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USER MANUAL

PRECISION SCALE

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PŘESNÁ VÁHA

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MANUAL DE INSTRUCCIONES

BALANZA DE PRECISIÓN

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SBS-LW-2000A

INHALT | CONTENT | TREŚĆ | OBSAH | CONTENU | CONTENUTO | CONTENIDO

▪ Deutsch	3
▪ English	10
▪ Polski	17
▪ Český	24
▪ Français	31
▪ Italiano	38
▪ Español	45

NAZWA PRODUKTU PRODUKTNAME PRODUCT NAME NOM DU PRODUIT NOME DEL PRODOTTO NOMBRE DEL PRODUCTO NÁZEV VÝROBKU	PRÄZISIONSWAAGEN PRECISION SCALE WAGA PRECYZYJNA PŘESNÁ VÁHA BALANCES DE PRÉCISION BILANCE DI PRECISIONE BALANZA DE PRECISIÓN
MODEL PRODUKTU MODELL PRODUCT MODEL MODÈLE MODELLO MODELO MODEL VÝROBKU	SBS-LW-2000A
HERSTELLER MANUFACTURER PRODUCENT VÝROBCE FABRICANT PRODUTTORE FABRICANTE	EXPONDO POLSKA SP. Z O.O. SP. K.
ANSCHRIFT DES HERSTELLERS MANUFACTURER ADDRESS ADRES PRODUCTENTA ADRESA VÝROBCE ADRESSE DU FABRICANT INDIRIZZO DEL PRODUTTORE DIRECCIÓN DEL FABRICANTE	UL. NOWY KISIELIN – INNOWACYJNA 7, 66-002 ZIELONA GÓRA POLAND, EU

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1. EINFÜHRUNG

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Sealed Lead Acid
Battery
Must be recycled
Properly

1. INTRODUCTION

Thank you for choosing our product. Each scale has a range of options and features that are ideal both for use in laboratory, classroom or for private use. We hope you enjoy working with our balance!

2. PACKAGE CONTENTS

- scale
- Weighing platform
- Power Cable
- Manual

3. BEFORE FIRST USE

INSERTING THE BATTERIES

Remove the battery box cover and insert the batteries (AA)

SCREEN AND KEYS

KEYS:

- ON/OFF - turn on / off of the scale
- UNITS – units change
- PCS - Counting function:
 1. per unit
 2. the scale can calculate a percentage
- ZERO – scale zeroing

SCREEN:

1. Low battery symbol.
2. Zero symbol – is displayed when the zero option is activated.
3. Tare symbol.
4. Stable symbol – is displayed when the result is stable.
5. Not used.
6. Not used.
7. Not used.
8. Not used.
9. Display showing the current weighing unit / pieces counting.
10. Not used.
11. Percentage counting mode

HHHHH – the weight is too big, the scale cannot weigh

PLACEMENT OF THE DEVICE AND PRECAUTIONARY MEASURES

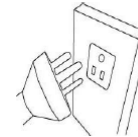
To keep your balance in a good condition, we recommend you follow these tips:


Avoid extreme temperatures. Do not place the unit in direct sunlight or near the air conditioning.

Make sure the scale is placed on a stable table and is not subject to vibrations.

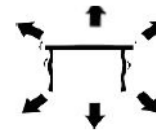Avoid unstable power sources. Do not use close to large electrical appliances such as welding machines or large motors. Do not leave batteries in the device if you do not use it for a long time.

Avoid vibrations. Do not place your scale near vibrating heavy machinery.

Avoid places where moisture can lead to condensation. Avoid direct contact with water.

Do not place the unit close to windows, doors or fans that may cause unstable results due to air currents.

Keep your machine clean. Do not store products on the scale when it is not used.

4. OPERATION

GENERAL INFORMATION

- Capacity / Size: 2000g / 0,01 g
- Display: LCD digital
- Backlight: blue
- Auto power save function: time off automatically programmable (2.5 or 8 min.)
- Units / Staging: g, ct, lb, oz, dr, GN, ozt, dwt, MM, tl.J, tl.T, tl.H, t (a total of 13 units)
- Counting function: 1. per unit 2. the scale can calculate a percentage
- Power: DC 9V adapter or AA batteries
- Weight: 550g

5. FUNCTIONS OF THE KEYS

ON / OFF

Press the ON / OFF button to start the unit. Hold the button pressed for 3 seconds to turn off the scale.

UNIT

The unit change is possible by pressing the UNIT button. Then press up to reach the desired unit.

PCS

Counting function: Press PCS to enter the counting function. The scales switches from the weighing mode to the counting mode. By pressing this button, you can switch from one mode to another.

ZERO

Press this button to subtract the weight of a container or any weight on the scale.

CALIBRATION AND PARAMETER SETTING

The calibration is performed in two stages. To calibrate the scale on 2kg, you need standards of 1 and 2 kg. Turn on the scale. Hold pressed the zero key for 3 seconds until the display shows „CALL“.

The identification number of the scale appears (SCL)

Press briefly again ON / OFF to display „CAL-1“. Wait until the scale displays „0“

Place the standard of 1 kg on the scale. It shows „PASS-1“ and „CAL-2“.

Then place the standard of 2 kg on the scale. It displays „PASS-2“

The scale is then calibrated.

FUNCTION OF TARE

- Place an empty container on the weighing platform until the balance indicates a stable value. Then press „ZERO“ Your balance now displays „0“.
- To exit the tare function, remove the container and repress the „ZERO“ button. Your scale is back to normal.

AUTO-ZERO

- Press „ZERO“ if the scale does not show „0“ when nothing is placed on it.
- The area of auto zero is $\pm 2\%$ of the total capacity of the scale.

DO NOT OVERLOAD THE BALANCE!

COUNTING MODE

- Press „PCS“ to activate the counting function. Your weight unit is maintained.
- Press „UNIT“, then choose between 100 and 1000 pieces and 10-1000 units by repeated pressure.
- Place a unit / object on the scale. Press ON / OFF for the scale to remind the weight of a unit.

PERCENTAGE MODE

- Press twice „PCS“ then press „UNIT“ to choose between 100.00 and 100.0
- After choosing the percentage of accuracy, place the weight on the scale and press „ON / OFF.“ This weight now accounts for 100%. You can now place a weight, the scale displays the percentage.

NOTE

To return to weighing mode, press „PCS“

ADVANCED SETUP:

	ACTION / STEP	Value which can be adjusted (bolded value is factory setting) Values can be changed by UNITS button	DESCRIPTION
1.	Scale must be turned off. Simultaneously press and hold ON/OFF and PCS buttons for ~5 sec. – display will show: [0] F01.	0/1	Units setting 0: without setting, move to next step 1: open or close some of the 13 units

2.	Press [ON/OFF] button – display will show: [0]F02	0/2/10/50	Zero returning range when the scale powers on: 2: If the weight exceeds 2 percent of the capacity before the scale powers on, the scale will not work after turning on
3.	Press [ON/OFF] button – display will show: [2.0]F03	2.0/0.0/0.3/0.5/0.7/1.0/1.5	Zero tracking range
4.	Press [ON/OFF] button – display will show: [0.5]F04	0.5/0.7/1.0/1.5/2.0/0.0/0.3	Dynamic tracking range
5.	Press [ON/OFF] button – display will show: [1.0]F05	1.0/2.0/3.0/0.5	Dynamic tracing time
6.	Press [ON/OFF] button – display will show: [0]F06	0/2/5/8	Auto-off after 2, 5 or 8 minutes of inactivity.
7.	Press [ON/OFF] button – display will show: [0] CAL Press [UNITS] button – display will show: [1] CAL	0/1	0 – after pressing the ON/OFF button, scale will exit the setup with saving changes 1 – after pressing the ON/OFF button scale will proceed with the remain of the setup
8.	Press [ON/OFF] button – display will show: [200000]WET	200000/300000/10000/20000/30000/50000/60000/100000/120000/150000	Capacity of the scale.
9.	Press [ON/OFF] button – display will show: [1]diV	1/2/5	Division 1, 2 or 5.
10.	Press [ON/OFF] button – display will show: [0.01]dEC	0.01/0.001/0.1	Changing the decimal place of the displayed weight value

11.	Press [ON/OFF] button – display will show: [1]LCP	1/0	Linear correction switch: 0 is off, 1 is on.
12.	Press [ON/OFF] button – display will show: [100000]WE1	100000/120000/150000/200000/300000/10000/20000/30000/50000/60000	Value of first standard weight for the calibration
13.	Press [ON/OFF] button – display will show: [200000]WE2	200000/300000/10000/20000/30000/50000/60000/100000/120000/150000	Value of second standard weight for the calibration
14.	Press [ON/OFF] button – display will show: [---CODE---]SCL	-	-
15.	Press [ON/OFF] button – display will show: CALL1 and [0]CAL	-	-
16.	Put on the scale 1 kg weight, display will show [---CODE---] and PASS-1 and CALL-2	-	-
17.	Very quickly take out the 1 kg weight from the scale and put 2 kg	-	-
18.	Display will show PASS – take out the weight from the scale	-	-
19.	End of setup	-	-

ATTENTION: All the parameters of above advanced functions have been thoroughly tested and set by the manufacturer. It is strongly recommended not to modify any of these parameters, otherwise the scale may not work properly. However positions no 6, 9 and 10 (auto –off, division and decimal point) can be modified according to the needs.

REGULAR CONTROL OF THE DEVICE

Control regularly that the device doesn't present any damage. If there is any damage, please stop using the device. Please contact your customer service to solve the problem.

What to do in case of a problem?

Please contact your customer service and prepare following information:

- Invoice number and serial number (the latter is to be found on the technical plate on the device).
- If relevant, a picture of the damaged, broken or defective part.
- The more detailed your information, the better the customer service will be able to solve your problem rapidly and efficiently!

CAUTION: Never open the device without the authorization of your customer service. This can lead to a loss of warranty!

6. RESPECT THE WEEE (ENVIRONMENTAL STANDARD)

According to the European standard 2002/95/EEC, electrical or electronic components and must be recycled or disposed so that they do not damage the environment and the health. The recycling of batteries is clearly regulated by laws of 2002 and 2005 and was updated in the text WEEE (en: Recycling of electrical and electronic equipment.)

Sealed Lead Acid
Battery
Must be recycled
Properly

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