USER MANUAL

Bedienungsanleitung
Instrukcja obsługi
Návod k použití
Manuel d'utilisation
Istruzioni per l'uso Manual
de instrucciones Használati
útmutató Brugsanvisning
Käyttöohje
Gebruiksaanwijzing
Bruksanvisning Instruções
de utilização Používateľská
príručka Ръководство за
употреба Οδηγίες χρήσης
Upute za uporabu
Naudojimo instrukcija
Manual de utilizare
Navodila za uporabo

PRECISION SCALE

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For products shipped in October 2025 or later

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OverLd – The weight is too big, the scale cannot weigh.

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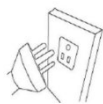
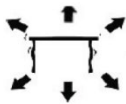
4.4. Device use

4.4.1. General information

- a) Capacity / Size: 2000g / 0.01g
- b) Display: LCD digital
- c) Backlight: blue
- d) Auto power save function: time off automatically programmed (0.2.5 or 8 min)
- e) Units / Staging: g, ct, lb, oz, dr, GN, ozt, dwt, MM, tl.J, tl.T, tl.H, t (a total of 13 units)
- f) Counting function:
 - 1) Per unit 2
 - 2) The scale can calculate a percentage
- g) Power: DC 8V adapter or AA batteries
- h) Weight: 550g

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4.4.6. Calibration and parameter setting

- 1) The calibration is performed in one stage. To calibrate the scale on 500g, you need standards of 500g.
- 2) Turn on the scale.
- 3) Press and hold the **ZERO** key for 3 seconds until the display shows "LoAd".
- 4) The identification number of the scale appears (WEIGHT NUMBER)
- 5) Then place the standard of 500g on the scale. It displays "PASS".
- 6) The scale is then calibrated.

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4.4.8. Auto-zero

- 1) Press **ZERO** if the scale does not show "0" when nothing is placed on it.
- 2) The area of auto zero is $\pm 4\%$ of the total capacity of the scale. **DO NOT OVERLOAD THE BALANCE!**

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4.4.11. Advanced setup

No.	Action	Value which can be adjusted (bolded value is factory setting) Values can be changed by PCS button	Description
1	Scale must be turned off. During the switch on process, press PCS 3 times + ZERO key. The display will show: [2E] -0..5	0.5 /1.0/1.5/2.0/2.5/3.0/3.5/ 4.0/4.5/5.0	Zero tracking range: Select by PCS Press ZERO to confirm and proceed to the next step
2	display will show: [HL-0.5]	0.5 /1.0/1.5/2.0/2.5/3.0/3.5/ 4.0/4.5/5.0/5.5/6.0/6.5/7.0/ 7.5/8.0/8.5/9.0	Zero returning range when the scale powers on: Select by PCS Press ZERO to confirm and proceed to the next step
3	Display will show: [Flt-3]	1/2/ 3 /4/5/6	Filter setting: Select by PCS Press ZERO to confirm and proceed to the next step
4	Display will show: [TrA-ON]	TrA-ON /trA-OF	Creep tracking switch: Select by PCS Press ZERO to confirm and proceed to the next step
5	Display will show: [CLS-0]	0 /2/5/8	Auto-off after 2, 5, or 8 minutes of inactivity: Select by PCS Press ZERO to confirm and return to weighing mode

ATTENTION: All the parameters of the above advanced functions have been thoroughly tested and set by the manufacturer. It is strongly recommended not to modify any of these parameters; otherwise, the scale may not work properly. However, No. 5 (auto-off) can be modified according to the needs.

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4.5.3. Disposing of used devices

Do not dispose of this device in municipal waste systems. Hand it over to an electric and electrical device recycling and collection point. Check the symbol on the product, instruction manual and packaging. The plastics used to construct the device can be recycled in accordance with their markings. By choosing to recycle you are making a significant contribution to the protection of our environment.

Contact local authorities for information on your local recycling facility.

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For products shipped in September 2025 or earlier

This User Manual has been translated using machine translation. We have made every effort to ensure the translation is accurate, but please note that automated translations are not perfect and are not meant to replace human translators. The official version of the User Manual is in English. Any differences between the translated version and the original English are not legally binding. If you have any questions about the accuracy of the translation, please refer to the English version, which is the official reference. More language versions are available upon request via info@expondo.com.

1. Technical data

Table 1: Technical data of the product

Parameter description	Parameter value
Product name	Precision scale
Model	SBS-LW-2000A LABORWAAGE / WAGA D
Max. load [g]	2000
Min. load [g]	0.05
Division [g]	0.01
Battery	6 x AA
Dimensions [width x depth x height; mm]	130 x 168 x 53
Weight [kg]	0.55

2. General description

The user manual is designed to assist in the safe and trouble-free use of the device. The product is designed and manufactured in accordance with strict technical guidelines, using state-of-the-art technologies and components. Additionally, it is produced in compliance with the most stringent quality standards.

DO NOT USE THE DEVICE UNLESS YOU HAVE THOROUGHLY READ AND UNDERSTOOD THIS USER MANUAL.

To increase the product life of the device and to ensure trouble-free operation, use it in accordance with this user manual and regularly perform maintenance tasks. The technical data and specifications in this user manual are up to date. The manufacturer reserves the right to make changes associated with quality improvement. The device is designed to reduce noise emission risks to a minimum, taking into account technological progress and noise reduction opportunities.

2.1. Legend

Icon	Description
	The product satisfies the relevant safety standards.

	Read instructions before use.
	The product must be recycled.
	WARNING! or CAUTION! or REMEMBER! Applicable to the given situation. (general warning sign)
	ATTENTION! Electric shock warning!

PLEASE NOTE! DRAWINGS IN THIS MANUAL ARE FOR ILLUSTRATION PURPOSES ONLY AND IN SOME DETAILS MAY DIFFER FROM THE ACTUAL PRODUCT.

3. Usage safety

ATTENTION! READ ALL SAFETY WARNINGS AND ALL INSTRUCTIONS. FAILURE TO FOLLOW THE WARNINGS AND INSTRUCTIONS MAY RESULT IN ELECTRIC SHOCK, FIRE AND/OR SERIOUS INJURY OR EVEN DEATH.

The terms "device" or "product" are used in the warnings and instructions to refer to:

Precision scale

3.1. Electrical safety

- a) Do not touch the device with wet or damp hands.
- b) **ATTENTION! DANGER TO LIFE!** While cleaning, never immerse the device in water or other liquids.
- c) Do not use in very humid environments or in the direct vicinity of water tanks.
- d) Prevent the device from getting wet. Risk of electric shock!

3.2. Safety in the workplace

- a) Make sure the workplace is clean and well lit. A messy or poorly lit workplace may lead to accidents. Try to think ahead, observe what is going on and use common sense when working with the device.
- b) If you discover damage or irregular operation, immediately switch the device off and report it to a supervisor without delay.

- c) If you are unsure about whether the product is operating correctly or if you find damage, please contact the manufacturer's service centre.
- d) Only the manufacturer's service centre may make repairs to the product. Do not attempt to make repairs yourself!
- e) In case of fire, use a powder or carbon dioxide (CO₂) fire extinguisher (one intended for use on live electrical devices) to put it out.
- f) Regularly inspect the condition of the safety labels. If the labels are illegible, they must be replaced.
- g) Please keep this manual available for future reference. If this device is passed on to a third party, the manual must be passed on with it.
- h) Keep packaging elements and small assembly parts in a place not available to children.

REMEMBER! WHEN USING THE DEVICE, PROTECT CHILDREN AND OTHER BYSTANDERS.

3.3. Personal safety

- a) Do not use the device when tired, ill or under the influence of alcohol, narcotics or medication which can significantly impair the ability to operate the device.
- b) The device is not designed to be handled by persons (including children) with limited mental and sensory functions or persons lacking relevant experience and/or knowledge unless they are supervised by a person responsible for their safety or they have received instruction on how to operate the device.

3.4. Safe device use

- a) Do not overload the device. Use the appropriate tools for the given task. A correctly-selected device will perform the task for which it was designed better and in a safer manner.
- b) Keep the device in perfect technical condition. Before each use check for general damage, especially check moving components for cracked parts or elements, and for any other conditions which may impact the safe operation of the device. If damage is discovered, hand over the device for repair before use.
- c) Device repair or maintenance should be carried out by qualified persons, only using original spare parts. This will ensure safe use.
- d) To ensure the operational integrity of the device, do not remove factory-fitted guards and do not loosen any screws.
- e) When transporting and handling the device between the warehouse and the destination, observe the occupational health and safety principles for manual transport operations which apply in the country where the device will be used.
- f) Clean the device regularly to prevent stubborn grime from accumulating.

- g) Ensure that the weight placed on the scales is stable.
- h) Check that all screws, rods, and brackets are secure.
- i) Do not weigh in case of extreme humidity and rain.
- j) Avoid working at high ambient temperatures.

ATTENTION! DESPITE THE SAFE DESIGN OF THE DEVICE AND ITS PROTECTIVE FEATURES, AND DESPITE THE USE OF ADDITIONAL ELEMENTS PROTECTING THE OPERATOR, THERE IS STILL A SLIGHT RISK OF ACCIDENT OR INJURY WHEN USING THE DEVICE. STAY ALERT AND USE COMMON SENSE WHEN USING THE DEVICE.

4. Use guidelines

Thank you for choosing our product. Each scale has a range of options and features that are ideal both for use in laboratory, classroom, or for private use.

We hope you enjoy working with our balance!

The user is liable for any damage resulting from unintended use of the device.

4.1. Product description

4.1.1. Keys

- ON/OFF - Turn on / off the scale
- UNITS – Units change
- PCS - Counting function:
 - 1) Per unit
 - 2) The scale can calculate a percentage
- ZERO – Scale zeroing

4.1.2. Screen

- 1) Low battery symbol.

- 2) Zero symbol – is displayed when the zero option is activated.
- 3) Tare symbol.
- 4) Stable symbol – is displayed when the result is stable.
- 5) Not used.
- 6) Not used.
- 7) Not used.
- 8) Not used.
- 9) Display showing the current weighing unit / pieces counting.
- 10) Not used.
- 11) Percentage counting mode

HHHHH – The weight is too big, the scale cannot weigh.

4.2. Package contents

- Scale
- Weighing platform
- Power Cable
- Manual

4.3. Preparing for use

4.3.1. Inserting the batteries

Remove the battery box cover and insert the batteries (AA)

4.3.2. Placement of the device and precautionary measures

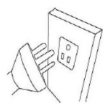
To keep your balance in good condition, we recommend you follow these tips:


Avoid extreme temperatures. Do not place the unit in direct sunlight or near the air conditioning.

Make sure the scale is placed on a stable table and is not subject to vibrations.

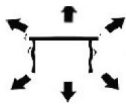Avoid unstable power sources. Do not use close to large electrical appliances such as welding machines or large motors. Do not leave batteries in the device if you do not use it for a long time.

Avoid vibrations. Do not place your scale near vibrating heavy machinery.

Avoid places where moisture can lead to condensation. Avoid direct contact with water.

Do not place the unit close to windows, doors or fans that may cause unstable results due to air currents.

Keep your machine clean. Do not store products on the scale when it is not used.

4.4. Device use

4.4.1. General information

- a) Capacity / Size: 2000g / 0.01 g
- b) Display: LCD digital
- c) Backlight: blue
- d) Auto power save function: time off automatically programmed (2.5 or 8 min.)
- e) Units / Staging: g, ct, lb, oz, dr, GN, ozt, dwt, MM, tl.J, tl.T, tl.H, t (a total of 13 units)
- f) Counting function:
 - 1) Per unit 2
 - 2) The scale can calculate a percentage
- g) Power: DC 9V adapter or AA batteries
- h) Weight: 550g

4.4.2. ON / OFF

Press the ON / OFF button to start the unit. Hold the button pressed for 3 seconds to turn off the scale.

4.4.3. UNIT

The unit change is possible by pressing the **UNIT** button. Then press up to reach the desired unit.

4.4.4. PCS

Counting function: Press **PCS** to enter the counting function. The scales switch from the weighing mode to the counting mode. By pressing this button, you can switch from one mode to another.

4.4.5. ZERO

Press this button to subtract the weight of a container or any weight on the scale.

4.4.6. Calibration and parameter setting

- 1) The calibration is performed in two stages. To calibrate the scale on 2 kg, you need standards of 1 and 2 kg.
- 2) Turn on the scale.
- 3) Hold the **ZERO** key for 3 seconds until the display shows "CALL".
- 4) The identification number of the scale appears (SCL).
- 5) Press briefly again **ON / OFF** to display "CAL-1". Wait until the scale displays "o".
- 6) Place the standard of 1 kg on the scale. It shows "PASS-1" and "CAL-2". Then place the standard of 2 kg on the scale. It displays "PASS-2".
- 7) The scale is then calibrated.

4.4.7. Function of tare

- 1) Place an empty container on the weighing platform until the balance indicates a stable value. Then press **ZERO**. Your balance now displays "0".
- 2) To exit the tare function, remove the container and repress the **ZERO** button. Your scale is back to normal.

4.4.8. Auto-zero

- 1) Press **ZERO** if the scale does not show "0" when nothing is placed on it.
- 2) The area of auto zero is $\pm 2\%$ of the total capacity of the scale. **DO NOT OVERLOAD THE BALANCE!**

4.4.9. Counting mode

- 1) Press **PCS** to activate the counting function. Your weight unit is maintained.
- 2) Press **UNIT**, then choose between 100 and 1000 pieces and 10-1000 units by repeated pressure.
- 3) Place a unit / object on the scale. Press **ON / OFF** for the scale to remind the weight of a unit.

4.4.10. Percentage mode

- 1) Press twice **PCS**, then press **UNIT** to choose between 100.00 and 100.0

- 2) After choosing the percentage of accuracy, place the weight on the scale and press ON / OFF . This weight now accounts for 100%. You can now place a weight; the scale displays the percentage.

Note: To return to weighing mode, press PCS .

4.4.11. Advanced setup

No.	Action	Value which can be adjusted (bolded value is factory setting) Values can be changed by UNIT button	Description
1	Scale must be turned off. Simultaneously press and hold ON / OFF and PCS buttons for ~5 sec. – display will show: [0] F01	0/1	Units setting 0: without setting, move to next step 1: open or close some of the 13 units
2	Press ON / OFF button – display will show: [0]F02	0/2/10/50	Zero returning range when the scale powers on: 2: If the weight exceeds 2 percent of the capacity before the scale powers on, the scale will not work after turning on
3	Press ON / OFF button – display will show: [2.0]F03	2.0/0.0/0.3/0.5/0.7/1.0/1.5	Zero tracking range
4	Press ON / OFF button – display will show: [0.5]F04	0.5/0.7/1.0/1.5/2.0/0.0/0.3	Dynamic tracking range
5	Press ON / OFF button – display will show: [1.0]F05	1.0/2.0/3.0/0.5	Dynamic tracing time
6	Press ON / OFF button – display will show: [0]F06	0/2/5/8	Auto-off after 2, 5, or 8 minutes of inactivity.
7	Press ON / OFF button – display will show: [0] CAL Press UNIT button – display will show: [1] CAL	0/1	0 – after pressing the ON / OFF button, the scale will exit the setup with saving changes 1 – after pressing the ON / OFF button, the scale will proceed with the remain of the setup
8	Press ON / OFF button – display will show: [200000]WET	200000/300000/10000/2000 0/30000/50000/60000/10000 0/120000/150000	Capacity of the scale.
9	Press ON / OFF button – display will show: [1]diV	1/2/5	Division 1, 2, or 5.

10	Press ON / OFF button – display will show: [0.01]dEC	0.01/0.001/0.1	Changing the decimal place of the displayed weight value
11	Press ON / OFF button – display will show: [1]LCP	1/0	Linear correction switch: 0 is off, 1 is on.
12	Press ON / OFF button – display will show: [100000]WE1	100000/120000/150000/200000/300000/10000/20000/30000/50000/60000	Value of the first standard weight for the calibration
13	Press ON / OFF button – display will show: [200000]WE2	200000/300000/10000/20000/30000/50000/60000/100000/120000/150000	Value of the second standard weight for the calibration
14	Press ON / OFF button – display will show: [---CODE---]SCL	-	-
15	Press ON / OFF button – display will show: CALL1 & [0]CAL	-	-
16	Put on the scale 1 kg weight, display will show [---CODE---] & PASS-1 & CALL-2	-	-
17	Quickly take out the 1 kg weight from the scale and put the 2 kg	-	-
18	Display will show PASS – take out the weight from the scale	-	-
19	End of setup	-	-

ATTENTION: All the parameters of the above advanced functions have been thoroughly tested and set by the manufacturer. It is strongly recommended not to modify any of these parameters; otherwise, the scale may not work properly. However, No. 6, 9, and 10 (auto-off, division, and decimal point) can be modified according to the needs.

4.4.12. Regular control of the device

Control regularly that the device doesn't present any damage. If there is any damage, please stop using the device. Please contact your customer service to solve the problem.

What to do in case of a problem?

Please contact your customer service and prepare the following information:

- Invoice number and serial number (the latter is to be found on the technical plate on the device).
- If relevant, a picture of the damaged, broken, or defective part.
- The more detailed your information, the better the customer service will be able to solve your problem rapidly and efficiently!

CAUTION: Never open the device without the authorization of your customer service. This can lead to a loss of warranty!

4.5. Cleaning and maintenance

4.5.1. General instructions

- a) Use only non-corrosive cleaners to clean the surface.
- b) After cleaning the device, all parts should be dried completely before using it again.
- c) Store the unit in a dry, cool place, free from moisture and direct exposure to sunlight.
- d) Do not spray the device with a water jet or submerge it in water.
- e) Do not allow water to get inside the device through vents in the housing of the device.
- f) The device must be regularly inspected to check its technical efficiency and spot any damage.
- g) Use a soft cloth for cleaning.
- h) Do not leave the battery in the device if it will not be used for a longer period of time.
- i) Do not use sharp and/or metal objects for cleaning (e.g. a wire brush or a metal spatula) because they may damage the surface material of the appliance.
- j) Do not clean the device with an acidic substance, agents of medical purposes, thinners, fuel, oils or other chemical substances because it may damage the device.

4.5.2. Safe removal of batteries and rechargeable batteries

Batteries are installed in the devices.

Remove used batteries from the device using the same procedure by which you installed them.

Recycle batteries with the appropriate organisation or company.

4.5.3. Respect the WEEE (environmental standard)

According to the European standard 2002/95/EEC, electrical or electronic components must be recycled or disposed, so that they do not damage the environment and health. The recycling of batteries is clearly regulated by the laws of 2002 and 2005 and was updated in the text WEEE (Recycling of electrical and electronic equipment.)

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