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USER MANUAL

ELECTRIC STIRRER

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AGITATEUR ELECTRIQUE

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MANUAL DE INSTRUCCIONES

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INHALT | CONTENU | CONTENT | CONTENUTO | CONTENIDO | TREŚĆ | OBSAH

Deutsch	3
English	5
Polski	7
Česky	9
Français	11
Italiano	13
Español	15

NAZWA PRODUKTU	MIESZADŁO ELEKTRYCZNE
PRODUKTNAME	ELEKTRISCHER RÜHRER
PRODUCT NAME	ELECTRIC STIRRER
NOM DU PRODUIT	AGITATEUR ELECTRIQUE
NOME DEL PRODOTTO	AGITATORE ELETTRICO
NOMBRE DEL PRODUCTO	AGITADOR ELÉCTRICO
NÁZEV VÝROBKU	ELEKTRICKÉ MÍCHADLO
MODEL PRODUKTU	SBS-ER-3000
MODELL	
PRODUCT MODEL	
MODÈLE	
MODELLO	
MODELO	EXPONDO POLSKA SP. Z O.O. SP. K.
MODEL VÝROBKU	
HERSTELLER	
MANUFACTURER	
PRODUCENT	
FABRICANT	UL. NOWY KISIELIN – INNOWACYJNA 7, 66-002 ZIELONA GÓRA POLAND, EU
PRODUTTORE	
FABRICANTE	
VÝROBCE	
ANSCHRIFT DES HERSTELLERS	
MANUFACTURER ADDRESS	
ADRES PRODUCENTA	
ADRESSE DU FABRICANT	
INDIRIZZO DEL PRODUTTORE	
DIRECCIÓN DEL FABRICANter	
ADRESA VÝROBCE	

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USER MANUAL

The operation manual must be reviewed.

Never dispose of electrical equipment together with household waste.

This machine conforms to the CE declarations.

For indoor use only.

HINWEIS! Drawings in this manual are for illustration purposes only and in some details it may differ from the actual machine.

The original operation manual is in German. Other language versions are translations from German.

USER MANUAL

SAFETY INSTRUCTIONS

General safety information for the use of electrical devices To avoid injury from fire or electric shock, please ensure the compliance with the safety instructions when using this device. Please read the instructions carefully and make sure that you have understood them well. Keep the manual near the equipment to be able to read it at any time. Always use current sources which are connected to the ground and which provide the necessary voltage (this info is indicated on the label on the device). If you have any doubts, let an electrician check that your outlet is properly grounded. Never use a damaged power cable. Do not open the unit in damp or wet environment, or if your hands or body are damped or wet. Protect the unit from solar radiation. Use the device in a protected location to avoid damaging the equipment or endangering others. Make sure the device is able to cool and avoid placing it too close to other devices that produce heat. Before cleaning the device, disconnect it from the power source. Use a soft damp cloth for cleaning. Avoid using detergents and make sure that no liquid enters the unit. No internal element of this device needs to be maintained by the user. Opening the device without our approval leads to a loss of warranty!

SAFETY GUIDELINES

- 1. Please read and understand these instructions carefully before operating the device! Please observe the safety guidelines carefully in order to prevent damages through improper use!
- 2. Please keep this manual available for future reference. If this device is passed on to third parties, the manual must be passed on with it.
- 3. Only use this device indoors and for its intended use.
- 4. We do not offer any warranty for damages resulting from improper use or incorrect operation.
- 5. Before the first use, check whether the main voltage type and current comply with the indicated data on the type plate.
- 6. This device is not intended to be used by persons (including children) with limited physical, sensory or mental aptitude, as well as lack of experience and/or knowledge, unless they are supervised by a person responsible for their safety or if they have received instructions as to how the device is used.
- 7. **ELECTRIC SHOCK HAZARD!** Do not attempt to repair the unit yourself. In case of device failures, repairs must be done by qualified experts.
- 8. Please check the main plug and the power cable regularly. If the power cable for this device is damaged, it has to be replaced by the manufacturer

or its customer service or another qualified person in order to prevent hazards.

- 9. Prevent damages to the power cable by avoiding squeezing or bending it or rubbing it on sharp edges. Also keep the power cable away from hot surfaces and open flames.
- 10. **ATTENTION! DANGER TO LIFE!** While cleaning, never immerse the device in water or other liquids.
- 11. Under no circumstances should the housing be opened.

TECHNICAL DATA

Product name	Electric Stirrer
Model	SBS-ER-3000
Rated voltage [V]	230~
Frequency [Hz]	50
Rated power [W]	100
Timer range (timer) [min]	0 – 120
Rotational speed range [rpm]	~3400
Weight [kg]	7,5

SCOPE OF APPLICATION

The electric agitator is a device designated for general use in laboratories, schools, universities and similar establishments. Used for contactless liquid agitation in glass laboratory vessels, such as flasks, beakers, etc). **ATTENTION!** The product is outside of the scope of MDD 93/42/EEC Medical Devices regulations. It is not a medical device and should not be used for medical research and tests in hospitals and other health facilities. **The user is liable for any damage resulting from nonintended use of the device.**

BEFORE THE FIRST USE

Upon receipt of the goods, check the packaging for integrity and open it. If the packaging is damaged, please contact your transport company and distributor within 3 days, and document the damages as detailed as possible. Do not turn the package upside down! When transporting the package, please ensure that it is kept horizontal and stable.

DISPOSAL OF THE PACKAGING

Please keep all packaging materials (cardboard, plastic tapes and Styrofoam), so that in case of a problem, the device can be sent back to the service centre in accurate condition!

INSTALLING THE APPLIANCE

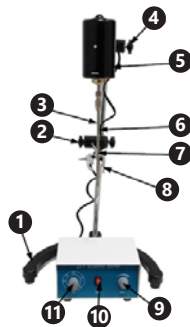
APPLIANCE LOCATION

The temperature of the environment must not exceed 45°C and the relative humidity should not exceed 85%. Ensure good ventilation in the room in which the device is used. The distance between the back of the chain toaster and the wall should be at least 20 cm, a minimum distance between the sides of the device and the walls should be at least 10 cm on each side and 35 cm between the top of the device and the ceiling. Keep the device away from hot surfaces. Operate the device on flat, stable, clean, refractory and dry surfaces out of the reach of children and people with mental disabilities.

Install the device making sure that constant access to the main plug is ensured. The power cord connected to the appliance must be properly grounded and correspond to the technical details. Before the first use, remove all elements and clean them along with the entire device.

PRINCIPLE OF OPERATION

1. base
2. clamp handle
3. foot
4. motor handle
5. motor
6. agitator arm
7. agitator
8. clamp
9. rotational speed adjuster
10. ON/OFF switch
11. timer

Before using the device, make sure the rotational speed adjustment knob is set to the minimum speed and that the timer is in the OFF position.

Place a container with the liquid in front of the agitator or suspend it in the clamp, adjust the height of the clamp and agitator. Use the ON/OFF switch to switch the device on. The switch will illuminate. Then use the timer to set the required time, or set the timer to „ON“ and use the revs controller to select an appropriate rotational speed. The agitator begins mixing. If a certain time was set on the timer, the agitator will stop after that time has elapsed. If the timer was set to „ON“ then the device will continue operating until the timer is set to „OFF“, the revs controller is set to its minimum setting or the device is switched off using the ON/OFF switch.

ATTENTION! If the device is in use for 2 h, a 30 min break is required!

ASSEMBLY/DISASSEMBLY OF THE DEVICE'S ELEMENTS

Lay the base on a flat surface. Screw two legs to the base. Slide a bracket onto each leg. Attach motor onto the upper bracket and tighten; attach and tighten the agitator to the motor shaft. Attach clamp onto the lower bracket and tighten securely. Insert lead into the back of the device.

TRANSPORTATION AND STORAGE

During transport, the machine should be protected from shaking, crashing and turning upside down. Store it in a properly ventilated place with dry air and without any corrosive gas.

CLEANING AND MAINTENANCE

- Always unplug the device before cleaning the equipment. Disconnect the device from the electricity when it is not in use. Let the device cool down before cleaning it.
- Use only cleaner without corrosive substances to clean the surface.
- After each cleaning, all the elements must be carefully dried and secured by greasing the heating plates.
- Store the unit in a dry, cool place, free from moisture and direct exposure to sunlight.

- Never clean the device with a pressurized water-jet, industrial cleaners, and brushes or any other type of cleaning utensil that could damage the surface of the appliance.

REGULAR CONTROL OF THE DEVICE

Check regularly the device to see if it presents any damages. If there is any damage, please stop using the device. Please contact your customer service to solve the problem.

What to do in case of a problem?

Please contact your customer service and prepare the following information:

- Invoice number and serial number (the latter is to be found on the technical plate on the device).
- If relevant, take a picture of the damaged, broken or defective part.
- It will be easier for your customer service clerk to determine the source of the problem if you give a detailed and precise description of the matter. The more detailed your information, the better the customer service will be able to solve your problem rapidly and efficiently!

CAUTION: Never open the device without the authorization of your customer service. This can lead to a loss of warranty!

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