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USER MANUAL

MAGNETIC STIRRER

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MANUAL DE INSTRUCCIONES

AGITADOR MAGNÉTICO

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INHALT | CONTENU | CONTENT | CONTENUTO | CONTENIDO | TREŚĆ | OBSAH

Deutsch	3
English	8
Polski	12
Česky	17
Français	21
Italiano	26
Español	31

NAZWA PRODUKTU	MIESZADŁO MAGNETYCZNE Z GRZANIEM	MIESZADŁO MAGNETYCZNE Z GRZANIEM I TIMEREM	MIESZADŁO MAGNETYCZNE Z GRZANIEM I TIMEREM
PRODUKTNAME	MAGNETRÜHRER 6FACH	MAGNETRÜHRER TIMER	MAGNETRÜHRER TIMER PROFESSIONAL
PRODUCT NAME	MAGNETIC STIRRER	MAGNETIC STIRRER TIMER	MAGNETIC STIRRER TIMER PROFESSIONAL
NOM DU PRODUIT	AGITATEUR MAGNÉTIQUE	AGITATEUR MAGNÉTIQUE AVEC MINUTERIE	AGITATEUR MAGNÉTIQUE PROFESSIONNEL AVEC MINUTERIE
NOME DEL PRODOTTO	AGITATORE MAGNETICO	AGITATORE MAGNETICO CON TIMER	AGITATORE MAGNETICO PROFESSIONALE CON TIMER
NOMBRE DEL PRODUCTO	AGITADOR MAGNÉTICO	AGITADOR MAGNÉTICO CON TEMPORIZADOR	AGITADOR MAGNÉTICO CON TEMPORIZADOR PROFESIONAL
NÁZEV VÝROBKU	MAGNETICKÉ MÍCHADLO	MAGNETICKÉ MÍCHADLO S ČASOVAČEM	MAGNETICKÉ MÍCHADLO S ČASOVAČEM PROFESIONÁLNÍ
MODEL PRODUKTU			
MODELL			
PRODUCT MODEL			
MODÈLE			
MODELLO			
MODELO			
MODEL VÝROBKU			
HERSTELLER			
MANUFACTURER			
PRODUCENT			
FABRICANT	EXPONDO POLSKA SP. Z O.O. SP. K.		
PRODUTTORE			
FABRICANTE			
VÝROBCE			
ANSCHRIFT DES HERSTELLERS			
MANUFACTURER ADDRESS			
ADRES PRODUCTENTA			
ADRESSE DU FABRICANT	UL. NOWY KISIELIN – INNOWACYJNA 7, 66-002 ZIELONA GÓRA POLAND, EU		
INDIRIZZO DEL PRODUTTORE			
DIRECCIÓN DEL FABRICANter			
ADRESA VÝROBCE			

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-  The operation manual must be reviewed.
-  Never dispose of electrical equipment together with household waste.
-  This machine conforms to CE declarations
-  Attention! Hot surface may cause burns.

PLEASE NOTE! Drawings in this manual are for illustration purposes only and in some details it may differ from the actual machine.

The original operation manual is in German. Other language versions are translations from German.

USER MANUAL
SAFETY INSTRUCTIONS

General safety information for the use of electrical devices: To avoid injury from fire or electric shock, please ensure compliance with safety instructions when using this device. Please read the instructions carefully and make sure that you have understood them well. Keep the manual near the equipment to be able to read it at any time. Always use current sources connected to the ground and providing the necessary voltage (indicated on the label on the device). If you have any doubt, let an electrician check that your outlet is properly grounded. Never use a damaged power cable. Do not open the unit in damp or wet environment, or if your hands or body are damp or wet. Protect the unit from solar radiation. Use the device in a protected location to avoid damaging the equipment or endangering others. Make sure the device is able to cool and avoid placing it too close to other devices that produce heat. Before cleaning, disconnect it from the power source. Use a soft damp cloth for cleaning. Avoid using detergents and make sure that no liquid enters the unit. No internal element of this device needs to be maintained by the user. Opening the device without our approval leads to a loss of warranty!

SAFETY GUIDELINES

1. Please read and understand these instructions carefully before operating the device! Please observe the safety guidelines carefully in order to prevent damages through improper use!
2. Please keep this manual available for future reference. If this device is passed on to third parties, the manual must be passed on with it.
3. Only use this device indoors and for its intended use.
4. We do not offer any warranty for damages resulting from improper use or incorrect operation.
5. Before the first use, please check whether the main voltage type and current comply with the indicated data on the type plate.
6. This device is not intended to be used by persons (including children) with limited physical, sensory or mental aptitude or lack of experience and/or knowledge, unless they are supervised by a person responsible for their safety or have received instruction from this person as to how the device is used.
7. **ELECTRIC SHOCK HAZARD!** Do not attempt to repair the unit yourself. In case of device failures, repairs must be done by qualified experts.
8. Please check main plug and power cable regularly. If the power cable for this device is damaged, it has to be replaced by the manufacturer or its customer service or another qualified person in order to prevent hazards.

9. Prevent damages to the power cable by avoiding squeezing or bending it or rubbing it on sharp edges. Also keep the power cable away from hot surfaces and open flames.
10. **ATTENTION! DANGER TO LIFE!** While cleaning, never immerse the device in water or other liquids.
11. Under no circumstances should the housing be opened.
12. The device is designed for internal use only.
13. Do not use the device at altitudes above 2000 m.
14. For use safety ensure the device housing is earthed.

TECHNICAL DATA

Product name	Magnetic Stirrer	Magnetic Stirrer Timer	Magnetic Stirrer Professional
Model	SBS-MR-1600/6	SBS-MR-1600/1T	SBS-MR-1600/1T PRO
Rated voltage [V~]/Frequency [Hz]	230/50		
Power [W]	870	310	660
Protection class	IP 30		
Temperature range [°C]	5-99	5 – 100	0-220
Maximum rotational speed [rev/min]	1600		
Volume of agitated substance [ml]	100 -1000	100 -5000	5000
Number of workstations	6	1	1
Work surface dimensions [mm]	φ120	161x161	187x187
Dimensions [mm]	301x963 x131	195x266 x168	310x209 x140
Maximum dimensions incl. mounted accessories [mm]	301x963 x390	195x266 x443	310x209 x525
Timer [min]	0-120	0-120	0-9999
Ambient temperature [°C]	15-25		
Weight [kg]	10,8	4,2	5,1

SCOPE OF APPLICATION

The magnetic agitator is a device designed for general use in laboratories, schools, universities and similar establishments. Used for contactless liquid agitation in glass laboratory vessels, such as flasks, beakers, etc.

The product is outside of the scope of MDD 93/42/EEC Medical Devices regulations.

It is not a medical device and should not be used for medical research and tests in hospitals and other health facilities.
The user is liable for any damage resulting from non-intended use of the device!

BEFORE THE FIRST USE

Upon receipt of the goods, check the packaging for integrity and open it. If the packaging is damaged, please contact your transport company and distributor within 3 days, and document the damages as detailed as possible. Do not turn the package upside down! When transporting the package, please ensure that it is kept horizontal and stable.

DISPOSING OF PACKAGING


Please keep all packaging materials (cardboard, plastic tapes and Styrofoam), so that in case of a problem, the device can be sent back to the service centre in accurate condition!

INSTALLING THE APPLIANCE
APPLIANCE LOCATION

Do not use in ambient temperatures outside of the 15-25°C range and if relative humidity exceeds 60%. Ensure good ventilation in the room in which the device is to be used. The distance between the back of the chain toaster and the wall should be at least 20 cm, minimum distance between the sides of the device and the walls should be at least 10 cm on each side and 35 cm between the top of the device and the ceiling. Keep the device away from hot surfaces. Operate the device on flat, stable, clean, refractory and dry surfaces out of the reach of children and people with mental disabilities. Install the device making sure that constant access to the main plug is ensured. The power cord connected to the appliance must be properly grounded and correspond to the technical details! Voltage fluctuation should not exceed ±10% of the rated value.

DEVICE OPERATION – BASIC PRINCIPLES

SBS-MR-1600/6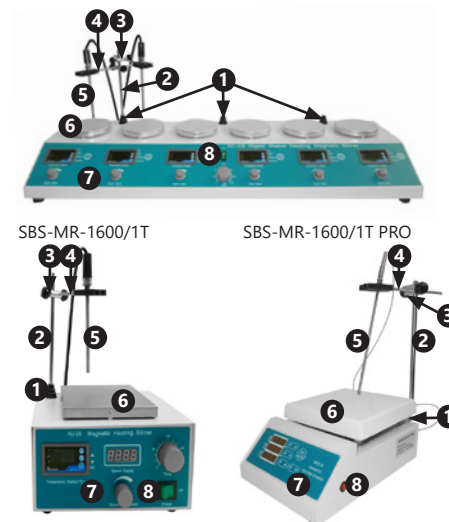


1. Stand base
2. Rod
3. Attachment
4. Clamp
5. Temperature sensor
6. Work surface
7. Control panel (see description below)
8. ON/OFF switch

User assembly of the temperature sensor stand is required.

To do that:

- Screw the threaded end of the rod (2) to base (1).
- Slide attachment (3) onto the rod (2) and tighten screw.
- Insert clamp (4) rod through the hole in the attachment and tighten screw.
- Insert temperature sensor into one of the holes in the clamp's plastic part.
- The device is ready for use.

Control panel of SBS-MR-1600/6 i SBS-MR-1600/1T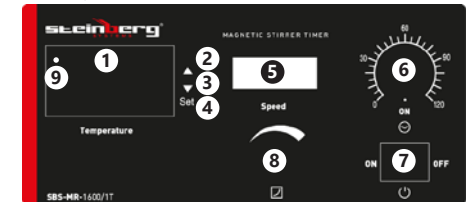


1. Temperature display
2. Increase value button
3. Decrease value button
4. „SET“ button
5. Rotational speed display
6. Timer
7. ON/OFF switch
8. Rotational speed adjustment knob
9. Diode

Temperature display abbreviations:

SV – Setting Value
PV – process variable / OUT – Output – process variable for PID controller (current temperature reading)

Control panel of SBS-MR-1600/1T PRO

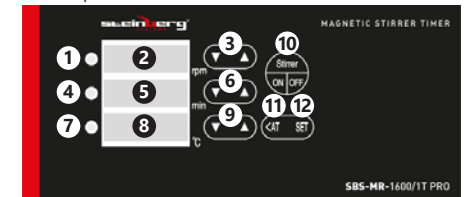1. Alarm diode
2. Rotational speed display
3. Rotational speed adjustment button
4. „OFF“ diode
5. Time display (timer)
6. Time setting adjustment button
7. „OUT“ diode
8. Temperature display
9. Temperature setting adjustment button
10. Heating, agitation and timer On/Off switch
11. Automatic adaptation function button
12. „SET“ button

OPERATION

Model SBS-MR-1600/6 | SBS-MR-1600/1T

1. Before use check state of the device
2. Position beaker (or another laboratory vessel) in the middle of the work surface.
3. Pour liquid into the vessel. Insert small magnetic rod (agitator) into the vessel with liquid.
4. Insert temperature sensor into the liquid (min. 2/3 of the sensor length should be submerged) by inserting it into a hole in the support. Adjust immersion height by changing the position of the attainment on the support.
5. Connect the device to a power source.
6. Use On/Off switch (7) to turn on; the switch will become illuminated.
7. „nHc“ will appear on the temperature display and after a while the current sensor temperature reading will be displayed.
8. To set the heating temperature:
 - Press the „SET“ (4) button, the temperature value will begin flashing.
 - Use Δ and ∇ (2; 3) buttons to set the temperature. ATTENTION: To change the temperature value up or down quickly press and hold the appropriate button.
 - Press the „SET“ (4) button again to confirm and return to operation mode.
 - The display will show the current temperature of the liquid with the submerged sensor.
 - The red diode on the bottom of the display will come on when heating is active.
 - Attention: The work surface heats up considerably!
9. Use knob (8) to adjust the agitator rotational speed. The rotational speed can only be adjusted once the timer (6) is set to „ON“ or countdown mode. The agitator will start moving.
10. The device features a PID controller to adjust the heating temperature. To enter controller parameter settings follow these steps:
 - press and hold the „SET“ (4) button for 3s. Ignore the „E00“ message on the display.
 - Press the „SET“ button twice, until a „P“ appears on the screen with a numerical value. The PID controller proportional term can be set in the 0.0 to 9.9 range. P=0“ is the threshold value and signifies no heating, P≠0 means that the liquid heating process is active. The output signal strength increases as the value of P increases, i.e. the liquid reaches the set temperature faster and the adjustment time decreases, and the maximum overshoot value increases.
 - press the „SET“ button to confirm the setting.
 - Additionally the user has an option to select the T adjustment time. In order to enter the adjustment time settings, proceed analogously to proportional term setting, however once the „E00“ message appears, press the „SET“ button three times. „T“ followed by a numerical value will appear on the display. The adjustment time can be set in the 1 – 99 range, with recommended settings between 20 and 60s.
 - To simplify use, controller parameters such as integration and differentiation times are hidden from the user and set to I=180 and D=45 by default.
11. Measurement error corrections If the user is certain that an incorrect temperature is shown on the display, it is possible to correct the reading in the – 4.9 to 4.9°C range.

To activate this option, press and hold the „SET“ button for 3s. „E00“ will appear on the display (threshold value – negligible parameter). Press the „SET“ button again, and then change the parameter using the Δ and ∇ buttons. Press the „SET“ button to confirm changes.

12. The device includes a timer. If it is set to „0“ then the device will heat up the liquid, but the agitation function will not be active. Set the knob to „ON“ for the device to operate without a time limit. The device may be operated in a time restricted mode with a maximum time setting of 120 min. The device switches off by itself once the set time has elapsed.
- For the SBS-MR-1600/6 model, the parameters described in points 7 – 11 can be set for each work station in use at the given time (6 in total). The device has one master On/Off switch and one timer for all work stations. It is not possible to set separate operation time values for given work stations.

Model SBS-MR-1600/1T PRO

1. Before use check state of the device
2. Position beaker (or another laboratory vessel) in the middle of the work surface.
3. Pour liquid into the vessel. Insert small magnetic rod (agitator) into the vessel with liquid.
4. Insert temperature sensor into the liquid (min. 2/3 of the sensor length should be submerged) by inserting it into a hole in the support. Adjust immersion height by changing the position of the attainment on the support.
5. Connect the device to a power source. Activate using the master On/Off switch on the side of the device.
6. Use Δ and ∇ (3 ; 6 ; 9) buttons to set the required rotational speed values, agitator operation time and heating temperature. Additionally the „<AT“ (11) button is used to switch between positions in the decimal numerical temperature value format. ATTENTION: To change the temperature value up or down quickly press and hold the appropriate button. The selected value will be automatically stored approx. 2 s after the button is released.
7. ATTENTION: The work surface heats up considerably! Press the „Stirrer On/Off“ (10) button, agitation and heating will commence and the timer will begin counting down. During heating element activation the „OUT“ (7) diode will be illuminated, and when the temperature is above the set temperature or the liquid is heated up to the required temperature the „OFF“ (4) diode will be on.
8. Once the time set by the user elapses, the timer alarm will sound, heating will switch off and the agitator will stop. The user may stop the device by pressing the „Stirrer On/Off“ (10) button.
9. To use the device without a time limit set 0000 as the time, Press the „Stirrer On/Off“ button to switch off the device.
10. The device features a PID controller to adjust the heating temperature. The user may set parameters shown in the table below. These parameters should only be set by a person with appropriate knowledge and qualifications. To enter the settings press and hold the „SET“ button for 3s. Use Δ and ∇ , buttons to set the values, use the „SET“ button to go to the next value. To confirm the selected parameter values press and hold the „SET“ button for approx. 3s.

Values will appear on the temperature display in the same order as in the table below. The „ALM“ (1) diode will come on if an alarm is triggered (a set alarm value is exceeded).

Parameter code	Parameter meaning	Settings range	Factory setting
LC	Lock 00 – No lock; 01 – Internal parameter changes locked 02 – Present temperature and internal parameter changes locked	---	01
H	Alarm value	±19°C	00.3
F	Alarm trip resistance zone	0-99.9	00.1
P	PID controller settings proportional term	0-99.9	02.0
I	PID controller constant integration time	0-999s	160
D	PID controller constant differentiation time	0-999s	040
T	Present temperature attainment period	1-200s	003
C	Sensor reading correction	±19°C	000
Y	No response of master output to changes zone (P=0)	0-99.9	00.1
E	Maximum overshoot limit	0-10.0	00.0
AT	Automatic PID value adaptation	OF/ON	OF

11. To activate automatic PID value adaptation option, press and hold the „<AT“ (11) button for 3s.

USE GUIDELINES

- Do not change the rotational speed violently, set it slowly by turning the knob. By setting the highest agitator speed tight at the outset might result in vibrations of the vessel with the liquid or even of the entire device.
- Do not activate heating under an empty vessel.
- To earth the housing, connect the earthing cable to a bolt located at the back of the device marked with the earth symbol.
- Monitor the liquid quantity in vessels. The liquid may evaporate.

- If the heating temperature is set to a value below ambient temperature the device will not heat up.
- The PID controller should only be configured by person with appropriate knowledge and qualifications. Configuration by a person with insufficient knowledge may result in faulty operation of the device.
- Always switch the device off after use.
- A user with appropriate knowledge and qualifications may replace the fuses at the back of the device housing. Change the fuse for an identical model with the same parameters.

TRANSPORTATION AND STORAGE

Shaking, crashing and turning upside down of the device should be prevented when transporting it. Store it in a properly ventilated location with dry air and without any corrosive gas.

CLEANING AND MAINTENANCE

- Always unplug the device before cleaning it.
- Use cleaner without corrosive substances to clean surface.
- Dry all parts well after cleaning before the device is used again.
- Store the unit in a dry, cool place, free from moisture and direct exposure to sunlight.
- Never spray the device with water jets.

REGULAR CONTROL OF THE DEVICE

Check regularly if the device is damaged. In case of damage, please stop using it immediately and contact customer service for solving the problem

What to do in case of a problem?

Please contact customer service and prepare the following information:

- Invoice number and serial number (the latter is to be found on the technical plate on the device).
- If necessary, take a picture of the damaged, broken or defective part.
- It will be easier for your customer service assistant to determine the source of the problem if you give a detailed and precise description of the problem. The more detailed your information, the better customer service will be able to solve your problem rapidly and efficiently!

CAUTION: Never open the device without the authorization of your customer service. This can lead to a loss of warranty!

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