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MIKROPIPETTE 100

USER MANUAL

MICROPIPETTE 100

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MIKROPIPETA 100

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MICROPIPETTE 100

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MICROPIPETTA 100

MANUAL DE INSTRUCCIONES

MICROPIPETA 100

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MIKROPIPETA 100

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INHALT | CONTENU | CONTENT | CONTENUTO | CONTENIDO | TREŚĆ | OBSAH

Deutsch	3
English	4
Polski	5
Česky	6
Français	7
Italiano	8
Español	9

NAZWA PRODUKTU	MIKROPIPETA 100
PRODUKTNAME	MIKROPIPETTE 100
PRODUCT NAME	MICROPIPETTE 100
NOM DU PRODUIT	MICROPIPETTE 100
NOME DEL PRODOTTO	MICROPIPETTA 100
NOMBRE DEL PRODUCTO	MICROPIPETA 100
NÁZEV VÝROBKU	MIKROPIPETA 100
MODEL PRODUKTU	
MODELL	
PRODUCT MODEL	
MODÈLE	SBS-PIP100
MODELLO	
MODELO	
MODEL VÝROBKU	
HERSTELLER	
MANUFACTURER	
PRODUCENT	
FABRICANT	EXPONDO POLSKA SP. Z O.O. SP. K.
PRODUTTORE	
FABRICANTE	
VÝROBCE	
ANSCHRIFT DES HERSTELLERS	
MANUFACTURER ADDRESS	
ADRES PRODUCENTA	
ADRESSE DU FABRICANT	UL. NOWY KISIELIN – INNOWACYJNA 7, 66-002 ZIELONA GÓRA POLAND, EU
INDIRIZZO DEL PRODUTTORE	
DIRECCIÓN DEL FABRICANER	
ADRESA VÝROBCE	

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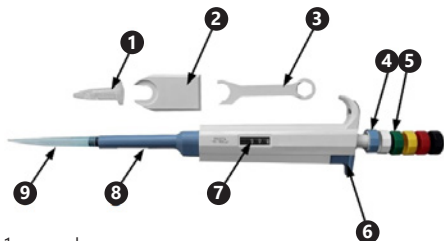
SCOPE OF APPLICATION

The micropipette is designed to accurately collect and dispense liquids.

The user is liable for any damage resulting from nonintended use of the device.

PRINCIPLE OF OPERATION

1. scale
2. 2hook
3. spanner/scale attachment
4. collect/dispense button, volume selector knob
5. removable sleeves for the collect/dispense button
6. pipette nozzle release / eject button
7. selected volume indicator
8. nozzle ejector
9. pipette nozzle

Attach nozzle to the pipette, use knob to select required volume (range: 100µl – 1000µl), immerse pipette nozzle in the liquid to a depth of about 1 cm and then press the collect button until resistance is discernible and release. Remove pipette nozzle from the liquid. To dispense the liquid from pipette nozzle, press the collect/dispense button all the way. To remove a nozzle, press the pipette nozzle release/eject button. To set the required liquid volume turn the volume selector knob anticlockwise or clockwise depending on whether more or less volume is required.

CHANGING THE SLEEVES

Grab a sleeve with your fingers and remove from the button, use fingers to insert new sleeve and press all the way.

CLEANING AND MAINTENANCE

Use a soft cloth dipped in ethanol or warm water to clean the nozzle ejector regularly.

PISTON MAINTENANCE

To access and lubricate the piston remove pipette nozzle by grabbing the pipette in your hand, press the ejector button all the way down, insert the spanner from the pipette ejector end and lever up the clamp locking the nozzle ejector to the release button. Gently remove the nozzle ejector whilst holding down the release button (8).

and release the pipette nozzle eject button (6) – ensure the spring does not fall out.

Insert the other end of the spanner onto the part uncovered by the removed nozzle ejector and turn the spanner anticlockwise to gently remove it.

Carefully remove the piston, paying attention to springs and the order in which the elements were removed, lubricate gently using silicone lubricant (ATTENTION! Excess lubricant may cause the piston to clog up), pay attention when re-inserting piston and screw the assembly together by turning the spanner clockwise. Then attach the release button, attach nozzle ejector, and assemble the device until the clamp locking the nozzle ejector to the release button locks.

CALIBRATION PROCEDURE

Once the piston has been greased the micropipette has to be calibrated. To do that insert the spanner under the collect button. Then gently turn the spanner anticlockwise (less volume) or clockwise (more volume) until the micropipette is calibrated correctly.

Remove spanner, the micropipette is ready for use.

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