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FEINWAAGE

## **USER MANUAL**

PRECISION SCALE

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BILANCIA DI PRECISIONE

## **MANUAL DE INSTRUCCIONES**

BALANZA DE PRECISIÓN

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PŘESNÁ VÁHA

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**SBS-LW-200N | SBS-LW-2000N | SBS-LW-3000N**

■ Deutsch	3
■ English	5
■ Polski	7
■ Česky	9
■ Français	11
■ Italiano	13
■ Español	15

NAZWA PRODUKTU PRODUKTNAME PRODUCT NAME NOM DU PRODUIT NOME DEL PRODOTTO NOMBRE DEL PRODUCTO NÁZEV VÝROBKU	WAGA PRECZYJNA FEINWAAGE PRECISION SCALE BALANCE DE PRÉCISION BILANCIA DI PRECISIONE BALANZA DE PRECISIÓN PŘESNÁ VÁHA
MODEL PRODUKTU MODELL PRODUCT MODEL MODÈLE MODELLO MODELO MODEL VÝROBKU	SBS-LW-200N SBS-LW-2000N SBS-LW-3000N
HERSTELLER MANUFACTURER PRODUCENT VÝROBCE FABRICANT PRODUTTORE FABRICANTE	EXPONDO POLSKA SP. Z O.O. SP. K.
ANSCHRIFT DES HERSTELLERS MANUFACTURER ADDRESS ADRES PRODUCTENTA ADRESA VÝROBCE ADRESSE DU FABRICANT INDIRIZZO DEL PRODUTTORE DIRECCIÓN DEL FABRICANTE	UL. NOWY KISIELIN-INNOWACYJNA 7, 66-002 ZIELONA GÓRA   POLAND, EU

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1. GENERAL DESCRIPTION

The user manual is designed to assist in the safe and trouble-free use of the device. The product is designed and manufactured in accordance with strict technical guidelines, using state-of-the-art technologies and components. Additionally, it is produced in compliance with the most stringent quality standards.

**DO NOT USE THE DEVICE UNLESS YOU HAVE THOROUGHLY READ AND UNDERSTOOD THIS USER MANUAL.**

To increase the product life of the device and to ensure a trouble-free operation, use it in accordance with this user manual and regularly perform the maintenance tasks. The technical data and specifications in this user manual are up to date. The manufacturer reserves the right to make changes associated with quality improvement. Taking into account technological progress and noise reduction opportunities, the device was designed to reduce noise emission risks to the minimum.

Legend

- The product satisfies the relevant safety standards.
- Read the instructions before use.
- The product must be recycled.

**PLEASE NOTE!** Drawings in this manual are for illustration purposes only and in some details may differ from the actual machine.

The original operation manual is in German. Other language versions are translations from German.

2. Usage safety

Electrical equipment:

**ATTENTION!** Read all safety warnings and all instructions. Failure to follow the warnings and instructions may result in an electric shock, fire and/or serious injury or even death.

Whenever „device“ or „product“ are used in the warnings and instructions, it shall mean Precision scale. Do not use in very humid environments or in the direct vicinity of water tanks. Prevent the device from getting wet. Risk of electric shock!

3.1 Electrical safety

- a) The plug has to fit the socket. Do not modify the plug in any way. Using original plugs and matching sockets reduces the risk of electric shock.
- b) Use the cable only in accordance with its designated use. Never use it to carry the device or to pull the plug out of a socket. Keep the cable away from heat sources, oil, sharp edges or moving parts. Damaged or tangled cables increase the risk of electric shock.

3.2 Safety at the workplace

- a) Upon discovering damage or irregular operation, immediately switch the device off and report it to a supervisor without delay.
- b) If there are any doubts as to the correct operation of the device, contact the manufacturer's support service.
- c) Only the manufacturer's service point may repair the device. Do not attempt any repairs independently!

3.3 Safe device use

- a) Device repair or maintenance should be carried out by qualified persons, only using original spare parts. This will ensure safe use.
- b) To ensure the operational integrity of the device, do not remove factory fitted guards and do not loosen any screws.
- c) When transporting and handling the device between the warehouse and the destination, take into account the occupational health and safety principles for manual transport operations which apply in the country where the device will be used.

4. Use guidelines

The precision scale is designed to weigh elements placed in the platter.

**The user is liable for any damage resulting from non-intended uses of the device!**

4.1 Device description

1. Weighing container
2. Platter
3. Tare/calibration button

4. Count button
5. Display

#### 4.2 Preparing for use

Place the device on a flat and level working surface, which will be at least as large as the device. The balance should be used in a place free from wind, corrosion, vibration, humidity and temperature variations.

#### 4.3 Device use

Use guidelines:

- Do not weigh items which are too heavy. If „OL” is displayed, this indicates that the item on the scale is too heavy; remove it at once to avoid damaging the scale.
- Error1 message appears on the display if the items on the scale are vibrating and the vibrations are interfering with the weighing.
- Error2 message appears if the load is unstable (not necessarily vibrating). Ensure the item to be weighed is stable on the scale.
- Do NOT leave load on the scale for long.
- Avoid extreme temperatures. Do not place the unit in direct sunlight or near the air conditioning.
- Avoid using strong electrical or magnetic fields during use.
- Make sure the scale is placed on a stable table and is not subject to vibrations.
- Do not place the unit close to windows, doors or fans that may cause unstable results due to air currents.
- Keep your machine clean. Do not store products on the scale when it is not used.

#### ON/OFF switch (I/O)

Use the main I/O switch on the back of the device to turn the scale on. Switching the device on initiates a self-test. The self-test begins with a display test. A correctly displayed series of digits (8.8.8.8.8.8.) has to appear on the display followed by the maximum acceptable weight (3000g/2000g/200g). Once the test is complete, the scale is ready to for use. The display will show „0”. To turn the scale off press the I/O button.

#### TARE

Put an empty container in the scale bowl and wait for the display to settle. Press the [TAR/CAL] button. Zero should appear on the display.

#### ZEROING

If the display is showing a value other than 0 when the scale bowl is empty, reset it by pressing the [TAR/CAL] button.

#### COUNTING PIECES

To enter count mode, press the [COU/AFF] button. Place a pre-counted number of items on the scale. Use the [TAR/CAL] button to select the number of items on the scale (10, 20, 50, 100, 250, 500). Press [COU/AFF] to confirm selection. The device is ready for use.

#### CALIBRATION

Switch the scale on with the empty platter. Wait for the scale to stabilise (up to approx. 15 min), press and hold the [TAR/CAL] button. A flashing „100.000” will appear on the display. Place a 100g weight on the scale. „-----”, and then „100.000” will appear on the display. Calibration is complete.

#### 4.4 Cleaning and maintenance

- Always unplug the device before cleaning it.
- Use cleaners without corrosive substances to clean each surface.
- It is prohibited to clean the scale with water, in order not to damage electronic components.
- Store the unit in a dry, cool place, free from moisture and direct exposure to sunlight

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