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USER MANUAL

OVERHEAD STIRRER

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MANUAL DE INSTRUCCIONES

AGITADOR DIGITAL DE VARILLA

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SBS-MR-2500

Deutsch	3
English	8
Polski	12
Česky	17
Français	21
Italiano	26
Español	31

PRODUKTNAME	DIGITALER ÜBERKOPFRÜHRER
PRODUCT NAME	OVERHEAD STIRRER
NAZWA PRODUKTU	MIESZADŁO LABORATORYJNE
NÁZEV VÝROBKU	HŘÍDELOVÉ MÍCHADLO
NOM DU PRODUIT	AGITATEUR À HÉLICE NUMÉRIQUE
NOME DEL PRODOTTO	AGITATORE DIGITALE AD ASTA
NOMBRE DEL PRODUCTO	AGITADOR DIGITAL DE VARILLA
MODELL	SBS-MR-2500
PRODUCT MODEL	
MODEL PRODUKTU	
MODEL VÝROBKU	
MODÈLE	
MODELLO	EXPONDO POLSKA SP. Z O.O. SP. K.
MODELO	
IMPORTEUR	
IMPORTER	
IMPORTER	
DOVOZCE	UL. NOWY KISIELIN-INNOWACYJNA 7, 66-002 ZIELONA GÓRA POLAND, EU
IMPORTATEUR	
IMPORTATORE	
IMPORTADOR	
ADRESSE VON IMPORTEUR	
IMPORTER ADDRESS	
ADRES IMPORTERA	
ADRESA DOVOZCE	
ADRESSE DE L'IMPORTATEUR	
INDIRIZZO DELL'IMPORTATORE	
DIRECCIÓN DEL IMPORTADOR	

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1. GENERAL DESCRIPTION

The user manual is designed to assist in the safe and trouble-free use of the device. The product is designed and manufactured in accordance with strict technical guidelines, using state-of-the-art technologies and components. Additionally, it is produced in compliance with the most stringent quality standards.

DO NOT USE THE DEVICE UNLESS YOU HAVE THOROUGHLY READ AND UNDERSTOOD THIS USER MANUAL

To increase the product life of the device and to ensure trouble-free operation, use it in accordance with this user manual and regularly perform maintenance tasks. The technical data and specifications in this user manual are up to date. The manufacturer reserves the right to make changes associated with quality improvement. The device is designed to reduce noise emission risks to a minimum, taking into account technological progress and noise reduction opportunities.

LEGEND

	The product satisfies the relevant safety standards.
	Read instructions before use.
	The product must be recycled.
	WARNING! or CAUTION! or REMEMBER! Applicable to the given situation (general warning sign).
	ATTENTION! Electric shock warning!
	ATTENTION! Hot surface, risk of burns!
	Connect the device to a grounded power supply.
	Do not use the device to experiment with flammable, explosive, poisonous and highly caustic substances.
	Only use indoors.

PLEASE NOTE! Drawings in this manual are for illustration purposes only and in some details may differ from the actual product. The original operation manual is written in German. Other language versions are translations from the German.

2. USAGE SAFETY

ATTENTION! Read all safety warnings and all instructions. Failure to follow the warnings and instructions may result in electric shock, fire and/or serious injury or even death.

The terms "device" or "product" are used in the warnings and instructions to refer to < Ceramic Magnetic Stirrer >. Do not use in very humid environments or in the direct vicinity of water tanks. Prevent the device from getting wet. Risk of electric shock! Do not cover the ventilation openings!

2.1. ELECTRICAL SAFETY

- The plug must fit the socket. Do not modify the plug in any way. Using original plugs and matching sockets reduces the risk of electric shock.
- Avoid touching earthed elements such as pipes, heaters, boilers and refrigerators. There is an increased risk of electric shock if the earthed device is exposed to rain, comes into direct contact with a wet surface or is operating in a damp environment. Water getting into the device increases the risk of damage to the device and of electric shock.
- Do not touch the device with wet or damp hands.
- Use the cable only for its designated use. Never use it to carry the device or to pull the plug out of a socket. Keep the cable away from heat sources, oil, sharp edges or moving parts. Damaged or tangled cables increase the risk of electric shock.
- If using the device in a damp environment cannot be avoided, a residual current device (RCD) should be applied. The use of an RCD reduces the risk of electric shock.
- Do not use the device if the power cord is damaged or shows obvious signs of wear. A damaged power cord should be replaced by a qualified electrician or the manufacturer's service centre.
- To avoid electric shock, do not immerse the cord, plug or device in water or other liquids. Do not use the device on wet surfaces.

2.2. SAFETY IN THE WORKPLACE

- Make sure the workplace is clean and well lit. A messy or poorly lit workplace may lead to accidents. Try to think ahead, observe what is going on and use common sense when working with the device.
- Do not use the device in a potentially explosive environment, for example in the presence of flammable liquids, gases or dust. The device generates sparks which may ignite dust or fumes.
- If you discover damage or irregular operation, immediately switch the device off and report it to a supervisor without delay.
- If there are any doubts as to the correct operation of the device, contact the manufacturer's support service.
- Only the manufacturer's service point may repair the device. Do not attempt any repairs independently!
- In case of fire, use a powder or carbon dioxide (CO₂) fire extinguisher (one intended for use on live electrical devices) to put it out.
- Please keep this manual available for future reference. If this device is passed on to a third party, the manual must be passed on with it.

- Keep packaging elements and small assembly parts in a place not available to children.
- Keep the device away from children and animals.

REMEMBER! When using the device, protect children and other bystanders.

2.3. PERSONAL SAFETY

- Do not use the device when tired, ill or under the influence of alcohol, narcotics or medication which can significantly impair the ability to operate the device.
- The device is not designed to be handled by persons (including children) with limited mental and sensory functions or persons lacking relevant experience and/or knowledge unless they are supervised by a person responsible for their safety or they have received instruction on how to operate the device.
- When working with the device, wear personal protective equipment that prevents:
 - splashing or contact with evaporating liquid,
 - contact with toxic or flammable gases.
- To prevent accidental starting, make sure the knobs are in off position before connecting to the power source.
- The device is not a toy. Children must be supervised to ensure that they do not play with the device.

2.4. SAFE DEVICE USE

- Do not use the device if the knobs or control buttons do not function properly. Devices that cannot be controlled by the controls are dangerous, may not be operated and must be repaired.
- Disconnect the device from the power supply before commencement of adjustment, cleaning and maintenance. Such a preventive measure reduces the risk of accidental activation.
- When not in use, store in a safe place, away from children and people not familiar with the device who have not read the user manual. The device may pose a hazard in the hands of inexperienced users.
- Keep the device in perfect technical condition. Before each use check for general damage and especially check for cracked parts or elements and for any other conditions which may impact the safe operation of the device. If damage is discovered, hand over the device for repair before use.
- Keep the device out of the reach of children.
- Device repair or maintenance should be carried out by qualified persons, only using original spare parts. This will ensure safe use.
- To ensure the operational integrity of the device, do not remove factory-fitted guards and do not loosen any screws.
- Do not move, adjust or rotate the device in the course of work.
- Do not leave this appliance unattended while it is in use.
- Clean the device regularly to prevent stubborn grime from accumulating.
- The device is not a toy. Cleaning and maintenance may not be carried out by children without supervision by an adult person.
- It is forbidden to interfere with the structure of the device in order to change its parameters or construction.
- Keep the device away from sources of fire and heat.
- Control the amount of liquid in containers. The liquid may evaporate.

- Spilling the flammable contents of the container onto the heating elements of the device may cause a fire.
- Special attention must be paid to the hazards arising from:
 - Overfilled containers
 - Containers made of brittle material or of low melting point
- Keep the device away from strong magnetic fields.
- When mixing viscous liquids reduce the rotational speed and mix a small volume of liquid only.
- The device can work continuously for 8 hours at moderate speed and 4 hours at high mixing speed

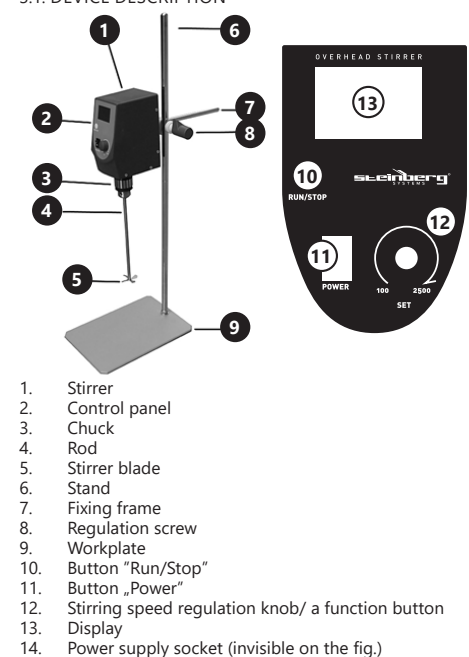
ATTENTION! Despite the safe design of the device and its protective features, and despite the use of additional elements protecting the operator, there is still a slight risk of accident or injury when using the device. Stay alert and use common sense when using the device.

3. USE GUIDELINES

The device is designed for use in laboratories, schools, universities, and similar institutions. It is used for non-contact stirring and for heating liquids in laboratory vessels. The product is outside of the scope of MDD 93/42/EEC Medical Devices regulations. It is not a medical device and should not be used for medical research and tests in hospitals and other health facilities.

The user is liable for any damage resulting from unintended use of the device.

3.1. DEVICE DESCRIPTION

3.1. PREPARING FOR USE

APPLIANCE LOCATION

Ensure good ventilation in the room in which the device is being used. There should be at least 10 cm distance between each side of the device and the wall or other objects.

The device should always be used when positioned on an even, stable, clean, fireproof and dry surface, and be out of the reach of children and persons with limited mental and sensory functions. Position the device such that you always have access to the power plug. The power cord connected to the appliance must be properly grounded and correspond to the technical details on the product label.

APPLIANCE LOCATION

- Unpack the device, check if the package contains the complete set of undamaged elements.
- Place the stirring plate on a stable and level surface.
- Assemble the two elements of the stand.
- Screw a ready stand to the stirring plate.
- Screw the fixing frame to a rear part of the main body of the stirrer.
- Assemble the regulation screw to the stand.
- Put the fixing frame into a regulation screw clamp.
- Place the stirrer rod in the chuck and tighten it.
- Connect a power cord to an adapter.
- Connect the adapter to the power supply and its socket.

3.2. DEVICE USE

- Place a vessel containing liquid under the main body of the overhead stirrer so that the stirrer blade is inside the vessel.
- Turn „Power“ button to „I“ position.
- Press „Run/Stop“ button to start stirring.
- Press and turn clockwise the regulation knob to increase stirrer blade revolutions.
- Press and turn anticlockwise the regulation knob to decrease stirrer blade revolutions.
- Press „Run/Stop“ button to stop the blade.
- Turn the device off by turning „Power“ button to „0“ position.

- 3.3.1. CAUTION: The device is equipped with a control panel for internal parameters presented in the table below. However, these settings should be changed only by a person with proper knowledge and qualifications.
- Press and hold for three seconds a function button [12] to operate the control panel. LC menu for selecting a function will be displayed. A user should comply with the table to change stirrer parameters.
 - Confirm a choosen code in LC menu by choosing the value with the speed regulation knob and pressing the function button.
 - In the menu, a user can change parameters by pressing the function button.
 - Parameters value is changed by turning the speed regulation knob.
 - Press and hold for three seconds the function button to back to an initial menu.

Message	Function name	Function description	(setting range) Factory value
Lc	Menu number	Selecting menu	3
Pd	Proportionation	Proportionation regulation	(1~100) 10
Id	Integration	Integration regulation	(1~100) 5

InT	Acceleration time	0-maximum speed time	(1~60) 10
dET	Decreasing speed time	maximum speed-0 time	(1~60) 10
SdL	revolutions minimum value	minimum speed	(20~6000) 20
SdH	revolutions maximum value	maximum speed	(20~6000) 2000
Message	Function name	Function description	(setting range) Factory value
Lc	Menu number	Selecting menu	9
Ear	Gearing	Gear/pinion diameter	(1,0~10,0) 1,0
PoL	Motor poles	Motor poles	(1~32) 4
InT	reference direction of the motor	0: clockwise revolutions 1: anticlockwise revolutions	(0~1) 0
FdS	Speed feedback	Velocity feedback	(1,0~10,0) 1,0
FdC	Amperage feedback		(1,0~10,0) 1,0
FrE	Carrier frequency	PWM carrier frequency	(5~15) 15
Po	Motor power	Motor power	(1~750) 40
CL	Excessive amperage multiples	Current multiples authorization	(1,0~10,0) 5,0
Fr	Motor revolutions direction	0: The motor works only forward 1: The motor work only backward 2: The motor works forward and backward	(0~2) 0
db	Speed dead zone	Speed dead zone	(0~100) 5
Message	Function name	Function description	(setting range) Factory value
Lc	Menu number	Selecting menu	101
Pd	Zero correction	Update a measurement error	(-99,9 ~99,9) 0,0
PK	Full correction	Update a measurement error PK= 1000x a measured actual value	(-999~999) 0

dot	Decimal display	0: Turn decimal display off 1: Turn decimal display on	(0~1) 1
Message	Function name	Function description	(setting range) Factory value
Lc	Menu number	Selecting menu	103
ET	Time display	0: Turn a time display off 1: Turn a time display on	(0~1) 0
Hn	Synchronization mode	0: minute-based measurement 1: hour-based measurement	(0~1) 0
rT	Time correction	Correction value=[current working time (a second) – actual time (a second)]* 10/ actual time (a minute)	(-999~999) 0

Message	Function name	Function description	(setting range) Factory value
Lc	Menu number	Selecting menu	110
rSET	Default settings	Restoring default settings	(0~1) 0

Warning: When an error occurs, Er (1-6) will be displayed, then the laboratory stirrer should be switched off with the "Power" button and then restarted.

3.3.2 TIME SETTING MENU

- After starting the time display (ET function) additional functions will appear on the rotation frequency change screen:
- Increase and decrease the function parameters by turning the speed control knob.

Message	Function name	Function description
ST	Counter	Setting time, after which the stirrer will turn off.
FT		Setting a displayed time value
PT		Stop a displayed time value
rT		Setting a reverse time value

3.4. CLEANING AND MAINTENANCE

- Unplug the mains plug and allow the device to cool completely before each cleaning, adjustment or replacement of accessories, or if the device is not being used.
- Use only non-corrosive cleaners to clean the surface.
- After cleaning the device, all parts should be dried completely before using it again.
- Store the unit in a dry, cool place, free from moisture and direct exposure to sunlight.
- Do not spray the device with a water jet or submerge it in water.
- Do not allow water to get inside the device through vents in the housing of the device.
- Clean the vents with a brush and compressed air.
- The device must be regularly inspected to check its technical efficiency and spot any damage.

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