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TISCHWAAGE

USER MANUAL

TABLE SCALE

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WAGA STOŁOWA

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STOLNÍ VÁHA

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BALANCE DE TABLE

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BILANCIA DA TAVOLO

MANUAL DE INSTRUCCIONES

BALANZA DE MESA

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DE | EN | PL | CZ | FR | IT | ES | HU

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1. GENERAL DESCRIPTION

The user manual is designed to assist in the safe and trouble-free use of the device. The product is designed and manufactured in accordance with strict technical guidelines, using state-of-the-art technologies and components. Additionally, it is produced in compliance with the most stringent quality standards.

DO NOT USE THE DEVICE UNLESS YOU HAVE THOROUGHLY READ AND UNDERSTOOD THIS USER MANUAL.

To increase the product life of the device and to ensure trouble-free operation, use it in accordance with this user manual and regularly perform maintenance tasks.

The technical data and specifications in this user manual are up to date. The manufacturer reserves the right to make changes associated with quality improvement.

LEGEND

	The product satisfies the relevant safety standards.
	Read instructions before use.
	The product must be recycled.
	WARNING! or CAUTION! or REMEMBER! Applicable to the given situation (general warning sign).
	Only use indoors.

PLEASE NOTE! Drawings in this manual are for illustration purposes only and in some details may differ from the actual machine.

The original operation manual is in German. Other language versions are translations from German.

2. USAGE SAFETY

ATTENTION! Read all safety warnings and all instructions. Failure to follow the warnings and instructions may result in an electric shock, fire and/or serious injury or even death.

The terms „device“ or „product“ are used in the warnings and instructions to refer to <Table scale>. Prevent the device from getting wet. Risk of electric shock! Do not cover the ventilation openings!

2.1. ELECTRICAL SAFETY

For battery charger

- a) The plug must fit the socket. Do not modify the plug in any way. Using original plugs and matching sockets reduces the risk of electric shock.
- b) Avoid touching earthed elements such as pipes, heaters, boilers and refrigerators. There is an increased risk of electric shock if the earthed device is exposed to rain, comes into direct contact with a wet surface or is operating in a damp environment. Water getting into the device increases the risk of damage to the device and of electric shock.
- c) Do not touch the device with wet or damp hands.
- d) Use the cable only for its designated use. Never use it to carry the device or to pull the plug out of a socket. Keep the cable away from heat sources, oil, sharp edges or moving parts. Damaged or tangled cables increase the risk of electric shock.
- e) If using the device in a damp environment cannot be avoided, a residual current device (RCD) should be applied. The use of an RCD reduces the risk of electric shock.
- f) Do not use the device if the power cord is damaged or shows obvious signs of wear. A damaged power cord should be replaced by a qualified electrician or the manufacturer's service centre.
- g) **ATTENTION! DANGER TO LIFE!** While cleaning, never immerse the device in water or other liquids.

2.2. SAFETY IN THE WORKPLACE

- a) If there are any doubts as to the correct operation of the device, contact the manufacturer's support service.

- b) Only the manufacturer's service point may repair the device. Do not attempt any repairs independently!
- c) In case of fire, use a powder or carbon dioxide (CO₂) fire extinguisher (one intended for use on live electrical devices) to put it out.
- d) Regularly inspect the condition of the safety labels. If the labels are illegible, they must be replaced.
- e) Please keep this manual available for future reference. If this device is passed on to a third party, the manual must be passed on with it.
- f) Keep packaging elements and small assembly parts in a place not available to children.
- g) Keep the device away from children and animals.

⚠ REMEMBER When using the device, protect children and other bystanders.

2.3. PERSONAL SAFETY

- a) The device is not designed to be handled by persons (including children) with limited mental and sensory functions or persons lacking relevant experience and/or knowledge unless they are supervised by a person responsible for their safety or they have received instruction on how to operate the device.
- b) The device is not a toy. Children must be supervised to ensure that they do not play with the device.

2.4. SAFE DEVICE USE

- a) Do not overload the device. A correctly-selected device will perform the task for which it was designed better and in a safer manner.
- b) Do not use the device if the ON/OFF switch does not function properly (does not switch the device on and off). Devices which cannot be switched on and off using the ON/OFF switch are hazardous, should not be operated and must be repaired.
- c) When not in use, store in a safe place, away from children and people not familiar with the device who have not read the user manual. The device may pose a hazard in the hands of inexperienced users.
- d) Keep the device in perfect technical condition. Before each use check for general damage and especially check for cracked parts or elements and for any other conditions which may impact the safe operation of the device. If damage is discovered, hand over the device for repair before use.
- e) Keep the device out of the reach of children.
- f) Device repair or maintenance should be carried out by qualified persons, only using original spare parts. This will ensure safe use.
- g) To ensure the operational integrity of the device, do not remove factory-fitted guards and do not loosen any screws.
- h) Avoid situations where the device stops working during use due to excessive loading.
- i) Do not move, adjust or rotate the device in the course of work.
- j) Clean the device regularly to prevent stubborn grime from accumulating.
- k) The device is not a toy. Cleaning and maintenance may not be carried out by children without supervision by an adult person.
- l) It is forbidden to interfere with the structure of the device in order to change its parameters or construction.
- m) Keep the device away from sources of fire and heat.
- n) Do not overload the scale. This can damage the load cells.

- o) Before making a measurement, make sure that the device is placed on a stable surface and is not exposed to vibrations and sudden temperature changes, which may adversely affect the measurement results.
- p) Do NOT leave load on the scale for long. This will decrease the scale's accuracy and shorten the load cell's life.
- q) Do NOT use the scale in thunder or rain.
- r) Avoid extreme temperatures. Do not place the unit in direct sunlight or near the air conditioning.
- s) Do not use the scale close to large electrical appliances such as welding machines or large motors. Do not leave batteries in the device if you do not use it for a long time.
- t) Avoid places where moisture can lead to condensation. Avoid direct contact with water.
- u) Do not place the unit close to windows, doors or fans that may cause unstable results due to air currents.
- v) Keep your machine clean. Do not store products on the scale when it is not used.

⚠ ATTENTION! Despite the safe design of the device and its protective features, and despite the use of additional elements protecting the operator, there is still a slight risk of accident or injury when using the device. Stay alert and use common sense when using the device.

3. USE GUIDELINES

The device is intended for measuring the weight of objects placed on the device's weighing platform.

The user is liable for any damage resulting from unintended use of the device.

3.1. DEVICE DESCRIPTION

1. Display
2. Weighing platform
3. Control panel
4. Main body
5. Foot
6. Charging indicator light
7. Power supply socket (not visible in the picture)
8. Spirit level bubble (not visible in the picture)
9. ON / OFF switch (not visible in the picture)
10. RS-232 port (not visible in the picture)

BUTTON	DESCRIPTION
MODE	Selecting function
SET	Setting limit values
↑	Entering a numeric value
CALI	Weight calibration
UNIT	Choice of measurement units

ZERO	Scale zeroing / backlight
ENTER	Confirming selection
TARE	Tare / Delete tare

3.2. PREPARING FOR USE APPLIANCE LOCATION

The temperature of environment must not be higher than 40°C and the relative humidity should be less than 85%. The device should always be used when positioned on an even, stable, clean, fireproof and dry surface, and be out of the reach of children and persons with limited mental and sensory functions. The power cord connected to the appliance must be properly grounded and correspond to the technical details on the product label.

3.3. DEVICE USE

NOTE: Before using the device, make sure that it is properly levelled.

3.3.1. USING THE DEVICE

NOTE: After starting the device, the scale will automatically go into weighing mode. „0“ and „kg“ will appear on the [1] display. To select a different operating mode, press the [MODE] button.

NOTE: To turn the display backlight on or off, press and hold the [ZERO / BL] button for a few seconds.

1. Weighing mode

Weighing an object.

- a) Change of units:
 - Press the [UNIT] button to change the unit of measurement kg > lb.
- b) Tare acquisition:
 - Place the container or weight on the [2] weighing platform.
 - Wait for a while until the device stabilizes.
 - Press the [TARE] button, the „1“ display will show the „0“ value and the „TARE“ message. After removing the weight, a minus value will appear on the [1] display.
 - Press the [TARE] button again to end tare acquisition.
- c) Reset:
 - Set the weight on the [2] weighing platform.
 - To zero the measurement value, press the [ZERO] button.

2. Counting mode

- Counting the pieces according to the single piece weight.
- a) Press the [SET] button to set the number of samples. The [1] display will show „SAP X“ (where X is the number of samples).
 - b) Press the [1] button to change the number of samples. Possible choices are 10, 20, 50, 100, 200.
 - c) Confirm the selection by pressing [ENTER].
 - d) „LOAD“ will appear on the [1] display. Place the selected number of samples, e.g. 10, on the [2] weighing platform and confirm by pressing [ENTER].
 - e) Place the total number of samples to be counted on the [2] weighing platform. The device will count the number of samples, the measurement result will appear on the [1] display.
 - f) Possible inaccuracy messages:
 - „-LAC-“ appears when the sample unit weight is less than 80% of the scale of the device. Counting the number of samples is possible in this case, but may be inaccurate.

- „-CSL-“ appears when the sample unit weight is insufficient to perform the counting operation.

3. Percentage mode

Counting the pieces according to the percentage value of the weighed objects.

- a) Press the [SET] button. „LOAD-P“ will appear on the [1] display. Put the samples on the [2] weighing platform. Press the [ENTER] button to make the measurement or
- b) press the [UNIT] button. „000000“ will appear on the [1] display. Press the [1] and the [ENTER] button to enter the weight.
- c) Possible inaccuracy messages:
 - „-CSL-“ appears when the sample unit weight is insufficient, i.e. it does not exceed 0.1% FS

4. Setting the weight alarm

- a) High alarm
 - Press the [SET] button, the display will show „-HH-“.
 - Confirm your selection by pressing [ENTER]
 - Press the [1] button to set the first digit.
 - Confirm with the [ENTER] button and repeat the steps until the total desired value is set.
- b) Low alarm
 - Press the [SET] button, the display will show „-LL-“.
 - Proceed in the same way as for high weight alarm.
- c) Signalling modes
 - The signalling mode can be modified using the [1] button.
 - „IN“ - the signal is activated when the measured value is within the accepted scale (between the „LL“ and „HH“ value).
 - „OUT“ - the alarm is activated when the measured value is outside the accepted scale (below the „LL“ or above the „HH“).
 - „NO“ - the alarm is deactivated.

3.3.2 MESSAGES

The [1] display shows the following messages:

1. „-OF-“ or „-ADC-“ and a continuous beep. This means that the load on the [2] weighing platform has exceeded the measuring scale of the device. Remove the load from the weighing platform.
2. „-Lo-“ means there is no load on the weighing platform. This means that the battery is almost flat. Connect the power supply.
3. „HHHH“ or „LLLL“ means the value exceeds the allowable measuring range during zeroing.
4. „UNSTA“ means the balance is not stabilized.
5. „-SYS-“ means the scale must be recalibrated or needs servicing.
6. „-SET-“ means incorrect alarm settings. The alarm must be reset.

3.3.3 RS-232 TRANSMISSION

1. To enter the RS-232 settings, press and hold [MODE] and [TARE] until a beep sounds.
 - a) Transmission method
 - Press the [SET] button to select the transmission method:
 - A. „Stb“ - when the balance is stabilized,
 - B. „Etb“ - after pressing the [ENTER] button,
 - C. „Ser“ - continuous transmission,
 - D. „CLOSE“ - transmission disabled,

- Press the [1] button to change the current transmission method. The method is selected in the following order:

- Press [ENTER] to confirm the selection and move to the next settings menu.

b) Transmission speed

- Press the [1] button to select the transmission speed. The selection is made in the following order:

- Press [ENTER] to confirm the selection.

2. RS-232 interface

- The balance has a standard RS-232 UART signal, 10bit.
- Data format:
 - 1: D0 ~ D7 - 0FFH
 - 2: D0 ~ D2 - decimal point
D3 ~ D4 - current mode: 00 (weighing), 01 (counting), 10 (percent)
D5 - 1 (negative weight), other (positive)
D6 - 1 (stable), other (unstable)
D7 - 2 (excessive weight), other (normal)
 - 3: D0 ~ D7 - BCD1 (LSB)
 - 4: D0 ~ D7 - BCD2 (MSB)
 - 5: D0 ~ D7 - BCD3 (HSB)
 - 6: D0 ~ D7 - weight units: kg, lb

3.3.4 DEVICE CALIBRATION

- Turn the the device on with the [9] switch.
- Press and hold the [CALI] button until „CALI” appears on the display.
 - Method I:
 - Prepare a weight equal to the maximum load of the scale.
 - Place the weight on the [2] weighing platform.
 - Confirm the displayed value by pressing [ENTER].
 - Method II:
 - Prepare the weight.
 - Press the [SET] button. The device will display the last calibration value.
 - Press the [1] button and the [ENTER] button to enter the desired value equal to the mass of the prepared weight.
 - Place the weight on the [2] weighing platform.
 - Confirm the displayed value by pressing [ENTER].

Note: The maximum and minimum weight values of the device are factory programmed. The value entered using the [1] button will not exceed the maximum weight value.

3.4. CLEANING AND MAINTENANCE

- Use only non-corrosive cleaners to clean the surface.
- After cleaning the device, all parts should be dried completely before using it again.
- Store the unit in a dry, cool place, free from moisture and direct exposure to sunlight.
- Do not allow water to get inside the device through vents in the housing of the device.

- The device must be regularly inspected to check its technical efficiency and spot any damage.
- Do not leave the battery in the device if it will not be used for a longer period of time.
- Use a soft, damp cloth for cleaning.

SAFE REMOVAL OF BATTERIES AND RECHARGEABLE BATTERIES

The devices are equipped with 9V / 300mA batteries. Recycle batteries with the appropriate organisation or company.

DISPOSING OF USED DEVICES

Do not dispose of this device in municipal waste systems. Hand it over to an electric and electrical device recycling and collection point. Check the symbol on the product, instruction manual and packaging. The plastics used to construct the device can be recycled in accordance with their markings. By choosing to recycle you are making a significant contribution to the protection of our environment. Contact local authorities for information on your local recycling facility.

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