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USER MANUAL

DIGITAL PRECISION SCALE

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PŘESNÁ VÁHA

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BILANCIA DI PRECISIONE

MANUAL DE INSTRUCCIONES

BALANZA DE PRECISIÓN

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SBS-LW-300001E | SBS-LW-600001E

INHALT | CONTENT | TREŠĆ | OBSAH | CONTENU | CONTENUTO | CONTENIDO

▪ Deutsch	3
▪ English	7
▪ Polski	11
▪ Český	15
▪ Français	19
▪ Italiano	23
▪ Español	27

PRODUKTNAME PRODUCT NAME NAZWA PRODUKTU NÁZEV VÝROBKU NOM DU PRODUIT NOME DEL PRODOTTO NOMBRE DEL PRODUCTO	PRÄZISIONSWAAGE DIGITAL PRECISION SCALE WAGA LABORATORYJNA PŘESNÁ VÁHA BALANCE DE PRÉCISION BILANCIA DI PRECISIONE BALANZA DE PRECISSION
MODELL PRODUCT MODEL MODEL PRODUKTU MODEL VÝROBKU MODÈLE MODELLO MODELO	SBS-LW-300001E SBS-LW-600001E
HERSTELLER MANUFACTURER PRODUCENT VÝROBCE FABRICANT PRODUTTORE FABRICANTE	EXPONDO POLSKA SP. Z O.O. SP. K.
ANSCHRIFT DES HERSTELLERS MANUFACTURER ADDRESS ADRES PRODUCENTA ADRESA VÝROBCE ADRESSE DU FABRICANT INDIRIZZO DEL PRODUTTORE DIRECCIÓN DEL FABRICANTE	UL. NOWY KISIELIN-INNOWACYJNA 7, 66-002 ZIELONA GÓRA POLAND, EU

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1.GENERAL DESCRIPTION

The user manual is designed to assist in the safe and trouble-free use of the device. The product is designed and manufactured in accordance with strict technical guidelines, using state-of-the-art technologies and components. Additionally, it is produced in compliance with the most stringent quality standards.

DO NOT USE THE DEVICE UNLESS YOU HAVE THOROUGHLY READ AND UNDERSTOOD THIS USER MANUAL.

To increase the product life of the device and to ensure trouble-free operation, use it in accordance with this user manual and regularly perform maintenance tasks. The technical data and specifications in this user manual are up to date. The manufacturer reserves the right to make changes associated with quality improvement.

LEGEND

- The product satisfies the relevant safety standards.
- Read instructions before use.

- The product must be recycled.
- WARNING! or CAUTION! or REMEMBER!** Applicable to the given situation. (general warning sign)
- ATTENTION! Electric shock warning!**

PLEASE NOTE! Drawings in this manual are for illustration purposes only and in some details may differ from the actual product. The original operation manual is written in German. Other language versions are translations from the German.

2. USAGE SAFETY

ATTENTION! Read all safety warnings and all instructions. Failure to follow the warnings and instructions may result in electric shock, fire and/or serious injury or even death.

The terms „device“ or „product“ are used in the warnings and instructions to refer to < Digital Precision Scale >. Do not use in very humid environments or in the direct vicinity of water tanks. Do not cover the ventilation openings!

2.1. ELECTRICAL SAFETY

- The plug must fit the socket. Do not modify the plug in any way. Using original plugs and matching sockets reduces the risk of electric shock.
- Avoid touching earthed elements such as pipes, heaters, boilers and refrigerators. There is an increased risk of electric shock if the earthed device is exposed to rain, comes into direct contact with a wet surface or is operating in a damp environment. Water getting into the device increases the risk of damage to the device and of electric shock.
- Do not touch the device with wet or damp hands.
- Use the cable only for its designated use. Never use it to carry the device or to pull the plug out of a socket. Keep the cable away from heat sources, oil, sharp edges or moving parts. Damaged or tangled cables increase the risk of electric shock.
- If using the device in a damp environment cannot be avoided, a residual current device (RCD) should be applied. The use of an RCD reduces the risk of electric shock.
- Do not use the device if the power cord is damaged or shows obvious signs of wear. A damaged power cord should be replaced by a qualified electrician or the manufacturer's service centre.
- To avoid electric shock, do not immerse the cord, plug or device in water or other liquids. Do not use the device on wet surfaces.

2.2. SAFETY IN THE WORKPLACE

- Do not use the device in a potentially explosive environment, for example in the presence of flammable liquids, gases or dust. The device generates sparks which may ignite dust or fumes.
- If you discover damage or irregular operation, immediately switch the device off and report it to a supervisor without delay.
- If there are any doubts as to the correct operation of the device, contact the manufacturer's support service.
- Only the manufacturer's service point may repair the device. Do not attempt any repairs independently!
- In case of fire, use a powder or carbon dioxide (CO₂) fire extinguisher (one intended for use on live electrical devices) to put it out.

- f) Regularly inspect the condition of the safety labels. If the labels are illegible, they must be replaced.
- g) Please keep this manual available for future reference. If this device is passed on to a third party, the manual must be passed on with it.
- h) Keep packaging elements and small assembly parts in a place not available to children.

REMEMBER! When using the device, protect children and other bystanders.

2.3. PERSONAL SAFETY

- a) Do not use the device when tired, ill or under the influence of alcohol, narcotics or medication which can significantly impair the ability to operate the device.
- b) The device is not designed to be handled by persons (including children) with limited mental and sensory functions or persons lacking relevant experience and/or knowledge unless they are supervised by a person responsible for their safety or they have received instruction on how to operate the device.
- c) To prevent the device from accidentally switching on, make sure the switch is on the OFF position before connecting to a power source.
- d) The device is not a toy. Children must be supervised to ensure that they do not play with the device.

2.4. SAFE DEVICE USE

- a) Do not use the device if the ON/OFF switch does not function properly (does not switch the device on and off). Devices which cannot be switched on and off using the ON/OFF switch are hazardous, should not be operated and must be repaired.
- b) Disconnect the device from the power supply before commencement of adjustment, cleaning and maintenance. Such a preventive measure reduces the risk of accidental activation.
- c) When not in use, store in a safe place, away from children and people not familiar with the device who have not read the user manual. The device may pose a hazard in the hands of inexperienced users.
- d) Keep the device out of the reach of children.
- e) Device repair or maintenance should be carried out by qualified persons, only using original spare parts. This will ensure safe use.
- f) To ensure the operational integrity of the device, do not remove factory-fitted guards and do not loosen any screws.
- g) Clean the device regularly to prevent stubborn grime from accumulating.
- h) The device is not a toy. Cleaning and maintenance may not be carried out by children without supervision by an adult person.
- i) It is forbidden to interfere with the structure of the device in order to change its parameters or construction.
- j) Keep the device away from sources of fire and heat.

ATTENTION! Despite the safe design of the device and its protective features, and despite the use of additional elements protecting the operator, there is still a slight risk of accident or injury when using the device. Stay alert and use common sense when using the device.

3. USE GUIDELINES

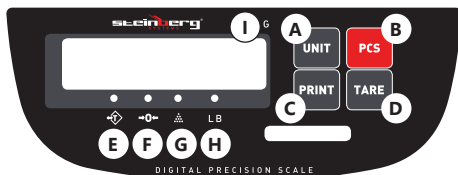
The device is intended to weigh items placed on the weighing pan. Warning! The device cannot be used for commercial purposes. **The user is liable for any damage resulting from unintended use of the device.**

The balance should be used in a space free of wind, corrosion, vibration, humidity and temperature variations. The scale should avoid vibration, violent air current, direct sunshine and high temperature.

3.1. DEVICE DESCRIPTION

- 1 Weighing pan
- 2 Control panel
- 3 RS-232 interface
- 4 I/O (ON/ OFF) button
- 5 Power supply socket
- 6 Bubble level

The keys:

- A. UNIT: Unit selection
- B. PRINT: Printing function.
- C. PCS: Counting function.
- D. TARE: Tare weight of container.

LEDs indicating function:

- E. Tare
- F. Zero status
- G. Counting
- H. Weight measurement unit: lb
- I. Weight measurement unit: g

3.2. PREPARING FOR USE LEVELLING THE DEVICE

The scale is equipped with a level indicator under the weighing pan and four adjustable legs. Before using the device, adjust the length of the legs until the bubble of the level indicator is in the middle position.

TURNING THE DEVICE ON

When the device is being turned on, the weighing pan should be empty. It is recommended to allow the device to warm up for about 15-30 minutes before further use. Perform calibration before first use.

3.3. DEVICE USE COUNTING FUNCTION

The weight can count the goods depending on their weight. This function can be used together with the tare function if needed. To do this, place the container on the pan before starting the counting and press the [TARE] button. While the display is showing zero, press the [PCS] button. The display will show the „—P 10—“ message. Press the [TARE] button to select the number of items to be sampled (10, 20, 50, 100, 200, 500). Place the same number of items on the weighing pan, then press the [PCS] button. Once the measurement has stabilized, the sampling procedure will end and you can use the counting function. The display will show „FAIN“ and the device will return to counting mode. Hint: It is better to use more samples to make the counting more accurate.

THE TARE FUNCTION

The tare is the weight of the container that can be subtracted to read the weight of the item being weighed. To use the tare function, place an empty container on the weighing pan. When the device stabilizes, press the [TARE] key. The display will show zero and the tare symbol will light up. The tare weight (and the weight of the item) must not exceed the maximum load of the device.

PRINT FUNCTION

During weighing, press the [PRINT] button. The measured value will be sent to the computer.

CALIBRATION

Press the [TARE] button for about 2 seconds. The display will show the weight that is to be placed on the device. Place a load of the indicated weight on the pan and press the [TARE] button after the device stabilizes. When the load value is displayed, the calibration procedure is over.

RS-232 TRANSMISSION METHOD

This function allows the measurement data to be sent to any terminal or serial port. The device uses a Simplex asynchronous serial communication (can be changed to Duplex if necessary) of the RS232 electrical interface. The data transmission is in ASCII code. Each data set has 10 bits. Information about the individual bits:

bit1	bit2	bit3	bit4	bit5	bit6	bit7	bit8	bit9	bit10
------	------	------	------	------	------	------	------	------	-------

- BIT1: start bit
- BIT2 - BIT9: data bit
- BIT10: stop bit
- Transmission speed: default „9600“ bps.

TRANSMISSION DATA FORMAT

- Bit 1 - 2: Operation mode:
 - » WT - Weighing
 - » CT - Counting
 - » PC - Percentage mode.
- Bit No. 3 - 4: Measurement status
 - » OL - Overload
 - » UL - Negative
 - » ST - Stable
 - » US - Unstable.
- Bit No. 5 - 12: The measurement value contained in 8 bits where one bit corresponds to the sign separating decimal from integer values (here „.“); if the data does not take up all 8 bits, a space will be used as the fill.

- The „+“ or „-“ symbol is 1 bit (spaces are used instead of the „+“ symbol).
- Bit 13 - 16: The current weight unit contained in 4 bits. If the unit does not take all the 4 bits in the weighing mode or in the percentage mode, the empty fields will be filled with spaces.
- Bit 17 - 18: Movement to the beginning of the current line and a new line are marked with „0DH“, „0AH“.

Exemplary data transmission recording:

Weighing, steady value, 38.25g:

W	T	S	T				3	8	.	2	5									g	0D	0A
---	---	---	---	--	--	--	---	---	---	---	---	--	--	--	--	--	--	--	--	---	----	----

Weighing, unsteady value, 300ct:

W	T	U	S							3	0	0								c	t	0D	0A
---	---	---	---	--	--	--	--	--	--	---	---	---	--	--	--	--	--	--	--	---	---	----	----

Counting mode, steady value, + 60:

C	T	S	T							6	0									p	c	s	0D	0A
---	---	---	---	--	--	--	--	--	--	---	---	--	--	--	--	--	--	--	--	---	---	---	----	----

Percentage mode, steady value + 25.4:

P	C	S	T							2	5	.	4							%	0D	0A
---	---	---	---	--	--	--	--	--	--	---	---	---	---	--	--	--	--	--	--	---	----	----

USE GUIDELINES

- Do NOT leave a load on the scale for long periods of time.
- Avoid extreme temperatures. Do not place the unit in direct sunlight or near air conditioning units.
- Avoid using strong electrical or magnetic fields during use.
- Avoid vibration. Do not place the device near heavy or vibrating machinery.
- Do not place the unit close to open windows, doors or fans that may cause unstable results due to air currents.
- Keep your machine clean. Do not store products on the scale when it is not being used.
- Avoid any strong shock or heavy weight. Don't put any article or give pressure over maximum capacity on the scale pan even the scale is not in use.
- Wrong wires connection of storage battery is forbidden. Right connection: Red wire should be connected with red end; Black wire should be connected with black end.

3.4. CLEANING AND MAINTENANCE

- a) Unplug the mains plug and allow the device to cool completely before each cleaning, adjustment or replacement of accessories, or if the device is not being used.
- b) Use only non-corrosive cleaners to clean the surface.
- c) After cleaning the device, all parts should be dried completely before using it again.
- d) Store the unit in a dry, cool place, free from moisture and direct exposure to sunlight.
- e) Do not spray the device with a water jet or submerge it in water.
- f) Do not allow water to get inside the device through vents in the housing of the device.
- g) The device must be regularly inspected to check its technical efficiency and spot any damage.
- h) Use a soft cloth for cleaning.
- i) Do not leave the battery in the device if it will not be used for a longer period of time.

3.4. CLEANING AND MAINTENANCE

The devices are equipped with 6V / 1.2Ah lead-acid batteries. Remove used batteries from the device using the same procedure by which you installed them. Recycle batteries with the appropriate organisation or company.

DISPOSING OF USED DEVICES

Do not dispose of this device in municipal waste systems. Hand it over to an electric and electrical device recycling and collection point. Check the symbol on the product, instruction manual and packaging. The plastics used to construct the device can be recycled in accordance with their markings. By choosing to recycle you are making a significant contribution to the protection of our environment. Contact local authorities for information on your local recycling facility.

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