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USER MANUAL

DIGITAL PRECISION SCALE

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PŘESNÁ VÁHA

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BILANCIA DI PRECISIONE

MANUAL DE INSTRUCCIONES

BALANZA DE PRECIÓN

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SBS-LW-2000E | SBS-LW-300001E

INHALT | CONTENT | TREŚĆ | OBSAH | CONTENU | CONTENUTO | CONTENIDO

- Deutsch 3
- English 7
- Polski 11
- Česky 15
- Français 19
- Italiano 23
- Español 27

PRODUKTNAME PRODUCT NAME NAZWA PRODUKTU NÁZEV VÝROBKU NOM DU PRODUIT NOME DEL PRODOTTO NOMBRE DEL PRODUCTO	PRÄZISIONSWAAGE DIGITAL PRECISION SCALE WAGA LABORATORYJNA PŘESNÁ VÁHA BALANCE DE PRÉCISION BILANCIA DI PRECISIONE BALANZA DE PRECISIÓN
MODELL PRODUCT MODEL MODEL PRODUKTU MODEL VÝROBKU MODÈLE MODELLO MODELO	SBS-LW-2000E SBS-LW-3000001E
IMPORTEUR IMPORTER IMPORTER DOVOZCE IMPORTATEUR IMPORTATORE IMPORTADOR	EXPONDO POLSKA SP. Z O.O. SP. K.
ADRESSE VON IMPORTEUR IMPORTER ADDRESS ADRES IMPORTERA ADRESA DOVOZCE ADRESSE DE L'IMPORTATEUR INDIRIZZO DELL'IMPORTATORE DIRECCIÓN DEL IMPORTADOR	UL. NOWY KISIELIN-INNOWACYJNA 7, 66-002 ZIELONA GÓRA POLAND, EU

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1. GENERAL DESCRIPTION

The user manual is designed to assist in the safe and trouble-free use of the device. The product is designed and manufactured in accordance with strict technical guidelines, using state-of-the-art technologies and components. Additionally, it is produced in compliance with the most stringent quality standards.

DO NOT USE THE DEVICE UNLESS YOU HAVE THOROUGHLY READ AND UNDERSTOOD THIS USER MANUAL.

To increase the product life of the device and to ensure trouble-free operation, use it in accordance with this user manual and regularly perform maintenance tasks. The technical data and specifications in this user manual are up to date. The manufacturer reserves the right to make changes associated with quality improvement.

LEGEND

- The product satisfies the relevant safety standards.
- Read instructions before use.
- The product must be recycled.

WARNING! or CAUTION! or REMEMBER!
Applicable to the given situation. (general warning sign).

ATTENTION! Electric shock warning!

Only use indoors.

PLEASE NOTE! Drawings in this manual are for illustration purposes only and in some details may differ from the actual machine.

The original operation manual is in German. Other language versions are translations from German.

2. USAGE SAFETY

ATTENTION! Read all safety warnings and all instructions. Failure to follow the warnings and instructions may result in an electric shock, fire and/or serious injury or even death.

The terms „device“ or „product“ are used in the warnings and instructions to refer to <Digital Precision Scale>. Do not use in very humid environments or in the direct vicinity of water tanks. Prevent the device from getting wet. Risk of electric shock! Do not cover air inlets / outlets! Do not cover the ventilation openings!

2.1. ELECTRICAL SAFETY

- The plug must fit the socket. Do not modify the plug in any way. Using original plugs and matching sockets reduces the risk of electric shock.
- Avoid touching earthed elements such as pipes, heaters, boilers and refrigerators. There is an increased risk of electric shock if the earthed device is exposed to rain, comes into direct contact with a wet surface or is operating in a damp environment. Water getting into the device increases the risk of damage to the device and of electric shock.
- Do not touch the device with wet or damp hands.
- Use the cable only for its designated use. Never use it to carry the device or to pull the plug out of a socket. Keep the cable away from heat sources, oil, sharp edges or moving parts. Damaged or tangled cables increase the risk of electric shock.
- If using the device in a damp environment cannot be avoided, a residual current device (RCD) should be applied. The use of an RCD reduces the risk of electric shock.
- Do not use the device if the power cord is damaged or shows obvious signs of wear. A damaged power cord should be replaced by a qualified electrician or the manufacturer's service centre
- To avoid electric shock, do not immerse the cord, plug or device in water or other liquids. Do not use the device on wet surfaces.

2.2. SAFETY IN THE WORKPLACE

- Do not use the device in a potentially explosive environment, for example in the presence of flammable liquids, gases or dust. The device generates sparks which may ignite dust or fumes.
- If you discover damage or irregular operation, immediately switch the device off and report it to a supervisor without delay.
- If there are any doubts as to the correct operation of the device, contact the manufacturer's support service.

- d) Only the manufacturer's service point may repair the device. Do not attempt any repairs independently!
- e) In case of fire, use a powder or carbon dioxide (CO₂) fire extinguisher (one intended for use on live electrical devices) to put it out.
- f) Regularly inspect the condition of the safety labels. If the labels are illegible, they must be replaced.
- g) Please keep this manual available for future reference. If this device is passed on to a third party, the manual must be passed on with it.
- h) Keep packaging elements and small assembly parts in a place not available to children.

REMEMBER When using the device, protect children and other bystanders.

2.3. PERSONAL SAFETY

- a) Do not use the device when tired, ill or under the influence of alcohol, narcotics or medication which can significantly impair the ability to operate the device.
- b) The device is not designed to be handled by persons (including children) with limited mental and sensory functions or persons lacking relevant experience and/or knowledge unless they are supervised by a person responsible for their safety or they have received instruction on how to operate the device.
- c) To prevent the device from accidentally switching on, make sure the switch is on the OFF position before connecting to a power source.
- d) The device is not a toy. Children must be supervised to ensure that they do not play with the device.

2.4. SAFE DEVICE USE

- a) Do not use the device if the ON/OFF switch does not function properly (does not switch the device on and off). Devices which cannot be switched on and off using the ON/OFF switch are hazardous, should not be operated and must be repaired.
- b) Disconnect the device from the power supply before commencement of adjustment, cleaning and maintenance. Such a preventive measure reduces the risk of accidental activation.
- c) When not in use, store in a safe place, away from children and people not familiar with the device who have not read the user manual. The device may pose a hazard in the hands of inexperienced users.
- d) Keep the device in perfect technical condition. Before each use check for general damage and especially check for cracked parts or elements and for any other conditions which may impact the safe operation of the device. If damage is discovered, hand over the device for repair before use.
- e) Keep the device out of the reach of children.
- f) Device repair or maintenance should be carried out by qualified persons, only using original spare parts. This will ensure safe use.
- g) To ensure the operational integrity of the device, do not remove factory-fitted guards and do not loosen any screws.
- h) Clean the device regularly to prevent stubborn grime from accumulating.
- i) The device is not a toy. Cleaning and maintenance may not be carried out by children without supervision by an adult person.
- j) It is forbidden to interfere with the structure of the device in order to change its parameters or construction.
- k) Keep the device away from sources of fire and heat.

ATTENTION! Despite the safe design of the device and its protective features, and despite the use of additional elements protecting the operator, there is still a slight risk of accident or injury when using the device. Stay alert and use common sense when using the device.

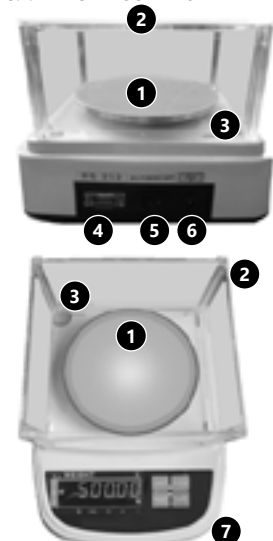
3. USE GUIDELINES

The device is intended to weigh items placed on the weighing pan. Warning! The device cannot be used for commercial purposes.

The user is liable for any damage resulting from unintended use of the device.

The balance should be used in a space free of wind, corrosion, vibration, humidity and temperature variations. The scale should avoid vibration, violent air current, direct sunshine and high temperature.

3.1. DEVICE DESCRIPTION

1. Weighing pan
2. Cover
3. Bubble level
4. RS-232 Interface
5. I/O (ON/ OFF)
6. Power socket
7. Control panel
 - A. UNIT button
 - B. PRINT button
 - C. PCS button
 - D. TARE button
 - E. Display

Display symbols

	Tare	Tare function active
	Steady reading	The device is stable
	Zero	Reset function active

	Piece counting	Piece counting function active
tola	Tola	Weighing unit: tola
lb	Pound	Weighing unit: pound
ct	Carat	Weighing unit: carat
g	Gram	Weighing unit: gram
AC	Power on	The device is connected to power supply

3.2. PREPARING FOR USE DEVICE USE

- ▶ **UNIT.** Changing weighing units:
In order to change the weighing unit, press the UNIT button, the unit symbol will change on the display.
- ▶ **TARE.** Tare weight:
Place an empty container on the weighing pan. When the reading is steady, press the TARE button and the display will zero. The display will show:
 - ZERO if the value is less than $\pm 5\%$ of the scale;
 - TARE if the value is more than $\pm 5\%$ of the scale;
- ▶ **PCS.** Piece counting
If required, place the container on the weighing pan and press „TARE“. While the zero function is active, press the PCS button. The „—SAP 10—“ message will be displayed. Press the UNIT button to select the sample size (10, 20, 50, 100, 200, 500, 1000 pieces). After selecting the number of samples, press the TARE button to confirm. The display will show „LOAD-C“. Place the previously selected number of samples on the weighing pan. After the steady reading has disappeared, the sampling procedure is completed, and you can start using the piece counting function.
NOTE: it is best to use more samples to make the count more accurate.
- ▶ **PRINT.** Printing function
Pressing the button during the weighing procedure will send the weight data to the computer.
- ▶ **CALIBRATION**
Press and hold the TARE button for 3 seconds, the „LOAD-2“ message will be displayed. By using the PCS key, you can select „LOAD-2“, „LOAD-3“ or „LOAD-4“. There are two ways to calibrate:
 - Constant load:
Press the TARE button again, a value will be displayed, place an equivalent load on the weighing pan. The display will show „ZERO“. Remove the load and put on a weight corresponding to the next value indicated. After this operation the calibration is completed.
 - Load selection:
Select „LOAD-1“ using the PCS key. Press the TARE button. While the display is flashing, press the „PCS“ button and enter the value. To change the digit position, press the TARE button. After moving to the last digit position, place the selected load on the weighing pan. When the reading stabilizes, the calibration procedure is completed.

- ▶ **RS-232 TRANSMISSION METHOD**
This function allows the measurement data to be sent to any terminal or serial port. The device uses a Simplex asynchronous serial communication (can be changed to Duplex if necessary) of the RS232 electrical interface. The data transmission is in ASCII code. Each data set has 10 bits. Information about the individual bits:

bit1	bit2	bit3	bit4	bit5	bit6	bit7	bit8	bit9	bit10
------	------	------	------	------	------	------	------	------	-------

BIT1: start bit
BIT2 - BIT9: data bit
BIT10: stop bit
Transmission speed: default „9600“ bps.

- ▶ **TRANSMISSION DATA FORMAT**
 - Bit 1 - 2: Operation mode:
 - » WT - Weighing
 - » CT - Counting
 - » PC - Percentage mode.
 - Bit No. 3 - 4: Measurement status
 - » OL - Overload
 - » UL - Negative
 - » ST - Stable
 - » US - Unstable.
 - Bit No. 5 - 12: The measurement value contained in 8 bits where one bit corresponds to the sign separating decimal from integer values (here „.“); if the data does not take up all 8 bits, a space will be used as the fill.
 - The „+“ or „-“ symbol is 1 bit (spaces are used instead of the „+“ symbol).
 - Bit 13 - 16: The current weight unit contained in 4 bits.
 - If the unit does not take all the 4 bits in the weighing mode or in the percentage mode, the empty fields will be filled with spaces.
 - Bit 17 - 18: Movement to the beginning of the current line and a new line are marked with „ODH“, „OAH“.

Exemplary data transmission recording:
Weighing, steady value, 38.25g:

W	T	S	T				3	8	.	2	5							g	OD	OAH
---	---	---	---	--	--	--	---	---	---	---	---	--	--	--	--	--	--	---	----	-----

Weighing, unsteady value, 300ct:

W	T	U	S							3	0	0							c	t	OD	OAH
---	---	---	---	--	--	--	--	--	--	---	---	---	--	--	--	--	--	--	---	---	----	-----

Counting mode, steady value, + 60:

C	T	S	T							6	0								p	c	s	OD	OAH
---	---	---	---	--	--	--	--	--	--	---	---	--	--	--	--	--	--	--	---	---	---	----	-----

Percentage mode, steady value + 25.4:

P	C	S	T							2	5	.	4									%	OD	OAH
---	---	---	---	--	--	--	--	--	--	---	---	---	---	--	--	--	--	--	--	--	--	---	----	-----

USE GUIDELINES

- Do NOT leave a load on the scale for long periods of time.
- Avoid extreme temperatures. Do not place the unit in direct sunlight or near air conditioning units.
- Avoid using strong electrical or magnetic fields during use.
- Avoid vibration. Do not place the device near heavy or vibrating machinery.
- Do not place the unit close to open windows, doors or fans that may cause unstable results due to air currents.
- Keep your machine clean. Do not store products on the scale when it is not being used.

- Avoid any strong shock or heavy weight. Don't put any article or give pressure over maximum capacity on the scale pan even the scale is not in use.
- Wrong wires connection of storage battery is forbidden. Right connection: Red wire should be connected with red end; Black wire should be connected with black end.

3.3. CLEANING AND MAINTENANCE

- Unplug the mains plug and allow the device to cool completely before each cleaning, adjustment or replacement of accessories, or if the device is not being used.
- Use only non-corrosive cleaners to clean the surface.
- again.
- Store the unit in a dry, cool place, free from moisture and direct exposure to sunlight.
- Do not spray the device with a water jet or submerge it in water.
- Do not allow water to get inside the device through vents in the housing of the device.
- The device must be regularly inspected to check its technical efficiency and spot any damage.
- Use a soft cloth for cleaning.
- Do not leave the battery in the device if it will not be used for a longer period of time.

SAFE REMOVAL OF BATTERIES AND RECHARGEABLE BATTERIES

The devices are equipped with 6V / 1.2Ah lead-acid batteries. Remove used batteries from the device using the same procedure by which you installed them. Recycle batteries with the appropriate organisation or company.

DISPOSING OF USED DEVICES

Do not dispose of this device in municipal waste systems. Hand it over to an electric and electrical device recycling and collection point. Check the symbol on the product, instruction manual and packaging. The plastics used to construct the device can be recycled in accordance with their markings. By choosing to recycle you are making a significant contribution to the protection of our environment. Contact local authorities for information on your local recycling facility.

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