Safety & Operating Instructions

DE	Sicherheits- und Betriebsanweisungen
PL	Instrukcje dotyczące bezpieczeństwa i obsługi
CZ	Bezpečnostní a provozní pokyny
FR	Consignes de sécurité et d'utilisation
IT	Istruzioni per la sicurezza e l'uso
ES	Instrucciones de seguridad y funcionamiento
HU	Biztonsági és használati útmutató
DK	Sikkerheds- og betjeningsvejledning
FI	Turvallisuus- ja käyttöohjeet
NO	Sikkerhets- og bruksanvisning
NL	Veiligheids- en bedieningsinstructies
SE	Säkerhets- och bruksanvisningar
PT	Instruções de segurança e de funcionamento
SK	Bezpečnostné a prevádzkové pokyny
BG	Инструкции за безопасност и експлоатация
GR	Οδηγίες ασφαλείας και λειτουργίας
HR	Sigurnosne i upute za rad
LT	Saugos ir naudojimo instrukcijos
RO	Instrucțiuni de siguranță și operare
SL	Varnost in navodila za uporabo

DIGITAL PRECISION SCALE

DE	Produktname	Digitale Präzisionswaage
EN	Product name	Digital Precision Scale
PL	Nazwa produktu	Cyfrowa waga precyzyjna
CZ	Název výrobku	Digitální přesná váha
FR	Nom du produit	Balance de précision numérique
IT	Nome del prodotto	Bilancia di precisione digitale
ES	Nombre del producto	Balanza de precisión digital
HU	Termék neve	Digitális pontos mérleg
DA	Produktnavn	Digital præcisionsvægt
FI	Tuotteen nimi	Digitaalinen tarkkuusvaaka
NL	Productnaam	Digitale precisiebalans
NO	Produktnavn	Digital presisjonsvekt
SE	Produktnamn	Digital precisionsväg
PT	Nome do produto	Balança de precisão digital
SK	Názov produktu	Digitálna presná váha
BG	Име на продукта	Цифрова прецизна везна
EL	Όνομα προϊόντος	Ψηφιακή ζυγαριά ακριβείας
HR	Naziv proizvoda	Digitalna precizna vaga
LT	Produkto pavadinimas	Skaitmeninės tikslios svarstyklės
RO	Numele produsului	Cântar digital de precizie
SL	Ime izdelka	Digitalna precizna tehtnica
DE Modell EN Product model PL Model produktu CZ Model výrobku FR Modèle IT Modello ES Modelo HU Modell DA Model FI Tuotteen malli NL Productmodel NO Produktmodell SE Produktmodell PT Modelo do produto SK Model BG Модел на продукт EL Μοντέλο προϊόντος HR Model proizvoda LT : Gaminio modelis RO : Model de produs SL : Model izdelka		SBS-LW-600E SBS-LW-300001E SBS-LW-600001E
DE Hersteller EN Manufacturer PL Producent CZ Výrobce FR Fabricant IT Produttore ES Fabricante HU Termelő DA Producent FI Valmistaja NL Producent NO Produsent SE Tillverkare PT Fabricante SK Výrobca BG Производител EL Κατασκευαστής HR Proizvođač LT Gamintojas RO Producător SL Proizvajalec		expondo Polska sp. z o.o. sp. k.
DE Anschrift des Herstellers EN Manufacturer Address PL Adres producenta CZ Adresa výrobce FR Adresse du fabricant IT Indirizzo del produttore ES Dirección del fabricante HU A gyártó címe DA Producentens adresse FI Valmistajan osoite NL Adres producent NO Produsentens adresse SE Tillverkarens adress PT Endereço do fabricante SK Adresa výrobcu BG Адрес на производителя EL : Διεύθυνση κατασκευαστή HR Adresa proizvođača LT Gamintojo adresas RO Adresa producătorului SL Naslov proizvajalca		ul. Nowy Kisielin – Innowacyjna 7, 66-002 Zielona Góra Poland, EU

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This User Manual has been translated using machine translation. We have made every effort to ensure the translation is accurate, but please note that automated translations are not perfect and are not meant to replace human translators. The official version of the User Manual is in English. Any differences between the translated version and the original English are not legally binding. If you have any questions about the accuracy of the translation, please refer to the English version, which is the official reference. More language versions are available upon request via info@expondo.com.

Technical data

Parameter description	Parameter value		
Product name	Digital Precision Scale		
Model	SBS-LW- 600E	SBS-LW- 300001E	SBS-LW- 600001E
Rated voltage [V~]/ Frequency [Hz]	230/50		
Working temperature/ Relative working humidity [°C/%RH]	0°C-40°C/<65%		
Storage temperature/ Relative storage humidity [°C/%RH]	10°C--50°C/<85%		
Storage battery capacity	6V/1,3Ah		
Power supply parameters	8V/500mA		
Division [g]	0.01	0.1	
Max capacity [g]	600	3000	6000
Min capacity [g]	0.05	0.5	
Battery life [h]	24		
Battery charging time [h]	8		
Units of measure- ment	g/lb		
Pan dimensions [mm]	Ø130	160x180	
Dimensions [mm]	265x196 x80	195x270x90	
Weight [kg]	1.4	1.7	

1. General description

The user manual is designed to assist in the safe and trouble-free use of the device. The product is designed and manufactured in accordance with strict technical guidelines, using state-of-the-art technologies and components. Additionally, it is produced in compliance with the most stringent quality standards.

DO NOT USE THE DEVICE UNLESS YOU HAVE THOROUGHLY READ AND UNDERSTOOD THIS USER MANUAL.

To increase the product life of the device and to ensure trouble-free operation, use it in accordance with this user manual and regularly perform maintenance tasks. The technical data and specifications in this user manual are up to date. The manufacturer reserves the right to make changes associated with quality improvement. The device is designed to reduce noise emission risks to a minimum, taking into account technological progress and noise reduction opportunities.

Legend

The product satisfies the relevant safety standards.

Read instructions before use.

The product must be recycled.

WARNING! or **CAUTION!** or **REMEMBER!** Applicable to the given situation.
(general warning sign)

ATTENTION! Electric shock warning!

PLEASE NOTE! Drawings in this manual are for illustration purposes only and in some details may differ from the actual product.

2. Usage safety

ATTENTION! Read all safety warnings and all instructions. Failure to follow the warnings and instructions may result in electric shock, fire and/or serious injury or even death.

The terms "device" or "product" are used in the warnings and instructions to refer to: Digital Precision Scale.

2.1. Electrical safety

- a) The plug must fit the socket. Do not modify the plug in any way. Using original plugs and matching sockets reduces the risk of electric shock.
- b) Avoid touching earthed elements such as pipes, heaters, boilers and refrigerators. There is an increased risk of electric shock if the earthed device is exposed to rain, comes into direct contact with a wet surface or is operating in a damp environment. Water getting into the device increases the risk of damage to the device and of electric shock.
- c) Do not touch the device with wet or damp hands.
- d) Use the cable only for its designated use. Never use it to carry the device or to pull the plug out of a socket. Keep the cable away from heat sources, oil, sharp edges or moving parts. Damaged or tangled cables increase the risk of electric shock.
- e) If using the device in a damp environment cannot be avoided, a residual current device (RCD) should be applied. The use of an RCD reduces the risk of electric shock.
- f) Do not use the device if the power cord is damaged or shows obvious signs of wear. A damaged power cord should be replaced by a qualified electrician or the manufacturer's service centre.
- g) To avoid electric shock, do not immerse the cord, plug or device in water or other liquids. Do not use the device on wet surfaces.

2.2. Safety in the workplace

- a) Do not use the device in a potentially explosive environment, for example in the presence of flammable liquids, gases or dust. The device generates sparks which may ignite dust or fumes.
- b) If you discover damage or irregular operation, immediately switch the device off and report it to a supervisor without delay.
- c) If you are unsure about whether the device is operating correctly or if you find damage, please contact the manufacturer's service centre.
- d) Only the manufacturer's service centre may make repairs to the device. Do not attempt to make repairs yourself!
- e) In case of fire, use a powder or carbon dioxide (CO₂) fire extinguisher (one intended for use on live electrical devices) to put it out.
- f) Regularly inspect the condition of the safety labels. If the labels are illegible, they must be replaced.
- g) Please keep this manual available for future reference. If this device is passed on to a third party, the manual must be passed on with it.
- h) Keep packaging elements and small assembly parts in a place not available to children.

Remember! When using the device, protect children and other bystanders.

2.3. Personal safety

- a) Do not use the device when tired, ill or under the influence of alcohol, narcotics or medication which can significantly impair the ability to operate the device.
- b) The device is not designed to be handled by persons (including children) with limited mental and sensory functions or persons lacking relevant experience and/or knowledge unless they are supervised by a person responsible for their safety or they have received instruction on how to operate the device.
- c) To prevent the device from accidentally switching on, make sure the switch is on the OFF position before connecting to a power source.
- d) The device is not a toy. Children must be supervised to ensure that they do not play with the device.

2.4. Safe device use

- a) Do not use the device if the "ON/OFF" switch does not function properly (does not switch the device on and off). Devices which cannot be switched on and off using the "ON/OFF" switch are hazardous, should not be operated and must be repaired.
- b) Disconnect the device from the power supply before commencement of adjustment, cleaning and maintenance. Such a preventive measure reduces the risk of accidental activation.
- c) When not in use, store in a safe place, away from children and people not familiar with the device who have not read the user manual. The device may pose a hazard in the hands of inexperienced users.
- d) Keep the device out of the reach of children.
- e) Device repair or maintenance should be carried out by qualified persons, only using original spare parts. This will ensure safe use.
- f) To ensure the operational integrity of the device, do not remove factory-fitted guards and do not loosen any screws.
- g) Clean the device regularly to prevent stubborn grime from accumulating.
- h) The device is not a toy. Cleaning and maintenance may not be carried out by children without supervision by an adult person.
- i) It is forbidden to interfere with the structure of the device in order to change its parameters or construction.
- j) Keep the device away from sources of fire and heat.

ATTENTION! Despite the safe design of the device and its protective features, and despite the use of additional elements protecting the operator, there is still a slight risk of accident or injury when using the device. Stay alert and use common sense when using the device.

3. Use guidelines

The device is intended to weigh items placed on the weighing pan. Warning! The device cannot be used for commercial purposes. The user is liable for any damage resulting from unintended use of the device.

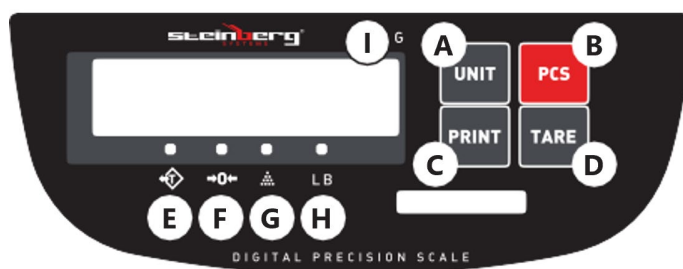
The user is liable for any damage resulting from unintended use of the device.

The balance should be used in a space free of wind, corrosion, vibration, humidity and temperature variations. The scale should avoid vibration, violent air current, direct sunshine and high temperature.

3.1. Device description

- 1- Weighing pan
- 2- Control panel
- 3- RS-232 interface
- 4- I/O (ON/OFF) button
- 5- Power supply socket
- 6- Bubble level

The keys:

- A. UNIT: Unit selection
- B. PRINT: Printing function.
- C. PCS: Counting function.
- D. TARE: Tare weight of container.

LEDs indicating function:

- E. Tare
- F. Zero status
- G. Counting
- H. Weight measurement unit: lb
- I. Weight measurement unit: g

3.2. Preparing for use

LEVELLING THE DEVICE

The scale is equipped with a level indicator under the weighing pan and four adjustable legs. Before using the device, adjust the length of the legs until the bubble of the level indicator is in the middle position.

TURNING THE DEVICE ON

When the device is being turned on, the weighing pan should be empty. It is recommended to allow the device to warm up for about 15-30 minutes before further use. Perform calibration before first use.

3.3. Device use

COUNTING FUNCTION

The weight can count the goods depending on their weight. This function can be used together with the tare function if needed. To do this, place the container on the pan before starting the counting and press the [TARE] button.

While the display is showing zero, press the [PCS] button. The display will show the "P 10" message. Press the [TARE] button to select the number of items to be sampled (10, 20, 50, 100, 200, 500). Place the same number of items on the weighing pan, then press the [PCS] button. Once the measurement has stabilized, the sampling procedure will end, and you can use the counting function. The display will show "FAIN," and the device will return to counting mode.

Hint: It is better to use more samples to make the counting more accurate.

THE TARE FUNCTION

The tare is the weight of the container that can be subtracted to read the weight of the item being weighed. To use the tare function, place an empty container on the weighing pan. When the device stabilizes, press the [TARE] key. The display will show zero, and the tare symbol will light up. The tare weight (and the weight of the item) must not exceed the maximum load of the device.

PRINT FUNCTION

During weighing, press the [PRINT] button. The measured value will be sent to the computer.

CALIBRATION

Press the **【TARE】** key for about 2 seconds, the display will show the "calibration weight value". Put the same weight on the pan, press the **【TARE】** key after the stable symbol appears. When the weight value is displayed, the calibration procedure is finished.

RS-232 TRANSMISSION METHOD

This function allows the measurement data to be sent to any terminal or serial port. The device uses a Simplex asynchronous serial communication (can be changed to Duplex if necessary) of the RS232 electrical interface.

The data transmission is in ASCII code.

Each data set has 10 bits. Information about the individual bits:

bit1	bit2	bit3	bit4	bit5	bit6	bit7	bit8	bit9	bit10
------	------	------	------	------	------	------	------	------	-------

- BIT1: start bit
- BIT2 - BIT9: data bit
- BIT10: stop bit
- Transmission speed: default "9600" bps.

TRANSMISSION DATA FORMAT

- Bit 1 - 2: Operation mode:
 - WT - Weighing
 - CT - Counting

- PC - Percentage mode
- Bit No. 3 - 4: Measurement status
 - OL - Overload
 - UL - Negative
 - ST - Stable
 - US - Unstable
- Bit No. 5 - 12: The measurement value contained in 8 bits, where one bit corresponds to the sign separating decimal from integer values (here "."); if the data does not take up all 8 bits, a space will be used as the fill.
- The "+" or "-" symbol is 1 bit (spaces are used instead of the "+" symbol).
- Bit 13 - 16: The current weight unit is contained in 4 bits. If the unit does not take all 4 bits in the weighing mode or in the percentage mode, the empty fields will be filled with spaces.
- Bit 17 - 18: Movement to the beginning of the current line and a new line are marked with "0DH", "0AH".

Exemplary data transmission recording:

Weighing, steady value, 38.25g:

W	T	S	T					3	8	.	2	5					g	0D	0A
---	---	---	---	--	--	--	--	---	---	---	---	---	--	--	--	--	---	----	----

Weighing, unsteady value, 300ct:

W	T	U	S								3	0	0					c	t	0D	0A
---	---	---	---	--	--	--	--	--	--	--	---	---	---	--	--	--	--	---	---	----	----

Counting mode, steady value, +60:

C	T	S	T									6	0					p	c	s	0D	0A
---	---	---	---	--	--	--	--	--	--	--	--	---	---	--	--	--	--	---	---	---	----	----

Percentage mode, steady value +25.4:

P	C	S	T									2	5	.	4					%	0D	0A
---	---	---	---	--	--	--	--	--	--	--	--	---	---	---	---	--	--	--	--	---	----	----

USE GUIDELINES

- Do NOT leave a load on the scale for long periods of time.
- Avoid extreme temperatures. Do not place the unit in direct sunlight or near air conditioning units.
- Avoid using strong electrical or magnetic fields during use.
- Avoid vibration. Do not place the device near heavy or vibrating machinery.
- Do not place the unit close to open windows, doors, or fans that may cause unstable results due to air currents.
- Keep your machine clean. Do not store products on the scale when it is not being used.

- Avoid any strong shock or heavy weight. Don't put any article or apply pressure over the maximum capacity on the scale pan even if the scale is not in use.
- Wrong wire connection of storage battery is forbidden.
- Right connection: Red wire should be connected with red end; Black wire should be connected with black end.

3.4. Cleaning and maintenance

- a) Unplug the mains plug and allow the device to cool completely before each cleaning, adjustment, or replacement of accessories, or if the device is not being used.
- b) Use only non-corrosive cleaners to clean the surface.
- c) After cleaning the device, all parts should be dried completely before using it again.
- d) Store the unit in a dry, cool place, free from moisture and direct exposure to sunlight.
- e) Do not spray the device with a water jet or submerge it in water.
- f) Do not allow water to get inside the device through vents in the housing of the device.
- g) The device must be regularly inspected to check its technical efficiency and spot any damage.
- h) Use a soft cloth for cleaning.
- i) Do not leave the battery in the device if it will not be used for a longer period of time.

SAFE REMOVAL OF BATTERIES AND RECHARGEABLE BATTERIES:

The devices are equipped with **6V / 1.2Ah lead-acid batteries**. Remove used batteries from the device using the same procedure by which you installed them. Recycle batteries with the appropriate organisation or company.

DISPOSING OF USED DEVICES:

Do not dispose of this device in municipal waste systems. Hand it over to an electric and electrical device recycling and collection point. Check the symbol on the product, instruction manual and packaging. The plastics used to construct the device can be recycled in accordance with their markings. By choosing to recycle you are making a significant contribution to the protection of our environment. Contact local authorities for information on your local recycling facility.

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Environmental and disposal information

Producer to customer

Dear Sir or Madam,

Let's take care of the environment! The product you own should not be disposed of with household waste. According to applicable regulations (Directive 2012/19/EU and other local regulations), used equipment must be taken to designated collection points for proper disposal and recycling.

How to Handle the Product at the End of Its Lifecycle?

- 1. Electrical and Electronic Equipment:** Products marked with the crossed-out wheeled bin symbol must not be disposed of with household waste. You can:
 - Take it to your local collection point for electrical and electronic equipment.
 - Use the “one-for-one” system (returning the old device when purchasing a new one, if available in your region).
- 2. Batteries and Accumulators:** If the product contains batteries, remove them before recycling and take them to designated battery collection points.
- 3. Packaging Materials:** Packaging, such as cardboard or plastic, should be sorted and placed in recycling bins.

Why is This Important?

Proper disposal of equipment and packaging materials:

- Protects the environment from harmful substances.
- Enables the recovery of valuable resources.
- Reduces the amount of waste sent to landfills.

Where to Find Information?

Information about waste collection points can be found on local authorities' websites or through your recycling service provider. If needed, contact our customer service team, who will gladly provide further details.

Thank you for your care for the environment and commitment to recycling!

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