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USER MANUAL

PRECISION SCALE

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PŘESNÁ VÁHA

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BILANCIA DI PRECISIONE

MANUAL DE INSTRUCCIONES

BALANZA DE PRECISIÓN

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**SBS-LW-300E | SBS-LW-600001 | SBS-LW-3000001
SBS-LW-300001 | SBS-LW-6000**

■ Deutsch	3
■ English	7
■ Polski	11
■ Česky	15
■ Français	19
■ Italiano	23
■ Español	27

PRODUKTNAME PRODUCT NAME NAZWA PRODUKTU NÁZEV VÝROBKU NOM DU PRODUIT NOME DEL PRODOTTO NOMBRE DEL PRODUCTO	PRÄZISIONSWAAGE PRECISION SCALE WAGA LABORATORYJNA PŘESNÁ VÁHA BALANCE DE PRÉCISION BILANCIA DI PRECISIONE BALANZA DE PRECISIÓN
MODELL PRODUCT MODEL MODEL PRODUKTU MODEL VÝROBKU MODÈLE MODELLO MODELO	SBS-LW-300E SBS-LW-600001 SBS-LW-3000001 SBS-LW-300001 SBS-LW-6000
IMPORTEUR IMPORTER IMPORTER DOVOZCE IMPORTATEUR IMPORTATORE IMPORTADOR	EXPONDO POLSKA SP. Z O.O. SP. K.
ADRESSE VON IMPORTEUR IMPORTER ADDRESS ADRES IMPORTERA ADRESA DOVOZCE ADRESSE DE L'IMPORTATEUR INDIRIZZO DELL'IMPORTATORE DIRECCIÓN DEL IMPORTADOR	UL. NOWY KISIELIN-INNOWACYJNA 7, 66-002 ZIELONA GÓRA POLAND, EU

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PLEASE NOTE! Drawings in this manual are for illustration purposes only and in some details may differ from the actual machine. The original operation manual is in German. Other language versions are translations from German.

2. USAGE SAFETY

ATTENTION! Read all safety warnings and all instructions. Failure to follow the warnings and instructions may result in an electric shock, fire and/or serious injury or even death.

The terms "device" or "product" are used in the warnings and instructions to refer to <Digital Precision Scale >. Do not use in very humid environments or in the direct vicinity of water tanks. Do not cover the ventilation openings!

2.1. ELECTRICAL SAFETY

- The plug must fit the socket. Do not modify the plug in any way. Using original plugs and matching sockets reduces the risk of electric shock.
- Avoid touching earthed elements such as pipes, heaters, boilers and refrigerators. There is an increased risk of electric shock if the earthed device is exposed to rain, comes into direct contact with a wet surface or is operating in a damp environment. Water getting into the device increases the risk of damage to the device and of electric shock.
- Do not touch the device with wet or damp hands.
- Use the cable only for its designated use. Never use it to carry the device or to pull the plug out of a socket. Keep the cable away from heat sources, oil, sharp edges or moving parts. Damaged or tangled cables increase the risk of electric shock.
- If using the device in a damp environment cannot be avoided, a residual current device (RCD) should be applied. The use of an RCD reduces the risk of electric shock.
- Do not use the device if the power cord is damaged or shows obvious signs of wear. A damaged power cord should be replaced by a qualified electrician or the manufacturer's service centre
- To avoid electric shock, do not immerse the cord, plug or device in water or other liquids. Do not use the device on wet surfaces.

2.2. SAFETY IN THE WORKPLACE

- Do not use the device in a potentially explosive environment, for example in the presence of flammable liquids, gases or dust. The device generates sparks which may ignite dust or fumes.
- If you discover damage or irregular operation, immediately switch the device off and report it to a supervisor without delay.
- If there are any doubts as to the correct operation of the device, contact the manufacturer's support service.
- Only the manufacturer's service point may repair the device. Do not attempt any repairs independently!
- In case of fire, use a powder or carbon dioxide (CO₂) fire extinguisher (one intended for use on live electrical devices) to put it out.
- Regularly inspect the condition of the safety labels. If the labels are illegible, they must be replaced.
- Please keep this manual available for future reference. If this device is passed on to a third party, the manual must be passed on with it.

TECHNICAL DATA

Parameter description	Parameter value				
Product name	DIGITAL PRECISION SCALE				
Model	SBS-LW-300E	SBS-LW-600 001	SBS-LW-300 0001	SBS-LW-300 001	SBS-LW-6000
Min. weight [g]	0.05			0.5	
Max. weight [g]	300	600	3,000	6,000	
Division [g]	0,01			0,1	
Operating/ storage temp [°C]	0-40/10				
Operating/ storage humidity [%]	50-85				
Battery type	6V/1.3AH				
Power supply	8V/500mA				
Plate dimensions [mm]	Ø130		160x180		
Weight [kg]	1.8				

1. GENERAL DESCRIPTION

The user manual is designed to assist in the safe and trouble-free use of the device. The product is designed and manufactured in accordance with strict technical guidelines, using state-of-the-art technologies and components. Additionally, it is produced in compliance with the most stringent quality standards.

DO NOT USE THE DEVICE UNLESS YOU HAVE THOROUGHLY READ AND UNDERSTOOD THIS USER MANUAL.

To increase the product life of the device and to ensure trouble-free operation, use it in accordance with this user manual and regularly perform maintenance tasks. The technical data and specifications in this user manual are up to date. The manufacturer reserves the right to make changes associated with quality improvement.

LEGEND

	The product satisfies the relevant safety standards.
	Read instructions before use.
	The product must be recycled.
	WARNING! or CAUTION! or REMEMBER! Applicable to the given situation (general warning sign).
	Only use indoors.

- h) Keep packaging elements and small assembly parts in a place not available to children.

REMEMBER When using the device, protect children and other bystanders.

2.3. PERSONAL SAFETY

- a) Do not use the device when tired, ill or under the influence of alcohol, narcotics or medication which can significantly impair the ability to operate the device.
- b) The device is not designed to be handled by persons (including children) with limited mental and sensory functions or persons lacking relevant experience and/or knowledge unless they are supervised by a person responsible for their safety or they have received instruction on how to operate the device.
- c) To prevent the device from accidentally switching on, make sure the switch is in the OFF position before connecting to a power source.
- d) The device is not a toy. Children must be supervised to ensure that they do not play with the device.

2.4. SAFE DEVICE USE

- a) Do not use the device if the ON/OFF switch does not function properly (does not switch the device on and off). Devices which cannot be switched on and off using the ON/OFF switch are hazardous, should not be operated and must be repaired.
- b) Disconnect the device from the power supply before commencement of adjustment, cleaning and maintenance. Such a preventive measure reduces the risk of accidental activation.
- c) When not in use, store in a safe place, away from children and people not familiar with the device who have not read the user manual. The device may pose a hazard in the hands of inexperienced users.
- d) Keep the device out of the reach of children.
- e) Device repair or maintenance should be carried out by qualified persons, only using original spare parts. This will ensure safe use.
- f) To ensure the operational integrity of the device, do not remove factory-fitted guards and do not loosen any screws.
- g) Clean the device regularly to prevent stubborn grime from accumulating.
- h) The device is not a toy. Cleaning and maintenance may not be carried out by children without supervision by an adult person.
- i) It is forbidden to interfere with the structure of the device in order to change its parameters or construction.
- j) Keep the device away from sources of fire and heat.

ATTENTION! Despite the safe design of the device and its protective features, and despite the use of additional elements protecting the operator, there is still a slight risk of accident or injury when using the device. Stay alert and use common sense when using the device.

3. USE GUIDELINES

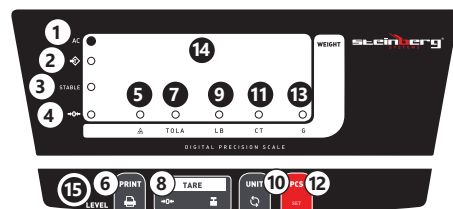
The device is intended to weigh items placed on the weighing pan. Warning! The device cannot be used for commercial purposes. The balance should be used in a space free of wind, corrosion, vibration, humidity and temperature variations. The scale should avoid vibration, violent air current, direct sunshine and high temperature

The user is liable for any damage resulting from unintended use of the device.

3.1. DEVICE DESCRIPTION

1. Cover
2. Cover hole
3. Weighing pan
4. Control panel
5. RS-232 socket
6. On / Off button
7. Power socket

1. Power indicator light
2. Tare lamp
3. Stability lamp
4. Reset lamp
5. Input sample lamp
6. Print button
7. Weight unit lamp: tola
8. Tare button
9. Weight unit lamp: pound
10. Unit of measure selection
11. Weight unit lamp: carat
12. Quantity of input sample
13. Weight unit lamp: gram
14. Display
15. Level indicator

3.2. PREPARING FOR USE APPLIANCE LOCATION

The temperature of environment must not be higher than 40°C and the relative humidity should be less than 85%. There should be at least 10 cm distance between each side of the device and the wall or other objects. The device should always be used when positioned on an even, stable, clean,

fireproof and dry surface, and be out of the reach of children and persons with limited mental and sensory functions. The power cord connected to the appliance must be properly grounded and correspond to the technical details on the product label.

3.3. DEVICE USE

NOTE: Make sure that the weighing pan is empty before you turn the device on.

IMPORTANT: Leave the device turned on for 15-30 minutes before use

- a) Select weight unit
 - Press the "UNIT" button to select the measurement unit.
- b) Tare the scale
 - Place a container or a weight on the pan.
 - Press the "TARE" button.
 - The balance will be zeroed, the tare lamp will come on.
- c) Print function
 - Connect the scale to the computer.
 - Press the "PRINT" button.
 - The weight data will be sent to the computer.
- d) Counting mode
 - If required, place the container on the weighing pan and press "TARE". While the zero function is active, press the PCS button. The "—SAP 10—" message will be displayed. Press the UNIT button to select the sample size (10, 20, 50, 100, 200, 500, 1000 pieces). After selecting the number of samples, press the TARE button to confirm. The display will show "LOAD-C". Place the previously selected number of samples on the weighing pan. After the steady reading has disappeared, the sampling procedure is completed and you can start using the piece counting function. Note: it is best to use more samples to make the count more accurate.
 - **IMPORTANT:** Calibrate the scale before first use.
- e) Calibration
 - Press and hold the TARE button for 3 seconds, the "LOAD-2" message will be displayed.
 - By using the PCS key, select "LOAD-2", "LOAD-3" or "LOAD-4".
 - » 1st calibration method: constant load range
 - 1. Press the TARE button again, a weight value will be displayed.
 - 2. Place the corresponding load on the weighing pan.
 - 3. Remove the load, the next weight value will be displayed.
 - 4. Place the corresponding weight. The calibration will be completed.
 - » 2nd calibration method: any load range
 - 1. Select "LOAD-1" using the PCS key.
 - 2. Press the TARE button to confirm.
 - 3. Select the proper value on the display.
 - 4. To enter the value on the display, press the "PCS" button and select the digit. Press "TARE" to move to the next digit.
 - 5. After entering the appropriate value, place the corresponding weight on the pan.
- f) RS-232 TRANSMISSION METHOD

This function allows the measurement data to be sent to any terminal or serial port. The device uses a Simplex asynchronous serial communication (can be changed to Duplex if necessary) of the RS232 electrical interface.

The data transmission is in ASCII code. Each data set has 10 bits. Information about the individual bits:

- BIT1: start bit
- BIT2 - BIT9: data bit
- BIT10: stop bit
- Transmission speed: default "9600" bps.

TRANSMISSION DATA FORMAT

- Bit 1 - 2: Operation mode:
 - » WT - Weighing
 - » CT - Counting
 - » PC - Percentage mode.
- Bit No. 3 - 4: Measurement status
 - » OL - Overload
 - » UL - Negative
 - » ST - Stable
 - » US - Unstable.
- Bit No. 5 - 12: The measurement value contained in 8 bits where one bit corresponds to the sign separating decimal from integer values (here "."); if the data does not take up all 8 bits, a space will be used as the fill.
- The "+" or "-" symbol is 1 bit (spaces are used instead of the "+" symbol).
- Bit 13 - 16: The current weight unit contained in 4 bits. If the unit does not take all the 4 bits in the weighing mode or in the percentage mode, the empty fields will be filled with spaces.
- Bit 17 - 18: Movement to the beginning of the current line and a new line are marked with "0DH", "0AH".

Exemplary data transmission recording:

Weighing, steady value, 38.25g:

```
W T S T | | | | 3 8 . | 2 5 | | | | g 0D 0A
```

Weighing, unsteady value, 300ct:

```
W T U S | | | | 3 0 0 | | | | c t 0D 0A
```

Counting mode, steady value, + 60:

```
C T S T | | | | 6 0 | | | | p c s 0D 0A
```

Percentage mode, steady value + 25.4:

```
P C S T | | | | 2 5 . 4 | | | | % 0D 0A
```

USE GUIDELINES

- Do NOT leave a load on the scale for long periods of time.
- Avoid extreme temperatures. Do not place the unit in direct sunlight or near air conditioning units.
- Avoid using strong electrical or magnetic fields during use.
- Avoid vibration. Do not place the device near heavy or vibrating machinery.
- Do not place the unit close to open windows, doors or fans that may cause unstable results due to air currents.
- Keep your machine clean. Do not store products on the scale when it is not being used.
- Avoid any strong shock or heavy weight. Don't put any article or give pressure over maximum capacity on the scale pan even the scale is not in use.
- Wrong wires connection of storage battery is forbidden. Right connection: Red wire should be connected with red end; Black wire should be connected with black end

3.4. CLEANING AND MAINTENANCE

- Use only non-corrosive cleaners to clean the surface.
- After cleaning the device, all parts should be dried completely before using it again.
- Store the unit in a dry, cool place, free from moisture and direct exposure to sunlight.
- Do not allow water to get inside the device through vents in the housing of the device.
- The device must be regularly inspected to check its technical efficiency and spot any damage.
- Do not leave the battery in the device if it will not be used for a longer period of time.
- Use a soft, damp cloth for cleaning.

SAFE REMOVAL OF BATTERIES AND RECHARGEABLE BATTERIES.

The devices are equipped with 6V / 1.2Ah lead-acid batteries. Recycle batteries with the appropriate organisation or company.

DISPOSING OF USED DEVICES

Do not dispose of this device in municipal waste systems. Hand it over to an electric and electrical device recycling and collection point. Check the symbol on the product, instruction manual and packaging. The plastics used to construct the device can be recycled in accordance with their markings. By choosing to recycle you are making a significant contribution to the protection of our environment. Contact local authorities for information on your local recycling facility.

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