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MIKROSKOP

## **USER MANUAL**

MICROSCOPE

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MIKROSKOP

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MIKROSKOP

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MICROSCOPE

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MICROSCOPIO

## **MANUAL DE INSTRUCCIONES**

MICROSCOPIO

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MIKROSZKÓP

**DE | EN | PL | CZ | FR | IT | ES | HU**

**SBS-MK-1**

## INHALT | CONTENT | TREŚĆ | OBSAH | CONTENU | CONTENUTO | CONTENIDO | INDHOLD

Deutsch	3
English	7
Polski	11
Česky	15
Français	19
Italiano	23
Español	27
Magyar	31

PRODUKTNAME	MIKROSKOP
PRODUCT NAME	MICROSCOPE
NAZWA PRODUKTU	MIKROSKOP
NÁZEV VÝROBKU	MIKROSKOP
NOM DU PRODUIT	MICROSCOPE
NOME DEL PRODOTTO	MICROSCOPIO
NOMBRE DEL PRODUCTO	MICROSCOPIO
TERMÉK NEVE	MIKROSKÓP
MODELL	SBS-MK-1
PRODUCT MODEL	
MODEL PRODUKTU	
MODEL VÝROBKU	
MODÈLE	
MODELLO	
MODELO	
TERMÉKMODELL	
HERSTELLER	EXPONDO POLSKA SP. Z O.O. SP. K.
MANUFACTURER	
PRODUCENT	
VÝROBCE	
FABRICANT	
PRODUTTORE	
FABRICANTE	
GYÁRTÓ	
ANSCHRIFT DES HERSTELLERS	UL. NOWY KISIELIN-INNOWACYJNA 7, 66-002 ZIELONA GÓRA   POLAND, EU
MANUFACTURER ADDRESS	
ADRES PRODUCENTA	
ADRESA VÝROBCE	
ADRESSE DU FABRICANT	
INDIRIZZO DEL PRODUTTORE	
DIRECCIÓN DEL FABRICANTE	
GYÁRTÓ CÍME	

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### 1.GENERAL DESCRIPTION

The user manual is designed to assist in the safe and trouble-free use of the device. The product is designed and manufactured in accordance with strict technical guidelines, using state-of-the-art technologies and components. Additionally, it is produced in compliance with the most stringent quality standards.

**DO NOT USE THE DEVICE UNLESS YOU HAVE THOROUGHLY READ AND UNDERSTOOD THIS USER MANUAL.**

To increase the product life of the device and to ensure trouble-free operation, use it in accordance with this user manual and regularly perform maintenance tasks. The technical data and specifications in this user manual are up to date. The manufacturer reserves the right to make changes associated with quality improvement.

### LEGEND

	The product satisfies the relevant safety standards.
	Read instructions before use.
	The product must be recycled.
	ATTENTION! Electric shock warning!
	Only use indoors.

**PLEASE NOTE!** Drawings in this manual are for illustration purposes only and in some details may differ from the actual product.

The original operation manual is written in German. Other language versions are translations from the German.

### 2. USAGE SAFETY

**ATTENTION!** Read all safety warnings and all instructions. Failure to follow the warnings and instructions may result in electric shock, fire and/or serious injury or even death.

The terms „device“ or „product“ are used in the warnings and instructions to refer to Microscope. Do not use in very humid environments or in the direct vicinity of water tanks. Prevent the device from getting wet. Risk of electric shock (when the device is powered from the mains) !

### 2.1. ELECTRICAL SAFETY

- a) The plug must fit the socket. Do not modify the plug in any way. Using original plugs and matching sockets reduces the risk of electric shock.
- b) Do not touch the device with wet or damp hands.
- c) Use the cable only for its designated use. Never use it to carry the device or to pull the plug out of a socket. Keep the cable away from heat sources, oil, sharp edges or moving parts. Damaged or tangled cables increase the risk of electric shock.
- d) If using the device in a damp environment cannot be avoided, a residual current device (RCD) should be applied. The use of an RCD reduces the risk of electric shock.
- e) Do not use the device if the power cord is damaged or shows obvious signs of wear. A damaged power cord should be replaced by a qualified electrician or the manufacturer's service centre.
- f) To avoid electric shock, do not immerse the cord, plug or device in water or other liquids. Do not use the device on wet surfaces.

### 2.2. SAFETY IN THE WORKPLACE

- a) Make sure the workplace is clean and well lit. A messy or poorly lit workplace may lead to accidents. Try to think ahead, observe what is going on and use common sense when working with the device.
- b) If you are unsure about whether the product is operating correctly or if you find damage, please contact the manufacturer's service centre.
- c) Only the manufacturer's service centre may make repairs to the product. Do not attempt to make repairs yourself!
- d) In case of fire, use a powder or carbon dioxide (CO<sub>2</sub>) fire extinguisher (one intended for use on live electrical devices) to put it out.
- e) Please keep this manual available for future reference. If this device is passed on to a third party, the manual must be passed on with it.
- f) Keep packaging elements and small assembly parts in a place not available to children.
- g) Keep the device away from children and animals.
- h) If this device is used together with another equipment, the remaining instructions for use shall also be followed.

### 2.3. PERSONAL SAFETY

- a) The device is not designed to be handled by persons (including children) with limited mental and sensory functions or persons lacking relevant experience and/or knowledge unless they are supervised by a person responsible for their safety or they have received instruction on how to operate the device.

### 2.4. SAFE DEVICE USE

- a) Do not overload the device. Use the appropriate tools for the given task. A correctly-selected device will perform the task for which it was designed better and in a safer manner.
- b) Disconnect the device from the power supply before commencement of adjustment, cleaning and maintenance. Such a preventive measure reduces the risk of accidental activation.
- c) When not in use, store in a safe place, away from children and people not familiar with the device who have not read the user manual. The device may pose a hazard in the hands of inexperienced users.

- d) Keep the device in perfect technical condition. If damage is discovered, hand over the device for repair before use.
- e) Keep the device out of the reach of children.
- f) Device repair or maintenance should be carried out by qualified persons, only using original spare parts. This will ensure safe use.
- g) To ensure the operational integrity of the device, do not remove factory-fitted guards and do not loosen any screws.
- h) Do not move, adjust or rotate the device in the course of work.
- i) Clean the device regularly to prevent stubborn grime from accumulating.
- j) The device is not a toy. Cleaning and maintenance may not be carried out by children without supervision by an adult person.
- k) It is forbidden to interfere with the structure of the device in order to change its parameters or construction.
- l) Keep the device away from sources of fire and heat.

**⚠ ATTENTION!** Despite the safe design of the device and its protective features, and despite the use of additional elements protecting the operator, there is still a slight risk of accident or injury when using the device. Stay alert and use common sense when using the device.

### 3. USE GUIDELINES

The microscope is used to observe small objects or details invisible to the unaided eye. The product is intended for home use only!

**The user is liable for any damage resulting from unintended use of the device.**

#### 3.1. DEVICE DESCRIPTION

**⚠ ATTENTION!** This product's view can be found on the last pages of the operating instructions (pp. 31).

- 1. WF 10x eyepiece
- 2. WF 16x eyepiece
- 3. Barlow lens
- 4. MicOcular
- 5. Tube
- 6. Stand
- 7. Screw
- 8. Revolver
- 9. Lens
- 10. Clips
- 11. Microscope table
- 12. LED lighting (transmitted light)
- 13. Base
- 14. Backlight selection dial
- 15. Power
- 16. Focus ring:
  - a) macrometric
  - b) micrometric
- 17. Colour filters
- 18. LED lighting (reflected light)
- 19. Disc with Ulead Photo Explorer SE software
- 20. 10 slides, 10 glass slips and 5 plastic specimen boxes
- 21. Specimens:
  - a) Yeast
  - b) „Gum Media“ adhesive
  - c) Sea salt
  - d) Shrimp eggs
- 22. Transformer with plug and power cable

- 23. Sample slicer
- 24. Shrimp hatchery
- 25. Tweezers
- 26. Pipette
- 27. Case

### 3.2. PREPARING FOR USE APPLIANCE LOCATION

The temperature of environment must not be higher than 40°C. Ensure good ventilation in the room in which the device is being used. There should be at least 10 cm distance between each side of the device and the wall or other objects. The device should always be used when positioned on an even, stable, clean, fireproof and dry surface. Position the device such that you always have access to the power plug. The power cord connected to the appliance must be properly grounded and correspond to the technical details on the product label.

### 3.3. DEVICE USE

#### 3.3.1 CORRECT MICROSCOPE ADJUSTMENT

- 1. Loosen the screw [7].
  - 2. Place the stand in a position convenient for the user.
  - 3. Adjust the microscope table to the lowest position.
  - 4. Turn the revolver until it clicks at the lowest magnification.
  - 5. Insert the 10x eyepiece into the tube.
  - 6. To double the image magnification, place a Barlow lens between the tube and eyepiece.
- NOTE: To avoid damage, always set the microscope table to the lowest position before changing the lens position.

#### 3.3.2. USING THE ELECTRIC LED LIGHTING

- 1. Connect the power cord to the socket on the back of the microscope.
- 2. Insert the plug into the power socket.

There are two diodes on the microscope. The first illuminates the sample from below, the second from above. Using the knob [14], you can select the lighting method depending on the type of object:

It.	Lighting type	Object type
I	from bottom (transmitted light)	transparent
II	from above (reflected light)	hard, non-transparent
II	from bottom and from above	half-transparent

### 3.3.3. USING COLOUR FILTERS

The colour filters [17] are located below the microscope table. This option is helpful when observing very bright objects. It makes it easier to recognize the components of colourless or transparent objects.

### 3.3.4. OBSERVING AN OBJECT

- 1. The slide should be inserted under the clips [10] on the microscope table [11] centrally under the lens [9] and above the lower LED lighting [12].

- 2. Specimen observation should be started at the lowest magnification. This will help centre the object and focus the image. The higher the magnification, the more light is needed to get a good image.
  - 3. In the next step, look into the eyepiece and gently turn the focus ring [16] until the image becomes bright and sharp. This procedure should be repeated after each change or adjustment of the eyepiece.
  - 4. To increase the magnification, install the 16x eyepiece and turn the revolver [8] to higher settings (10x or 40x).
- NOTE: Please note that the highest magnification is not the best for all specimen.

### 3.3.5. SPECIMEN PREPARATION

Place the object on the slide and use a pipette to put a few drops of distilled water on it. Place the glass slip vertically on the edge of the water droplet and slowly lower it.

### 3.3.6. SAMPLE PREPARATION - SLICES

NOTE: Samples should be prepared by adults or under their supervision. Heat a little wax or paraffin and immerse in it the object from which the sample will be prepared. Leave the object to harden. Using a sample slicer [23] (alternately a knife or a scalpel), cut very thin slices from the waxed object. The slice should be placed on a slide and covered with a glass slip.

### 3.3.7. MICROOCULAR INSTALLATION

NOTE: after the MicOcular is inserted, the zoom settings change and must be readjusted.

- 1. Remove the Barlow lens [3] and the eyepiece [1/2].
- 2. Join the MicOcular [4] with the eyepiece [1/2] and insert in the tube.
- 3. Install software and drivers.
- 4. Reboot the computer after downloading the system and connect the MicOcular to a USB port.

### 3.3.8. INSTALLING THE ULEAD PHOTOEXPLORER 7.0 SE SOFTWARE

After assembling the microscope, do not connect the MicOcular to the computer. Work with the microscope should be started from software installation.

- 1. Insert the CD into the CD drive, the menu will open automatically.
- 2. Click „Install PhotoExplorer 7.0 SE“.

- 3. Select your preferred language from the list and click „OK“.
- 4. On the welcome screen, click „Continue“.
- 5. A License Agreement will appear, which you must accept by clicking „Yes“.
- 6. Write the name and serial number from the CD packaging; the „Company Name“ field can be left blank. Click „Continue“.
- 7. In the next 4 windows select „Target Path“, „Create Icons“, „Select Program Folder“ and „TV system“ and accept the information by clicking „Continue“.
- 8. The „Setup Status“ window will appear, the program files are then copied to your computer, which may take several minutes.
- 9. When the a/m window disappears, another one will appear after a while. If the program is used in the future to display and edit photos, click „Continue“, if not select „Do not Change Current File Allocation“, then select „Continue“.
- 10. If the Internet Explorer opens, close it.
- 11. When the „Install Shield Wizard Complete“ window appears, you can click „No, I will restart computer later“ and press „Finish“. The software has been installed.

### 3.3.9 DRIVER INSTALLATION

Make sure that MicOcular is not connected to the computer.

- 1. Click „Install Bresser PC-Ocular Driver“.
- 2. Select your preferred language from the list and click „OK“.
- 3. On the welcome screen, click „Continue“.
- 4. Select „Complete“ and click „Continue“.
- 5. In the „Start Copying“ window, select „Continue“ again.
- 6. The „Setup Status“ window will appear, the program files are then copied to the computer, which may take several minutes.
- 7. When the „InstallShield Wizard Complete“ window appears press „Finish“. Drivers have been installed. Click „Exit“ in the menu.

### 3.3.10 MICROOCULAR - DISPLAYING AND SAVING PHOTOS

- 1. Turn on Ulead PhotoExplorer 7.0 SE.
  - 2. Select „Video“, then „Record Picture“ from the menu.
  - 3. The „Record Picture from Video Unit“ window will appear, click „Live View“.
  - 4. The image from the MicOcular should be visible on the screen (if it is not, click „Setup“ and select „Options“; the „Recording Unit“ window will appear. In „Select Camera“, select „Microocular“ and accept by clicking „Ok“).
  - 5. Focus again. If the image is clear click „Record Now“.
  - 6. The image has been saved to your computer.
  - 7. Click „Close“ and the photo will appear in Ulead PhotoExplorer.
- NOTE: the detailed program manual can be found on the CD in the „Document / Manual“ folder. If the folder does not include the language of the user, a message can be sent to Ulead asking for the availability of the user's language.

### 3.4. CLEANING AND MAINTENANCE

- a) Unplug the mains plug before each cleaning, adjustment or replacement of accessories, or if the device is not being used.
- b) Use only non-corrosive cleaners to clean the surface.

01.08.2022

- c) After cleaning the device, all parts should be dried completely before using it again.
- d) Do not spray the device with a water jet or submerge it in water.
- e) The device must be regularly inspected to check its technical efficiency and spot any damage.
- f) Use a soft cloth for cleaning.
- g) Do not leave the battery in the device if it will not be used for a longer period of time.
- h) Do not use sharp and/or metal objects for cleaning (e.g. a wire brush or a metal spatula) because they may damage the surface material of the appliance.
- i) Unplug the mains plug and remove the batteries before each cleaning, changing accessories and also when the device is not being used.
- j) The lens should be cleaned with a soft, smooth (lint-free) cloth.
- k) Excessive pressure during cleaning may scratch the lens.
- l) Heavily soiled lens pieces should be cleaned with a special liquid for glasses.
- m) Store the device in a special bag or the original packaging.
- n) Protect the device against dust and moisture.

#### SAFE REMOVAL OF BATTERIES AND RECHARGEABLE BATTERIES

Remove used batteries from the device using the same procedure by which you installed them. Recycle batteries with the appropriate organisation or company.

#### DISPOSING OF USED DEVICES

Do not dispose of this device in municipal waste systems. Hand it over to an electric and electrical device recycling and collection point. Check the symbol on the product, instruction manual and packaging. The plastics used to construct the device can be recycled in accordance with their markings. By choosing to recycle you are making a significant contribution to the protection of our environment. Contact local authorities for information on your local recycling facility.

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