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TASCHENWAAGE

USER MANUAL

POCKET SCALE

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KAPESNÍ VÁHA

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BALANCE DE POCHE

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BILANCIA TASCABILE

MANUAL DE INSTRUCCIONES

BALANZA DE BOLSILLO

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**SBS-DW-500HA | SBS-DW-1000HA | SBS-DW-500VAS
SBS-DW-500VA | SBS-DW-3000VA**

▪ Deutsch	3
▪ English	6
▪ Polski	9
▪ Český	12
▪ Français	15
▪ Italiano	18
▪ Español	21

PRODUKTNAME	TASCHENWAAGE
PRODUCT NAME	POCKET SCALE
NAZWA PRODUKTU	WAGA KIESZONKOWA
NÁZEV VÝROBKU	KAPESNÍ VÁHA
NOM DU PRODUIT	BALANCE DE POCHE
NOME DEL PRODOTTO	BILANCIA TASCABILE
NOMBRE DEL PRODUCTO	BALANZA DE BOLSILLO
MODELL	
PRODUCT MODEL	
MODEL PRODUKTU	SBS-DW-500HA
MODEL VÝROBKU	SBS-DW-1000HA
MODÈLE	SBS-DW-500VAS
MODELLO	SBS-DW-500VA
MODELO	SBS-DW-3000VA
IMPORTEUR	
IMPORTER	
IMPORTER	
DOVOZCE	EXPONDO POLSKA SP. Z O.O. SP. K.
IMPORTATEUR	
IMPORTATORE	
IMPORTADOR	
ADRESSE VON IMPORTEUR	
IMPORTER ADDRESS	
ADRES IMPORTERA	
ADRESA DOVOZCE	UL. NOWY KISIELIN-INNOWACYJNA 7, 66-002 ZIELONA GÓRA POLAND, EU
ADRESSE DE L'IMPORTATEUR	
INDIRIZZO DELL'IMPORTATORE	
DIRECCIÓN DEL IMPORTADOR	

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Parameter	Werte		
Produktname	TASCHENWAAGE		
Modell	SBS-DW-500HA	SBS-DW-1000HA	SBS-DW-500VAS
Abmessungen [mm]	127x79x20		115x64x17,5
Gewicht [kg]	0,120		0,100
Mindestlast [g]	0,05	0,5	0,05
Maximallast [kg]	500	1000	500
Genauigkeit [g]	0,01	0,1	0,01
Batterieparameter	1,5V 2xAAA		
Automatische Abschaltzeit [s]	300		
Produktname	TASCHENWAAGE		
Modell	SBS-DW-500VA	SBS-DW-3000VA	
Abmessungen [mm]	165x105x23	165x105x22,5	
Gewicht [kg]	0,270		
Mindestlast [g]	0,05		
Maximallast [kg]	500	3000	
Genauigkeit [g]	0,01		
Batterieparameter	1,5V 2xAAA		
Automatische Abschaltzeit [s]	300		

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1. GENERAL DESCRIPTION

The user manual is designed to assist in the safe and trouble-free use of the device. The product is designed and manufactured in accordance with strict technical guidelines, using state-of-the-art technologies and components. Additionally, it is produced in compliance with the most stringent quality standards.

DO NOT USE THE DEVICE UNLESS YOU HAVE THOROUGHLY READ AND UNDERSTOOD THIS USER MANUAL.

To increase the product life of the device and to ensure a trouble-free operation, use it in accordance with this user manual and regularly perform the maintenance tasks. The technical data and specifications in this user manual are up-to-date. The manufacturer reserves the right to make changes associated with quality improvement.

LEGEND

The product satisfies the relevant safety standards.

Read instructions before use.

- The product must be recycled.
- WARNING!** or **CAUTION!** or **REMEMBER!** Applicable to the given situation (general warning sign).
- Only use indoors.

PLEASE NOTE! Drawings in this manual are for illustration purposes only and in some details may differ from the actual machine. The original operation manual is in German. Other language versions are translations from German.

2. USAGE SAFETY

ATTENTION! Read all safety warnings and all instructions. Failure to follow the warnings and instructions may result in an electric shock, fire and/or serious injury or even death.

The terms „device“ or „product“ are used in the warnings and instructions to refer to POCKET SCALE.

2.1. SAFETY IN THE WORKPLACE

- a) Make sure the workplace is clean and well lit. A messy or poorly lit workplace may lead to accidents. Try to think ahead, observe what is going on and use common sense when working with the device.
- b) If you are unsure about whether the product is operating correctly or if you find damage, please contact the manufacturer's service centre.
- c) Only the manufacturer's service centre may make repairs to the product. Do not attempt to make repairs yourself!
- d) Please keep this manual available for future reference. If this device is passed on to a third party, the manual must be passed on with it.
- e) Keep packaging elements and small assembly parts in a place not available to children.

2.2. PERSONAL SAFETY

- a) The device is not designed to be handled by persons (including children) with limited mental and sensory functions or persons lacking relevant experience and/or knowledge unless they are supervised by a person responsible for their safety or they have received instruction on how to operate the device.
- b) The device can be handled only by physically fit persons who are capable of handling it, properly trained, familiar with this manual and trained within the scope of occupational health and safety.
- c) The device is not a toy. Children must be supervised to ensure that they do not play with the device.

2.3. SAFE DEVICE USE

- a) When not in use, store in a safe place, away from children and people not familiar with the device who have not read the user manual.
- b) Keep the device out of the reach of children.
- c) Device repair or maintenance should be carried out by qualified persons, only using original spare parts. This will ensure safe use.
- d) To ensure the operational integrity of the device, do not remove factory-fitted guards and do not loosen any screws.
- e) Clean the device regularly to prevent stubborn grime from accumulating.

- f) The device is not a toy. Cleaning and maintenance may not be carried out by children without supervision by an adult person.
- g) It is forbidden to interfere with the structure of the device in order to change its parameters or construction.
- h) Keep the device away from sources of fire and heat.
- i) Do not overload the device.

ATTENTION! Despite the safe design of the device and its protective features, and despite the use of additional elements protecting the operator, there is still a slight risk of accident or injury when using the device. Stay alert and use common sense when using the device.

3. USE GUIDELINES

The device is used for precise weight measurement. **The user is liable for any damage resulting from improper use of the device.**

3.1. DEVICE DESCRIPTION

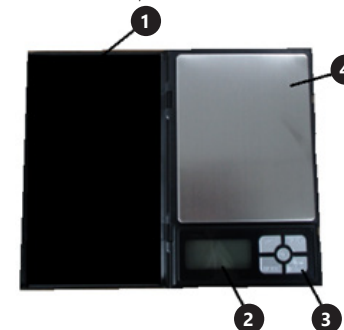
SBS-DW-500HA, SBS-DW-1000HA

SBS-DW-500VAS

SBS-DW-500VA, SBS-DW-3000VA

02.09.2020

1. Cover
2. Display
3. Control panel
4. Scale pan

3.2. DEVICE USE

3.2.1. CONTROL PANEL

Button name	Marking on the panel			Function
	SBS-DW-500HA, SBS-DW-1000HA	SBS-DW-500VAS	SBS-DW-500VA, SBS-DW-3000VA	
ON/OFF button	ON/OFF			used to turn the device on or off
The button of counting	PCS	PCS	PCS	used for counting the number of weighed items
The button of taring	TARE	TARE	TARE	used to tare; resets the counter regardless of a product placed on the pan
The button of unit changing	MODE	UNIT	MODE	used to change the measurement unit
The button of reading lock	-	-	HOLD	used to lock the reading value

CAUTION: To use the counting function, first press and hold the PCS button and then press the UNIT button to select the appropriate number of samples (25/50/75/100); place the desired number of samples on the pan and press PCS to confirm; place all objects on the pan and the display will show the quantity).

3.2.2. WEIGHTING

1. Insert the batteries into the device.
 2. Place the scale on an even, stable, dry and hard surface, away from any hot surfaces and out of the reach of children.
 3. Remove the cover.
 4. Turn the scale on.
 5. Wait until the display shows „0“.
 6. Put an object on the pan.
 7. Read the value after a few seconds when the measurement is stabilized (press UNIT/M to change the unit).
- CAUTION:** If the device is not used, it will turn off automatically after 5 minutes.

3.2.3. CALIBRATION

SBS-DW-500HA, SBS-DW-1000HA, SBS-DW-500VA, SBS-DW-3000VA

- Place the scale on an even, stable and hard surface.
- Remove the cover
- Turn the scale on.
- Press and hold the MODE unit button (until the display shows CAL).
- Briefly press the unit change button again and the display will show the number corresponding to the weight of loading in grams.
- Put appropriate weights on the pan one by one:

SBS-DW-500HA	SBS-DW-1000HA	SBS-DW-500VA	SBS-DW-3000VA
1. 300g 2. 500g	1. 1000g	1. 200g 2. 500g	1. 3000g

- When the (last) reading is stabilized, the display will show „PASS”.
CAUTION: If the weight becomes unstable or the calibration is interrupted, the display will show “Err” and will turn off automatically.
- Turn the scale off.

SBS-DW-500VAS

- Place the scale on an even, stable and hard surface.
- Remove the cover
- Turn the scale on.
- Press and hold the UNIT unit button.
- The display will show „500” (this is the weight of the loading that should be placed on the pan for calibration).
CAUTION: It is possible to change the weight by pressing the TARE button. Then the display will show „200”.
- Place a 200 g or 500 g weight on the pan according to your choice.
- When the (last) reading is stabilized, the display will show „PASS”.
CAUTION: If the weight becomes unstable or the calibration is interrupted, the display will show “Err” and will turn off automatically.
- Turn the scale off.

3.3. CLEANING AND MAINTENANCE

CAUTION: Skin contact with the pan may cause corrosion of the device, therefore it is recommended to put objects to be weighed on the pan with gloves or, if possible, with tweezers.

- Always unplug the device before cleaning it.
- Store the unit in a dry, cool place, free from moisture and direct exposure to sunlight.
- Do not spray the device with a water jet or submerge it in water.
- The device must be regularly inspected to check its technical efficiency and spot any damage.
- Use a soft cloth for cleaning.
- Do not leave the battery in the device if it will not be used for a longer period of time.
- Do not use sharp and/or metal objects for cleaning (e.g. a wire brush or a metal spatula) because they may damage the surface material of the appliance.

SAFE REMOVAL OF BATTERIES AND RECHARGEABLE BATTERIES

1.5V AAA batteries are installed in the devices. Remove used batteries from the device using the same procedure by which you installed them. Recycle batteries with the appropriate organisation or company.

DISPOSING OF USED DEVICES

Do not dispose of this device in municipal waste systems. Hand it over to an electric and electrical device recycling and collection point. Check the symbol on the product, instruction manual and packaging. The plastics used to construct the device can be recycled in accordance with their markings. By choosing to recycle you are making a significant contribution to the protection of our environment. Contact local authorities for information on your local recycling facility.

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