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LABORZENTRIFUGE

USER MANUAL

ELECTRIC CENTRIFUGE

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LABORATORNÍ CENTRIFUGA

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CENTRIFUGA DA LABORATORIO

MANUAL DE INSTRUCCIONES

CENTRÍFUGA PARA LABORATORIO

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SBS-LZ-2000SLS

Deutsch	3
English	8
Polski	13
Česky	18
Français	23
Italiano	28
Español	33

PRODUKTNAME	LABORZENTRIFUGE
PRODUCT NAME	ELECTRIC CENTRIFUGE
NAZWA PRODUKTU	WIRÓWKA LABORATORYJNA
NÁZEV VÝROBKU	LABORATORNÍ CENTRIFUGA
NOM DU PRODUIT	CENTRIFUGEUSE DE LABORATOIRE
NOME DEL PRODOTTO	CENTRIFUGA DA LABORATORIO
NOMBRE DEL PRODUCTO	CENTRÍFUGA PARA LABORATORIO
MODELL	SBS-LZ-2000SLS
PRODUCT MODEL	
MODEL PRODUKTU	
MODEL VÝROBKU	
MODÈLE	
MODELLO	
MODELO	
HERSTELLER	EXPONDO POLSKA SP. Z O.O. SP. K.
MANUFACTURER	
PRODUCENT	
VÝROBCE	
FABRICANT	
PRODUTTORE	
FABRICANTE	
ANSCHRIFT DES HERSTELLERS	UL. NOWY KISIELIN-INNOWACYJNA 7, 66-002 ZIELONA GÓRA POLAND, EU
MANUFACTURER ADDRESS	
ADRES PRODUCENTA	
ADRESA VÝROBCE	
ADRESSE DU FABRICANT	
INDIRIZZO DEL PRODUTTORE	
DIRECCIÓN DEL FABRICANTE	

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1. GENERAL DESCRIPTION

The user manual is designed to assist in the safe and trouble-free use of the device. The product is designed and manufactured in accordance with strict technical guidelines, using state-of-the-art technologies and components. Additionally, it is produced in compliance with the most stringent quality standards.

DO NOT USE THE DEVICE UNLESS YOU HAVE THOROUGHLY READ AND UNDERSTOOD THIS USER MANUAL.

To increase the product life of the device and to ensure trouble-free operation, use it in accordance with this user manual and regularly perform maintenance tasks. The technical data and specifications in this user manual are up to date. The manufacturer reserves the right to make changes associated with quality improvement. The device is designed to reduce noise emission risks to a minimum, taking into account technological progress and noise reduction opportunities.

LEGEND

	The product satisfies the relevant safety standards.
	Read instructions before use.
	The product must be recycled.
	WARNING! or CAUTION! or REMEMBER! Applicable to the given situation. (general warning sign).
	Wear protective goggles.
	Wear protective gloves.

Wear a protective apron.

ATTENTION! Electric shock warning!

ATTENTION! Rotating parts, entanglement hazard!

ATTENTION! Fire hazard - flammable materials!

WARNING! Toxic substances, danger of poisoning!

Only use indoors.

PLEASE NOTE! Drawings in this manual are for illustration purposes only and in some details may differ from the actual product.

The original operation manual is in German. Other language versions are translations from German.

2. USAGE SAFETY

ATTENTION! Read all safety warnings and all instructions. Failure to follow the warnings and instructions may result in an electric shock, fire and/or serious injury or even death.

The terms „device“ or „product“ are used in the warnings and instructions to refer to ELECTRIC CENTRIFUGE.

2.1. ELECTRICAL SAFETY

- The plug must fit the socket. Do not modify the plug in any way. Using original plugs and matching sockets reduces the risk of electric shock.
- Do not touch the device with wet or damp hands.
- Use the cable only for its designated use. Never use it to carry the device or to pull the plug out of a socket. Keep the cable away from heat sources, oil, sharp edges or moving parts. Damaged or tangled cables increase the risk of electric shock.
- If using the device in a damp environment cannot be avoided, a residual current device (RCD) should be applied. The use of an RCD reduces the risk of electric shock.
- Do not use the device if the power cord is damaged or shows obvious signs of wear. A damaged power cord should be replaced by a qualified electrician or the manufacturer's service centre.
- To avoid electric shock, do not immerse the cord, plug or device in water or other liquids. Do not use the device on wet surfaces.
- ATTENTION! DANGER TO LIFE! While cleaning, never immerse the device in water or other liquids.
- Do not use in very humid environments or in the direct vicinity of water tanks.
- Prevent the device from getting wet. Risk of electric shock!

2.2. SAFETY IN THE WORKPLACE

- Make sure the workplace is clean and well lit. A messy or poorly lit workplace may lead to accidents. Try to think ahead, observe what is going on and use common sense when working with the device.
- If you discover damage or irregular operation, immediately switch the device off and report it to a supervisor without delay.
- If there are any doubts as to the correct operation of the device, contact the manufacturer's support service.

- Only the manufacturer's service point may repair the device. Do not attempt any repairs independently!
- In case of fire, use a powder or carbon dioxide (CO₂) fire extinguisher (one intended for use on live electrical devices) to put it out.
- Children or unauthorised persons are forbidden to enter a work station.
- Use the device in a well-ventilated space.
- Please keep this manual available for future reference. If this device is passed on to a third party, the manual must be passed on with it.
- Keep packaging elements and small assembly parts in a place not available to children.
- Keep the device away from children and animals.
- If this device is used together with another equipment, the remaining instructions for use shall also be followed.

2.3. PERSONAL SAFETY

- Do not use the device when tired, ill or under the influence of alcohol, narcotics or medication which can significantly impair the ability to operate the device.
- The device is not designed to be handled by persons (including children) with limited mental and sensory functions or persons lacking relevant experience and/or knowledge unless they are supervised by a person responsible for their safety or they have received instruction on how to operate the device.
- The device can be handled only by physically fit persons who are capable of handling it, properly trained, familiar with this manual and trained within the scope of occupational health and safety.
- When working with the device, use common sense and stay alert. Temporary loss of concentration while using the device may lead to serious injuries.
- Use personal protective equipment as required for working with the device, specified in section 1 (Legend). The use of correct and approved personal protective equipment reduces the risk of injury.
- The device is not a toy. Children must be supervised to ensure that they do not play with the device.
- Do not put your hands or other items inside the device while it is in use!

2.4. SAFE DEVICE USE

- Do not overload the device. Use the appropriate tools for the given task. A correctly-selected device will perform the task for which it was designed better and in a safer manner.
- Do not use the device if the ON/OFF switch does not function properly (does not switch the device on and off). Devices which cannot be switched on and off using the ON/OFF switch are hazardous, should not be operated and must be repaired.
- Make sure the plug is disconnected from the socket before attempting any adjustments, accessory replacements or before putting the device aside. Such precautions will reduce the risk of accidentally activating the device.
- Disconnect the device from the power supply before commencement of adjustment, cleaning and maintenance. Such a preventive measure reduces the risk of accidental activation.

- When not in use, store in a safe place, away from children and people not familiar with the device who have not read the user manual. The device may pose a hazard in the hands of inexperienced users.
- Keep the device in perfect technical condition. Before each use check for general damage and especially check for cracked parts or elements and for any other conditions which may impact the safe operation of the device. If damage is discovered, hand over the device for repair before use.
- Keep the device out of the reach of children.
- Device repair or maintenance should be carried out by qualified persons, only using original spare parts. This will ensure safe use.
- To ensure the operational integrity of the device, do not remove factory-fitted guards and do not loosen any screws.
- When transporting and handling the device between the warehouse and the destination, observe the occupational health and safety principles for manual transport operations which apply in the country where the device will be used.
- Avoid situations where the device stops working during use due to excessive loading. This may result in overheating of the drive elements and damage to the device.
- Do not touch articulated parts or accessories unless the device has been disconnected from the power source.
- Do not move, adjust or rotate the device in the course of work.
- Do not leave this appliance unattended while it is in use.
- Clean the device regularly to prevent stubborn grime from accumulating.
- The device is not a toy. Cleaning and maintenance may not be carried out by children without supervision by an adult person.
- Do not run the device when empty.
- It is forbidden to interfere with the structure of the device in order to change its parameters or construction.
- Keep the device away from sources of fire and heat.
- Do not overload the device.
- Do not cover the ventilation openings!
- Keep 30 cm of free space around the device.
- Do not use force to open the lid.
- Appropriate test tubes must be used for the device.
- Tubes must not be filled up to more than 75%.

ATTENTION! Despite the safe design of the device and its protective features, and despite the use of additional elements protecting the operator, there is still a slight risk of accident or injury when using the device. Stay alert and use common sense when using the device.

3. USE GUIDELINES


The device is used to divide various mixtures, suspensions and body fluids into components of different densities. The product is outside of the scope of MDD 93/42/EEC Medical Devices regulations. It is not a medical device and should not be used for medical research and tests in hospitals and other health facilities.

The user is liable for any damage resulting from unintended use of the device.

3.1. DEVICE DESCRIPTION

Control panel

No.	Name	Function
1.	PROGRAM AND DIAL BUTTON (P / R) -	Centrifuge program and number selection
2.	SELECTION BUTTON	Speed, RCF and time parameters selection
3.	INCREASE BUTTON	Parameters increase
4.	DECREASE BUTTON	Parameters decrease
5.	STOP	Centrifugation stop
6.	START	Device activation
7.	OPERATION INDICATOR	Operation status
8.	STOP LIGHT	Operation status stop
9.	SPEED/RCF DISPLAY	Speed and RCF displaying
10.	TIME DISPLAY	Time displaying
11.	ACC./DEC.	Up and down speed control

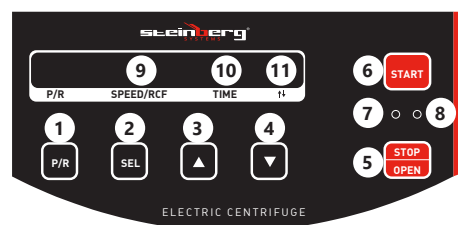3.2. PREPARING FOR USE

3.2.1. APPLIANCE LOCATION

The temperature of environment must not be higher than 40°C and the relative humidity should be less than 85%. Ensure good ventilation in the room in which the device is being used. There should be at least 10 cm distance between each side of the device and the wall or other objects. Keep the product away from hot surfaces. The device should always be used when positioned on an even, stable, clean, fireproof and dry surface, and be out of the reach of children and persons with limited mental and sensory functions. Position the device such that you always have access to the power plug. The power cord connected to the appliance must be properly grounded and correspond to the technical details on the product label.

ASSEMBLING THE DEVICE

CAUTION!

If there is the concern that the device is contaminated with toxic, pathogenic, radioactive substances or infectious test tubes with blood that are harmful to health promptly sterilize the equipment in accordance to valid laboratory methods. The device should be placed on a stable worktable, so that the four rubber feet of the device adhere well to the even surface of the table. Then, insert one power plug into the power socket at the back of the device. Connect the other end of the plug to the power socket. CAUTION! The power socket must be grounded. This will reduce the risk of electric shock.

3.3. DEVICE USE

Step 1:

- Set the power switch (in the lower right corner of the control panel) to the „I“ position, the device turns on. When the power switch is in the „O“ position, the device is turned off.

Step 2:

- Check if the round nut in the impeller is loose. If so, tighten it.

Step 3:

- Prepare test tubes. Place the test fluids in the tubes. Remember that they should be on the same level.

Step 4:

- Place the tubes symmetrically on the tube rack. Tubes must be arranged symmetrically, otherwise it will cause strong vibrations in the centrifugation process which may lead to a serious accident.

Step 5:

- Close the lid and make sure it is locked.

Step 6:

- Set the parameters.

CAUTION! The device is not designed to work with materials that can generate flammable or explosive vapors. Such substances should not be centrifuged or kept within the device.

3.3.1. IMPELLER CHAMBER

Do not pour any solutions such as water, detergent or disinfectant directly into the impeller chamber. This will damage the drive unit bearing. If the impeller chamber is wet, remove the moisture with a soft cloth or sponge. Any dirt in the impeller chamber should be cleaned with a damp cloth or a sponge using a diluted, neutral detergent. Leave the device open until it is completely dry.

3.3.2. DRIVE SHAFT (CROWN)

At least once a month, clean the inside of the impeller's drive slot (crown) and the surface of drive shaft. The crown transmits the driving force to the impeller. If the slot or the drive shaft is dirty, foreign objects should be removed immediately with a slightly damp cloth. If left unclean, the impeller may not be fitted correctly and may fall off during operation.

3.3.3. HOUSING

Always keep the housing and the worktop clean to prevent dust and other materials from entering the impeller chamber. For cleaning, use a soft cloth or sponge moistened with a diluted neutral detergent. If any harmful solution enters the equipment, take the necessary measures in accordance to appropriate laboratory methods.

3.3.4. IMPELLER

To prevent corrosion, after finishing work, remove the impeller from the chamber and remove the impeller lid to dry the inside. If any test tube is spilled inside the impeller, wash and dry it thoroughly, and then lightly apply engine grease to the impeller. Apply grease to the threaded parts of the impeller regularly.

3.3.5. PARAMETER SETTING

- Speed setting**
Press the selection button [2], „A1 4000“ will flash on the display [9], then press the increase [3] or decrease [4] button to adjust the speed. The range is from 0 to 4,000 rpm.
- Time setting**
Press the selection button [2], „10:00“ will flash on the display [9], then press the increase [3] or decrease [4] button to set the time. The range is from 0 to 99.
- Brake setting (acceleration and deceleration)**
Press the selection button [2] to pause time. BX will appear in the time display [10]. Set the brake speed using the increase [3] or decrease [4] buttons. The speed range is from b0 to b9. The speed is set at b5 at the factory.
- Starting centrifugation**
Press the START button [6], the operation indicator [7] will brighten. The device will start working. The speed display will show RPM and the time display will begin to count down the remaining time. The device takes approximately 1 minute to reach maximum speed.
- Stopping centrifugation**
When centrifugating is finished or the STOP button [5] is pressed, the time display will show „0“, the stop indicator [8] will light up and the speed display [9] will show „0“. You can open the lid and remove test tubes.

3.4. CLEANING AND MAINTENANCE

- Unplug the mains plug and allow the device to cool completely before each cleaning, adjustment or replacement of accessories, or if the device is not being used. Wait for the rotating elements to stop.
- Use only non-corrosive cleaners to clean the surface.
- After cleaning the device, all parts should be dried completely before using it again.
- Store the unit in a dry, cool place, free from moisture and direct exposure to sunlight.
- Do not spray the device with a water jet or submerge it in water.
- Do not allow water to get inside the device through vents in the housing of the device.
- The device must be regularly inspected to check its technical efficiency and spot any damage.
- Use a soft cloth for cleaning.
- Do not use sharp and/or metal objects for cleaning (e.g. a wire brush or a metal spatula) because they may damage the surface material of the appliance.

- Avoid places exposed to UV rays when working and storing the device. If installation in such a place is not possible, after finishing work, cover the device with a cloth to protect it.

SAFE REMOVAL OF BATTERIES AND RECHARGEABLE BATTERIES

1.5V AAA batteries are installed in the devices. Remove used batteries from the device using the same procedure by which you installed them. Recycle batteries with the appropriate organisation or company.

DISPOSING OF USED DEVICES

Do not dispose of this device in municipal waste systems. Hand it over to an electric and electrical device recycling and collection point. Check the symbol on the product, instruction manual and packaging. The plastics used to construct the device can be recycled in accordance with their markings. By choosing to recycle you are making a significant contribution to the protection of our environment. Contact local authorities for information on your local recycling facility.

ERROR CODE FORM

E1	Lid unlocked
E2	Device failure
E3	Parameters set incorrectly
E4	Sending error
E5	Receiving error
E6	ECC error
E7	Exceeding the speed limit

TROUBLESHOOTING

CAUTION! The device may be repaired by qualified personnel.

Problem	Possible cause	Solution
The display window does not work when the AC adapter is connected	No 220V power	Check the power
	Defective fuse	Check and replace a defective fuse
Impeller does not rotate when START button is pressed	The connector is not connected properly	Check the connection
	The power transformer was damaged	The transformer must be replaced
	The motor, although connected to the power supply, does not work. Damaged motor	Replace the motor

Centrifuge malfunction – incorrect vibrations	The pipe in the impeller has not been placed symmetrically	Check the position and arrange it properly
	Broken test tube	Check a test tube and replace it with the correct one
	The impeller has not been turned hard enough	Check the rotation of the impeller
	Damaged shock absorber	The shock absorber needs to be replaced

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