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LABORZENTRIFUGE

USER MANUAL

ELECTRIC CENTRIFUGE

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LABORATORNÍ CENTRIFUGA

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CENTRIFUGA DA LABORATORIO

MANUAL DE INSTRUCCIONES

CENTRÍFUGA PARA LABORATORIO

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SBS-LZ-5000LS | SBS-LZ-6000HS

Deutsch	3
English	8
Polski	12
Česky	17
Français	21
Italiano	26
Español	31

PRODUKTNAME	LABORZENTRIFUGE
PRODUCT NAME	ELECTRIC CENTRIFUGE
NAZWA PRODUKTU	WIRÓWKA LABORATORYJNA
NÁZEV VÝROBKU	LABORATORNÍ CENTRIFUGA
NOM DU PRODUIT	CENTRIFUGEUSE DE LABORATOIRE
NOME DEL PRODOTTO	CENTRIFUGA DA LABORATORIO
NOMBRE DEL PRODUCTO	CENTRÍFUGA PARA LABORATORIO
MODELL	
PRODUCT MODEL	
MODEL PRODUKTU	
MODEL VÝROBKU	SBS-LZ-5000LS
MODÈLE	SBS-LZ-6000HS
MODELLO	
MODELO	
HERSTELLER	
MANUFACTURER	
PRODUCENT	
VÝROBCE	EXPONDO POLSKA SP. Z O.O. SP. K.
FABRICANT	
PRODUTTORE	
FABRICANTE	
ANSCHRIFT DES HERSTELLERS	
MANUFACTURER ADDRESS	
ADRES PRODUCENTA	
ADRESA VÝROBCE	UL. NOWY KISIELIN-INNOWACYJNA 7, 66-002 ZIELONA GÓRA POLAND, EU
ADRESSE DU FABRICANT	
INDIRIZZO DEL PRODUTTORE	
DIRECCIÓN DEL FABRICANTE	

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1. GENERAL DESCRIPTION

The user manual is designed to assist in the safe and trouble-free use of the device. The product is designed and manufactured in accordance with strict technical guidelines, using state-of-the-art technologies and components. Additionally, it is produced in compliance with the most stringent quality standards.

DO NOT USE THE DEVICE UNLESS YOU HAVE THOROUGHLY READ AND UNDERSTOOD THIS USER MANUAL.

To increase the product life of the device and to ensure trouble-free operation, use it in accordance with this user manual and regularly perform maintenance tasks. The technical data and specifications in this user manual are up to date. The manufacturer reserves the right to make changes associated with quality improvement. The device is designed to reduce noise emission risks to a minimum, taking into account technological progress and noise reduction opportunities.

LEGEND

	The product satisfies the relevant safety standards.
	Read instructions before use.
	The product must be recycled.
	WARNING! or CAUTION! or REMEMBER! Applicable to the given situation. (general warning sign).
	Wear protective goggles.
	Wear protective gloves.

Wear a protective apron.

ATTENTION! Electric shock warning!

ATTENTION! Rotating parts, entanglement hazard!

ATTENTION! Fire hazard - flammable materials!

WARNING! Toxic substances, danger of poisoning!

Only use indoors.

PLEASE NOTE! Drawings in this manual are for illustration purposes only and in some details may differ from the actual product.

The original operation manual is in German. Other language versions are translations from German.

2. USAGE SAFETY

ATTENTION! Read all safety warnings and all instructions. Failure to follow the warnings and instructions may result in an electric shock, fire and/or serious injury or even death.

The terms „device“ or „product“ are used in the warnings and instructions to refer to ELECTRIC CENTRIFUGE.

2.1. ELECTRICAL SAFETY

- The plug must fit the socket. Do not modify the plug in any way. Using original plugs and matching sockets reduces the risk of electric shock.
- Avoid touching earthed elements such as pipes, heaters, boilers and refrigerators. There is an increased risk of electric shock if the earthed device is exposed to rain, comes into direct contact with a wet surface or is operating in a damp environment. Water getting into the device increases the risk of damage to the device and of electric shock.
- Do not touch the device with wet or damp hands.
- Use the cable only for its designated use. Never use it to carry the device or to pull the plug out of a socket. Keep the cable away from heat sources, oil, sharp edges or moving parts. Damaged or tangled cables increase the risk of electric shock.
- If using the device in a damp environment cannot be avoided, a residual current device (RCD) should be applied. The use of an RCD reduces the risk of electric shock.
- Do not use the device if the power cord is damaged or shows obvious signs of wear. A damaged power cord should be replaced by a qualified electrician or the manufacturer's service centre.
- To avoid electric shock, do not immerse the cord, plug or device in water or other liquids. Do not use the device on wet surfaces.
- ATTENTION! DANGER TO LIFE! While cleaning, never immerse the device in water or other liquids.
- Do not use in very humid environments or in the direct vicinity of water tanks.
- Prevent the device from getting wet. Risk of electric shock!

2.2. SAFETY IN THE WORKPLACE

- Make sure the workplace is clean and well lit. A messy or poorly lit workplace may lead to accidents. Try to think ahead, observe what is going on and use common sense when working with the device.

- If you discover damage or irregular operation, immediately switch the device off and report it to a supervisor without delay.
- If there are any doubts as to the correct operation of the device, contact the manufacturer's support service.
- Only the manufacturer's service point may repair the device. Do not attempt any repairs independently!
- In case of fire, use a powder or carbon dioxide (CO₂) fire extinguisher (one intended for use on live electrical devices) to put it out.
- Children or unauthorised persons are forbidden to enter a work station. (A distraction may result in loss of control over the device).
- Regularly inspect the condition of the safety labels. If the labels are illegible, they must be replaced.
- Please keep this manual available for future reference. If this device is passed on to a third party, the manual must be passed on with it.
- Keep packaging elements and small assembly parts in a place not available to children.
- Keep the device away from children and animals.

REMEMBER! When using the device, protect children and other bystanders.

2.3. PERSONAL SAFETY

- Do not use the device when tired, ill or under the influence of alcohol, narcotics or medication which can significantly impair the ability to operate the device.
- The device is not designed to be handled by persons (including children) with limited mental and sensory functions or persons lacking relevant experience and/or knowledge unless they are supervised by a person responsible for their safety or they have received instruction on how to operate the device.
- The device can be handled only by physically fit persons who are capable of handling it, properly trained, familiar with this manual and trained within the scope of occupational health and safety.
- When working with the device, use common sense and stay alert. Temporary loss of concentration while using the device may lead to serious injuries.
- Use personal protective equipment as required for working with the device, specified in section 1 (Legend). The use of correct and approved personal protective equipment reduces the risk of injury.
- To prevent the device from accidentally switching on, make sure the switch is on the OFF position before connecting to a power source.
- Do not wear loose clothing or jewellery. Keep hair, clothes and gloves away from moving parts. Loose clothing, jewellery or long hair may get caught in moving parts.
- The device is not a toy. Children must be supervised to ensure that they do not play with the device.
- Do not put your hands or other items inside the device while it is in use!

2.4. SAFE DEVICE USE

- Do not use the device if the ON/OFF switch does not function properly (does not switch the device on and off). Devices which cannot be switched on and off using the ON/OFF switch are hazardous, should not be operated and must be repaired.

- Disconnect the device from the power supply before commencement of adjustment, cleaning and maintenance. Such a preventive measure reduces the risk of accidental activation.
- When not in use, store in a safe place, away from children and people not familiar with the device who have not read the user manual. The device may pose a hazard in the hands of inexperienced users.
- Keep the device in perfect technical condition. Before each use check for general damage and especially check for cracked parts or elements and for any other conditions which may impact the safe operation of the device. If damage is discovered, hand over the device for repair before use.
- Keep the device out of the reach of children.
- Device repair or maintenance should be carried out by qualified persons, only using original spare parts. This will ensure safe use.
- To ensure the operational integrity of the device, do not remove factory-fitted guards and do not loosen any screws.
- When transporting and handling the device between the warehouse and the destination, observe the occupational health and safety principles for manual transport operations which apply in the country where the device will be used.
- Avoid situations where the device stops working during use due to excessive loading. This may result in overheating of the drive elements and damage to the device.
- Do not touch articulated parts or accessories unless the device has been disconnected from the power source.
- Do not move, adjust or rotate the device in the course of work.
- Do not leave this appliance unattended while it is in use.
- Clean the device regularly to prevent stubborn grime from accumulating.
- The device is not a toy. Cleaning and maintenance may not be carried out by children without supervision by an adult person.
- It is forbidden to interfere with the structure of the device in order to change its parameters or construction.
- Keep the device away from sources of fire and heat.
- Do not overload the device.
- Do not cover the ventilation openings!
- A flammable, explosive or corrosive test sample may damage or shorten the life of the device. Any damage caused by the above-mentioned type test samples are not covered by the warranty.
- The maximum rotor speed must not be exceeded.
- The condition of the rotor should be frequently checked (for cracks, corrosion, damaged seals, scratches, etc.). Do not use the device if the rotor is not in good technical condition.
- You must never open the device cover while it is working! The lid can be opened with the device switched off and the rotor at a standstill.
- Uneven alignment or use of damaged test tubes can cause severe vibrations and damage the device.
- For the user's safety, observe min. 30cm distance between the operating device and other objects and devices or people.

- y) Regularly clean the device's chamber as recommended. Dirt on the drive shaft can cause the rotor to fall and get damaged.
- z) Remove the rotor before moving the machine to avoid damaging the drive shaft.

⚠ ATTENTION! Despite the safe design of the device and its protective features, and despite the use of additional elements protecting the operator, there is still a slight risk of accident or injury when using the device. Stay alert and use common sense when using the device.

3. USE GUIDELINES

The device is used to divide various mixtures, suspensions and body fluids into components of different densities.

The product is outside of the scope of MDD 93/42/EEC Medical Devices regulations. It is not a medical device and should not be used for medical research and tests in hospitals and other health facilities.

The user is liable for any damage resulting from unintended use of the device.

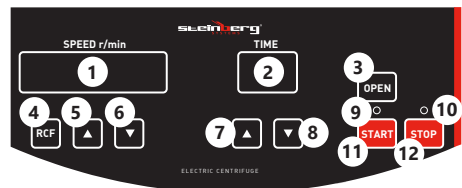
3.1. DEVICE DESCRIPTION

1. Lid
2. Rotor
3. On / Off button
4. Feet
5. Vents
6. Control panel

NOTE: The power socket is located on the back of the device (not shown in the figure above).

CONTROL PANEL DESCRIPTION

1. Display
2. Timer display
3. Lid release button
4. RCF button
5. „Up“ button for speed settings
6. „Down“ button for speed setting
7. „Up“ button for run time setting
8. „Down“ button for run time setting
9. Device operation LED
10. Device stop LED
11. START button
12. STOP button

3.2. PREPARING FOR USE

APPLIANCE LOCATION

The ambient temperature must not exceed 45°C and the relative humidity should not exceed 85%. Only use the device in well-ventilated areas. Do not obstruct the air outlet of the device. Keep the device away from any hot surfaces. The device should always be used when positioned on an even, stable, clean, fireproof and dry surface, and be out of the reach of children and persons with limited mental and sensory functions. The valve that disables the compressed air supply should be immediately accessible from the area where the device is being used. Make sure that the pressure of the air supplying the device does not exceed the recommended range!

ASSEMBLING THE DEVICE

Place the device on a flat, stable surface so that all four rubber feet are firmly set on the surface. Connect the power cord to the device and its other end to a 230VAC socket. Please note: The power socket must be grounded. This will reduce the risk of electric shock. Do not use excessive force when installing the rotor. Do it with care, so as not to damage the rotor or the drive shaft.

3.3. DEVICE USE

1. Check that the nut on the rotor has not loosened, if so, tighten it.
2. Get the test tubes with a preparation ready for centrifugation and place them in the device so that they meet the following premises:
 - the test tube should be filled to a maximum of 70% of its total volume,
 - place the test tubes symmetrically in the rotor.
 - The sample-filled tubes placed opposite each other must be of equal weight.
 - Always verify the condition of the test tube used. Do not insert cracked or deformed test tubes in the device!
3. Close the device lid.
4. Switch the device connected to the power supply on, using the ON/OFF button at the back of the device.
5. Set the device parameters.
 - Speed setting
 Set the speed value using buttons (5) and (6), the current value will be shown on the display (1). When the displayed value flashes twice, the setting is saved.

The maximum speed must not be exceeded.

Please note: If the density of the sample intended for centrifugation exceeds 1.2 g/ml, the maximum speed value should be recalculated acc. to the formula below:

$$N = n \sqrt{(1.2c / S)}$$

Where,

N - new maximum rotation speed

n- original maximum rotation speed

S- density of the tested sample

- Setting the maximum relative centrifugal force (RCF) Press the RCF button (4), the display (1) shows the value of the centrifugal force. Change the value using buttons (5) and (6). When the displayed value flashes twice, the setting is saved.
 - Rotor selection Press button (4) until the display shows the rotor number. The last two digits of the number are displayed, i.e. for the rotor no. 31277, the displayed setting is 77. When replacing the rotor with a different model, the number on the device must be changed.
 - Time setting Set the timer using buttons (7) and (8), the current value will be displayed on the display (2). When the displayed value flashes twice, the setting is saved.
 - Set the acceleration and deceleration values To enter this option, the device must not be working. Press and hold the STOP button (12) until the device beeps and the time display (2) shows a value such as „b8“. Set the value using buttons (7) and (8). The device has 10 speed settings (range from 0 to 9). The higher the value is selected, the faster the rate of deceleration and acceleration will be. After setting the parameter, wait approx. 5 seconds, the value will be saved automatically. Note: The device is factory set to either „b4“ or „b5“.
6. The speed value is shown as standard on the display (1). By pressing the RCF button (4), the following parameter values will appear in turn on the display: the rotor number, rotational speed and relative centrifugal force.
 7. To start the process, press the START button (11), and the power LED (9) lights up green. The display will show the current speed value (the set value will be reached in about 1-2 minutes), and the countdown time to the end of spinning.
 8. To stop the device, press the STOP (12) button or wait until the timer counts down to the value 0. The LED (10) will turn red. The device will not be fully stopped until the actual speed value is 0 on the display. The device will beep after it has stopped completely.
 9. Release the lid by pressing the button (3). You will hear a humming sound when the lid is unlocked, then you can open it.
 10. Remove the test tubes with centrifuged samples.
 11. Switch the device off using the On/Off button. Even after the device is turned off, it is still live. Disconnect the device from the power source when not in use.
 12. After finishing work, check that there is no condensation in the chamber of the device. Condensation should be removed as soon as possible with a sponge or a hose to prevent it from entering the drive unit.
 13. Leave the lid open until the chamber is completely dry.

3.4. CLEANING AND MAINTENANCE

- a) Always unplug the device before cleaning it.
- b) Use only non-corrosive cleaners to clean the surface.
- c) After cleaning the device, all parts should be dried completely before using it again.
- d) Store the unit in a dry, cool place, free from moisture and direct exposure to sunlight.
- e) Do not spray the device with a water jet or submerge it in water.
- f) Do not allow water to get inside the device through vents in the housing of the device.
- g) Clean the vents with a brush and compressed air.
- h) The device must be regularly inspected to check its technical efficiency and spot any damage.
- i) Use a soft, damp cloth for cleaning.
- j) Do not use sharp and/or metal objects for cleaning (e.g. a wire brush or a metal spatula) because they may damage the surface material of the appliance.
- k) Do not clean the device with an acidic substance, agents of medical purposes, thinners, fuel, oils or other chemical substances because it may damage the device.
- l) If there is the concern that the device is contaminated with toxic, pathogenic, radioactive substances or infectious test tubes with blood that are harmful to health promptly sterilize the equipment in accordance to valid laboratory methods.
- m) Periodically lubricate the central part of the rotor.
- n) Avoid exposing the device to UV rays during operation or storage. UV rays can cause discoloration of the casing and the peeling of the surface. It is recommended to cover the device during storage.
- o) The device chamber, including the rotor drive, should be cleaned once a month.
- p) Never pour water, detergents, disinfectants, etc. into the chamber of the device. This can damage the device and cause corrosion.

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