USER MANUAL

Bedienungsanleitung
Instrukcja obsługi
Návod k použití
Manuel d'utilisation
Istruzioni per l'uso Manual
de instrucciones Használati
útmutató Brugsanvisning
Käyttöohje
Gebruiksaanwijzing
Bruksanvisning Instruções
de utilização Používateľská
príručka Ръководство за
употреба Οδηγίες χρήσης
Upute za uporabu
Naudojimo instrukcija
Manual de utilizare
Navodila za uporabo

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1. GENERAL DESCRIPTION

The user manual is designed to assist in the safe and trouble-free use of the device. The product is designed and manufactured in accordance with strict technical guidelines, using state-of-the-art technologies and components. Additionally, it is produced in compliance with the most stringent quality standards.

DO NOT USE THE DEVICE UNLESS YOU HAVE THOROUGHLY READ AND UNDERSTOOD THIS USER MANUAL.

To increase the product life of the device and to ensure trouble-free operation, use it in accordance with this user manual and regularly perform maintenance tasks. The technical data and specifications in this user manual are up to date. The manufacturer reserves the right to make changes associated with quality improvement. The device is designed to reduce noise emission risks to a minimum, taking into account technological progress and noise reduction opportunities.

LEGEND

	The product satisfies the relevant safety standards.
	Read instructions before use.
	The product must be recycled.
	WARNING! or CAUTION! or REMEMBER! Applicable to the given situation. (general warning sign)
	Wear protective gloves.
	ATTENTION! Electric shock warning!
	ATTENTION! Hot surface, risk of burns!
	Only use indoors.

PLEASE NOTE! Drawings in this manual are for illustration purposes only and in some details may differ from the actual product.

The original operation manual is written in German. Other language versions are translations from the German.

2. USAGE SAFETY

ATTENTION! Read all safety warnings and all instructions. Failure to follow the warnings and instructions may result in electric shock, fire and/or serious injury or even death.

The terms „device“ or „product“ are used in the warnings and instructions to refer to ELECTRO THERMOSTATIC AIR DRYING OVEN.

2.1. ELECTRICAL SAFETY

- The plug must fit the socket. Do not modify the plug in any way. Using original plugs and matching sockets reduces the risk of electric shock.
- Avoid touching earthed elements such as pipes, heaters, boilers and refrigerators. There is an increased risk of electric shock if the earthed device is exposed to rain, comes into direct contact with a wet surface or is operating in a damp environment. Water getting into the device increases the risk of damage to the device and of electric shock.
- Do not touch the device with wet or damp hands.
- Use the cable only for its designated use. Never use it to carry the device or to pull the plug out of a socket. Keep the cable away from heat sources, oil, sharp edges or moving parts. Damaged or tangled cables increase the risk of electric shock.
- If using the device in a damp environment cannot be avoided, a residual current device (RCD) should be applied. The use of an RCD reduces the risk of electric shock.
- Do not use the device if the power cord is damaged or shows obvious signs of wear. A damaged power cord should be replaced by a qualified electrician or the manufacturer's service centre.
- To avoid electric shock, do not immerse the cord, plug or device in water or other liquids. Do not use the device on wet surfaces.
- ATTENTION! DANGER TO LIFE! While cleaning, never immerse the device in water or other liquids.
- Prevent the device from getting wet. Risk of electric shock!

2.2. SAFETY IN THE WORKPLACE

- Make sure the workplace is clean and well lit. A messy or poorly lit workplace may lead to accidents. Try to think ahead, observe what is going on and use common sense when working with the device.
- Do not use the device in a potentially explosive environment, for example in the presence of flammable liquids, gases or dust. The device generates sparks which may ignite dust or fumes.
- If you discover damage or irregular operation, immediately switch the device off and report it to a supervisor without delay.
- If there are any doubts as to the correct operation of the device, contact the manufacturer's support service.
- Only the manufacturer's service point may repair the device. Do not attempt any repairs independently!
- In case of fire, use a powder or carbon dioxide (CO2) fire extinguisher (one intended for use on live electrical devices) to put it out.

- Children or unauthorised persons are forbidden to enter a work station. (A distraction may result in loss of control over the device).
- Please keep this manual available for future reference. If this device is passed on to a third party, the manual must be passed on with it.
- Keep packaging elements and small assembly parts in a place not available to children.
- Keep the device away from children and animals.
- If this device is used together with another equipment, the remaining instructions for use shall also be followed.

REMEMBER! When using the device, protect children and other bystanders.

2.3. PERSONAL SAFETY

- Do not use the device when tired, ill or under the influence of alcohol, narcotics or medication which can significantly impair the ability to operate the device.
- The device is not designed to be handled by persons (including children) with limited mental and sensory functions or persons lacking relevant experience and/or knowledge unless they are supervised by a person responsible for their safety or they have received instruction on how to operate the device.
- The device can be handled only by physically fit persons who are capable of handling it, properly trained, familiar with this manual and trained within the scope of occupational health and safety.
- When working with the device, use common sense and stay alert. Temporary loss of concentration while using the device may lead to serious injuries.
- Use personal protective equipment as required for working with the device, specified in section 1 (Legend). The use of correct and approved personal protective equipment reduces the risk of injury.
- To prevent the device from accidentally switching on, make sure the switch is on the OFF position before connecting to a power source.
- The device is not a toy. Children must be supervised to ensure that they do not play with the device.

2.4. SAFE DEVICE USE

- Do not use the device if the ON/OFF switch does not function properly (does not switch the device on and off). Devices which cannot be switched on and off using the ON/OFF switch are hazardous, should not be operated and must be repaired.
- Make sure the plug is disconnected from the socket before attempting any adjustments, accessory replacements or before putting the device aside. Such precautions will reduce the risk of accidentally activating the device.
- When not in use, store in a safe place, away from children and people not familiar with the device who have not read the user manual. The device may pose a hazard in the hands of inexperienced users.
- Keep the device in perfect technical condition. Before each use check for general damage and especially check for cracked parts or elements and for any other conditions which may impact the safe operation of the device. If damage is discovered, hand over the device for repair before use.

- Keep the device out of the reach of children.
- Device repair or maintenance should be carried out by qualified persons, only using original spare parts. This will ensure safe use.
- To ensure the operational integrity of the device, do not remove factory-fitted guards and do not loosen any screws.
- When transporting and handling the device between the warehouse and the destination, observe the occupational health and safety principles for manual transport operations which apply in the country where the device will be used.
- Do not move, adjust or rotate the device in the course of work.
- Do not leave this appliance unattended while it is in use.
- Clean the device regularly to prevent stubborn grime from accumulating.
- The device is not a toy. Cleaning and maintenance may not be carried out by children without supervision by an adult person.
- It is forbidden to interfere with the structure of the device in order to change its parameters or construction.
- Keep the device away from sources of fire and heat.
- Do not overload the device.
- Do not cover the ventilation openings! Flammable objects must not be put into the device.
- Objects inside the device should not be placed too close to each other. Adequate space should be left to facilitate circulation of the hot air.

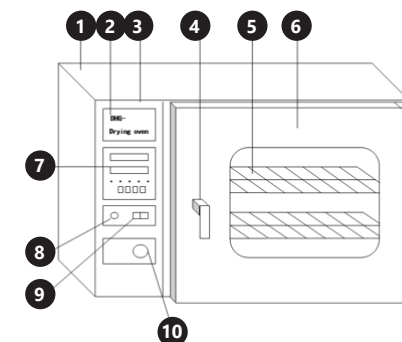
ATTENTION! Despite the safe design of the device and its protective features, and despite the use of additional elements protecting the operator, there is still a slight risk of accident or injury when using the device. Stay alert and use common sense when using the device.

3. USE GUIDELINES

The SBS-ADO-1000, SBS-ADO-2000 drying oven is a device used for gentle and thorough drying of various materials. The SBS-ADO-1000, SBS-ADO-2000 drying oven are products outside the scope of MDD 93/42 / EEC regulations for medical devices.


The user is liable for any damage resulting from unintended use of the device.

3.1. DEVICE DESCRIPTION

1. Housing
2. Nameplate
3. Control panel
4. Door handle
5. Shelf
6. Door
7. Temperature regulation
8. Power light
9. ON/OFF switch
10. Air vacuum valve

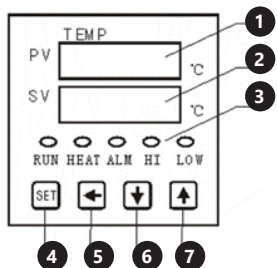Temperature control panel:

1. Temperature inside the device
2. Set temperature
3. Indicators:
 - RUN – operation indicator
 - HEAT – heating indicator
 - ALM – alarm indicator
 - HI – high speed indicator
 - LOW – low speed indicator
4. SET function button
5. SHIFT button
6. „Down“ „-“ setting button
7. „Up“ „+“ setting button

Devices have:

- precise temperature control: the temperature can be set every 10°C,
- a display showing the set temperature and the temperature inside the device,
- an audible and visual alarm to indicate a higher temperature inside the device than that set on the control panel. The device will automatically turn off the heater and activate an alarm,
- an air vacuum valve (with an adjustment knob you can regulate the inflow and outflow of air from inside the device).

PLEASE NOTE! Device is equipped with temperature safety switch – it turns off power when temperature inside device exceeds temperature value set on the switch. In each case switch should be set to a temperature value of 10°C higher than temperature set on main control panel.

3.2. PREPARING FOR USE

APPLIANCE LOCATION

The temperature of environment must not be higher than 35°C and the relative humidity should be less than 85%. Ensure good ventilation in the room in which the device is being used. A minimum distance of 20 cm to the rear wall, 15 cm to the sides and 35 cm upwards must be maintained. The device should always be used when positioned on an even, stable, clean, fireproof and dry surface, and be out of the reach of children and persons with limited mental and sensory functions. Position the device such that you always have access to the power plug. The power cord connected to the appliance must be properly grounded and correspond to the technical details on the product label. Disassemble the device and all its components and clean them before the first use.

3.3. DEVICE USE

BASIC PRINCIPLES

1. Place the items to be dried inside the device and close the door securely.
2. Set the power switch to the ON position. The power indicator light will illuminate and the numbers will appear on the temperature controller.
3. Set the temperature. The display will show the temperature set by the user and the temperature inside the device. The heating ON indicator light will come on. This means that the device is heating and raising the temperature. After a while, when the set temperature is close to the temperature inside the device, the indicator light starts blinking. Under normal conditions, after heating for 90 minutes, the device will only maintain the set temperature.
4. If the required operating temperature is low, e.g. 80°C, the two-step temperature setting method should be used. This method will prevent the temperature from dropping and allow the device to reach a constant temperature faster. For example, if the desired temperature is 80°C, set the temperature to 70°C first. After heating the device to the set temperature, set the target temperature, i.e. 80°C.
5. The drying time should be adjusted according to the moisture content of different items. If the items to be dried are very humid, turn the air volume control to the „≡“ position. Then the humid air inside the device will be exhausted.
6. After completing the drying process, leave the air supply open, set the power switch to the OFF position, and then open the door. Be careful not to burn yourself! If the user does not want to take the items out after the drying process is finished, just shut off the air supply and items inside the device will be kept in safe conditions.

PROCEDURE SHOWING THE COURSE BETWEEN THE FUNCTIONS

ATTENTION! This product's view can be found on the last pages of the operating instructions (pp. 39).

CONTROLLER OPERATION

When the temperature controller is on, the PV display shows the value of measured temperature, and the SV display shows the preset temperature value.

BASIC SETTINGS:

Temperature setting: Press the SET button, the PV display shows „SP“ the SV displays shows the adjustable temperature value. When last digit blinks, set its value with the ▲ and ▼ buttons. Use the button ◀ to switch to the next digit and set its value using ▲ and ▼. After setting the temperature, press the SET button to set the timer.

Timer setting: Press the SET button, the PV display shows „SI“ the SV displays shows the adjustable timer value. When last digit blinks, set its value with the ▲ and ▼ buttons. Use the ◀ button to switch to the next digit and set its value using ▲ and ▼. After setting the time, press the SET button for the device to start operation with the defined parameters.

ADVANCED SETTINGS:

CAUTION: All the controllers and parameters have been thoroughly tested and set by the manufacturer. If the functional parameters of the vacuum dryer meet your expectations and if the vacuum dryer runs properly, there is no need to modify these parameters.

How to open the advanced settings: Press and hold ▲, then press SET. The display will show „LCK“. Release the SET button (while pressing ▲) and set the value of the blinking digit to „7“. Release ▲ and press SET. The display will show „AL1“.

- AL1 – temperature deviation alarm – set the required value with the ▲ and ▼ buttons. Adjustable values: range between -99,9°C and 999,9°C. The device will trigger an alarm when the sum of the set temperature and the set temperature deviation exceeds the actual temperature value. After setting the temperature deviation value, press SET to enter the high/low speed „FAN“ setting.
- FAN – high/low speed control. Adjustable values: range between -99,9°C and 999,9°C. After setting, press the SET button to go to the settings for correction of the measured temperature value „SC“. After setting, press SET in order to adjust the measured temperature value „SC“.
- SC – measured temperature value adjustment – the PV display shows the adjusted temperature value. Available values: -99,9°C to +100°C. After setting the measured temperature adjustment value, press SET in order to adjust the PID controller self-regulation – the display shows „ATU“.
- ATU – enable or disable self-regulation of the PID controller – available values: „0“ – PID self-regulation disabled; „1“ – PID self-regulation enabled. When this parameter is set to „1“, the device can be started in self-regulation mode. In order to start the self-regulation mode, leave the advanced settings and, in the normal operation mode, press and hold ▼ for about 5 seconds. When this mode is on, the SV displays blinks and shows „ATU“. During this process the actual temperature may differ from the preset one – this is normal. The duration of the self-regulation function is about 3 to 5 minutes. After this time the PID controller will adjust its settings, in order to work more precisely. After setting, press SET in order to

adjust the proportionality range “P”.

- P – proportionality range – heating control adjustment. Available values: from 0 to 100.0 After setting, press SET in order to adjust the integration time constant “P”.
- I – integration time constant. Available values: from 1 to 4320 After setting, press SET in order to adjust the differentiation time constant “d”.
- d – differentiation time constant. Available values: from 0 to 1200 After setting, press SET in order to adjust the heating period “P”.
- T – heating time. Available values: from 1 to 60 After setting, press SET in order to adjust the thermal accuracy control range “HL”.
- HL – thermal accuracy control range. Available values: from 500 to 1.500 After setting, press the SET button to enter the „OUT“ setting of the heater output power.
- OUT – output power of the heating body. The values are adjustable from 0 up to 100. After setting, press and hold SET for about 5 seconds in order to leave the advanced settings and save the changes.

3.4. CLEANING AND MAINTENANCE

- a) Unplug the mains plug and allow the device to cool completely before each cleaning, adjustment or replacement of accessories, or if the device is not being used.
- b) Use only mild, food-safe detergents to wash the device.
- c) After cleaning the device, all parts should be dried completely before using it again.
- d) Store the unit in a dry, cool place, free from moisture and direct exposure to sunlight.
- e) Do not spray the device with a water jet or submerge it in water.
- f) Do not allow water to get inside the device through vents in the housing of the device.
- g) Clean the vents with a brush and compressed air.
- h) The device must be regularly inspected to check its technical efficiency and spot any damage.
- i) Use a soft, damp cloth for cleaning.
- j) Do not use sharp and/or metal objects for cleaning (e.g. a wire brush or a metal spatula) because they may damage the surface material of the appliance.
- k) Do not clean the device with an acidic substance, agents of medical purposes, thinners, fuel, oils or other chemical substances because it may damage the device.

DISPOSING OF USED DEVICES

Do not dispose of this device in municipal waste systems. Hand it over to an electric and electrical device recycling and collection point. Check the symbol on the product, instruction manual and packaging. The plastics used to construct the device can be recycled in accordance with their markings. By choosing to recycle you are making a significant contribution to the protection of our environment. Contact local authorities for information on your local recycling facility.

TROUBLESHOOTING

Problem	Possible cause	Solution
No power	The plug is inserted incorrectly or the cord is disconnected.	Insert the plug/ connect the cord.
	Blown fuse.	Replace the fuse.
The temperature in the device does not rise	The set temperature is too low.	Adjust set temperature.
	The electric heater is not working.	Replace the electric heater.
	The temperature controller is not working.	Replace the temperature controller.
The set temperature differs significantly from the temperature in the device	Cycle fan does not work.	Replace the fan.
	The temperature sensor is not working.	Replace the temperature sensor.
The set temperature differs significantly from the temperature in the device	Fine tuning potentiometer for setting the temperature is not well adjusted.	Adjust the potentiometer.
	The set temperature is too low.	Adjust set temperature.
Overtemperature alarm is invalid.	The temperature controller is not working.	Replace the temperature controller.

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