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USER MANUAL

VACUUM PUMP

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POMPE À VIDE

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POMPA A VUOTO

MANUAL DE INSTRUCCIONES

BOMBA DE VACÍO

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SBS-VP-200

DE	Produktname:	VAKUUMPUMPUMPE
EN	Product name:	VACUUM PUMP
PL	Nazwa produktu:	POMPA PRÓŻNIOWA
CZ	Název výrobku	VAKUOVÁ PUMPA
FR	Nom du produit:	POMPE À VIDE
IT	Nome del prodotto:	POMPA A VUOTO
ES	Nombre del producto:	BOMBA DE VACÍO
HU	Termék neve	VÁKUUMSZIVATTYÚ
DA	Produktnavn	VAKUUMPUMPE
DE	Modell:	SBS-VP-200
EN	Product model:	
PL	Model produktu:	
CZ	Model výrobku	
FR	Modèle:	
IT	Modello:	
ES	Modelo:	
HU	Modell	
DA	Model	
DE	Hersteller	expondo Polska sp. z o.o. sp. k.
EN	Manufacturer	
PL	Producent	
CZ	Výrobce	
FR	Fabricant	
IT	Produttore	
ES	Fabricante	
HU	Termelő	
DA	Producent	
DE	Anschrift des Herstellers	ul. Nowy Kisielin – Innowacyjna 7, 66-002 Zielona Góra Poland, EU
EN	Manufacturer Address	
PL	Adres producenta	
CZ	Adresa výrobce	
FR	Adresse du fabricant	
IT	Indirizzo del produttore	
ES	Dirección del fabricante	
HU	A gyártó címe	
DA	Producentens adresse	

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1. General Description

The instruction manual is intended to assist in safe and reliable use. The product is designed and manufactured strictly according to technical specifications using the latest technology and components and maintaining the highest quality standards.

**PLEASE CAREFULLY READ
AND UNDERSTAND THIS INSTRUCTION MANUAL BEFORE
OPERATION,**

To ensure long and reliable operation of the device, make sure to operate and maintain it properly in accordance with the guidelines in this instruction manual. The technical data and specifications contained in this instruction manual are up to date. The manufacturer reserves the right to make changes in order to improve the quality. Taking the technical progress and the possibility of reducing noise into account, the device is designed and built in such a way so that risks resulting from noise emissions are reduced to the lowest possible level.

Explanation of symbols

	<p>The product complies with applicable safety standards.</p>
	<p>Please read the instructions before use.</p>
	<p>Recyclable product.</p>
	<p>CAUTION! or WARNING! or REMINDER! describing a situation. (general warning sign).</p>
	<p>CAUTION! Warning of electric shock!</p>

CAUTION! The illustrations in this instruction manual are for reference only and may differ from the actual product in some details.

The original instruction manual is in the German language version. Other language versions are translations from German.

2. Safety of use

 **CAUTION!** Read all safety warnings and all instructions. Failure to follow the warnings and instructions may result in electric shock, fire and/or severe personal injury or death.

The term "device" or "product" in the warnings and in the description of the instructions refers to the VACUUM PUMP

Electrical safety

- a) The plug of this device must fit into the outlet. Do not modify the plug in any way. Original plugs and matching outlets reduce the risk of electric shock.

-
- b) Avoid touching grounded parts, such as pipes, heaters, ovens, and refrigerators. There is an increased risk of electric shock if your body is grounded and touches the device exposed to direct rain, wet pavement, and operation in a humid environment. If water enters the device, there is an increased risk of damage to the device and electric shock.
 - c) Do not touch the device with wet or damp hands.
 - d) Do not use the cord in an unintended manner. Never use it to carry the device or to pull the plug out of the socket. Keep the cord away from heat sources, oil, sharp edges or moving parts. Damaged or tangled cords increase the risk of electric shock.
 - e) If you cannot avoid using the device in a wet environment, use a residual current device (RCD). Using an RCD reduces the risk of electric shock.
 - f) Do not use the device if the power cord is damaged or shows signs of wear. A damaged power cord should be replaced by a qualified electrician or the manufacturer's service department.
 - g) To avoid electric shock, do not immerse the cable, plug, or device itself in water or other liquid. Do not use the device on wet surfaces.

Safety in the workplace

- a) Keep the work area tidy and well lit. Disorder or poor lighting can lead to accidents. Be foresighted, watch what you are doing and use common sense when using the device.
- b) Do not use the appliance in an explosive area, for example in the presence of flammable liquids, gases or dust.
- c) If you find any damage or irregularities in the operation of the device, immediately turn it off and report it to an authorized person.
- d) • If you have any doubts as to whether the product is working properly or if it is damaged, contact the manufacturer's service department.
- e) Only the manufacturer's service department can repair the device. Do not carry out repairs yourself!
- f) In case of open flames or fire, use only dry powder or snow (CO₂) fire extinguishers to extinguish the live equipment.
- g) Keep these instructions for use for future reference. If the device is to be passed on to a third party, the operating instructions must also be handed over together with the device.
- h) Keep the packaging and small assembly parts out of the reach of children.
- i) Keep the device away from children and animals.
- j) When using this device together with other devices, also follow the other instructions for use.

Personal safety

- a) Do not operate this product if you are tired, ill or under the influence of alcohol, drugs or medication that could impair your ability to operate the product.
- b) The device is not intended to be used by persons (including children) with reduced mental, sensory or intellectual functions or persons who lack experience and/or knowledge unless they are supervised or have been instructed by a person responsible for their safety on how to operate the device.
- c) Use caution and common sense when operating this device. A moment's inattention during operation may result in serious personal injury.
- d) To prevent accidental start-up, make sure the switch is in the off position before connecting to a power source.
- e) The device is not a toy. Children should be watched to ensure that they do not play with the device.
- f) Do not place your hands or any objects inside the running device!

Safe use of the device

- a) Do not overload the device. Use tools that are suitable for the application. A correctly selected device will do a better and safer job for which it was designed.
- b) Do not use the device if the ON/OFF switch does not function properly (does not turn on and off). Devices that cannot be controlled by the switch are unsafe, cannot operate, and must be repaired.
- c) Unplug the device before making adjustments, changing accessories, or putting it away. This precaution reduces the risk of accidental start-up.
- d) Keep unused equipment out of the reach of children and out of the reach of anyone unfamiliar with the device or this instruction manual. These devices are dangerous in the hands of inexperienced users.
- e) Keep the device in good working condition. Check before each use for general damage or damage to moving parts (cracks in parts and components or any other condition that may affect the safe operation of the device). If damaged, have the device repaired before use.
- f) Keep the device out of the reach of children.
- g) Repairs and maintenance should be carried out by qualified personnel using only original spare parts. This will ensure the safety of use.
- h) To ensure the designed operational integrity of the device, do not remove factory-installed covers or loosen screws.

-
- i) When transporting or moving the device from storage to the place of use, observe the health and safety rules for manual handling applicable in the country where the device is used.
 - j) Do not move, shift, or rotate the machine while in operation.
 - k) Do not leave the device switched on unattended.
 - l) Clean the device regularly to prevent permanent dirt build-up.
 - m) The device is not a toy. Cleaning and maintenance must not be performed by children without adult supervision.
 - n) Do not tamper with the device to alter its performance or design.
 - o) Keep the device away from sources of fire and heat.
 - p) Do not block the ventilation openings of the device!
 - q) Do not use the machine without water
 - r) The water level should be between the minimum and maximum level.
 - s) Do not allow water to freeze in the device.

CAUTION! Although the product has been designed to be safe, with adequate safeguards, and despite the additional safety features provided to the user, there is still a slight risk of accident or injury when handling the device. You are advised to use caution and common sense when using this product.

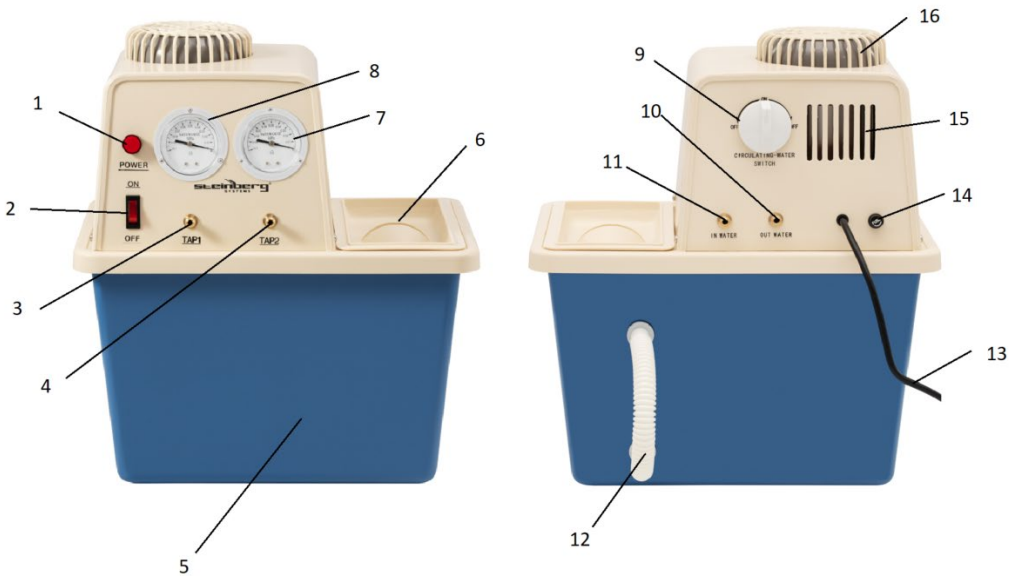
3. Rules of use

The device is designed to create a vacuum, which is necessary in processes such as: evaporation, distillation, crystallization, drying, degassing.

The product is intended only for use in university laboratories or industrial plants (e.g. in chemical or food processing industry).

The user is responsible for any damage resulting from misuse.

3.1. Description

1. Power light
2. ON/OFF switch
3. Vacuum connection I
4. Vacuum connection II
5. Water tank
6. Water tank cover
7. Manometer II
8. Manometer I
9. Water circulation switch
10. Water outlet
11. Water inlet
12. Overflow pipe
13. Power cord
14. Fuse
15. Ventilation hole
16. Motor cover

3.2. Preparation for operation

POSITIONING OF THE device

The ambient temperature must not exceed 40°C and ambient humidity should not exceed 85%. Place the device in a way that ensures good air circulation. Maintain a

minimum clearance of 100 cm from any wall of the device. Keep the device away from any hot surfaces. Always operate the device on a level, stable, clean, fireproof and dry surface and out of the reach of children and persons of impaired mental, sensory and intellectual functions. Place the device in such a way that the mains plug can be reached at any time. Ensure that the power supply to the device corresponds to that specified on the identification plate!

3.3. Working with the device

1. Before using for the first time, make sure that the overflow pipe is properly seated on the connections (on the back of the device).
2. Open the cover of the water tank.

Pour clean, cool water into the tank.

The water level should be between the minimum and maximum level.

3. Connect the vacuum connection. The machine works with one, the other or both connections connected.

Connect the hose on one side to the vacuum connection and on the other side to the device in which you want to create a vacuum.

4. Connect the plug to a power outlet.
5. Make sure that the circulation switch is in the OFF position.
6. Set the ON/OFF switch to the ON position. The ON/OFF switch will illuminate.

The pressure gauge will indicate the appropriate amount of vacuum generated.

7. When finished, turn the ON/OFF switch to the OFF position.

All the lights will go out. Pour the water out of the tank and drain it.

CAUTION! If the device stops working during operation, replace the water in the tank.

Pour the water out of the tank, clean the tank of any debris, then pour the water back in.

Make sure that each time you fill the water tank, the water is clean and cool

and that the water level is between the minimum and maximum levels specified.

CAUTION! If the water level in the tank is too high, the pressure gauge values may not be accurate. The water will automatically flow out through the overflow pipe.

WATER CIRCULATION FUNCTION

To start the cold water circulation function, you must:

1. Connect the water inlet and outlet of the device with the inlet and outlet of the laboratory device needing cooling.
2. Set the ON/OFF switch to the ON position. The ON/OFF switch will illuminate.
3. Set the water circulation switch to the ON position.
4. When finished, turn the ON/OFF switch to the OFF position. All the lights will go out. Pour the water out of the tank and drain it.

3.4. Cleaning and maintenance

- a) Cleaning and maintenance
Pull the mains plug and let the device cool down completely before cleaning, adjusting or replacing accessories and when the device is not in use.
 - Wait until the rotating parts stop.
- b) Pull out the mains plug before each cleaning and when the device is not in use.
- c) Use only non-corrosive cleaning agents for cleaning the surfaces.
- d) After each cleaning, all the parts should be dried well before the device is used again.
- e) Store the device in a dry and cool place protected from moisture and direct sunlight.
- f) Do not spray the device with a stream of water or immerse it in water.
- g) Make sure that no water enters through the ventilation openings in the casing.
- h) Clean the ventilation openings with a brush and compressed air.
- i) Perform regular inspections of the device checking technical fitness and any damages.
- j) Use a soft, damp cloth for cleaning.
- k) Do not clean the device with acidic substances, medical products, diluent, fuel, oil or other chemicals as this may cause damage to the device.

DISPOSAL OF USED UNITS.

At the end of its useful life, this product should not be disposed of with normal household waste but should be taken to a collection point for the recycling of electrical and electronic equipment. This is indicated by the symbol on the product, operating instructions or packaging. The materials used in this device are recyclable according to their marking. You will be making an important contribution to protecting our environment by reusing, recycling or otherwise disposing of used devices.

Your local administration will provide you with information about the appropriate disposal point for used devices.

TROUBLESHOOTING

Problem	Possible cause	Action
No vacuum or inaccurate gauge reading	Motor speed is not sufficient or the voltage is too low	Check that the voltage is stable
	Clogged water inlet, filters	Remove accumulated dirt inside. Clean the device thoroughly.
	Rotor failure	Replace the impeller
The motor does not work	Protective pipe is burnt through	Replace the protective tube
	The bearing is broken	Replace the bearing
	Motor winding burnt	Replace the device with a new one
Warm motor, too much noise	Bearing worn out	Replace the bearing
	Wear, deformation of the rotor	Replace the impeller
	Voltage instability or voltage is too low	Check if the line is overloaded and if low voltage occurs
Water tank	Water level is too low	Replenish water in the tank
Leakage	Pass-through water nozzle in the water tank is loose	Tighten the pass-through water nozzle
	Old seal	Replace the sealing ring
	Water tank cracked	Replace the water tank

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