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LABORDESTILLATIONSGERÄT

## **USER MANUAL**

SHORT PATH DISTILLATION FLASKS

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# Technical data

Parameter value	Parameter value
Name of the product	SHORT PATH DISTILLATION FLASK
Model	SBS-DK-1000
Supply voltage [V~] / Frequency [Hz]	230/50
Rated power [W]	1100
Maximum heating temperature [°C]	300
Temperature precision [°C]	+/-1
Maximum mixing speed [rpm]	1800
Jacket volume [l]	5
IP Code	IPX0
Dimensions [Width x depth x height; cm]	100x55x50
Weight [kg]	32

## 1. General description

These instructions are intended to assist you in safe and reliable use. The product is designed and manufactured strictly according to technical specifications using the latest technology and components and maintaining the highest quality standards.

**READ THE INSTRUCTIONS CAREFULLY AND  
UNDERSTAND THEM BEFORE USE.**

To ensure long and reliable operation of the device, care should be taken to operate and maintain the device correctly in accordance with the instructions. The technical data and specifications contained in these instructions are up-to-date. The manufacturer reserves the right to make changes for the purpose of quality improvement. Taking technological progress and the possibility to reduce noise into account, the device is designed and built so that risks resulting from noise emissions are reduced to the lowest possible level.

### Explanation of symbols

	The product meets the requirements of relevant safety standards.
---	--

	Read the instructions before use.
	Recyclable product.
	<b>NOTE!</b> or <b>WARNING!</b> or <b>REMEMBER!</b> describing a given situation (general warning sign).
	NOTE! Risk of electric shock!
	Caution! Touching hot surfaces may lead to burns!
	For indoor use only.
	The device shall be connected to grounded power supply sources.
	It is forbidden to use this equipment for experiments involving flammable, explosive, toxic and highly corrosive substances.

**NOTE!** The illustrations in this manual are for illustrative purposes only and may differ in some details from the actual product appearance.

The original version of the instructions is the German version. Other language versions are translations from German.

## 2. Use safety

**NOTE!** Read all safety warnings and all instructions. Failure to follow the warnings and instructions may result in electric shock, fire and/or serious injury or death.

---

The term "device" or "product" in the warnings and instructions refers to the SHORT PATH DISTILLATION APPARATUS.

## **2.1. Electrical safety**

- a) The device plug must fit into the socket. Do not modify the plug in any way. Original plugs and matching sockets reduce the risk of electric shock.
- b) Avoid touching earthed items such as tubes, radiators, heaters and refrigerators. There is an increased risk of electric shock if your body is earthed and touches the device exposed to direct rain, wet floor and humid environment. If water penetrates into the device, there is an increased risk of damage to the device and electric shock.
- c) Do not touch the device with wet or damp hands.
- d) Do not use the cable in a wrong way. Never use it to carry the device or to pull the plug out of the socket. Keep the cable away from heat sources, oil, sharp edges or moving parts. Damaged or tangled cables increase the risk of electric shock.
- e) If you cannot avoid using the device in a wet environment, use a residual current device (RCD). Using an RCD reduces the risk of electric shock.
- f) Do not use the device if the power cable is damaged or shows clear signs of wear. A damaged power cable should be replaced by a qualified electrician or the manufacturer's service.
- g) To avoid electric shock, do not immerse the cable, plug or device itself in water or other liquid. Do not use the device on wet surfaces.

## **2.2. Safety in the work area**

- a) Keep the work area tidy and well lit. Disorder or poor lighting can lead to accidents. Be careful, watch what you are doing and use common sense when working with the device.
- b) Do not use the device in an area with a risk of explosion, for example in the presence of flammable liquids, gases or dust. The device produces sparks which can ignite dust or vapours.
- c) In the event of damage or malfunction of the device, it must be switched off immediately and the incident must be reported to an authorised person.
- d) If in doubt as to whether the device is working properly, contact the manufacturer's service.
- e) Repairs to the device may only be carried out by the manufacturer's service. Do not carry out repairs yourself!
- f) In the event of fire or fire, use only powder or snow (CO<sub>2</sub>) extinguishers to extinguish fire in the device where there is voltage.

- 
- g) Keep these instructions for future reference. If the device is to be passed on to third parties, the instructions for use must be handed over together with the device.
  - h) Keep packaging parts and small assembly parts out of the reach of children.
  - i) Keep the device away from children and animals.

**Remember!** Keep children and other bystanders safe while operating the device.

## **2.3. Personal safety**

- a) Do not operate the device if you are tired, ill or under the influence of alcohol, drugs or medication which might impair your ability to operate the device.
- b) The device is not intended to be used by persons (including children) with reduced mental, sensory or intellectual functions or lack of experience and/or knowledge, unless they are supervised by a person responsible for their safety or have been given instructions by them on how to operate the device.
- c) When operating the device use personal protection equipment preventing:
  - splashing or contact with evaporating liquid,
  - contact with toxic or flammable gases.
- d) To prevent accidental start-up, make sure the switch is in the OFF position before connecting the device to the power source.
- e) The device is not a toy. Children should be supervised to ensure that they do not play with the device.

## **2.4. Safe use of the device**

- a) Do not use the device if the knobs or control buttons do not operate in efficient manner. Devices that cannot be controlled by control elements are hazardous, must not be operated and must be repaired.
- b) The device must be disconnected from the power supply before adjustment, cleaning and maintenance. This preventive measure reduces the risk of accidental activation.
- c) Keep unused devices out of the reach of children and anyone not familiar with the device or these instructions. Devices are dangerous in the hands of inexperienced users.
- d) Keep the device in good working order. Before each use, check for general damage or damage to moving parts (cracks in parts and components or any other condition that may affect the safe operation of the device). If damaged, have the device repaired before use.
- e) Keep the device out of the reach of children.

- 
- f) Repair and maintenance of the device should be carried out by qualified persons using only original spare parts. This will ensure a safe use of the device.
  - g) To ensure the designed operational integrity of the device, do not remove factory-installed covers or loosen any screws.
  - h) Do not move, shift or rotate the device during operation.
  - i) Do not leave the switched on device unattended.
  - j) Clean the device regularly to prevent the permanent accumulation of dirt.
  - k) The device is not a toy. Cleaning and maintenance must not be performed by children without adult supervision.
  - l) It is forbidden to tamper with the device in order to change its parameters or construction.
  - m) Keep the device away from sources of fire and heat.
  - n) NOTE! During active heating function, the jacket heats up to a high temperature. Do not touch hot elements of the apparatus, because it poses a hazard of burns.
  - o) When handling flammable materials pay special care to the set heating temperature to be not too high.
  - p) Make sure that the container with the liquid can withstand the set heating temperature. A molten container may cause fire.
  - q) Never heat an empty container.
  - r) Do not use the device without a vessel/sample placed in the jacket.
  - s) Control the amount of liquid in the vessels. Liquid may evaporate.
  - t) Spilling a flammable contents of a vessel on the heating elements of the device may lead to a fire.
  - u) The set heating temperature must be lower at least by 25°C from the ignition temperature of the mixed substance.
  - v) Pay special attention to the hazards resulting from:
    - flammable materials and substances with low boiling point,
    - overful vessels,
    - vessels made of brittle material or material with a low melting temperature.
  - w) Pathogenic materials can be only mixed in closed vessels.
  - x) Do not cover the device when it is operating.
  - y) Make sure that the supply cable does not come into contact with hot elements of the device.
  - z) Keep the device away from strong magnetic field.
  - aa) Use caution not to damage the glass vessels.
  - bb) Wash and dry glass vessels after each use.
  - cc) It is recommended to first switch off heating, and only after the device has cooled down to turn off mixing.

**NOTE!** Although the device has been designed to be safe and has been equipped with adequate safeguards, and despite the use of additional

---

**safety measures, there is still a small risk of accident or injury when operating the device. Caution and common sense are advised when using the device.**


### **3. Rules of use**

The device is intended for use in laboratories, schools, universities and similar facilities. It is intended for contactless mixing and heating of liquids in laboratory vessels and for controlled liquid evaporation.

**The user is responsible for any damage caused by improper use.**

#### **3.1. Description of the device**


**Description of the entire set**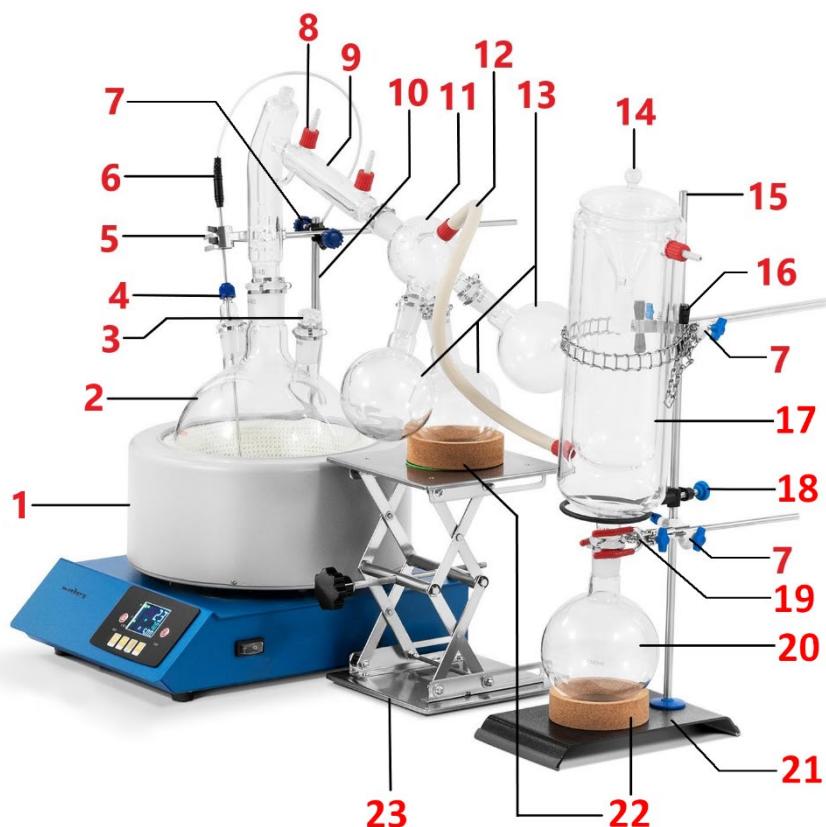

- 1 - Heating jacket
- 2 - Flask with 3 necks
- 3 - Plug
- 4 - Vacuum port
- 5 - Boom with handle
- 6 - External thermocouple
- 7 - Cross connector
- 8 - Connection adapter
- 9 - Distillation head
- 10 - Stand #1

- 
- 11 - 3-way adapter for receiving flasks
  - 12 - Connection hose
  - 13 - Receiving flask (500 ml)
  - 14 - Chiller lid
  - 15 - Stand #2
  - 16 - Boom with chain
  - 17 - Chiller
  - 18 - Ring support
  - 19 - Boom with handle
  - 20 - Receiving flask (1000 ml)
  - 21 - Chiller base
  - 22 - Cork ring
  - 23 - Regulated base

**Description of heating jacket with magnetic agitator**

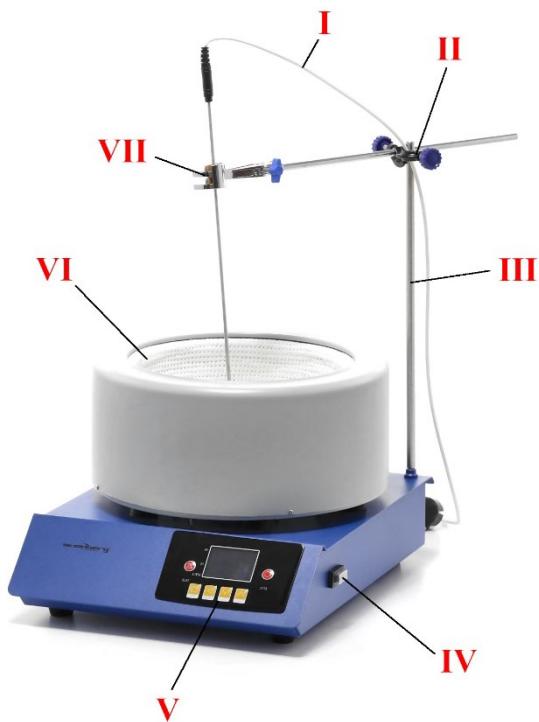I - external thermocouple wire

II - cross connector

III - stand

IV - I/O switch (on/off)

V - control panel

VI - chamber with heating jacket

VII - handle on a boom

View of the heating jacket housing (rear):

VIII - external/internal thermocouple socket

IX - supply cable

X - fuse

Control panel view:

A - PV - measured (current) temperature value

B - SV - set (target) temperature value

C - Heating activation button

D - R/MIN - Rotational speed of the agitator (presented in rpm)

E - "SET" button - enables setting the temperature and mixing speed values

F - "R/AT" value - automatic calibration / change of the location of digit edition on the display

G - “-” button - reducing the mixing speed or temperature

H - “+” button - increasing the mixing speed or temperature

I - Mixing activation button

## 3.2. Preparation for work

### LOCATION OF THE DEVICE

Position the device in a way that allows for good air circulation. Maintain a minimum distance of 10 cm from any wall. Keep the device away from any hot surfaces. Always operate the device on a level, stable, clean, fireproof and dry surface and out of the reach of children and persons with reduced mental, sensory and intellectual abilities. Position the device in such a way that the power supply plug can be reached at any time. Make sure that the device's power supply corresponds to the data on the rating plate!

### SET ASSEMBLY

1. Place the flask with 3 necks [2] in the jacket [1].

- 
2. Screw in the stand [10] in the corner of the housing of the jacket base [1].
  3. Apply the cross connector [7] on the stand [10] and screw on the clamping knob.

*Cross connector*

4. Place the plug [3] into one of the flask necks [2] and secure it with a clip.

5. Insert vacuum port [4] to one of the smaller flask necks [2].

*Vacuum port*

6. Insert the external thermocouple rod [6] into the vacuum port opening [4], so that the tip of the rod is located several centimetres above the flask bottom [2] (but it does not touch the bottom).
7. Insert the boom with handle [5] into the cross connector [7] and grasp the external thermocouple rod [6]. Secure parts in this position by turning the cross connector and boom with handle knobs.

8. Withdraw the internal thermocouple (if connected) from the socket in the base of the heating jacket.

*Internal thermocouple*

9. Connect the external thermocouple plug [6] into the socket in the base of the heating jacket.

10. Connect the distillation head [9] to the central neck of the flask inside the heating jacket [1] and secure the connection with a clip.

*Clip*


11. Connect the 3-way adapter for receiving flasks [11] to the free connection of the distillation head [9].
12. Connect one flask (500 ml) [13] to each neck of the 3-way receiving flasks adapter [11] and secure the connections with clips.
13. Install two connection adapters [8] on the glass ports of the distillation head [9] and 3-way adapter for receiving flasks [11].

14. Place the regulated base [23] for the lowest receiving flask (500 ml) [13].
15. Place the green washer and cork ring [22] on the regulated base [23].
16. Regulate the base height [23] so that the receiving flask (500 ml) [13] supports its weight lightly on the cork ring [22].

17. Place the chiller base [21] on a flat surface.
18. Screw in stand [15] into the chiller base opening [21].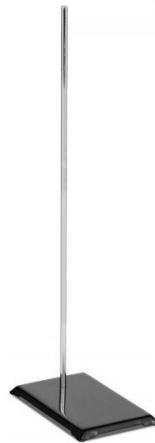


19. Install cross connector [7] on the stand [15].
20. Install the boom with chain [16] in the cross connector [7] and screw in the clamping knob.
21. Place the cork ring [22] on the base [21].
22. Place the receiving flask (1000 ml) [20] on the cork ring [22].

23. Install the ring support [18] on the stand [15]. Screw in the ring support screw on the stand.
24. Install the chiller neck [17] into the opening of the receiving flask (1000 ml) [20].
25. Wrap the chiller [17] using the chain on the boom [16]. Attach the chain on the hook on the other side of the boom [16].

26. Secure the connection between the chiller [17] and receiving flask (1000 ml) [20] with a clip.
27. Install the cross connector [7] on the stand [15].
28. Install the boom with handle [19] in the cross connector [7]. Insert the neck of the receiving flask (1000 ml) [20] into the handle [19] and screw in the screws of the handle and cross connector.

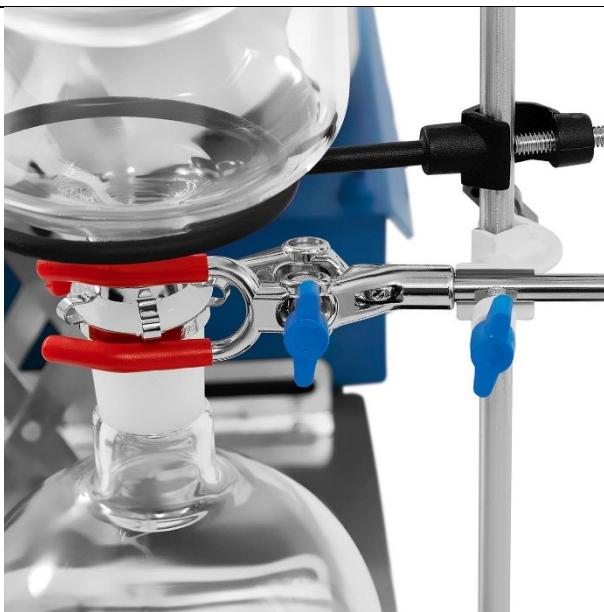29. Install two connection adapters [8] on glass ports of the chiller [17].
30. Connect the chiller port [17] with the port of the 3-way adapter for receiving flasks [11].

### 3.3. Working with the device

**NOTE!** The magnetic stirring heating mantles can operate as a separate device independent of the rest of the set. Operation of the control panel remains unchanged, but the following procedures refer to using the jacket with magnetic agitator as a separate device.

#### Heating jacket operation:

- Pour the liquid to the appropriate vessel and insert the magnetic agitator to the vessel.

*Magnetic agitator*

- Place the vessel with liquid inside the chamber with heating jacket and dip the tip of the temperature sensor into the liquid.

- 
- c) Connect the power supply cable to the power socket.
  - d) Switch on the device with I/O button.
  - e) Start mixing and set the appropriate speed following the “Mixing speed regulation” chapter.
  - e) Start heating and set the appropriate temperature following the “Temperature setting” chapter.
  - f) After completing operation of the device, switch off heating and mixing, switch off the device with I/O button, and disconnect the device from the power source.

### Liquid temperature measurement

- a) Remove the internal thermocouple from the socket located in rear portion of the device.

*Internal thermocouple*

*External thermocouple*


- b) Connect the external thermocouple to the socket in the rear portion of the device. The display will indicate the ambient temperature PV and set temperature SV.
- c) Insert the metal part of thermocouple into the liquid for which temperature will be measured. The display will indicate the measured temperature PV and set temperature SV.
- d) The safeguard will be activated when the external thermocouple is disconnected. In such a case the display will show “hhhh” message, the green indicator light switches off and the device stops working. When the above situation occurs, check the technical condition of the device.
- e) If it is not necessary to use external thermocouple, disconnect it and reconnect the internal thermocouple.

### Assembly of external thermocouple

Assembly of external thermocouple should be performed when it is necessary to heat the liquid and its temperature must be monitored.

- a) Screw in stand into the opening in the corner of the device housing.
- b) Apply the cross connector on the stand and screw on the clamping knob.
- c) Install the boom with handle in the cross connector and screw in the clamping knob.

- 
- d) Remove the metal part of external thermocouple. The sensor should be directed with the wire up.
  - e) Connect the external thermocouple wire to the socket in the rear portion of the device.
  - f) The device is ready for use.

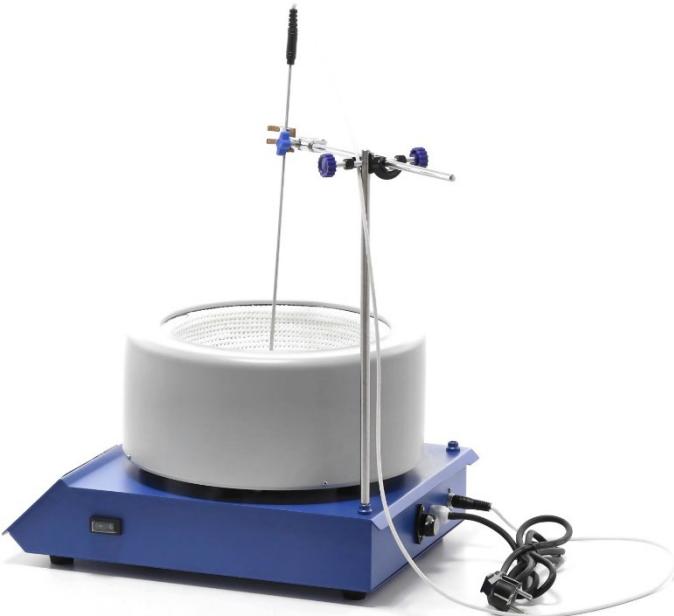### **Temperature setting**

- a) Switch the device on with I/O switch [IV].
- b) Press the heating activation button [C], the display will show the current the temperature PV and last set target temperature SV.  
Caution: The device will heat up to the last set temperature automatically, if the user does not set a new value.
- c) Press the "SET" button [E]. The last digit on the right in the SV field will flash.
- d) Press "+" or "-" [H or G], to reduce or increase the target temperature. In order to alternate between the edited digits, press "R/AT" [F].
- e) To complete temperature setting, press "SET" [E].

### **Automatic calibration**

- a) Switching the device off with I/O [IV], if the measured temperature PV increases above the set temperature SV.
- b) Wait until the device cools to room temperature.
- c) Press and hold "R/AT" [F] for 8 seconds.

- 
- d) Automatic calibration will end when the SV field stops flashing and the measured temperature PV is equal to the set temperature SV.

### **Mixing speed regulation**

**Caution:** Although the maximum mixing speed that can be set in the device is greater than 1800 rpm, it is not recommended to exceed this value!

The method of regulating the mixing speed differs depending on whether the heating function is active or inactive.

When the device is operated with heating function active:

- a) Press mixing activation button [I]. Field R/MIN [D] will show the rotating speed of the agitator.
- b) Press "SET" [E]. Editing of the last digit in the SV field will be activated (the digit starts flashing).
- c) Press the "SET" button [E] again. The R/MIN [D] field will display the set mixing speed and it will be possible to edit the last digit of this field (the digit will be flashing).
- d) Press "+" or "-" [H or G], to increase or reduce mixing speed. In order to alternate between the edited digits, press "R/AT" [F].
- e) Press "SET" [E], to complete mixing speed regulation. R/MIN field [D] will display the current mixing speed.

When the device is operated without heating function active:

- a) Switch the device on with I/O [IV].
- b) Press mixing activation button [I]. R/MIN field [D] will display the current mixing speed.
- c) Press "SET" [E]. The R/MIN [D] field will display the set mixing speed and it will be possible to edit the last digit of this field (the digit will be flashing).
- d) Press "+" or "-" [H or G], to increase or reduce mixing speed. In order to alternate between the edited digits, press "R/AT" [F].
- e) Press "SET" [E], to complete mixing speed regulation. R/MIN field [D] will display the current mixing speed.

### **Overload protection**

The device is equipped with a fuse (X) protecting against overload. In the event of the safeguard is activated (the device switches off), find and remove the cause. Subsequently, replace the fuse with a new one.

---

### **3.4. Cleaning and maintenance**

- a) Before cleaning, adjustment or replacement of accessories, and when the device is not in use, pull the power supply plug and let the device cool down completely.
- b) Use only non-corrosive agents for cleaning the surface.
- c) After each cleaning, all components should be well dried before the device is used again.
- d) Store the device in a dry and cool place protected from moisture and direct sunlight.
- e) Do not spray the device with a stream of water and do not immerse it in water.
- f) Ensure that no water enters through the ventilation openings in the housing.
- g) Clean the ventilation openings with a brush and compressed air.
- h) Regularly inspect the device for technical failure and damage.
- i) Use a soft cloth for cleaning.
- j) When selecting cleaning agents, take into consideration type of substance previously used in the device.

#### **DISPOSAL OF USED DEVICES.**

At the end of its service life, this product must not be disposed of with normal household waste, but should be taken to a collection point for the recycling of electrical and electronic devices. This is shown by the symbol placed on the product, operating manual or the packaging. The materials used in the device can be reused according to their intended use. By reusing materials or conducting other forms of using used devices, you make a significant contribution to the protection of our environment.

Local administration will provide you with information about the appropriate point for disposal of used devices.

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