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ROTATIONSVERDAMPFER SET

USER MANUAL

ROTARY EVAPORATOR SET

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EVAPORATORE ROTANTE - SET

MANUAL DE INSTRUCCIONES

SET DE EVAPORADOR ROTATORIO

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SBS-RV-5000 | SBS-RV-2000

DE	Produktname:	ROTATIONSVERDAMPFER SET
EN	Product name:	ROTARY EVAPORATOR SET
PL	Nazwa produktu:	ROTACYJNY DESTYLATOR PRÓŻNIOWY
CZ	Název výrobku	ROTAČNÍ VÝPARNÍK – SADA
FR	Nom du produit:	ÉVAPORATEUR ROTATIF
IT	Nome del prodotto:	EVAPORATORE ROTANTE - SET
ES	Nombre del producto:	SET DE EVAPORADOR ROTATORIO
HU	Termék neve	ROTÁCIÓS VÁKUUMDESZTILLÁTOR
DA	Produktnavn	ROTERENDE VAKUUMDESTILLER
DE	Modell:	SBS-RV-5000 SBS-RV-2000
EN	Product model:	
PL	Model produktu:	
CZ	Model výrobku	
FR	Modèle:	
IT	Modello:	
ES	Modelo:	
HU	Modell	
DA	Model	
DE	Hersteller	expondo Polska sp. z o.o. sp. k.
EN	Manufacturer	
PL	Producent	
CZ	Výrobce	
FR	Fabricant	
IT	Produttore	
ES	Fabricante	
HU	Termelő	
DA	Producent	
DE	Anschrift des Herstellers	ul. Nowy Kisielin – Innowacyjna 7, 66-002 Zielona Góra Poland, EU
EN	Manufacturer Address	
PL	Adres producenta	
CZ	Adresa výrobce	
FR	Adresse du fabricant	
IT	Indirizzo del produttore	
ES	Dirección del fabricante	
HU	A gyártó címe	
DA	Producentens adresse	

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Technical Data

Description of the parameter	Value of the parameter	
Product name	ROTARY EVAPORATOR SET	
Model	SBS-RV-5000	SBS-RV-2000
Supply voltage [V~] / Frequency [Hz]	230/50	
Rated power [W].	2000	1500
Motor power [W].	40	40
Rotational speed [RPM]	0-120	0-120
Heating temperature[°C]	0-99	0-99
Heating tub capacity [l]	6	5
Height adjustment range of the water bath bowl [mm]	120	120
Temperature accuracy [°]	±1	±1
Vacuum pressure [Mpa]	0.098	0.098
Receiving flask [l]	2	1
Rotating flask [l]	5	5
IP code	IP20	IP20
Dimensions [Width x Depth x Height; mm]	800x400x1000	730x330x1000
Weight [kg]	33	25

1. General Description

The instruction manual is intended to assist in safe and reliable use. The product is designed and manufactured strictly according to technical specifications using the latest technology and components and maintaining the highest quality standards.

**PLEASE CAREFULLY READ
AND UNDERSTAND THIS INSTRUCTION MANUAL BEFORE
OPERATION,**

To ensure long and reliable operation of the unit, make sure to operate and maintain it properly in accordance with the guidelines in this instruction manual. The technical data and specifications contained in this instruction manual are up to date. The manufacturer reserves the right to make changes in order to improve the quality. Taking the technical progress and the possibility of reducing noise into account, the unit

is designed and built in such a way so that risks resulting from noise emissions are reduced to the lowest possible level.

Explanation of symbols

	The product complies with applicable safety standards.
	Please read the instructions before use.
	Recyclable product.
	CAUTION! or WARNING! or REMINDER! describing a situation. (general warning sign).
	CAUTION! Warning of electric shock!
	CAUTION! Spinning elements!
	Caution! Hot surface can cause burns!
	For indoor use only.

CAUTION! The illustrations in this instruction manual are for reference only and may differ from the actual product in some details.

The original instruction manual is in the German language version. Other language versions are translations from German.

2. Safety of use

CAUTION! Read all safety warnings and all instructions. Failure to follow the warnings and instructions may result in electric shock, fire and/or severe personal injury or death.

The term "unit" or "product" in the warnings and in the description of the instructions refers to the ROTARY EVAPORATOR SET.

2.1. Electrical safety

- a) The plug of this unit must fit into the outlet. Do not modify the plug in any way. Original plugs and matching outlets reduce the risk of electric shock.
- b) Avoid touching grounded parts, such as pipes, heaters, ovens, and refrigerators. There is an increased risk of electric shock if your body is grounded and touches the unit exposed to direct rain, wet pavement, and operation in a humid environment. If water enters the unit, there is an increased risk of damage to the unit and electric shock.
- c) Do not touch the unit with wet or damp hands.
- d) Do not use the cord in an unintended manner. Never use it to carry the unit or to pull the plug out of the socket. Keep the cord away from heat sources, oil, sharp edges or moving parts. Damaged or tangled cords increase the risk of electric shock.
- e) If you cannot avoid using the unit in a wet environment, use a residual current unit (RCD). Using an RCD reduces the risk of electric shock.
- f) Do not use the unit if the power cord is damaged or shows signs of wear. A damaged power cord should be replaced by a qualified electrician or the manufacturer's service department.
- g) To avoid electric shock, do not immerse the cable, plug, or unit itself in water or other liquid. Do not use the unit on wet surfaces.

2.2. Safety in the workplace

- a) Keep the work area tidy and well lit. Disorder or poor lighting can lead to accidents. Be foresighted, watch what you are doing and use common sense when using the unit.
- b) Do not use the unit in an explosive area, for example in the presence of flammable liquids, gases or dust.
- c) If you find any damage or irregularities in the operation of the unit, immediately turn it off and report it to an authorized person.

-
- d) • If you have any doubts as to whether the product is working properly or if it is damaged, contact the manufacturer's service department.
 - e) Only the manufacturer's service department can repair the unit. Do not carry out repairs yourself!
 - f) In case of open flames or fire, use only dry powder or snow (CO₂) fire extinguishers to extinguish the live equipment.
 - g) Keep these instructions for use for future reference. If the unit is to be passed on to a third party, the operating instructions must also be handed over together with the unit.
 - h) Keep the packaging and small assembly parts out of the reach of children.
 - i) Keep the unit away from children and animals.
 - j) When using this machine together with other machines, also follow the other instructions for use.

2.3. Personal safety

- a) Do not operate this product if you are tired, ill or under the influence of alcohol, drugs or medication that could impair your ability to operate the product.
- b) The unit is not intended to be used by persons (including children) with reduced mental, sensory or intellectual functions or persons who lack experience and/or knowledge unless they are supervised or have been instructed by a person responsible for their safety on how to operate the unit.
- c) Use caution and common sense when operating this unit. A moment's inattention during operation may result in serious personal injury.
- d) To prevent accidental start-up, make sure the switch is in the off position before connecting to a power source.
- e) Do not wear loose clothing or jewelry. Keep hair, clothing, and gloves away from moving parts. Loose clothing, jewelry, or long hair can be caught in moving parts.
- f) Before switching the unit on, remove any regulating tools or keys. Any tools or keys left in the rotating part of the unit may cause injury.
- g) The unit is not a toy. Children should be watched to ensure that they do not play with the unit.
- h) Do not place your hands or any objects inside the running unit!

2.4. Safe use of the unit

- a) Do not overload the unit. Use tools that are suitable for the application. A correctly selected unit will do a better and safer job for which it was designed.

-
- b) Do not use the unit if the ON/OFF switch does not function properly (does not turn on and off). Units that cannot be controlled by the switch are unsafe, cannot operate, and must be repaired.
 - c) Unplug the unit before making adjustments, changing accessories, or putting it away. This precaution reduces the risk of accidental start-up.
 - d) Keep unused equipment out of the reach of children and out of the reach of anyone unfamiliar with the unit or this instruction manual. These units are dangerous in the hands of inexperienced users.
 - e) Keep the unit in good working condition. Check before each use for general damage or damage to moving parts (cracks in parts and components or any other condition that may affect the safe operation of the unit). If damaged, have the unit repaired before use.
 - f) Keep the unit out of the reach of children.
 - g) Repairs and maintenance should be carried out by qualified personnel using only original spare parts. This will ensure the safety of use.
 - h) To ensure the designed operational integrity of the unit, do not remove factory-installed covers or loosen screws.
 - i) When transporting or moving the unit from storage to the place of use, observe the health and safety rules for manual handling applicable in the country where the unit is used.
 - j) Do not move, shift, or rotate the machine while in operation.
 - k) Do not leave the unit switched on unattended.
 - l) Clean the unit regularly to prevent permanent dirt build-up.
 - m) The unit is not a toy. Cleaning and maintenance must not be performed by children without adult supervision.
 - n) Do not tamper with the unit to alter its performance or design.
 - o) Keep the unit away from sources of fire and heat.
 - p) Do not block the ventilation openings of the unit!
 - q) Use care not to damage the glassware.
 - r) Wash and dry the glassware after each use.
 - s) All round glassware joints should be lubricated with sealing grease before connecting.
 - t) The water bath container must be filled with water before connecting to the power supply. Heating without water is prohibited.
 - u) To obtain an accurate temperature, use a thermometer to measure.
 - v) When you have finished working with the unit, turn it off and pull the plug out from the electrical outlet.
 - w) Never use explosive or flammable substances in the unit.
 - x) Some materials during the evaporation process may explode, produce poisonous or flammable gases. Use caution!
 - y) If hazardous material is spilled, it is the user's responsibility to properly disinfect or dispose of the material.

-
- z) The room in which the unit is used must have an efficient ventilation system.
 - aa) Ensure that vapors are extracted at the place of use.

CAUTION! Although the product has been designed to be safe, with adequate safeguards, and despite the additional safety features provided to the user, there is still a slight risk of accident or injury when handling the unit. You are advised to use caution and common sense when using this product.

3. Rules of use

The unit is designed for controlled evaporation of liquids.

The user is responsible for any damage resulting from misuse.

3.1. Description

Model SBS-RV-5000

Model SBS-RV-2000

- 1 - cooler
- 2 - speed controller
- 3 - safety clamp
- 4 - vacuum meter
- 5 - motor
- 6 - rotating flask
- 7 - water bath
- 8 - knob for height adjustment of water bath bowl
- 9 - temperature controller
- 10 - filling valve
- 11 - safety clip
- 12 - receiving flask
- 13 - circular support
- 14 - drain valve for receiving flask
- 15 - rack
- 16 - rack base

Speed controller description (2):

- A - speed display
- B - rotational speed adjustment knob
- C - Speed controller ON/OFF switch
- D - Speed controller power cable
- E - rack mounting hole
- F - control cable

Temperature Controller Description (9):

Display:

PV - current measured temperature value (process value)

SV - target value of water bath temperature (set value)

Buttons:

Set - enter target temperature setting mode/approve settings

Swift - switching between display digits

+ - increase value

- decreasing the value

I/O - water bath on/off switch (on the SBS-RV-2000 this button is located on the back of the temperature controller casing)

Indicator lights:

AT - automatic settings

OUT - start heating

ALM1 - alarm 1

ALM2 - alarm 2

3.2. Preparation for operation

POSITIONING OF THE UNIT

The ambient temperature must not exceed 40°C and ambient humidity should not exceed 85%. Place the unit in a way that ensures good air circulation. Maintain a minimum clearance of 10 cm from any wall of the unit. Keep the unit away from any hot surfaces. Always operate the unit on a level, stable, clean, fireproof and dry surface and out of the reach of children and persons of impaired mental,

sensory and intellectual functions. Place the unit in such a way that the mains plug can be reached at any time. Ensure that the power supply to the unit corresponds to that specified on the identification plate!

ASSEMBLY OF THE UNIT

CAUTION! All round glassware joints should be lubricated with sealing grease before connecting.

Model SBS-RV-5000

- 1) Position the base (16) with the rack (15) next to the water bath (7).
- 2) Install the motor (5) to the rightmost tube of the rack (15) (closest to the water bath). To install the motor to the rack, insert the motor pin into the mounting hole. Position the motor so that it is facing the water bath at a height of approximately 40 cm and tilted 25° from horizontal. Tighten the clamping screw to secure the connection.
- 3) Mount the speed controller (2) to the end of the right rack tube (15), secure the connection with the clamp screw. Plug the 5-pin control cable connector routed from the motor (5) into the controller's lower socket.
- 4) Connect the right cooler connector (1) to the motor connector (5) and tighten the lock nut. Secure the position of the cooler (1) with a clamp (3) attached to the left rack tube (15).
- 5) Insert the fill valve (10) into the flexible tube and place the connected parts into the left cooler connector (1).
- 6) Connect the receiving flask (12) to the lower radiator connector (1) and secure the connection with the clip (11), then tighten the clip screw.
- 7) Attach the round support (13) to the lower bracket of the left rack tube (15) so that it holds the receiving flask (12) in place. Secure the connection by tightening the clamping screw.
- 8) Install the rotary flask (6) to the motor connector (5) and tighten the locking collar.
- 9) Place the water bath (7) under the rotary flask (6). Pour clean water into the bowl of the water bath. The water level should not exceed 2/3 of the bowl height.
- 10) Connect the vacuum hose to the upper vacuum gauge connection (4), then connect the vacuum pump or laboratory vacuum switch to the lower vacuum gauge connection.

Model SBS-RV-2000

-
- 1) Place the rack (15) in the opening of the base (16) and tighten the clamp screw to secure the connection.
 - 2) Position the base (16) with the rack (15) next to the water bath (7).
 - 3) Install the motor (5) into the center tube holder of the rack (15). To install the motor to the rack, insert the motor pin into the mounting hole. Position the motor so that it is facing the water bath at a height of approximately 40 cm and tilted 25° from horizontal. Tighten the clamping screw to secure the connection.
 - 4) Mount the speed controller (2) to the upper rack bracket (15), secure the connection with the clamp screw. Plug the 5-pin control cable connector routed from the motor (5) into the controller's lower socket.
 - 5) Connect the right cooler connector (1) to the motor connector (5) and tighten the lock nut.
 - 6) Insert the fill valve (10) into the flexible tube and place the connected parts into the left cooler connector (1).
 - 7) Connect the receiving flask (12) to the lower radiator connector (1) and secure the connection with the clip (11), then tighten the clip screw.
 - 8) Attach the round support (13) to the lower rack bracket (15) so that it holds the receiving flask (12) in place. Secure the connection by tightening the clamping screw.
 - 9) Install the rotary flask (6) to the motor connector (5) and tighten the locking collar.
 - 10) Place the water bath (7) under the rotary flask (6). Pour clean water into the bowl of the water bath. The water level should not exceed 2/3 of the bowl height.
 - 11) Connect the vacuum hose to the upper vacuum gauge connection (4), then connect the vacuum pump or laboratory vacuum switch to the lower vacuum gauge connection.

Connection of the cooling water circuit

For both models, connect the water circuit as shown in the figure below:

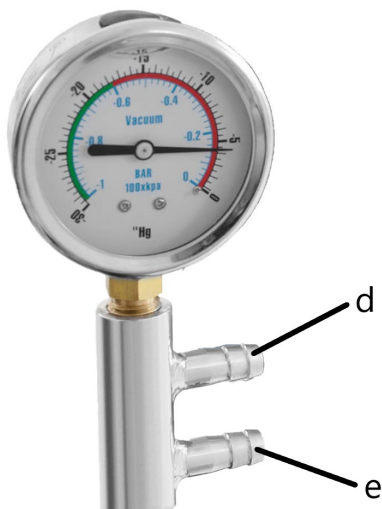
The connections are located at the rear of the cooler at the appropriately marked locations:

a - cooling water outlet

b - vacuum hose connection (should connect to upper vacuum gauge connection)

c - cooling water inlet

Vacuum pump connection

d - pressure hose connection (should connect with cooler)

e - vacuum pump connection

3.3. Working with the unit

The evaporation process should be as follows:

1. Make sure that an adequate amount of water has been poured into the bath bowl (7), then connect the unit to the power supply (by performing these steps in the order given, you reduce the risk of electrical shock).
Start the water bath and set the appropriate heating temperature using the controller (9).
2. Start the vacuum pump.
3. Close the filling valve (10).
4. Start rotation of the flask (6):
 - a) make sure that the speed controller (2) is connected to the power supply
 - b) Switch the speed controller (2) on at the ON/OFF switch (C)
 - c) set the selected speed using the adjustment knob (B)Caution: Avoid rocking, vibrating or shaking the water bath and motor during operation
5. Make sure the water bath (7) has reached the proper temperature and raise the water bowl to the proper level using the knob (8).

-
6. Allow the solvent to evaporate.
 7. Perform steps 1-5 in reverse order to complete the process. These include:
 - a) lower the water bath bowl
 - b) stop rotation of the flask
 - c) open the fill valve to equalize the pressure
 - d) stop the vacuum pump
 - e) turn off the water bath
 8. Recover the solvent from the receiving flask (12) and the non-volatile substance from the rotating flask (6). Depending on model:
 - Model SBS-RV-5000: Open the drain valve (14) to allow the solvent to flow out.
 - Model SBS-RV-2000: Loosen the screw of the clip (11) securing the receiving flask (12), then while holding the flask, disengage the clip and remove the flask and its contents.

Vacuum

If no vacuum is generated in the system despite the operation of the vacuum pump, check that all connections are sealed and that the vacuum pump has no leaks or damage.

Refilling the rotary flask

Liquids should be added to the flask through a flexible tube connected to the fill valve (10) using vacuum systems.

The rotary flask (6) can be filled to a maximum of half its capacity.

The substance in the rotary flask can be refilled while the instrument is in operation, but in this case the following steps must be taken:

- a) turn off the pump,
- b) stop heating,
- c) when evaporation is complete, slowly open the delivery valve to prevent backflow.

Heating

- The temperature control range for the water bath is 0-99°C.
- Due to thermal inertia, the actual water temperature is 3 degrees higher than the set temperature. It is possible to change the value setting: for example, if the water temperature needed is 40 degrees, set it to 38 degrees.
- To determine the exact actual water temperature, measure directly with a thermometer.

-
- The rotating flask (6) should be submerged in water to a depth of 1/3 to 1/2 its size.

Evaporation rate

The rate of evaporation depends on three factors:

1. The rate of evaporation the higher the degree of vacuum.
2. The higher the water bath temperature, the higher the evaporation rate. The water bath temperature should be selected according to the requirements of the substance.
3. The lower the cooling water temperature, the higher the evaporation rate.

3.4. Cleaning and maintenance

- a) Pull the mains plug and let the unit cool down completely before cleaning, adjusting or replacing accessories and when the unit is not in use.
 - Wait until the rotating parts stop.
- b) Use only non-corrosive cleaning agents for cleaning the surfaces.
- c) After each cleaning, all the parts should be dried well before the unit is used again.
- d) Store the unit in a dry and cool place protected from moisture and direct sunlight.
- e) Do not spray the unit with a stream of water or immerse it in water.
- f) Make sure that no water enters through the ventilation openings in the casing.
- g) Clean the ventilation openings with a brush and compressed air.
- h) Perform regular inspections of the unit checking technical fitness and any damages.
- i) Use a soft cloth for cleaning.
- j) When selecting cleaning agents, take the type of substance that was previously used in the unit into account.

DISPOSAL OF USED UNITS.

At the end of its useful life, this product should not be disposed of with normal household waste but should be taken to a collection point for the recycling of electrical and electronic equipment. This is indicated by the symbol on the product, operating instructions or packaging. The materials used in this unit are recyclable according to their marking. You will be making an important contribution to protecting our environment by reusing, recycling or otherwise disposing of used units.

Your local administration will provide you with information about the appropriate disposal point for used units.

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