USER MANUAL

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Mode d'emploi

Manuale d'uso

Manual de instrucciones

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Käyttöohje

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Bruksanvisning

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Návod na použitie

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1. Technical Data

Description of the parameter	Value of the parameter	
Product name	Lower temp liquid circulating chiller	
Model	SBS-LCC-3000	SBS-LCC-5000
Supply voltage [V~] / Frequency [Hz]	230/50	
Rated power [W].	660	900
Current intensity [A]	2.87	3.6
Accuracy [°C]	±2	±2
Type of temperature sensor	Pt100	
Tank capacity [L]	3	5
Circulation pump flow rate [L/min]	20	20
Circulation pump head [m]	5-7	5-7
Tank dimensions [mm]	ø180x130	ø250x130
Hose length [m]	3.6	-
Working temperature range[°C]	-20 ÷ 20 (max. 25)	-20 ÷ 20 (max. 25)
Ambient relative humidity during operation [%].	≤60	≤60
Compressor refrigerant	R404A	R404A
IP code	IPX4	
Dimensions [Width x Depth x Height; mm]	420x230x500	480x410x1000
Weight [kg]	27	40

2. General Description

The instruction manual is intended to assist in safe and reliable use. The product is designed and manufactured strictly according to technical specifications using the latest technology and components and maintaining the highest quality standards.

**PLEASE CAREFULLY READ
AND UNDERSTAND THIS INSTRUCTION MANUAL BEFORE
OPERATION,**

To ensure long and reliable operation of the unit, make sure to operate and maintain it properly in accordance with the guidelines in this instruction manual. The technical data and specifications contained in this instruction manual are up to date. The manufacturer reserves the right to make changes in order to improve the quality. Taking the technical progress and the possibility of reducing noise into account, the unit is designed and built in such a way so that risks resulting from noise emissions are reduced to the lowest possible level.

Explanation of symbols

	<p>The product complies with applicable safety standards.</p>
	<p>Please read the instructions before use.</p>
	<p>Recyclable product.</p>
	<p>CAUTION! or WARNING! or REMINDER! describing a situation. (general warning sign).</p>
	<p>CAUTION! Warning of electric shock!</p>
	<p>CAUTION! Spinning elements!</p>
	<p>Caution! Hot surface can cause burns!</p>
	<p>For indoor use only.</p>

CAUTION!

The illustrations in this instruction manual are for reference only and may differ from the actual product in some details.

The original instruction manual is in the German language version. Other language versions are translations from German.

3. Safety of use

CAUTION! Read all safety warnings and all instructions. Failure to follow the warnings and instructions may result in electric shock, fire and/or severe personal injury or death.

The term "unit" or "product" in the warnings and in the description of the instructions refers to the Lower temp liquid circulating chiller.

2.1. Electrical safety

- a) The plug of this unit must fit into the outlet. Do not modify the plug in any way. Original plugs and matching outlets reduce the risk of electric shock.
- b) Avoid touching grounded parts, such as pipes, heaters, ovens, and refrigerators. There is an increased risk of electric shock if your body is grounded and touches the unit exposed to direct rain, wet pavement, and operation in a humid environment. If water enters the unit, there is an increased risk of damage to the unit and electric shock.
- c) Do not touch the unit with wet or damp hands.
- d) Do not use the cord in an unintended manner. Never use it to carry the unit or to pull the plug out of the socket. Keep the cord away from heat sources, oil, sharp edges or moving parts. Damaged or tangled cords increase the risk of electric shock.
- e) If you cannot avoid using the unit in a wet environment, use a residual current unit (RCD). Using an RCD reduces the risk of electric shock.
- f) Do not use the unit if the power cord is damaged or shows signs of wear. A damaged power cord should be replaced by a qualified electrician or the manufacturer's service department.
- g) To avoid electric shock, do not immerse the cable, plug, or device itself in water or other liquid. Do not use the device on wet surfaces.

2.2. Safety in the workplace

- a) Keep the work area tidy and well lit. Disorder or poor lighting can lead to accidents. Be foresighted, watch what you are doing and use common sense when using the unit.

-
- b) Do not use the unit in an explosive area, for example in the presence of flammable liquids, gases or dust.
 - c) If you find any damage or irregularities in the operation of the unit, immediately turn it off and report it to an authorized person.
 - d) If you have any doubts as to whether the unit is working properly or if it is damaged, contact the manufacturer's service department.
 - e) Only the manufacturer's service department can repair the unit. Do not carry out repairs yourself!
 - f) In case of open flames or fire, use only dry powder or snow (CO₂) fire extinguishers to extinguish the live equipment.
 - g) No children or unauthorized persons are allowed in the work area. (Inattention may result in loss of control of the unit.)
 - h) Use the unit in a well-ventilated area.
 - i) Check the condition of the safety stickers regularly. Replace them if they are illegible.
 - j) Keep these instructions for use for future reference. If the unit is to be passed on to a third party, the operating instructions must also be handed over together with the unit.
 - k) Keep the packaging and small assembly parts out of the reach of children.
 - l) Keep the unit away from children and animals.
 - m) When using this machine together with other machines, also follow the other instructions for use.

Please note! Keep children and other bystanders safe while operating the unit.

2.3. Personal safety

- a) Do not operate this product if you are tired, ill or under the influence of alcohol, drugs or medication that could impair your ability to operate the product.
- b) The unit may be operated by persons who are physically fit, capable of operating it and appropriately trained, and who have read this instruction manual and have been trained in occupational safety and health.
- c) The machine is not intended to be used by persons (including children) with reduced mental, sensory or intellectual functions or persons who lack experience and/or knowledge unless they are supervised or have been instructed by a person responsible for their safety on how to operate the machine.
- d) To prevent accidental start-up, make sure the switch is in the off position before connecting to a power source.

-
- e) The unit is not a toy. Children should be watched to ensure that they do not play with the unit.
 - f) Do not place your hands or any objects inside the running unit!

2.4. Safe use of the unit

- a) Do not overload the unit. Use tools that are suitable for the application. A correctly selected unit will do a better and safer job for which it was designed.
- b) Do not use the unit if the ON/OFF switch does not function properly (does not turn on and off). Devices that cannot be controlled by the switch are unsafe, cannot operate, and must be repaired.
- c) Disconnect the unit from the power supply before adjusting, cleaning, or servicing. This precaution reduces the risk of accidental start-up.
- d) Keep unused equipment out of the reach of children and out of the reach of anyone unfamiliar with the unit or this instruction manual. These units is dangerous in the hands of inexperienced users.
- e) Keep the unit in good working condition. Check before each use for general damage or damage to moving parts (cracks in parts and components or any other condition that may affect the safe operation of the unit). If damaged, have the unit repaired before use.
- f) Keep the unit out of the reach of children.
- g) Repairs and maintenance should be carried out by qualified personnel using only original spare parts. This will ensure the safety of use.
- h) To ensure the designed operational integrity of the unit, do not remove factory-installed covers or loosen screws.
- i) When transporting or moving the unit from storage to the place of use, observe the health and safety rules for manual handling applicable in the country where the unit is used.
- j) Do not move, shift, or rotate the machine while in operation.
- k) Do not leave the unit switched on unattended.
- l) Clean the unit regularly to prevent permanent dirt build-up.
- m) The unit is not a toy. Cleaning and maintenance must not be performed by children without adult supervision.
- n) Do not start up an empty machine.
- o) Do not tamper with the unit to alter its performance or design.
- p) Keep the unit away from sources of fire and heat.
- q) Do not block the ventilation openings of the unit!
 - r) The liquid level should be between minimum and maximum.
- s) Do not allow water to freeze in the unit.
- t) Replace the coolant if the unit's performance has deteriorated or if the refrigerant has become viscous.

-
- u) Only qualified persons should perform maintenance on the unit.
High-pressure gas is present in the unit.
 - v) Do not operate the unit's pump with an empty tank, when the liquid level is too low or when the liquid is frozen.
 - w) Always turn off the unit before opening the cover.
 - x) Do not pour corrosive, explosive or other substances into the unit.
 - y) Do not start cooling or circulation if the unit has an empty tank or the liquid level is too low.

CAUTION! Although the product has been designed to be safe, with adequate safeguards, and despite the additional safety features provided to the user, there is still a slight risk of accident or injury when handling the unit. You are advised to use caution and common sense when using this product.

3. Rules of use

The unit is designed to control and maintain a specific temperature by circulating fluid through the system.



The unit can be connected to electron microscopes, lasers, distillers, and other laboratory equipment requiring appropriate operating conditions.

The user is responsible for any damage resulting from misuse.

3.1. Description


SBS-LCC-3000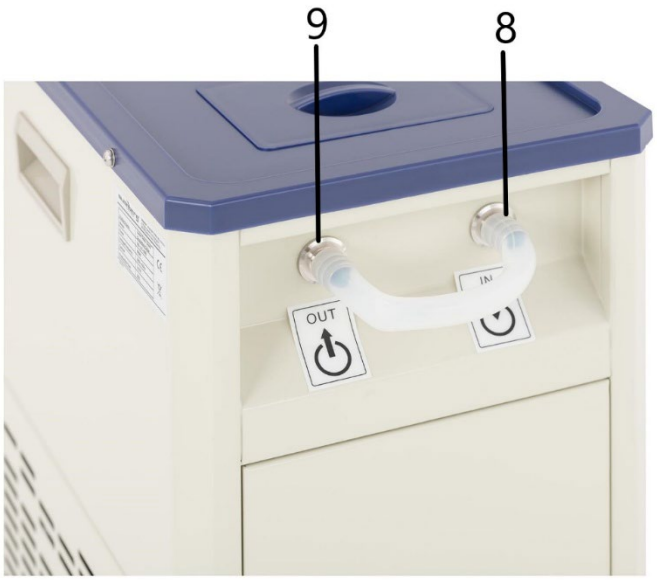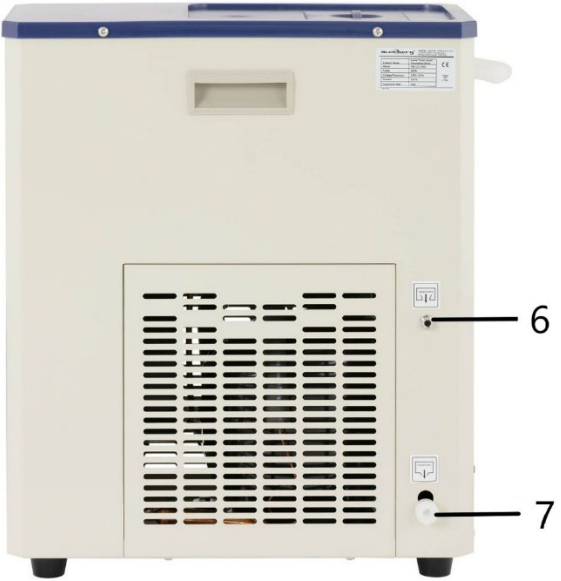

-
1. Cover
 2. Control panel
 3. Feet
 4. Fuse
 5. Handle
 6. Overflow
 7. Drain hole
 8. Input connector (inlet)
 9. Output connector (outlet)

Control panel

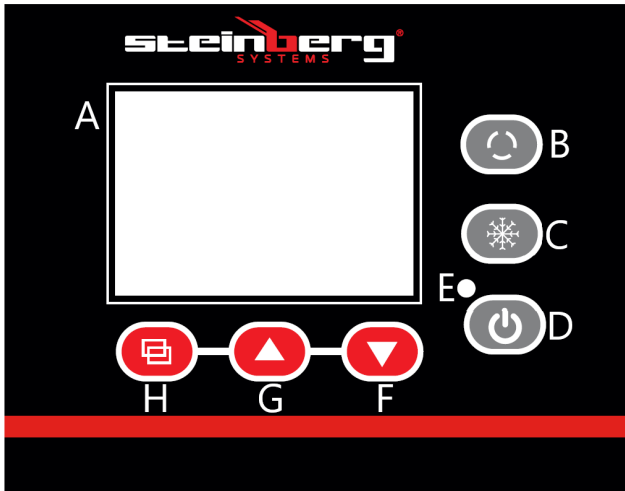- A - Display
- B - Cycle switch
- C - Cooling switch
- D - On/Off button
- E - Power light
- F - Decrease cooling button
- G - Increase cooling button
- H - SET button

SBS-LCC-5000

1. Cover
2. Control panel
3. Feet
4. Rod
5. Input connector (inlet)
6. Output valve (outlet)

Control panel

- A - Display
- B - Cycle switch
- C - Cooling switch
- D - On/Off button
- E - Power light
- F - Decrease cooling button
- G - Increase cooling button
- H - SET button

3.2. Preparation for operation

POSITIONING OF THE UNIT

The ambient temperature must not exceed 40°C and ambient humidity should not exceed 60%. Place the unit in a way that ensures good air circulation. Maintain a minimum clearance of 30 cm from any wall of the unit. Keep the unit away from any hot surfaces. Always operate the unit on a level, stable, clean, fireproof and dry surface and out of the reach of children and persons of impaired mental, sensory and intellectual functions. Place the unit in such a way that the mains plug can be reached at any time. Ensure that the power supply to the unit corresponds to that specified on the identification plate!

3.3. Working with the unit

3.3.1. Description of the buttons on the control panel

1. Cycle switch (B) - pressing the button will open/close the outlet
2. Cooling switch (C) - pressing the button turns the cooling mode on/off.
3. On/Off button (D) - pressing the button turns the control panel on/off.
Note: turning the display off does not cut off power to the unit.
4. Decrease value button (F) - in set-up mode pressing the button decreases the value and longer holding the button decreases the value in continuous mode.
5. Increase button (G) - in set-up mode pressing the button increases the value and holding the button longer increases the value in continuous mode.
In normal operation mode the button allows you to turn on/off the backlight.
6. Settings button (H) - pressing the button allows you to display temperature values and related parameters and to edit them.

3.3.2. Operation

Model SBS-LCC-3000

1. Open the cover (1) and fill with an appropriate amount of liquid (the type of liquid, e.g. demineralized water, glycol, etc., should be selected according to the temperature to be maintained by the unit so that it will not freeze).
The liquid level should be approx. 2 cm below the tank edges.
Examples of liquid mixtures:

-
- a) Ethylene glycol and demineralized water in proportion 55:45 - freezing point -40°C
 - b) Glycerol and demineralized water in proportion 70:30 - freezing point -30°C
2. Connect the inlet and outlet of the unit with the inlet and outlet of the laboratory instrument needing temperature control.

For the SBS-LCC-5000, you can also connect the inlet and outlet with each other when cooling a component placed in a tank. The item to be cooled (e.g., a laboratory beaker - not included) should be supported by, for example, a bracket attached to a rod (10).

Caution: Use an insulated hose for the connection.

Caution: The liquid at the inlet (coming out of the external device) should not exceed 40 °C for water, and 10 °C for solutions containing alcohol (e.g. ethanol).

3. Connect the device plug to a power source.
4. Switch on the display with the On/Off key (D). The display switches to the normal operating state after approx. 3s. The upper line of the display shows the current temperature (PV), and the lower line shows the desired temperature (SV).
5. Set the desired temperature by following the steps below:
 - enter the settings by pressing the settings button (H), the upper line of the display will show "SP", the lower line will show the temperature value to be edited using the buttons (F) and (G).
 - To save the settings and return to the main screen press the setting button (H).

Caution: After entering the settings, if the user does not press any button for about 1 min, the unit will automatically return to the main screen.

-
6. Press the cooling switch (C) to start the process of lowering the system temperature. The compressor will start after approx. 3 minutes. When the unit has reached the set temperature inside its system, open the output from the unit by pressing button (B). On the SBS-LCC-5000 model additionally remember to open the output valve (12).

 7. When you have finished using the unit, turn off the cooling by pressing button (C) and the display power by pressing button (D).
Disconnect the unit from the power supply.

Alarms:

- sensor failure / out of range temperature alarm - the top line of the display will show a flashing "E0" instead of the current temperature value

- Too low temperature alarm - the top line of the display will show a flashing message "E2" instead of the current temperature value

The beeping sound produced by the unit during an alarm can be silenced by pressing any button on the display.

Overload protection

The device is equipped with a fuse (4) for overload protection.

If the protection is triggered (the device switches off), find the cause and remove it.

Then replace the fuse with a new one.

3.3.3. Advanced settings

The device allows you to edit internal parameters of temperature measurement.

To enter the internal parameters setting mode press and hold the settings button (H) for over 3s.

The display screen will show "LC", which is a protection (password) against uncontrolled change of parameters.

Use the buttons (G) and (F) to edit the parameter value, and confirm your choice by pressing the setting button (H).

If the password is incorrect the display will return to the main screen, if correct it will go on to edit the parameter settings described in the table below.

Press the settings button (H) to move between the parameters, edit the parameter values with the buttons (F) and (G).

Press and hold the settings button (H) for more than 3 seconds to save the settings and return to the main screen.

Symbol	Name	Function description	Factory value (range)
Lc-	Password	"Lc=3" - parameter values can be viewed and modified	0
cL-	Deviation	If the difference between the measured value and the setpoint is greater than or equal to cL the compressor will be started, if the difference is less than the cL value the	2.0 (0.0~50.0°C)

		compressor will remain off.	
dL-	Low temperature alarm	If the measured value is less than the set value minus dL value, E2 alarm will appear.	2.0 (model SBS-LCC-3000) 5.0 (model SBS-LCC-5000) (0.0~50.0°C)
cb-	Zero adjustment (1)	Measured value <0, sensor measurement error (low temperature) $cb = x - y^{**}$	0.0 (-20.0~20.0°C)
ck-	Full adjustment (1)	Measured value <0, sensor measurement error (high temperature) $ck = 1000 * (x - y) / y^{**}$	0 (-999~999)
Pb-	Zero adjustment (2)	Measured value >0, correction of sensor measurement error (low temperature) $Pb = x - y^{**}$	0.0 (-20.0~20.0°C)
PK-	Full adjustment (2)	Measured value >0, correction for sensor measurement error (high temperature) $PK = 1000 * (x - y) / y^{**}$	0 (-999~999)

CT-	Compressor operation delay	Time between starting the cooling from the control level and the actual start of the compressor	3 (0~10)Min
-----	----------------------------	---	----------------

**where x - actual temperature value, y - measured value

3.4. Cleaning and maintenance

- a) Pull out the mains plug before each cleaning and when the unit is not in use.
- b) Use only non-corrosive cleaning agents for cleaning the surfaces.
- c) After each cleaning, all the parts should be dried well before the unit is used again.
- d) Store the unit in a dry and cool place protected from moisture and direct sunlight.
- e) Do not spray the unit with a stream of water or immerse it in water.
- f) Make sure that no water enters through the ventilation openings in the housing.
- g) Clean the ventilation openings with a brush and compressed air.
- h) Perform regular inspections of the unit checking technical fitness and any damages.
- i) Do not use sharp and/or metal objects (e.g. a wire brush or metal spatula) for cleaning, as these may damage the surface of the material from which the unit is made.
- j) Do not clean the unit with acidic substances, medical products, diluent, fuel, oil or other chemicals as this may cause damage to the unit.
- k) When not in use for an extended period of time or before storing, empty the unit's tank of any coolant poured in, following the instructions below.

Dispose of the collected liquid in accordance with local ordinances.

Model SBS-LCC-3000

Pull the hose extending from the drain hole (7). Place a container for the liquid, and then open the cap. Drain the liquid, replace the cap, and insert the hose.

Model SBS-LCC-5000

Place a liquid container under the outlet valve (12). Open the valve, drain liquid, and close the valve.

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