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LABORMISCHER

USER MANUAL

LABORATORY MIXER

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LABORATORNÍ MÍCHAČKA

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MISCELATORE DA LABORATORIO

MANUAL DE INSTRUCCIONES

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LABORATÓRIUMI KEVERŐ

BRUGSANVISNING

DREJESKIVE TIL MIKRORØR

DE | EN | PL | CZ | FR | IT | ES | HU | DA

SBS-LBM-200

Deutsch	3
English	6
Polski	9
Česky	12
Français	15
Italiano	18
Español	21
Magyar	24
Dansk	27

PRODUKTNAME	LABORMISCHER
PRODUCT NAME	LABORATORY MIXER
NAZWA PRODUKTU	MIESZADŁO LABORATORYJNE
NÁZEV VÝROBKU	LABORATORNÍ MÍCHAČKA
NOM DU PRODUIT	MÉLANGEUR DE LABORATOIRE
NOME DEL PRODOTTO	MISCELATORE DA LABORATORIO
NOMBRE DEL PRODUCTO	MEZCLADOR DE LABORATORIO
TERMÉK NEVE	LABORATÓRIUMI KEVERŐ
PRODUKTNÁVN	DREJESKIVE TIL MIKRORØR
MODELL	SBS-LBM-200
PRODUCT MODEL	
MODEL PRODUKTU	
MODEL VÝROBKU	
MODÈLE	
MODELLO	
MODELO	
MODELL	EXPONDO POLSKA SP. Z O.O. SP. K.
MODEL	
HERSTELLER	
MANUFACTURER	
PRODUCENT	
VÝROBCE	
FABRICANT	
PRODUTTORE	
FABRICANTE	
TERMELŐ	
PRODUCENT	
ANSCHRIFT DES HERSTELLERS	UL. NOWY KISIELIN-INNOWACYJNA 7, 66-002 ZIELONA GÓRA POLAND, EU
MANUFACTURER ADDRESS	
ADRES PRODUCENTA	
ADRESA VÝROBCE	
ADRESSE DU FABRICANT	
INDIRIZZO DEL PRODUTTORE	
DIRECCIÓN DEL FABRICANTE	
A GYÁRTÓ CÍME	
PRODUCENTENS ADRESSE	

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TECHNICAL DATA

Description of the parameter	Value of the parameter
Product name	LABORATORY MIXER
Model	SBS-LBM-200
Type of movement	orbital
Dimensions of the working platform surface [mm]	220
Number of mixing handles	15
Supply voltage [V~]/ Frequency [Hz]	230/50
Rated power [W].	15
Unit dimensions (Width x Depth x Height) [mm]	220x210x315
Weight [kg]	3,4
Rotational speed [RPM]	0-40
Permissible ambient temperature [°C]	5-40

1. GENERAL DESCRIPTION

The user manual is designed to assist in the safe and trouble-free use of the device. The product is designed and manufactured in accordance with strict technical guidelines, using state-of-the-art technologies and components. Additionally, it is produced in compliance with the most stringent quality standards.

DO NOT USE THE DEVICE UNLESS YOU HAVE THOROUGHLY READ AND UNDERSTOOD THIS USER MANUAL.

To increase the product life of the device and to ensure trouble-free operation, use it in accordance with this user manual and regularly perform maintenance tasks. The technical data and specifications in this user manual are up to date. The manufacturer reserves the right to make changes associated with quality improvement. The device is designed to reduce noise emission risks to a minimum, taking into account technological progress and noise reduction opportunities.

LEGEND

	The product satisfies the relevant safety standards.
	Read instructions before use.
	The product must be recycled.
	WARNING! or CAUTION! or REMEMBER! Applicable to the given situation (general warning sign).
	ATTENTION! Electric shock warning!

PLEASE NOTE! Drawings in this manual are for illustration purposes only and in some details may differ from the actual machine.

The original operation manual is in German. Other language versions are translations from German.

2. USAGE SAFETY

ATTENTION! Read all safety warnings and all instructions. Failure to follow the warnings and instructions may result in an electric shock, fire and/or serious injury or even death.

The terms „device“ or „product“ are used in the warnings and instructions to refer to LABORATORY MIXER.

2.1. ELECTRICAL SAFETY

- The plug must fit the socket. Do not modify the plug in any way. Using original plugs and matching sockets reduces the risk of electric shock.
- Avoid touching earthed elements such as pipes, heaters, boilers and refrigerators. There is an increased risk of electric shock if the earthed device is exposed to rain, comes into direct contact with a wet surface or is operating in a damp environment. Water getting into the device increases the risk of damage to the device and of electric shock.
- Do not touch the device with wet or damp hands.
- Use the cable only for its designated use. Never use it to carry the device or to pull the plug out of a socket. Keep the cable away from heat sources, oil, sharp edges or moving parts. Damaged or tangled cables increase the risk of electric shock.
- If using the device in a damp environment cannot be avoided, a residual current device (RCD) should be applied. The use of an RCD reduces the risk of electric shock.
- Do not use the device if the power cord is damaged or shows obvious signs of wear. A damaged power cord should be replaced by a qualified electrician or the manufacturer's service centre.
- To avoid electric shock, do not immerse the cord, plug or device in water or other liquids. Do not use the device on wet surfaces.
- Use only grounded power sources.

2.2. SAFETY IN THE WORKPLACE

- Make sure the workplace is clean and well lit. A messy or poorly lit workplace may lead to accidents. Try to think ahead, observe what is going on and use common sense when working with the device.
- If you discover damage or irregular operation, immediately switch the device off and report it to a supervisor without delay.
- If there are any doubts as to the correct operation of the device, contact the manufacturer's support service.
- Only the manufacturer's service point may repair the device. Do not attempt any repairs independently!
- In case of fire, use a powder or carbon dioxide (CO₂) fire extinguisher (one intended for use on live electrical devices) to put it out.
- Please keep this manual available for future reference. If this device is passed on to a third party, the manual must be passed on with it.
- Keep packaging elements and small assembly parts in a place not available to children.

- Keep the device away from children and animals.
- If this device is used together with another equipment, the remaining instructions for use shall also be followed.

REMEMBER! When using the device, protect children and other bystanders.

2.3. PERSONAL SAFETY

- Do not use the device when tired, ill or under the influence of alcohol, narcotics or medication which can significantly impair the ability to operate the device.
- The device is not designed to be handled by persons (including children) with limited mental and sensory functions or persons lacking relevant experience and/or knowledge unless they are supervised by a person responsible for their safety or they have received instruction on how to operate the device.
- To prevent the device from accidentally switching on, make sure the switch is on the OFF position before connecting to a power source.
- Do not wear loose clothing or jewellery. Keep hair, clothes and gloves away from moving parts. Loose clothing, jewellery or long hair may get caught in moving parts.
- The device is not a toy. Children must be supervised to ensure that they do not play with the device.

2.4. SAFE DEVICE USE

- Do not use the device if the knobs or control buttons do not function properly. Devices that cannot be controlled by the controls are dangerous, may not be operated and must be repaired.
- Make sure the plug is disconnected from the socket before attempting any adjustments, accessory replacements or before putting the device aside. Such precautions will reduce the risk of accidentally activating the device.
- When not in use, store in a safe place, away from children and people not familiar with the device who have not read the user manual. The device may pose a hazard in the hands of inexperienced users.
- Keep the device in perfect technical condition. Before each use check for general damage and especially check for cracked parts or elements and for any other conditions which may impact the safe operation of the device. If damage is discovered, hand over the device for repair before use.
- Keep the device out of the reach of children.
- Device repair or maintenance should be carried out by qualified persons, only using original spare parts. This will ensure safe use.
- To ensure the operational integrity of the device, do not remove factory-fitted guards and do not loosen any screws.
- Avoid situations where the device stops working during use due to excessive loading. This may result in overheating of the drive elements and damage to the device.
- Do not touch articulated parts or accessories unless the device has been disconnected from the power source.
- Do not move, adjust or rotate the device in the course of work.
- Do not leave this appliance unattended while it is in use.

- Clean the device regularly to prevent stubborn grime from accumulating.
- The device is not a toy. Cleaning and maintenance may not be carried out by children without supervision by an adult person.
- It is forbidden to interfere with the structure of the device in order to change its parameters or construction.
- Keep the device away from sources of fire and heat.
- Do not put your hands between the rollers of the device and do not put any foreign objects between them, otherwise the device may be damaged.

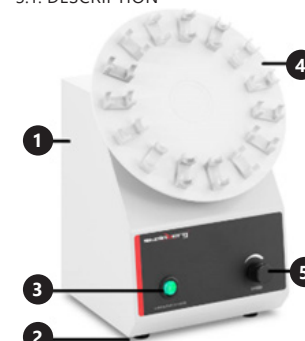
ATTENTION! Despite the safe design of the device and its protective features, and despite the use of additional elements protecting the operator, there is still a slight risk of accident or injury when using the device. Stay alert and use common sense when using the device.

3. USE GUIDELINES

The unit with a special design of the inclined circular stirrer is intended for mixing biological liquids characterized by high density and high effective properties.

The user is liable for any damage resulting from unintended use of the device.

3.1. DESCRIPTION


- Unit housing
- Anti-slip base
- Power switch
- Circular agitator platform
- Speed regulation knob

3.2. PREPARATION FOR OPERATION POSITIONING OF THE UNIT

The ambient temperature must not exceed 40°C. Place the unit in a way that ensures good air circulation. Maintain a minimum clearance of 10 cm from any wall of the unit. Keep the unit away from any hot surfaces. Always operate the unit on a level, stable, clean, fireproof and dry surface and out of the reach of children and persons of impaired mental, sensory and intellectual functions. Place the unit in such a way that the mains plug can be reached at any time. Ensure that the power supply to the unit corresponds to that specified on the identification plate!

3.3. WORKING WITH THE UNIT

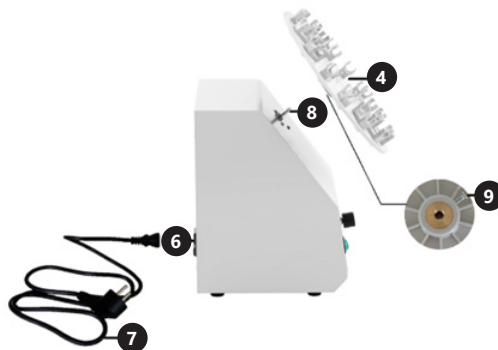CAUTION! Before working on the unit, make sure that all components are complete and as expected! In case of discrepancies, contact the supplier immediately!

Slide the circular agitator platform (4) onto the stem of the main unit (8). Secure from the bottom with the assembly screw (9). Check for stability, tighten if necessary. Connect the power cord (7) to the unit's power output (6). After completing these installation steps, you can start working with the unit.

3.3. WORKING WITH THE UNIT

- Plug the machine into a power outlet.
- Start up the unit by setting the power switch to the "I" position. The power indicator light will turn on. Setting the switch to the "0" position will turn off the unit.
- When the unit is turned on, the mixing process will begin automatically.
- Set the desired speed value using the dial. Turn the knob clockwise to increase the rpm, Turn the knob to the left to decrease rpm.
- If the process time is over, turn the unit off by reducing the speed and setting the power switch to the "0" position.

3.4. CLEANING AND MAINTENANCE

- Pull out the mains plug before each cleaning and when the unit is not in use.
- Use only non-corrosive cleaning agents for cleaning the surfaces.
- After each cleaning, all the parts should be dried well before the unit is used again.
- Store the unit in a dry and cool place protected from moisture and direct sunlight.
- Do not spray the unit with a stream of water or immerse it in water.
- Make sure that no water enters through the ventilation openings in the casing.
- Clean the ventilation openings with a brush and compressed air.
- Perform regular inspections of the unit checking technical fitness and any damages.
- Use a soft cloth for cleaning.

- Do not use sharp and/or metal objects (e.g. a wire brush or metal spatula) for cleaning, as these may damage the surface of the material from which the unit is made.
- Do not clean the unit with acidic substances, medical products, diluent, fuel, oil or other chemicals as this may cause damage to the unit.

DISPOSAL OF USED UNITS.

At the end of its useful life, this product should not be disposed of with normal household waste but should be taken to a collection point for the recycling of electrical and electronic equipment. This is indicated by the symbol on the product, operating instructions or packaging. The materials used in this unit are recyclable according to their marking. You will be making an important contribution to protecting our environment by reusing, recycling or otherwise disposing of used units. Your local administration will provide you with information about the appropriate disposal point for used units.

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