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USER MANUAL

MAGNETIC STIRRER

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MANUAL DE INSTRUCCIONES

AGITADOR MAGNÉTICO

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BRUGSANVISNING

MAGNETOMRØRER

DE | EN | PL | CZ | FR | IT | ES | HU | DA

SBS-MR-I500/H | SBS-MR-I500

Deutsch	3
English	7
Polski	10
Česky	14
Français	17
Italiano	21
Español	25
Magyar	30
Dansk	32

PRODUKTNAME	MAGNETRÜHRER
PRODUCT NAME	MAGNETIC STIRRER
NAZWA PRODUKTU	MIESZADŁO MAGNETYCZNE
NÁZEV VÝROBKU	MAGNETICKÉ MÍCHADLO
NOM DU PRODUIT	AGITATEUR MAGNÉTIQUE
NOME DEL PRODOTTO	AGITATORE MAGNETICO
NOMBRE DEL PRODUCTO	AGITADOR MAGNÉTICO
TERMÉK NEVE	MÉGNESES KEVERŐ
PRODUKTNÁVN	MAGNETOMRØRER
MODELL	SBS-MR-1500/H SBS-MR-1500
PRODUCT MODEL	
MODEL PRODUKTU	
MODEL VÝROBKU	
MODÈLE	
MODELLO	
MODELO	
MODELL	
MODEL	
HERSTELLER	EXPONDO POLSKA SP. Z O.O. SP. K.
MANUFACTURER	
PRODUCENT	
VÝROBCE	
FABRICANT	
PRODUTTORE	
FABRICANTE	
TERMELŐ	
PRODUCENT	
ANSCHRIFT DES HERSTELLERS	UL. NOWY KISIELIN-INNOWACYJNA 7, 66-002 ZIELONA GÓRA POLAND, EU
MANUFACTURER ADDRESS	
ADRES PRODUCENTA	
ADRESA VÝROBCE	
ADRESSE DU FABRICANT	
INDIRIZZO DEL PRODUTTORE	
DIRECCIÓN DEL FABRICANTE	
A GYÁRTÓ CÍME	
PRODUCENTENS ADRESSE	

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1. GENERAL DESCRIPTION

The user manual is designed to assist in the safe and trouble-free use of the device. The product is designed and manufactured in accordance with strict technical guidelines, using state-of-the-art technologies and components. Additionally, it is produced in compliance with the most stringent quality standards.

DO NOT USE THE DEVICE UNLESS YOU HAVE THOROUGHLY READ AND UNDERSTOOD THIS USER MANUAL.

To increase the product life of the device and to ensure trouble-free operation, use it in accordance with this user manual and regularly perform maintenance tasks. The technical data and specifications in this user manual are up to date. The manufacturer reserves the right to make changes associated with quality improvement. The device is designed to reduce noise emission risks to a minimum, taking into account technological progress and noise reduction opportunities.

EXPLANATION OF SYMBOLS

	The product satisfies the relevant safety standards.
	Read instructions before use.
	Recyclable product.
	CAUTION! or WARNING! or REMINDER! describing a situation (general warning sign).
	CAUTION! Warning of electric shock!
	PLEASE NOTE! The illustrations in this instruction manual are for reference only and may differ from the actual product in some details.

2. SAFETY OF USE

⚠ ATTENTION! Read all safety warnings and all instructions. Failure to follow the warnings and instructions may result in electric shock, fire and/or severe personal injury or death.

The term „appliance“ or „product“ in the warnings and in the description of the instructions refers to the MAGNETIC MIXER.

⚠ CAUTION! Do not use the device in rooms with very high humidity / in the immediate vicinity of water tanks! Do not allow the machine to get wet. Risk of electric shock!

2.1. ELECTRICAL SAFETY

- The plug of this unit must fit into the outlet. Do not modify the plug in any way. Original plugs and matching outlets reduce the risk of electric shock.
- Avoid touching grounded parts, such as pipes, heaters, ovens, and refrigerators. There is an increased risk of electric shock if your body is grounded and touches the device exposed to direct rain, wet pavement, and operation in a humid environment. If water enters the device, there is an increased risk of damage to the device and electric shock.
- Do not touch the unit with wet or damp hands.
- Do not use the cord in an unintended manner. Never use it to carry the unit or to pull the plug out of the socket. Keep the cord away from heat sources, oil, sharp edges or moving parts. Damaged or tangled cords increase the risk of electric shock.
- If you cannot avoid using the unit in a wet environment, use a residual current unit (RCD). Using an RCD reduces the risk of electric shock.
- Do not use the unit if the power cord is damaged or shows signs of wear. A damaged power cord should be replaced by a qualified electrician or the manufacturer's service department.
- To avoid electric shock, do not immerse the cable, plug, or unit itself in water or other liquid. Do not use the unit on wet surfaces.

2.2. SAFETY IN THE WORKPLACE

- Keep the work area tidy and well lit. Disorder or poor lighting can lead to accidents. Be foresighted, watch what you are doing and use common sense when using the unit.
- Do not use the appliance in an explosive area, for example in the presence of flammable liquids, gases or dust. The appliance produces sparks that can ignite dust or fumes.
- If you find any damage or irregularities in the operation of the unit, immediately turn it off and report it to an authorized person.
- If you have any doubts as to whether the unit is working properly or if it is damaged, contact the manufacturer's service department.
- Only the manufacturer's service department can repair the unit. Do not carry out repairs yourself!
- In case of open flames or fire, use only dry powder or snow (CO₂) fire extinguishers to extinguish the live equipment.
- Keep these instructions for use for future reference. If the unit is to be passed on to a third party, the operating instructions must also be handed over together with the unit.
- Keep the packaging and small assembly parts out of the reach of children.
- Keep the unit away from children and animals.

2.3. PERSONAL SAFETY

- Do not operate this unit if you are tired, ill or under the influence of alcohol, drugs or medication that could impair your ability to operate the unit.
- The unit is not intended to be used by persons (including children) with reduced mental, sensory or intellectual functions or persons who lack experience and/or knowledge unless they are supervised or have been instructed by a person responsible for their safety on how to operate the unit.
- To prevent accidental start-up, make sure the switch is in the off position before connecting to a power source.
- The unit is not a toy. Children should be watched to ensure that they do not play with the unit.

2.4. SAFE USE OF THE UNIT

- Do not use the unit if the knobs or controls do not operate smoothly. Units that cannot be controlled by the switch are unsafe, cannot operate, and must be repaired.
- Unplug the unit before making adjustments, changing accessories, or putting it away. This precaution reduces the risk of accidental start-up.
- Disconnect the unit from the power supply before adjusting, cleaning, or servicing. This precaution reduces the risk of accidental start-up.
- Keep unused equipment out of the reach of children and out of the reach of anyone unfamiliar with the unit or this instruction manual. These units are dangerous in the hands of inexperienced users.
- Keep the unit in good working condition. Check before each use for general damage or damage to moving parts (cracks in parts and components or any other condition that may affect the safe operation of the unit). If damaged, have the unit repaired before use.
- Keep the unit out of the reach of children.
- Repairs and maintenance should be carried out by qualified personnel using only original spare parts. This will ensure the safety of use.
- To ensure the designed operational integrity of the unit, do not remove factory-installed covers or loosen screws.
- Do not move, shift, or rotate the machine while in operation.
- Do not leave the unit switched on unattended.
- Clean the unit regularly to prevent permanent dirt build-up.
- The unit is not a toy. Cleaning and maintenance must not be performed by children without adult supervision.
- Do not tamper with the unit to alter its performance or design.
- Keep the unit away from sources of fire and heat.
- Keep the unit away from strong magnetic fields.
- Do not switch on the machine under an empty vessel.
- If the machine is turned on at the highest speed, the contents of the vessel placed on the mixing plate may splash or the vessel may be thrown off balance and overturned.
- Mixing flammable, explosive, corrosive or similar substances with this equipment is prohibited.

ATTENTION! Despite the safe design of the device and its protective features, and despite the use of additional elements protecting the operator, there is still a slight risk of accident or injury when using the device. Stay alert and use common sense when using the device.

3. USE GUIDELINES

The device is used for non-contact mixing of liquids of water density. The product is designed for use in laboratory conditions.

The user is liable for any damage resulting from unintended use of the device.

3.1. DEVICE DESCRIPTION

1. Work stand surface
2. Speed regulation knob
3. LCD display
4. Casing

CONTROL PANEL

3.2. PREPARATION FOR OPERATION POSITIONING OF THE UNIT

The ambient temperature must not exceed 40°C and ambient humidity should not exceed 80%. Place the unit in a way that ensures good air circulation. Maintain a minimum clearance of 10 cm from any wall of the unit. Keep the unit away from any hot surfaces. Always operate the unit on a level, stable, clean, fireproof and dry surface and out of the reach of children and persons of impaired mental, sensory and intellectual functions. Place the unit in such a way that the mains plug can be reached at any time. Ensure that the power supply to the unit corresponds to that specified on the identification plate!

3.3. WORKING WITH THE UNIT

CAUTION! Before working on the device, make sure that all components are complete and as expected! In case of discrepancies, contact the supplier immediately!

DEFINING THE OPERATION OF THE DEVICE

- Plug the machine into a power outlet.
- Turn on the unit by setting the power switch to the „I“ position:
SBS-MR-1500/H - rear of the unit,
SBS-MR-1500 - side of the device.

- Set the desired speed of the device with the knob. Turn the knob clockwise to increase the rpm, Turn the knob to the left to decrease rpm.
- The selected device speed will be displayed at the bottom of the LCD, with the message „SET“ (1 - see figure).
- Once defined, the value of the device speed will be saved.
- To start the appliance, press the speed control knob.
- The current device speed value is shown in the upper part of the LCD display, with the message „STIR“ (2 - see figure). This value gradually increases until the defined value is reached.

3.4. CLEANING AND MAINTENANCE

- Cleaning and maintenance**
Pull the mains plug and let the unit cool down completely before cleaning, adjusting or replacing accessories and when the unit is not in use. Wait until the rotating parts stop.
- Pull out the mains plug before each cleaning and when the device is not in use.
- Use only non-corrosive cleaning agents for cleaning the surfaces.
- After each cleaning, all the parts should be dried well before the unit is used again.
- Store the unit in a dry and cool place protected from moisture and direct sunlight.

- Do not spray the unit with a stream of water or immerse it in water.
- Make sure that no water enters through the ventilation openings in the casing.
- Clean the ventilation openings with a brush and compressed air.
- Perform regular inspections of the unit checking technical fitness and any damages.
- Use a soft, damp cloth for cleaning.
- Do not use sharp and/or metal objects (e.g. a wire brush or metal spatula) for cleaning, as these may damage the surface of the material from which the unit is made.
- Do not clean the unit with acidic substances, medical products, diluent, fuel, oil or other chemicals as this may cause damage to the unit.

DISPOSAL OF USED UNITS.

At the end of its useful life, this product should not be disposed of with normal household waste but should be taken to a collection point for the recycling of electrical and electronic equipment. This is indicated by the symbol on the product, operating instructions or packaging. The materials used in this unit are recyclable according to their marking. You will be making an important contribution to protecting our environment by reusing, recycling or otherwise disposing of used units. Your local administration will provide you with information about the appropriate disposal point for used units.

TROUBLESHOOTING

Problem	Possible cause	Action
The unit does not turn on	1. The power cord is not connected to the electrical outlet 2. The fuse is damaged	1. Plug the power cord into an electrical outlet 2. Replace the fuse

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