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LABORZENTRIFUGE

USER MANUAL

ELECTRIC CENTRIFUGE

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LABORATORNÍ CENTRIFUGA

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CENTRIFUGA DA LABORATORIO

MANUAL DE INSTRUCCIONES

CENTRÍFUGA PARA LABORATORIO

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ASZTALI CENTRIFUGA

BRUGSANVISNING

LABORATORIECENTRIFUGE

DE | EN | PL | CZ | FR | IT | ES | HU | DA

SBS-LZ-400B

Deutsch	3
English	7
Polski	11
Česky	15
Français	19
Italiano	23
Español	27
Magyar	31
Dansk	35

PRODUKTNAME PRODUCT NAME NAZWA PRODUKTU NÁZEV VÝROBKU NOM DU PRODUIT NOME DEL PRODOTTO NOMBRE DEL PRODUCTO TERMÉK NÉV PRODUKTNÁVN	LABORZENTRIFUGE ELECTRIC CENTRIFUGE WIRÓWKA LABORATORYJNA LABORATORNÍ CENTRIFUGA CENTRIFUGEUSE DE LABORATOIRE CENTRIFUGA DA LABORATORIO CENTRÍFUGA PARA LABORATORIO LABOR CENTRIFUGA LABORATORIECENTRIFUGE
MODELL PRODUCT MODEL MODEL PRODUKTU MODEL VÝROBKU MODÈLE MODELLO MODELO TÍPUS MODEL	SBS-LZ-400B
HERSTELLER MANUFACTURER PRODUCENT VÝROBCE FABRICANT PRODUTTORE FABRICANTE GYÁRTÓ PRODUCENT	EXPONDO POLSKA SP. Z O.O. SP. K.
ANSCHRIFT DES HERSTELLERS MANUFACTURER ADDRESS ADRES PRODUCENTA ADRESA VÝROBCE ADRESSE DU FABRICANT INDIRIZZO DEL PRODUTTORE DIRECCIÓN DEL FABRICANTE GYÁRTÓ CÍME PRODUCENTENS ADRESSE	UL. NOWY KISIELIN-INNOWACYJNA 7, 66-002 ZIELONA GÓRA POLAND, EU

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1. GENERAL DESCRIPTION

The user manual is designed to assist in the safe and trouble-free use of the device. The product is designed and manufactured in accordance with strict technical guidelines, using state-of-the-art technologies and components. Additionally, it is produced in compliance with the most stringent quality standards.

DO NOT USE THE DEVICE UNLESS YOU HAVE THOROUGHLY READ AND UNDERSTOOD THIS USER MANUAL.

To increase the product life of the device and to ensure trouble-free operation, use it in accordance with this user manual and regularly perform maintenance tasks. The technical data and specifications in this user manual are up to date. The manufacturer reserves the right to make changes associated with quality improvement. The device is designed to reduce noise emission risks to a minimum, taking into account technological progress and noise reduction opportunities.

LEGEND

	The product satisfies the relevant safety standards.
	Read instructions before use.
	The product must be recycled.
	WARNING! or CAUTION! or REMEMBER! Applicable to the given situation. (general warning sign).
	Wear protective goggles.
	Wear protective gloves.

	Wear protective gloves.
	Wear a protective apron.
	CAUTION! Warning of electric shock!
	CAUTION! Spinning elements!

PLEASE NOTE! Drawings in this manual are for illustration purposes only and in some details may differ from the actual product.

The original operation manual is in German. Other language versions are translations from German.

2. USAGE SAFETY

ATTENTION! Read all safety warnings and all instructions. Failure to follow the warnings and instructions may result in an electric shock, fire and/or serious injury or even death.

The terms „device“ or „product“ are used in the warnings and instructions to refer to ELECTRIC CENTRIFUGE.

2.1. ELECTRICAL SAFETY

- a) The plug must fit the socket. Do not modify the plug in any way. Using original plugs and matching sockets reduces the risk of electric shock.
- b) Avoid touching earthed elements such as pipes, heaters, boilers and refrigerators. There is an increased risk of electric shock if the earthed device is exposed to rain, comes into direct contact with a wet surface or is operating in a damp environment. Water getting into the device increases the risk of damage to the device and of electric shock.
- c) Do not touch the device with wet or damp hands.
- d) Use the cable only for its designated use. Never use it to carry the device or to pull the plug out of a socket. Keep the cable away from heat sources, oil, sharp edges or moving parts. Damaged or tangled cables increase the risk of electric shock.
- e) If using the device in a damp environment cannot be avoided, a residual current device (RCD) should be applied. The use of an RCD reduces the risk of electric shock.
- f) Do not use the device if the power cord is damaged or shows obvious signs of wear. A damaged power cord should be replaced by a qualified electrician or the manufacturer's service centre.
- g) To avoid electric shock, do not immerse the cord, plug or device in water or other liquids. Do not use the device on wet surfaces.
- h) ATTENTION! DANGER TO LIFE! While cleaning, never immerse the device in water or other liquids.
- i) Do not use in very humid environments or in the direct vicinity of water tanks.
- j) Prevent the device from getting wet. Risk of electric shock!

2.2. SAFETY IN THE WORKPLACE

- a) Make sure the workplace is clean and well lit. A messy or poorly lit workplace may lead to accidents. Try to think ahead, observe what is going on and use common sense when working with the device.
- b) If you discover damage or irregular operation, immediately switch the device off and report it to a supervisor without delay.

- c) If there are any doubts as to the correct operation of the device, contact the manufacturer's support service.
- d) Only the manufacturer's service point may repair the device. Do not attempt any repairs independently!
- e) In case of fire, use a powder or carbon dioxide (CO₂) fire extinguisher (one intended for use on live electrical devices) to put it out.
- f) Children or unauthorised persons are forbidden to enter a work station. (A distraction may result in loss of control over the device).
- g) Regularly inspect the condition of the safety labels. If the labels are illegible, they must be replaced.
- h) Please keep this manual available for future reference. If this device is passed on to a third party, the manual must be passed on with it.
- i) Keep packaging elements and small assembly parts in a place not available to children.
- j) Keep the device away from children and animals.

REMEMBER! When using the device, protect children and other bystanders.

2.3. PERSONAL SAFETY

- a) Do not use the device when tired, ill or under the influence of alcohol, narcotics or medication which can significantly impair the ability to operate the device.
- b) The device is not designed to be handled by persons (including children) with limited mental and sensory functions or persons lacking relevant experience and/or knowledge unless they are supervised by a person responsible for their safety or they have received instruction on how to operate the device.
- c) The device can be handled only by physically fit persons who are capable of handling it, properly trained, familiar with this manual and trained within the scope of occupational health and safety.
- d) When working with the device, use common sense and stay alert. Temporary loss of concentration while using the device may lead to serious injuries.
- e) Use personal protective equipment as required for working with the device, specified in section 1 (Legend). The use of correct and approved personal protective equipment reduces the risk of injury.
- f) To prevent the device from accidentally switching on, make sure the switch is on the OFF position before connecting to a power source.
- g) Do not wear loose clothing or jewellery. Keep hair, clothes and gloves away from moving parts. Loose clothing, jewellery or long hair may get caught in moving parts.
- h) The device is not a toy. Children must be supervised to ensure that they do not play with the device.
- i) Do not put your hands or other items inside the device while it is in use!

2.4. SAFE DEVICE USE

- a) Do not use the device if the ON/OFF switch does not function properly (does not switch the device on and off). Devices which cannot be switched on and off using the ON/OFF switch are hazardous, should not be operated and must be repaired.
- b) Disconnect the device from the power supply before commencement of adjustment, cleaning and maintenance. Such a preventive measure reduces the risk of accidental activation.

- c) When not in use, store in a safe place, away from children and people not familiar with the device who have not read the user manual. The device may pose a hazard in the hands of inexperienced users.
- d) Keep the device in perfect technical condition. Before each use check for general damage and especially check for cracked parts or elements and for any other conditions which may impact the safe operation of the device. If damage is discovered, hand over the device for repair before use.
- e) Keep the device out of the reach of children.
- f) Device repair or maintenance should be carried out by qualified persons, only using original spare parts. This will ensure safe use.
- g) To ensure the operational integrity of the device, do not remove factory-fitted guards and do not loosen any screws.
- h) Avoid situations where the device stops working during use due to excessive loading. This may result in overheating of the drive elements and damage to the device.
- i) Do not touch articulated parts or accessories unless the device has been disconnected from the power source.
- j) Do not move, adjust or rotate the device in the course of work.
- k) Do not leave this appliance unattended while it is in use.
- l) Clean the device regularly to prevent stubborn grime from accumulating.
- m) The device is not a toy. Cleaning and maintenance may not be carried out by children without supervision by an adult person.
- n) It is forbidden to interfere with the structure of the device in order to change its parameters or construction.
- o) Keep the device away from sources of fire and heat.
- p) Do not overload the device.
- q) Do not cover the ventilation openings!
- r) A flammable, explosive or corrosive test sample may damage or shorten the life of the device. Any damage caused by the above-mentioned type test samples are not covered by the warranty.
- s) The maximum rotor speed must not be exceeded.
- t) The condition of the rotor should be frequently checked (for cracks, corrosion, damaged seals, scratches, etc.). Do not use the device if the rotor is not in good technical condition.
- u) You must never open the device cover while it is working! The lid can be opened with the device switched off and the rotor at a standstill.
- v) Uneven alignment or use of damaged test tubes can cause severe vibrations and damage the device.

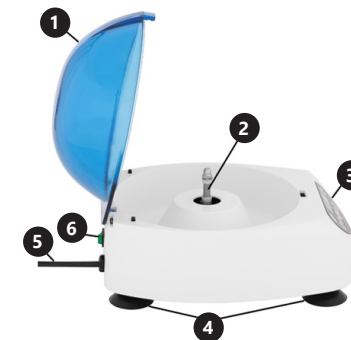
ATTENTION! Despite the safe design of the device and its protective features, and despite the use of additional elements protecting the operator, there is still a slight risk of accident or injury when using the device. Stay alert and use common sense when using the device.

3. USE GUIDELINES

The unit is used to separate various mixtures, suspensions and body fluids into components of different densities. The product is outside the scope of the MDD 93/42/ EEC regulations for medical units. It is not a medical unit and may not be used for examinations and tests in hospitals and medical centers.

The user is liable for any damage resulting from unintended use of the device.

3.1. DEVICE DESCRIPTION


1. Cover
2. Stem
3. Control panel
4. Feet
5. Power cord
6. On/Off button

CONTROL PANEL DESCRIPTION

- A. Program selection button
- B. Display
- C. Cover opening button
- D. Button for parameter selection: rotational speed (RMP) or relative centrifugal force (RCF)
- E. Up button
- F. SET button
- G. Down button
- H. Start/Stop button

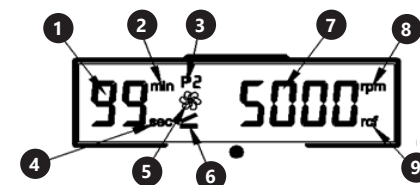Description of the displayed parameters (B)

1. Remaining time to the end of the program (timer)
2. Time unit - minutes
3. Number of the currently selected program
4. Time unit - seconds
5. Symbol - unit in operation
6. Symbol - cover open
7. Speed or centrifugal force value
8. Type of displayed value - speed
9. Type of displayed value - relative centrifugal force

3.2. PREPARATION FOR OPERATION

POSITIONING OF THE UNIT

The ambient temperature must not exceed 40°C and ambient humidity should not exceed 85%. Place the unit in a way that ensures good air circulation. Maintain a minimum clearance of 30 cm from any wall of the unit. Keep the unit away from any hot surfaces. Always operate the unit on a level, stable, clean, fireproof and dry surface and out of the reach of children and persons of impaired mental, sensory and intellectual functions. Place the unit in such a way that the mains plug can be reached at any time. Ensure that the power supply to the unit corresponds to that specified on the identification plate!

ASSEMBLY OF THE UNIT

Place the unit on a flat, stable surface so that all four rubber feet touch the surface of the ground. Caution! The power outlet must be grounded. This will reduce the risk of electric shock.

INSTALLING THE ROTOR

Unscrew the nut on the stem using the supplied wrench. Put on the rotor. Tighten the nut.

3.3. WORKING WITH THE UNIT

1. Check if rotor nut is loose, if so tighten it.
2. Prepare test tubes with the preparation for centrifugation and place them in the unit so that they meet the following requirements:
 - the tube should be filled to a maximum of 75% of its total volume,
 - place the tubes symmetrically in the rotor.
 - The sample tubes facing each other must be of equal weight.
 - Always verify the condition of the tubes being used. Do not place cracked or deformed tubes in the unit!

3. Close the unit cover.
4. Switch on the unit connected to the power supply using the I/O button on the back of the unit. The unit beeps and the display starts up.
5. Set the unit parameters or select one of the stored programs (select your preferred program by pressing button (A)):
 - by pressing the setting button „SET” (F) select which parameter is to be edited. Possible parameters to edit are: speed value, maximum relative centrifugal force value, timer and acceleration and deceleration values.
 - Use the up (E) and down (G) buttons to set the preferred parameter values.
6. Press button (H) to start the unit. The current value of the speed and the time remaining to the end of centrifuging will appear on the display.
7. To stop the machine before the set time has elapsed press button (H).
8. To open the unit cover press the (C) button. Caution: Opening the cover is possible only after the unit has come to a complete stop.
9. Remove the tubes with the centrifuged samples.
10. Turn off the unit by pressing the I/O button. Disconnect the unit from the power source when not in use.

3.4. CLEANING AND MAINTENANCE

- a) Pull out the mains plug before each cleaning and when the unit is not in use.
- b) Use only non-corrosive cleaning agents for cleaning the surfaces.
- c) After each cleaning, all the parts should be dried well before the unit is used again.
- d) Store the unit in a dry and cool place protected from moisture and direct sunlight.
- e) Do not spray the unit with a stream of water or immerse it in water.
- f) Make sure that no water enters through the ventilation openings in the casing.
- g) Clean the ventilation openings with a brush and compressed air.
- h) Perform regular inspections of the unit checking technical fitness and any damages.
- i) Use a soft, damp cloth for cleaning.
- j) Do not use sharp and/or metal objects (e.g. a wire brush or metal spatula) for cleaning, as these may damage the surface of the material from which the unit is made.
- k) Do not clean the unit with acidic substances, medical products, diluent, fuel, oil or other chemicals as this may cause damage to the unit.
- l) CAUTION! If there is concern that the equipment is contaminated with toxic, pathogenic, radioactive, or infectious blood samples that are harmful to health. Immediately sterilize the equipment according to current laboratory methods.
- m) Periodically lubricate the central part of the rotor.

TROUBLESHOOTING

Error code	Meaning
E1	Cover not closed
E2	Motor drive does not work
E6	Parameter setting error
E7	
E8	Speed exceeded
E9	Unbalanced load distribution

DISPOSAL OF USED UNITS.

At the end of its useful life, this product should not be disposed of with normal household waste but should be taken to a collection point for the recycling of electrical and electronic equipment. This is indicated by the symbol on the product, operating instructions or packaging. The materials used in this unit are recyclable according to their marking. You will be making an important contribution to protecting our environment by reusing, recycling or otherwise disposing of used units.

Your local administration will provide you with information about the appropriate disposal point for used units.

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