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PH-MESSGERÄT

USER MANUAL

PH METER

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MIERNIK PH

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MĚŘIČ PH

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PH-MÈTRE

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MANUAL DE INSTRUCCIONES

MEDIDOR DE PH

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PH MÉRŐ

BRUGSANVISNING

PH-MÅLER

DE | EN | PL | CZ | FR | IT | ES | HU | DA

SBS-PH-140

DE	Produktname:	PH-MESSGERÄT
EN	Product name:	PH METER
PL	Nazwa produktu:	MIERNIK PH
CZ	Název výrobku	MĚŘIČ PH
FR	Nom du produit:	PH-MÈTRE
IT	Nome del prodotto:	MISURATORE DI PH
ES	Nombre del producto:	MEDIDOR DE PH
HU	Termék neve	PH MÉRŐ
DA	Produktnavn	PH-MÅLER
DE	Modell:	SBS-PH-140
EN	Product model:	
PL	Model produktu:	
CZ	Model výrobku	
FR	Modèle:	
IT	Modello:	
ES	Modelo:	
HU	Modell	
DA	Model	
DE	Hersteller	expondo Polska sp. z o.o. sp. k.
EN	Manufacturer	
PL	Producent	
CZ	Výrobce	
FR	Fabricant	
IT	Produttore	
ES	Fabricante	
HU	Termelő	
DA	Producent	
DE	Anschrift des Herstellers	ul. Nowy Kisielin – Innowacyjna 7, 66-002 Zielona Góra Poland, EU
EN	Manufacturer Address	
PL	Adres producenta	
CZ	Adresa výrobce	
FR	Adresse du fabricant	
IT	Indirizzo del produttore	
ES	Dirección del fabricante	
HU	A gyártó címe	
DA	Producentens adresse	

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This User Manual has been translated for your convenience using machine translation. Reasonable efforts have been made to provide an accurate translation; however, no automated translation is perfect nor is it intended to replace human translators. The official User Manual is the English version. Any discrepancies or differences created in the translation are not binding and have no legal effect for compliance or enforcement purposes. If any questions arise related to the accuracy of the information contained in the User Manual, please refer to the English version of those contents which is the official version.

Technical data

Parameter description	Parameter value
Product name	pH METER
Model	SBS-PH-140
Power supply input parameters	AC 230-240V ~ / 50 Hz; 0.1A
Power supply output parameters	DC 6V; 250mA
Battery capacity [mAh]	110
Time of fully charging the power battery [h]	2-3
pH range [pH]	0-14
Temperature range [° C]	-50 - ~70
EC range [EC]	0 - ~19.99
CF range [CF]	0 - ~199.9
TDS range [ppm]	10 ~19990
Resolution	pH: 0.01 temperature: 0.1 EC: 0.1 CF: 0.1 TDS: 10
Accuracy of measurement display	pH: ± 0.1 temperature: ± 1 EC / CF / TDS: $\pm 2\%$ F. S
Dimensions (Width x Depth x Height) [mm]	152 x 79 x 26.4
Weight [g]	295


1. General overview

This manual is intended to assist you in safe and reliable use. The product is designed and manufactured strictly according to technical specifications using the latest technology and components and maintaining the highest quality standards.

**CAREFULLY READ AND UNDERSTAND THIS MANUAL
BEFORE STARTING THE WORK.**

To ensure long and reliable operation of the device, make sure to operate and maintain it properly in accordance with the guidelines in this instruction manual. The technical data and specifications in this manual are up-to-date. The manufacturer reserves the right to modifications for the purpose of quality improvement.

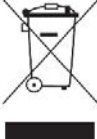Explanation of symbols

	The product complies with applicable safety standards.
	Read the manual before use.
	Recyclable product.
	CAUTION! or WARNING! or REMEMBER! General warning sign.
	CAUTION! Risk of electric shock!

CAUTION! The figures in this manual are illustrative only and may vary in some details from the actual appearance of the product.

2. Safety of use

CAUTION! Read all safety warnings and all instructions. Failure to follow the warnings and instructions may result in electric shock, fire and/or serious injury or death.

The term "device" or "product" in the warnings and in the description of the manual refers to the Ph Meter.

2.1. Electrical safety

- a) The charger plug on the device must fit into the outlet. Do not modify the plug in any way. Original plugs and matching outlets reduce the risk of electric shock.
- b) Do not use the charger cord improperly. Never use it to carry the device or to pull the plug out of the socket. Keep the cord away from heat sources, oil, sharp edges or moving parts. Damaged or tangled cords increase the risk of electric shock.
- c) Do not use the device if the charger cable is damaged or shows signs of wear. A damaged cord should be replaced by a qualified electrician or the manufacturer's service
- d) To avoid electric shock, do not immerse the charger cord, plug or the unit itself in water or any other liquid. Do not use the appliance on wet surfaces.
- e) CAUTION – DANGER TO LIFE! When cleaning or using the appliance, never immerse it in water or other liquids.
- f) Do not get the entire device wet. Risk of electric shock!

2.2. Safety in the workplace

- a) Do not touch the device with wet or damp hands.
- a) Keep the work area tidy and well lit. Disorder or poor lighting can lead to accidents. Be foresighted, watch what you are doing and use common sense when using the unit.
- b) Do not use the unit in an explosive area, for example in the presence of flammable liquids, gases or dust. The unit produces sparks that can ignite dust or fumes.
- c) If you have any doubts as to whether the product is working properly or if it is damaged, contact the manufacturer's service department.
- d) Repairs to the device may only be carried out by the manufacturer's service. Do not attempt to repair the product on your own!
- e) In case of open flames or fire, use only dry powder or snow (CO₂) fire extinguishers to extinguish the live equipment.
- f) Check the condition of the safety stickers regularly. Replace them if they are illegible.
- g) Keep these instructions for use for future reference. If the product is to be handed over to a third party, hand it over with this user manual.
- h) Keep packaging components and small installation parts out of the reach of children.
- i) Keep the device away from children and animals.

Remember! Keep children and other bystanders safe while operating the equipment.

2.3. Personal safety

- a) Do not operate this device if you are tired, ill or under the influence of alcohol, drugs or medication that could impair your ability to operate the device.
- b) The device is not intended to be used by persons (including children) with reduced mental, sensory or intellectual functions or persons who lack experience and/or knowledge unless they are supervised or have been instructed by a person responsible for their safety on how to operate the device.
- c) Do not overestimate your capabilities. Maintain body balance and equilibrium at all times during operation. This allows for better control of the machine in unexpected situations.
- d) The product is not a toy. Children should be watched to ensure that they do not play with the appliance.

2.4. Safe use of the device

- g) Do not use the device if the ON/OFF switch does not function properly (does not turn on and off). Units that cannot be controlled by the switch are unsafe, cannot operate, and must be repaired.
- h) Disconnect the device from the power supply before adjusting, cleaning, or servicing. This precaution reduces the risk of accidental start-up.
- i) Keep unused product out of the reach of children and anyone unfamiliar with the device or this manual. Products are dangerous when used by inexperienced users.
- j) Keep the product in good working order. Check before each use for general damage or damage to moving parts (cracks in parts and components or any other condition that may affect the safe operation of the device). If damaged, have the device repaired before use.
- k) Keep the product out of the reach of children.
- l) Repairs and maintenance should be carried out by qualified personnel using only original spare parts. This will ensure the safety of use.
- m) To ensure the designed operational integrity of the device, do not remove factory-installed covers or loosen screws.
- n) Clean the device regularly to prevent permanent dirt build-up.
- o) The product is not a toy. Cleaning and maintenance must not be performed by children without adult supervision.
- p) Do not tamper with the device to alter its performance or design.
- q) Keep the unit away from sources of fire and heat.


CAUTION! Although the product has been designed to be safe, with adequate safeguards, and despite the additional safety features provided to the user, there is still a slight risk of accident or injury when handling the product. Caution and common sense are advised when using the product.

3. Instructions for use

The device is designed to carry out measurements for solutions regarding: temperature, pH, EC/CF/TDS.

The user is responsible for any damage resulting from misuse.

3.1. Product overview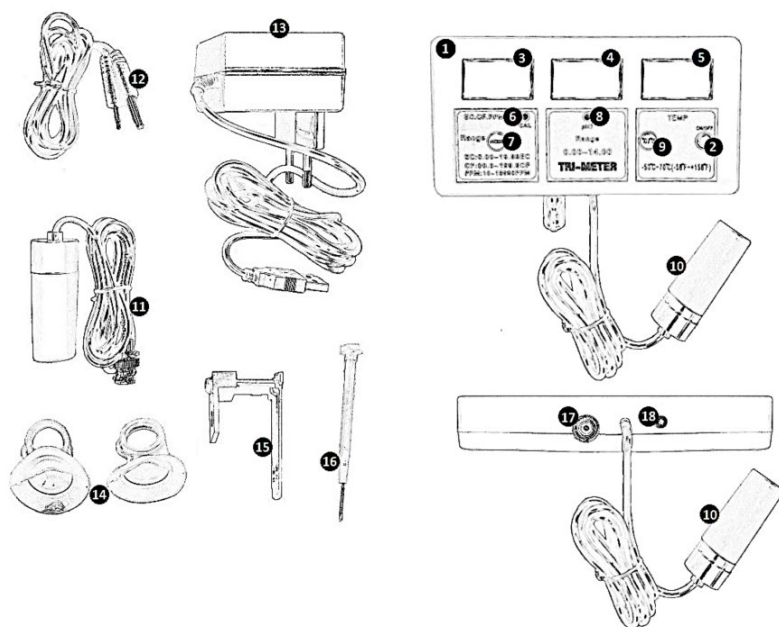

- 1- Control panel
- 2 - "ON / OFF" switch (on/off)
- 3 - LCD display for conductivity measurements
- 4 - LCD display for pH measurements

- 9 - °C/°F unit change button
- 10 - Conductivity measurement probe
- 11 - pH measurement probe
- 12 - Temperature measurement probe
- 13 - Power supply
- 14 - Electrode suction cup x2

5 - LCD display for temperature measurements	15 - Mounting bracket
6 - Conductivity calibration timer	16 - Screwdriver
7 - MODE button	17 - Inlet for connecting the pH probe
8 - pH calibration trimmer	18 - Inlet for connecting the temperature probe


3.2. Preparation for operation

POSITIONING OF THE UNIT

Place the unit in a way that ensures good air circulation. Keep the unit away from any hot surfaces. Always operate the unit on a level, stable, clean, fireproof and dry surface and out of the reach of children and persons of impaired mental, sensory and intellectual functions. Place the unit in such a way that the mains plug can be reached at any time. Ensure that the power supply to the unit corresponds to that specified on the identification plate!

ASSEMBLY OF THE UNIT

The device is designed to be installed on a working surface. If there is a need to mount the device in a vertical position when working with it, install the mounting bracket (15) or mount the work surface directly into the dedicated mounting indentations of the rear housing of the unit. When installing the mounting bracket (15), attach the longer arm to the dedicated hooks on the rear housing of the unit, and attach the shorter arm to the operating surface. Mounting the device onto the work surface should be performed in accordance with the standards and regulations prevailing in a given country.

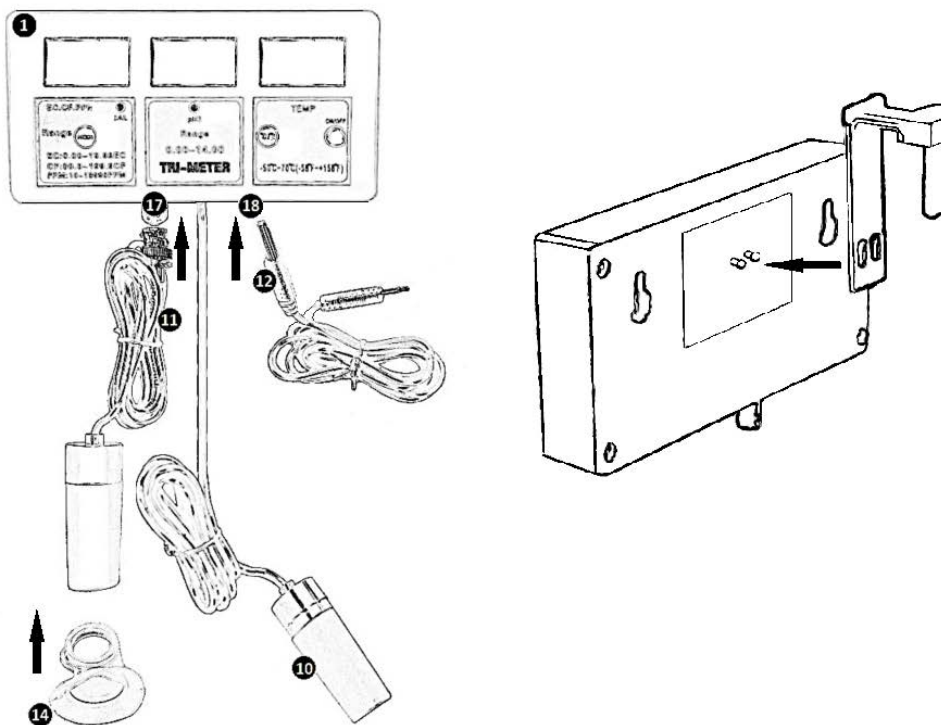3.3. Working with the unit

When the probe tip is immersed in the test solution, air bubbles can settle on its surface, leading to errors in readings. Therefore, when measuring and calibrating the device, it is recommended to slowly move the probe in a circular motion to minimize the occurrence of this issue.

If too large measurement deviations occur with a properly performed calibration, the measurement results do not reset, etc., clean the electrode with ethyl alcohol.

CONNECTING THE PROBES / CHARGING THE BATTERY

- Each time before working with the unit, fully charge the unit's battery in accordance with the guidelines in the technical data table. The inlet for connecting to the power supply is located in the side housing of the device.
NOTE: If during operation of the device the backlight of the LCD display darkens or the device displays measurements with too large deviations, recharge the battery of the device.
- The device has three dedicated measurement probes. The conductivity measurement probe for EC/CF/TDS is integrated into the device. Before each use of the device, check the correct installation of the measuring probes.

-
- c) Insert the pH probe connection output (11) into the pH probe connection input (17). Lock the connection by turning the probe connection output clockwise. The disassembly is done in reverse order to its assembly.
 - d) Connect the temperature probe connection output (12) to the temperature probe connection input (18).
 - e) In order to facilitate working with the device, the conductivity and pH probes can be installed in dedicated suction cups and attached to the operating surface.

CHANGING THE TEMPERATURE UNIT

To adjust the temperature unit (°C/°F), select the unit change button (9). The LCD will display the current temperature unit configuration.

PERFORMING MEASUREMENTS

- a) Remove the probe covers.
- b) Before and after each measurement, clean the probes with distilled water. Dip it in the solution and then dry with a cotton cloth.
- c) Turn on the device with the "ON/OFF" switch (2). Pressing the switch again will turn off the device.
- d) Immerse the tips of the probes in the test solution.
- e) Wait until the measurement is stabilized on the LCD displays (3, 4, 5), and then read the correct values.
NOTE: The conductivity values are displayed on the LCD in a specific order EC/CF/TDS. Switching between the measurement results may be done by pressing the [MODE] key.
- f) After taking the measurements, turn off the device using the power switch.
- g) Clean the probe.
- h) Protect the probe with a cover.

DEVICE CALIBRATION

The calibration process should be performed after the following occur:

- a) The first use of the device.
- b) Device was inactive for a long time.
- c) Measurement under specific conditions (strong solution).
- d) Replacing the probe with a new one

PH CALIBRATION PROCEDURE

- a) Prepare calibration solutions with pH: 6.86; 4.0; 9.18 and distilled water.
- b) Before and after each measurement, clean the probe with distilled water. Immerse the probe tip in distilled water, remove it and dry with a cotton cloth.
- c) Turn on the device with the power switch (2).
- d) Immerse the probe tip in a pH 6.86 solution at 25°C.
- e) Wait until the measurement stabilizes on the LCD display.

-
- f) Adjust the measurement reading with the pH calibration trimmer (8) using a screwdriver (16). The value displayed on the LCD screen should be equivalent to the tested calibration solution.
 - g) Clean the probe.
 - h) Check the correctness of the performed calibration. To do this, immerse the probe tip in a solution with a pH of 9.18 or 4.0. Wait until the measurement stabilizes on the LCD display. If the measurement is within the permissible error tolerance, the calibration should be considered correct. Otherwise, the calibration steps must be repeated.

CONDUCTIVITY CALIBRATION PROCEDURE

- a) Prepare calibration solutions with EC: 12.88 and 1.41 and distilled water.
- b) Before and after each measurement, clean the probe with distilled water. Immerse the probe tip in distilled water, remove it and dry with a cotton cloth.
- c) Turn on the device with the power switch (2).
- d) Immerse the probe tip in the dedicated 12.88 EC calibration solution at 25°C.
- e) Wait until the measurement stabilizes on the LCD display (3).
NOTE: When taking the measurement, remember to set the display properly to read the EC value.
- i) Adjust the measurement reading with the calibration trimmer (6) using a screwdriver (16). The value displayed on the LCD screen should be equivalent to the tested calibration solution.
- f) Clean the probe.
- g) Check the correctness of the performed calibration. To do this, dip the probe tip into the 1.41EC solution. Wait until the measurement stabilizes on the LCD display. If the measurement is within the permissible error tolerance, the calibration should be considered correct. Otherwise, the calibration steps must be repeated.

3.4. Cleaning and maintenance

- a) Store the appliance in a dry and cool place protected from moisture and direct sunlight.
- b) Perform regular inspections of the unit checking technical fitness and any damages.
- c) Do not leave the batteries in the unit when it will not be used for a long time.
- d) Avoid touching the probe directly with your hand, this can damage it!
- e) Always put the probe cover back on when the device will not be used for a long time.

DISPOSAL OF WASTE APPLIANCES

At the end of its useful life, this product should not be disposed of with normal household waste but should be taken to a collection point for the recycling of electrical and electronic equipment. This is indicated by the symbol on the

product, operating instructions or packaging. The materials used in this appliance are recyclable according to their marking. By reuse, recycling or applying other forms of use of waste machines, you make a significant contribution to the protection of our environment.

Your local administration will provide you with information about the appropriate disposal point for used appliances.

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