## USER MANUAL

Bedienungsanleitung  
Instrukcja obsługi  
Návod k použití  
Manuel d'utilisation  
Istruzioni per l'uso  
Manual de instrucciones  
Használati útmutató  
Brugsanvisning  
Käyttöohje  
Gebruiksaanwijzing  
Bruksanvisning  
Instruções de utilização  
Používateľská príručka  
Ръководство за употреба  
Οδηγίες χρήσης  
Upute za uporabu  
Naudojimo instrukcija  
Manual de utilizare  
Navodila za uporabo

---

---

# LAMINAR FLOW CABINET (VERTICAL TYPE)

<b>DE</b>	Produktname	LAMINAR-STRÖMUNG-SCHRANK (VERTIKALER TYP)
<b>EN</b>	Product name	LAMINAR FLOW CABINET (VERTICAL TYPE)
<b>PL</b>	Nazwa produktu	KOMORA LAMINARNA (TYP PIONOWY)
<b>CZ</b>	Název výrobku	LAMINÁRNÍ PROUDOVÁ SKŘÍŇ (VERTIKÁLNÍ TYP)
<b>FR</b>	Nom du produit	HOTTE A FLUX LAMINAIRE (TYPE VERTICAL)
<b>IT</b>	Nome del prodotto	CABINA A FLUSSO LAMINARE (TIPO VERTICALE)
<b>ES</b>	Nombre del producto	CABINA DE FLUJO LAMINAR (TIPO VERTICAL)
<b>HU</b>	Termék neve	LAMINÁRIS ÁRAMLÁSÚ FŰLKE (FÜGGŐLEGES TÍPUS)
<b>DA</b>	Produktnavn	LAMINAR FLOW KABINET (LODRET TYPE)
<b>FI</b>	Tuotteen nimi	LAMINAARIVIRTAUSKAAPPI (PYSTYTYYPPI)
<b>NL</b>	Productnaam	LAMINAIRE STROMINGSKAST (VERTICAAL TYPE)
<b>NO</b>	Produktnavn	LAMINARSTRØMSKABINETT (VERTIKAL TYPE)
<b>SE</b>	Produktnamn	LAMINÄRT FLÖDESKÅP (VERTIKAL TYP)
<b>PT</b>	Nome do produto	GABINETE DE FLUXO LAMINAR (TIPO VERTICAL)
<b>SK</b>	Názov produktu	LAMINÁRNY PRIETOKOVÝ KABINET (VERTIKÁLNY TYP)
<b>BG</b>	Име на продукта	ШКАФ ЛАМИНАРНОГО ПОТОКА (ВЕРТИКАЛЬНЫЙ ТИП)
<b>EL</b>	Όνομα προϊόντος	ΝΤΟΥΛΑΠΙ ΡΟΗΣ ΑΕΡΑ (ΚΑΤΑΚΟΡΥΦΟΣ ΤΥΠΟΣ)
<b>HR</b>	Naziv proizvoda	ORMARIĆ S LAMINARNIM PROTOKOM (VERTIKALNI TIP)
<b>LT</b>	Produkto pavadinimas	LAMINARIUMO SRAUTO SPINTA (VERTIKALUS TIPAS)
<b>RO</b>	Numele produsului	DULAP DE FLUX LAMINAR (TIP VERTICAL)
<b>SL</b>	Ime izdelka	LAMINARNA PRETOČNA OMARICA (VERTIKALNI TIP)
<b>DE</b> Modell   <b>EN</b> Product model   <b>PL</b> Model produktu   <b>CZ</b> Model výrobku   <b>FR</b> Modèle   <b>IT</b> Modello   <b>ES</b> Modelo   <b>HU</b> Modell   <b>DA</b> Model   <b>FI</b> Tuotteen malli   <b>NL</b> Productmodel   <b>NO</b> Produktmodell   <b>SE</b> Produktmodell   <b>PT</b> Modelo do produto   <b>SK</b> Model   <b>BG</b> Модел на продукт   <b>EL</b> Μοντέλο προϊόντος   <b>HR</b> Model proizvoda   <b>LT</b> : Gaminio modelis   <b>RO</b> : Model de produs   <b>SL</b> : Model izdelka		<b>SBS-LFC-100</b> <b>SBS-LFC-200</b>
<b>DE</b> Hersteller   <b>EN</b> Manufacturer   <b>PL</b> Producent   <b>CZ</b> Výrobce   <b>FR</b> Fabricant   <b>IT</b> Produttore   <b>ES</b> Fabricante   <b>HU</b> Termelő   <b>DA</b> Producent   <b>FI</b> Valmistaja   <b>NL</b> Producent   <b>NO</b> Produsent   <b>SE</b> Tillverkare   <b>PT</b> Fabricante   <b>SK</b> Výrobca   <b>BG</b> Производител   <b>EL</b> Κατασκευαστής   <b>HR</b> Proizvođač   <b>LT</b> Gamintojas   <b>RO</b> Producător   <b>SL</b> Proizvajalec		expondo Polska sp. z o.o. sp. k.
<b>DE</b> Anschrift des Herstellers   <b>EN</b> Manufacturer Address   <b>PL</b> Adres producenta   <b>CZ</b> Adresa výrobce   <b>FR</b> Adresse du fabricant   <b>IT</b> Indirizzo del produttore   <b>ES</b> Dirección del fabricante   <b>HU</b> A gyártó címe   <b>DA</b> Producentens adresse   <b>FI</b> Valmistajan osoite   <b>NL</b> Adres producent   <b>NO</b> Produsentens adresse   <b>SE</b> Tillverkarens adress   <b>PT</b> Endereço do fabricante   <b>SK</b> Adresa výrobcu   <b>BG</b> Адрес на производителя   <b>EL</b> : Διεύθυνση κατασκευαστή   <b>HR</b> Adresa proizvođača   <b>LT</b> Gamintojo adresas   <b>RO</b> Adresa producătorului   <b>SL</b> Naslov proizvajalca		ul. Nowy Kisielin – Innowacyjna 7, 66-002 Zielona Góra   Poland, EU

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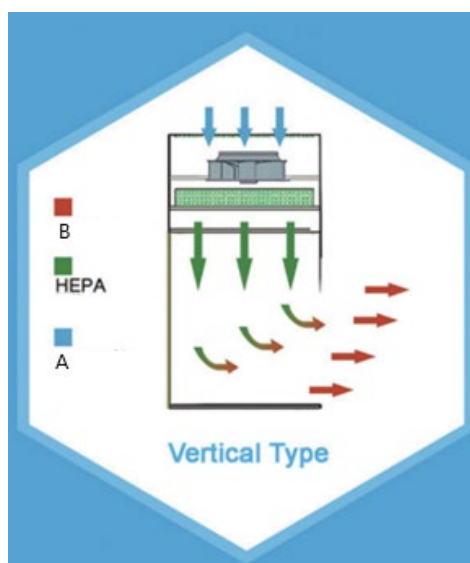
This User Manual has been translated using machine translation. We have made every effort to ensure the translation is accurate, but please note that automated translations are not perfect and are not meant to replace human translators. The official version of the User Manual is in English. Any differences between the translated version and the original English are not legally binding. If you have any questions about the accuracy of the translation, please refer to the English version, which is the official reference. More language versions are available upon request via [info@expondo.com](mailto:info@expondo.com)

## Technical data


Parameter description	Parameter value	
Product name	Laminar flow cabinet (vertical type)	
Model	SBS-LFC-100	SBS-LFC-200
Rated voltage [V~] / frequency [Hz]	230 / 50	
Rated power [W]	240	440
Max power of socket	800	800
UV lamp power [W]	20	30
LED lamp power [W]	16 (x2)	24 (x2)
Protection class	I	
Dimensions [width x depth x height; mm]	1060 x 795 x 1830	1460 x 792 x 1850
Weight [kg]	142.5	171
Inside chamber dimensions [mm]	920 x 690 x 650	1320 x 690 x 650
Duty cycle	S1	
Filter type	HEPA 99.999% efficiency at 0.3 $\mu\text{m}$	
Emitted noise [dB(A)]	60	
Max load capacity [kg]	50	
Air velocity measurement [m/s]	0.2~0.6	
Remote controller battery type	CR2025	

This appliance is a type of workstation designed to create a clean, particle-free working environment by directing filtered air vertically over the workspace. It operates by drawing air from the room, which is then filtered through a High-Efficiency Particulate Air (HEPA) filter before being directed downwards in a laminar (unidirectional) flow. It is crucial in environments requiring high standards of cleanliness, such as in life sciences, certain manufacturing processes etc.

**This product is outside the scope of REGULATION (EU) 2017/745 for medical devices. It has not been designed as a medical device.**

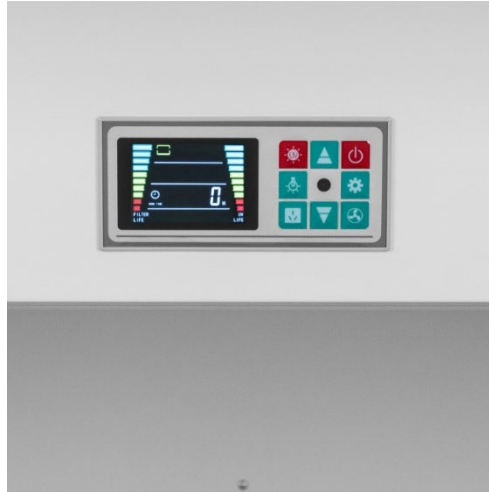- A. Room air
- B. Cleaned air

## Installation

- Do not tilt the appliance over 45° while transportation or operation.
- The appliance should be installed in the clean laboratory or factory building, away from dust and quiver source. Unpack it carefully and after setting it, please revolve the leg to make it exert itself, and then position the worktable horizontal.
- The earth terminal of the casing must be connected to a reliable grounding (see the universal ground symbol on the housing).
- Connect the power cord to a grounded socket.
- The air flow speed of the worktable should be adjusted to the best state while dispatching from the factory. It should be adjusted to the low-grade location in the initial stage, and generally improve in half a year.

## Device use

- If the worktable is newly installed or has not been used for a long time, clean it and the surrounding environment with dust catcher or the tool that will not produce fiber before using, and then use medicine sterilization or the ultraviolet sterilization method to deal with sterilization.
- Start the appliance 5 minutes in advance and turn on the UV-light.
- Shut off the sterilisation light after 30 minutes, and then turn on the air blower.
- The operation can be started 20 minutes later, turn on the fluorescent lamp while operating.

## Maintenance

- Unplug the mains plug and allow the device to cool completely before each cleaning, adjustment or replacement of accessories, or if the device is not being used.
  - Wait for the rotating elements to stop.
- Store the unit in a dry, cool place, free from moisture and direct exposure to sunlight.
- Do not spray the device with a water jet or submerge it in water.
- Do not allow water to get inside the device through vents in the housing of the device.
- Clean the vents with a brush and compressed air.
- Use a soft cloth for cleaning.
- Do not use sharp and/or metal objects for cleaning (e.g. a wire brush or a metal spatula) because they may damage the surface material of the appliance.
- HEPA-filter lifetime:
 

When the fan had been running for 2500 hours, the indicator light and gear of the fan will flash when the fan is running, indicating that the user needs to change the filter. While changing the high-efficient filter, open the top cover, pay attention to the arrow sign on the filter, the direction of arrow indicates the laminar flow. Check carefully while changing the filter, make sure the frame is sealed well all around; otherwise, the filter result will be influenced. After changing the filter, press the "Fan" button for 5 seconds until the four lights in the middle of the indicator bar are lit at the same time. At this time, the "Fan" button can be released to restore the normal working state.
- UV-sterilising lamp lifetime:

When the UV-sterilising lamp had been on for 1000 hours, the sterilisation indicator will flash when the this lamp is on, indicating that the user needs to replace it. After finishing changing the UV lamp, press the "sterilization" button for 5 seconds until the four lights in the middle of the indicator bar are lit at the same time. At this time, the "sterilisation" button can be released to restore the normal working state.

- j) According to the pollution degree of the environment, disassemble the filter cloth for wash or change (usually every 2-3 months).
- k) Sterilise the environment periodically (usual once a week), Use the gauze with alcohol for sterilisation. At the same time, keep the surface clean. Otherwise it will influence sterilisation result.

## Troubleshooting

Issue	Possible cause	Solution
No power supply.	1. The plug and socket are poorly connected. 2. The fuse is burnt.	1. Adjust the plug and socket, or check the circuit. 2. Change the fuse.
The air fan does not work.	The control does not work or the air fan is broken down.	Change the control panel or check the air fan.
The light does not work	The light socket is broken down. The light tube is burnt.	Adjust the lamp socket or change the light tube.
Cannot reach required air flow speed	The high-efficient filter wears out.	Change the filter.

## Disposing of used devices

Do not dispose of this device in municipal waste systems. Hand it over to an electric and electrical device recycling and collection point. Check the symbol on the product, instruction manual and packaging. The plastics used to construct the device can be recycled in accordance with their markings. By choosing to recycle you are making a significant contribution to the protection of our environment.

Contact local authorities for information on your local recycling facility.

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