USER MANUAL

Bedienungsanleitung
Instrukcja obsługi
Návod k použití
Manuel d'utilisation
Istruzioni per l'uso
Manual de instrucciones
Használati útmutató
Brugsanvisning
Käyttöohje
Gebruiksaanwijzing
Bruksanvisning
Instruções de utilização
Používateľská príručka
Ръководство за употреба
Οδηγίες χρήσης
Upute za uporabu
Naudojimo instrukcija
Manual de utilizare
Navodila za uporabo

MINI LAMINAR FLOW CABINET

DE	Produktname	MINI-LAMINAR-FLOW-SCHRANK
EN	Product name	MINI LAMINAR FLOW CABINET
PL	Nazwa produktu	MINI KOMORA LAMINARNA
CZ	Název výrobku	MINI LAMINÁRNÍ PROUDOVÁ SKŘÍŇ
FR	Nom du produit	MINI HOTTE A FLUX LAMINAIRE
IT	Nome del prodotto	MINI CABINA A FLUSSO LAMINARE
ES	Nombre del producto	MINI CABINA DE FLUJO LAMINAR
HU	Termék neve	MINI LAMINÁRIS ÁRAMLÁSÚ FÜLKE
DA	Produktnavn	MINI LAMINAR FLOW KABINET
FI	Tuotteen nimi	MINI LAMINAARIVIRTAUSKAAPPI
NL	Productnaam	MINI LAMINAIRE STROMINGSKAST
NO	Produktnavn	MINI LAMINARSTRØMSKABINETT
SE	Produktnamn	MINI LAMINÄRT FLÖDESKÅP
PT	Nome do produto	MINI GABINETE DE FLUXO LAMINAR
SK	Názov produktu	MINI LAMINÁRNY PRIETOKOVÝ KABINET
BG	Име на продукта	МИНИ ШКАФ ЛАМИНАРНОГО ПОТОКА
EL	Όνομα προϊόντος	ΜΙΝΙ ΝΤΟΥΛΑΠΙ ΡΟΗΣ ΑΕΡΑ
HR	Naziv proizvoda	MINI ORMARIĆ S LAMINARNIM PROTOKOM
LT	Produkto pavadinimas	MINI LAMINARIUMO SRAUTO SPINTA
RO	Numele produsului	MINI DULAP DE FLUX LAMINAR
SL	Ime izdelka	MINI LAMINARNA PRETOČNA OMARICA
DE Modell EN Product model PL Model produktu CZ Model výrobku FR Modèle IT Modello ES Modelo HU Modell DA Model FI Tuotteen malli NL Productmodel NO Produktmodell SE Produktmodell PT Modelo do produto SK Model BG Модел на продукт EL Μοντέλο προϊόντος HR Model proizvoda LT : Gaminio modelis RO : Model de produs SL : Model izdelka		SBS-LFC-10
DE Hersteller EN Manufacturer PL Producent CZ Výrobce FR Fabricant IT Produttore ES Fabricante HU Termelő DA Producent FI Valmistaja NL Producent NO Produsent SE Tillverkare PT Fabricante SK Výrobca BG Производител EL Κατασκευαστής HR Proizvođač LT Gamintojas RO Producător SL Proizvajalec		expondo Polska sp. z o.o. sp. k.
DE Anschrift des Herstellers EN Manufacturer Address PL Adres producenta CZ Adresa výrobce FR Adresse du fabricant IT Indirizzo del produttore ES Dirección del fabricante HU A gyártó címe DA Producentens adresse FI Valmistajan osoite NL Adres producent NO Produsentens adresse SE Tillverkarens adress PT Endereço do fabricante SK Adresa výrobcu BG Адрес на производителя EL : Διεύθυνση κατασκευαστή HR Adresa proizvođača LT Gamintojo adresas RO Adresa producătorului SL Naslov proizvajalca		ul. Nowy Kisielin – Innowacyjna 7, 66-002 Zielona Góra Poland, EU

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This User Manual has been translated using machine translation. We have made every effort to ensure the translation is accurate, but please note that automated translations are not perfect and are not meant to replace human translators. The official version of the User Manual is in English. Any differences between the translated version and the original English are not legally binding. If you have any questions about the accuracy of the translation, please refer to the English version, which is the official reference. More language versions are available upon request via info@expondo.com

Technical data

Parameter description	Parameter value
Product name	Mini laminar flow cabinet
Model	SBS-LFC-10
Rated voltage [V~] / frequency [Hz]	230 / 50
Rated power [W]	110
UV lamp power [W]	8
LED lamp power [W]	8
Protection class	I
Dimensions [width x depth x height; mm]	550 x 460 x 713
Weight [kg]	35.65
Inside chamber dimensions [mm]	480 x 340 x 370
Duty cycle	S1
Filter type	HEPA 99.999% efficiency at 0.3 µm
Emitted noise [dB(A)]	60
Max load capacity [kg]	10
Air velocity measurement [m/s]	0.2~0.6

This appliance is a type of workstation designed to create a clean, particle-free working environment by directing filtered air vertically over the workspace. It operates by drawing air from the room, which is then filtered through a High-Efficiency Particulate Air (HEPA) filter before being directed downwards in a laminar (unidirectional) flow. It is crucial in environments requiring high standards of cleanliness, such as in life sciences, certain manufacturing processes etc.

This product is outside the scope of REGULATION (EU) 2017/745 for medical devices. It has not been designed as a medical device.

Installation

- Do not tilt the appliance over 45° while transportation or operation.
- The appliance should be installed in the clean laboratory or factory building, away from dust and quiver source. Unpack it carefully and after setting it, please revolve the leg to make it exert itself, and then position the worktable horizontal.

- The earth terminal of the casing must be connected to a reliable grounding (see the universal ground symbol on the housing).
- Connect the power cord to a grounded socket.
- The air flow speed of the worktable should be adjusted to the best state while dispatching from the factory. It should be adjusted to the low-grade location in the initial stage, and generally improve in half a year.

Device use

- If the worktable is newly installed or has not been used for a long time, clean it and the surrounding environment with dust catcher or the tool that will not produce fiber before using, and then use medicine sterilization or the ultraviolet sterilisation method to deal with sterilization.
- Start the appliance 5 minutes in advance and turn on the UV-light.
- Shut off the sterilisation light after 30 minutes, and then turn on the air blower.
- The operation can be started 20 minutes later, turn on the fluorescent lamp while operating.

Maintenance

- Unplug the mains plug and allow the device to cool completely before each cleaning, adjustment or replacement of accessories, or if the device is not being used.
 - Wait for the rotating elements to stop.
- Store the unit in a dry, cool place, free from moisture and direct exposure to sunlight.
- Do not spray the device with a water jet or submerge it in water.
- Do not allow water to get inside the device through vents in the housing of the device.
- Clean the vents with a brush and compressed air.
- Use a soft cloth for cleaning.
- Do not use sharp and/or metal objects for cleaning (e.g. a wire brush or a metal spatula) because they may damage the surface material of the appliance.
- HEPA-filter lifetime:
When the fan had been running less efficiently and not reaching the air flow speed it is indicating that the user needs to change the filter. While changing the high-efficient filter, open the top cover, pay attention to the arrow sign on the filter, the direction of arrow indicates the laminar flow. Check carefully while changing the filter, make sure the frame is sealed well all around; otherwise, the filter result will be influenced.
- UV-sterilising lamp lifetime:
When the UV-sterilising lamp had been on for about 1000 hours it needs to be replaced.
- Sterilise the environment periodically (usual once a week), Use the gauze with alcohol for sterilisation. At the same time, keep the surface clean. Otherwise it will influence sterilisation result.

Troubleshooting

Issue	Possible cause	Solution
No power supply.	1. The plug and socket are poorly connected. 2. The fuse is burnt.	1. Adjust the plug and socket, or check the circuit. 2. Change the fuse.
The air fan does not work.	The control does not work or the air fan is broken down.	Change the control panel or check the air fan.
The light does not work	The light socket is broken down. The light tube is burnt.	Adjust the lamp socket or change the light tube.
Cannot reach required air flow speed	The high-efficient filter wears out.	Change the filter.

Disposing of used devices

Do not dispose of this device in municipal waste systems. Hand it over to an electric and electrical device recycling and collection point. Check the symbol on the product, instruction manual and packaging. The plastics used to

construct the device can be recycled in accordance with their markings. By choosing to recycle you are making a significant contribution to the protection of our environment.
Contact local authorities for information on your local recycling facility.

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