USER MANUAL

Bedienungsanleitung
Instrukcja obsługi
Návod k použití
Manuel d'utilisation
Istruzioni per l'uso
Manual de instrucciones
Használati útmutató
Brugsanvisning
Käyttöohje
Gebruiksaanwijzing
Bruksanvisning
Instruções de utilização
Používateľská príručka

THERMOSTATIC WATER BATH

DE	Produktname	THERMOSTATISCHES WASSERBAD
EN	Product name	THERMOSTATIC WATER BATH
PL	Nazwa produktu	LABORATORYJNA ŁAŹNIA WODNA
CZ	Název výrobku	TERMOSTATICKÁ VODNÍ LÁZEŇ
FR	Nom du produit	BAINS THERMOSTATÉS
IT	Nome del prodotto	BAGNO TERMOSTATICO
ES	Nombre del producto	BAÑO TERMOSTÁTICO
HU	Termék neve	TERMO SZTATIKUS VÍZFÜRDŐ
DA	Produktnavn	TERMOSTATISK VANDBAD
FI	Tuotteen nimi	TERMOSTAATTINEN VESIHAUDE
NL	Productnaam	THERMOSTATISCH WATERBAD
NO	Produktnavn	TERMOSTATISK VANNBAD
SE	Produktnamn	TERMOSTATERAT VATTENBAD
PT	Nome do produto	BANHO-MARIA DE LABORATÓRIO
SK	Názov produktu	TERMOSTATICKÝ VODNÝ KÚPEĽ
DE	Modell	SBS-TWB-400 SBS-TWB-600
EN	Product model	
PL	Model produktu	
CZ	Model výrobku	
FR	Modèle	
IT	Modello	
ES	Modelo	
HU	Modell	
DA	Model	
FI	Tuotteen malli	
NL	Productmodel	
NO	Produktmodell	
SE	Produktmodell	
PT	Modelo do produto	
SK	Model	
DE	Hersteller	expondo Polska sp. z o.o. sp. k.
EN	Manufacturer	
PL	Producent	
CZ	Výrobce	
FR	Fabricant	
IT	Produttore	
ES	Fabricante	
HU	Termelő	
DA	Producent	
FI	Valmistaja	
NL	Producent	
NO	Produsent	
SE	Tillverkare	
PT	Produtor	
SK	Výrobca	
DE	Anschrift des Herstellers	ul. Nowy Kisielin – Innowacyjna 7, 66-002 Zielona Góra Poland, EU
EN	Manufacturer Address	
PL	Adres producenta	
CZ	Adresa výrobce	
FR	Adresse du fabricant	
IT	Indirizzo del produttore	
ES	Dirección del fabricante	
HU	A gyártó címe	
DA	Producentens adresse	
FI	Valmistajan osoite	
NL	Adres producent	
NO	Produsentens adresse	
SE	Tillverkarens adress	
PT	Endereço do produtor	
SK	Adresa výrobcu	

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This User Manual has been translated for your convenience using machine translation. Reasonable efforts have been made to provide an accurate translation; however, no automated translation is perfect nor is it intended to replace human translators. The official User Manual is the English version. Any discrepancies or differences created in the translation are not binding and have no legal effect for compliance or enforcement purposes. If any questions arise related to the accuracy of the information contained in the User Manual, please refer to the English version of those contents which is the official version.

Technical data

Parameter description	Parameter value	
Product name	Thermostatic water bath	
Model	SBS-TWB-400	SBS-TWB-600
Supply voltage [V~] / frequency [Hz]	230/50	
Rated power [W]	1000	2200
Capacity [l]	11	36
Temperature range [°C]	5-100	
Dimensions (width x depth x height) [mm]	590 x 230 x 290	820 x 350 x 335
Weight [kg]	9	16.5

1. General Description

The manual is intended to assist in safe and reliable use. The product is designed and manufactured strictly according to technical specifications using the latest technology and components and maintaining the highest quality standards.

CAREFULLY READ AND UNDERSTAND THIS MANUAL BEFORE STARTING THE WORK.

To ensure the long and reliable operation of the device, make sure to operate and maintain it properly following the guidelines in this instruction manual. The technical data and specifications in this manual are up-to-date. The manufacturer reserves the right to make changes to improve the quality. Taking the technical progress and the possibility of reducing noise into account, the unit is designed and built in such a way that risks resulting from noise emissions are reduced to the lowest possible level.

Explanation of symbols

The product complies with applicable safety standards.

Read the manual before use.

Recyclable product.

CAUTION! or **WARNING!** or **REMEMBER!** describing a given situation (general warning sign).

Wear protective goggles.

Wear protective gloves.

Wear a protective apron.

CAUTION! Warning of electric shock!

Caution! Hot surface can cause burns!

For indoor use only.

CAUTION! The figures in this manual are illustrative only and may vary in some details from the actual appearance of the product.

2. Safety of use

CAUTION! Read all safety warnings and instructions. Failure to follow the warnings and instructions may result in electric shock, fire, and/or serious injury or death.

The term "device" or "product" in the warnings and the description of the instructions refers to ***Thermostatic water bath***

2.1. Electrical safety

- a) The plug of this device must fit into the outlet. Do not modify the plug in any way. Original plugs and matching outlets reduce the risk of electric shock.
- b) Avoid touching grounded parts, such as pipes, heaters, ovens, and refrigerators. There is an increased risk of electric shock if your body is grounded and touches the device while exposed to direct rain, wet pavement, or while working in a damp environment. If water enters the device, there is an increased risk of damage to the unit and electric shock.
- c) Do not touch the device with wet or damp hands.
- d) Do not use the cord in an unintended manner. Never use it to carry the device or to pull the plug out of the socket. Keep the cord away from heat sources, oil, sharp edges, or moving parts. Damaged or tangled cords increase the risk of electric shock.
- e) If you cannot avoid using the product in a wet environment, use a residual current device (RCD) to connect it to electrical mains. Using an RCD reduces the risk of electric shock.
- f) Do not use the device if the power cord is damaged or shows signs of wear. A damaged power cord should be replaced by a qualified electrician or the manufacturer's service department.
- g) To avoid electric shock, do not immerse the cable, plug, or the device itself in water or other liquid. Do not use the product on wet surfaces.
- h) Use only power sources that are grounded.

2.2. Safety in the workplace

- a) Keep the work area tidy and well-lit. Disorder or poor lighting can lead to accidents. Be foresighted, watch what you are doing and use common sense when using the device.
- b) If you find any damage or irregularities in the operation of the product, immediately turn it off and report it to an authorized person.
- c) Repairs to the device may only be carried out by the manufacturer's service. Do not attempt to repair the product on your own!
- d) In case of open flames or fire, use only dry powder or snow (CO₂) fire extinguishers to extinguish the live equipment.
- e) No children or unauthorized persons are allowed in the work area. (Inattention may result in loss of control of the unit.)
- f) Use the unit in a well-ventilated area.
- g) Check the condition of the safety stickers regularly. Replace them if they are illegible.
- h) Keep these instructions for use for future reference. If the product is to be handed over to a third party, hand it over with this user manual.
- i) Keep packaging components and small installation parts out of the reach of children.

- j) When using this product together with other devices, also follow the other instructions for use.

Remember! Keep children and other bystanders safe while operating the appliance.

2.3. Personal safety

- a) Do not operate this device if you are tired, ill, or under the influence of alcohol, drugs, or medication that could impair your ability to operate the device.
- b) The device is not intended to be used by persons (including children) with reduced mental, sensory, or intellectual functions or persons who lack experience and/or knowledge unless they are supervised or have been instructed by a person responsible for their safety on how to operate the device.
- c) The unit may be operated by persons who are physically fit, capable of operating it, and appropriately trained, and who have read this instruction manual and have been trained in occupational safety and health.
- d) Use caution and common sense when operating this unit. A moment's inattention during the operation may result in serious personal injury.
- e) Use personal protective equipment as required when operating the unit as specified in Section 1 of the explanation of symbols.
The use of appropriate, approved personal protective equipment reduces the risk of injury.
- f) To prevent accidental start-up, make sure the switch is in the off position before connecting to a power source.
- g) Do not wear loose clothing or jewelry. Keep hair, clothing, and gloves away from moving parts. Loose clothing, jewelry, or long hair can be caught in moving parts.
- h) The product is not a toy. Children should be watched to ensure that they do not play with the appliance.

2.4. Safe use of the device

- a) Do not overload the device. Use tools that are suitable for the application. A correctly selected product will do a better and safer job for which it was designed.
- b) Do not use the unit if the ON/OFF (I/O) switch does not function properly (does not turn on and off). Units that cannot be controlled by the switch are unsafe, cannot operate, and must be repaired.
- c) Disconnect the device from the power supply before adjusting, cleaning, or servicing. This precaution reduces the risk of accidental start-up.
- d) Keep unused product out of the reach of children and anyone unfamiliar with the device or this manual. Products are dangerous when used by inexperienced users.
- e) Keep the product in good working order. Check before each use for general damage or damage to moving parts (cracks in parts and components or any other condition that may affect the safe operation of the device). If damaged, return the device for repair before use.
- f) Repairs and maintenance should be carried out by qualified personnel using only original spare parts. This will ensure the safety of use.
- g) To ensure the designed operational integrity of the device, do not remove factory-installed covers or loosen screws.
- h) When transporting or moving the unit from storage to the place of use, observe the health and safety rules for manual handling applicable in the country where the unit is used.
- i) Do not move, shift, or rotate the machine while in operation.
- j) Do not leave the device switched on unattended.
- k) Clean the device regularly to prevent permanent dirt build-up.
- l) The product is not a toy. Cleaning and maintenance must not be performed by children without adult supervision.
- m) Do not tamper with the device to alter its performance or design.
- n) Keep the unit away from sources of fire and heat.
- o) It is forbidden to start an empty device! The minimum water level in the unit is 50 mm above the build plate.

CAUTION! Although the product has been designed to be safe, with adequate safeguards, and despite the additional safety features provided to the user, there is still a slight risk of accident or injury when handling the unit. Caution and common sense are advised when using the product.

3. Instructions for use

The device acts as an intermediary in heating or cooling laboratory vessels with their contents. The device is a vessel filled with a suitable substance that can be heated to transfer heat through the walls of the laboratory vessel to the substance being heated.

The product is intended for home use only!

The user is responsible for any damage resulting from misuse.

3.1. Product overview

1 - Main switch ON/OFF

2 - Water drain

3 - Cover

4 - Control panel:

1 - "PV" temperature display

- 2 - Setpoint temperature display/time display "SV"
- 3 - Indicators RUN (unit in operation) / OUT (heating is on) / AT (AUTO function) / ALM (alarm)
- 4 - SET button
- 5 - SHIFT function button
- 6 - Buttons for decreasing / increasing the entered values

3.2. Preparation for operation

POSITIONING OF THE UNIT

The ambient temperature must not exceed 40°C and ambient humidity should not exceed 85%. Place the unit in a way that ensures good air circulation. Maintain a minimum clearance of 10 cm from any wall of the unit. Keep the unit away from any hot surfaces. Always operate the unit on a level, stable, clean, fireproof, and dry surface and out of the reach of children and persons with impaired mental, sensory, and intellectual functions. Place the unit in such a way that the main plug can be reached at any time. Ensure that the power supply to the unit corresponds to that specified on the identification plate!

Before using the unit for the first time, disassemble all parts and wash them as well as the entire unit.

3.3. Working with the device

- Position the device as described above.
- Open the lid and add clean or distilled water to the water tank. The water level must be higher than the heating element and the temperature sensor. CAUTION! If the error "Er-1" appears on the display, add non-distilled water.
- Connect the appropriate power supply.
- The upper unit display (PV) shows the actual temperature and the lower display (SV) shows the set temperature or time.
- Press the SET button to go to the temperature setting. Set the desired temperature value using the input decrease/increase buttons. Confirm by pressing the SET button again.
- Set the desired operating time:
 - Press the SET button twice.
 - Use the input decrease/increase buttons to set the required time value.
 - Confirm by pressing the SET button.
 - The unit will start counting down when the set temperature is reached.

CAUTION! If only the temperature value is set (time value = 0), the unit will operate in continuous mode.

The AT function is designed to eliminate the possibility of control errors based on the settings of the parameter group (proportional, integral, derivative):

- Hold down the SHIFT function button for a few seconds. The "AT" indicator light will start flashing.
- When finished, the "AT" indicator light will turn off and the unit will return to normal operation.
- To stop the process, hold down the SHIFT function button again.

3.4. Cleaning and maintenance

- a) Pull the mains plug and let the unit cool down completely before cleaning, adjusting or replacing accessories and when the unit is not in use.
- b) Use only non-corrosive cleaning agents for cleaning the surfaces.
- c) After each cleaning, all parts should be thoroughly dried before the device is reused.
- d) Store the unit in a dry and cool place protected from moisture and direct sunlight.
- e) Do not spray the unit with a stream of water or immerse it in water.
- f) Make sure that no water enters through the ventilation openings in the casing.
- g) Clean the ventilation openings with a brush and compressed air.
- h) Perform regular inspections of the unit checking technical fitness and any damages.

- i) Use a soft, damp cloth for cleaning.
- j) Do not use sharp and/or metal objects (e.g. wire brush or metal spatula) as they may damage the surface of the material from which the appliance is made.
- k) Do not clean the unit with acidic substances, medical agents, thinners, fuel, oils, or other chemicals. It may cause damage to the device.

DISPOSAL OF USED UNITS.

At the end of its useful life, this product should not be disposed of with normal household waste but should be taken to a collection point for the recycling of electrical and electronic equipment. This is indicated by the symbol on the product, operating instructions, or packaging. The materials used in this appliance are recyclable according to their marking. By reusing, recycling, or applying other forms of use of waste machines, you make a significant contribution to the protection of our environment.

Your local administration will provide you with information about the appropriate disposal point for used appliances.

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