USER MANUAL

Bedienungsanleitung
Instrukcja obsługi
Návod k použití
Manuel d'utilisation
Istruzioni per l'uso
Manual de instrucciones
Használati útmutató
Brugsanvisning
Käyttöohje
Gebruiksaanwijzing
Bruksanvisning
Instruções de utilização
Používateľská príručka

THERMOSTATIC WATER BATHS

DE	Produktname:	Thermostatisches wasserbad
EN	Product name:	Thermostatic water baths
PL	Nazwa produktu:	Laboratoryjna łaźnia wodna
CZ	Název výrobku	Termostatická vodní lázeň
FR	Nom du produit:	Bains thermostatés
IT	Nome del prodotto:	Bagno termostatico
ES	Nombre del producto:	Baño termostático
HU	Termék neve	Termosztatikus vízfürdő
DA	Produktnavn	Termostatisk vandbad
DE	Modell:	SBS-TWB-2000 SBS-TWB-2002 SBS-TWB-2004 SBS-TWB-2006
EN	Product model:	
PL	Model produktu:	
CZ	Model výrobku	
FR	Modèle:	
IT	Modello:	
ES	Modelo:	
HU	Modell	
DA	Model	
DE	Hersteller	expondo Polska sp. z o.o. sp. k.
EN	Manufacturer	
PL	Producent	
CZ	Výrobce	
FR	Fabricant	
IT	Produttore	
ES	Fabricante	
HU	Termelő	
DA	Producent	
DE	Anschrift des Herstellers	ul. Nowy Kisielin – Innowacyjna 7, 66-002 Zielona Góra Poland, EU
EN	Manufacturer Address	
PL	Adres producenta	
CZ	Adresa výrobce	
FR	Adresse du fabricant	
IT	Indirizzo del produttore	
ES	Dirección del fabricante	
HU	A gyártó címe	
DA	Producentens adresse	

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This User Manual has been translated for your convenience using machine translation. Reasonable efforts have been made to provide an accurate translation; however, no automated translation is perfect nor is it intended to replace human translators. The official User Manual is the English version. Any discrepancies or differences created in the translation are not binding and have no legal effect for compliance or enforcement purposes. If any questions arise related to the accuracy of the information contained in the User Manual, please refer to the English version of those contents which is the official version.

Technical data

Parameter description	Parameter value			
Product name	Thermostatic water bath			
Model	SBS-TWB-2000	SBS-TWB-2002	SBS-TWB-2004	SBS-TWB-2006
Rated voltage [V~] / frequency [Hz]	230V~/50Hz			
Rated power [W]	400	500	1000	1500
Capacity [l]	3.4	6.1	14.6	22.5
Temperature range [°C]	RT+5~100			
Dimensions [width x depth x height; mm]	170 x 150 x 220	330 x 170 x 220	330 x 220 x 350	535 x 320 x 230
Weight [kg]	2.5	3.9	6	9.34

1. General description

The user manual is designed to assist in the safe and trouble-free use of the device. The product is designed and manufactured in accordance with strict technical guidelines, using state-of-the-art technologies and components. Additionally, it is produced in compliance with the most stringent quality standards.

DO NOT USE THE DEVICE UNLESS YOU HAVE THOROUGHLY READ AND UNDERSTOOD THIS USER MANUAL.

To increase the product life of the device and to ensure trouble-free operation, use it in accordance with this user manual and regularly perform maintenance tasks. The technical data and specifications in this user manual are up to date. The manufacturer reserves the right to make changes associated with quality improvement. The device is designed to reduce noise emission risks to a minimum, taking into account technological progress and noise reduction opportunities.

Legend

	The product satisfies the relevant safety standards.
	Read instructions before use.
	The product must be recycled.
	WARNING! or CAUTION! or REMEMBER! Applicable to the given situation. (general warning sign)
	Wear protective goggles.
	Wear protective gloves.

	Wear a protective apron.
	ATTENTION! Electric shock warning!
	ATTENTION! Hot surface, risk of burns!
	Only use indoors.

PLEASE NOTE! Drawings in this manual are for illustration purposes only and in some details may differ from the actual product.

2. Usage safety

ATTENTION! Read all safety warnings and all instructions. Failure to follow the warnings and instructions may result in electric shock, fire and/or serious injury or even death.

The terms "device" or "product" are used in the warnings and instructions to refer to:
Thermostatic water bath

2.1. Electrical safety

- a) The plug must fit the socket. Do not modify the plug in any way. Using original plugs and matching sockets reduces the risk of electric shock.
- b) Avoid touching earthed elements such as pipes, heaters, boilers and refrigerators. There is an increased risk of electric shock if the earthed device is exposed to rain, comes into direct contact with a wet surface or is operating in a damp environment. Water getting into the device increases the risk of damage to the device and of electric shock.
- c) Do not touch the device with wet or damp hands.
- d) Use the cable only for its designated use. Never use it to carry the device or to pull the plug out of a socket. Keep the cable away from heat sources, oil, sharp edges or moving parts. Damaged or tangled cables increase the risk of electric shock.
- e) If using the device in a damp environment cannot be avoided, a residual current device (RCD) should be applied. The use of an RCD reduces the risk of electric shock.
- f) Do not use the device if the power cord is damaged or shows obvious signs of wear. A damaged power cord should be replaced by a qualified electrician or the manufacturer's service centre.
- g) To avoid electric shock, do not immerse the cord, plug or device in water or other liquids. Do not use the device on wet surfaces.
- h) Only use grounded power sources.

2.2. Safety in the workplace

- a) Make sure the workplace is clean and well lit. A messy or poorly lit workplace may lead to accidents. Try to think ahead, observe what is going on and use common sense when working with the device.
- b) If you discover damage or irregular operation, immediately switch the device off and report it to a supervisor without delay.
- c) Only the manufacturer's service point may repair the device. Do not attempt any repairs independently!
- d) In case of fire, use a powder or carbon dioxide (CO₂) fire extinguisher (one intended for use on live electrical devices) to put it out.
- e) Children or unauthorised persons are forbidden to enter a work station. (A distraction may result in loss of control over the device).
- f) Use the device in a well-ventilated space.

- g) Regularly inspect the condition of the safety labels. If the labels are illegible, they must be replaced.
- h) Please keep this manual available for future reference. If this device is passed on to a third party, the manual must be passed on with it.
- i) Keep packaging elements and small assembly parts in a place not available to children.
- j) If this device is used together with another equipment, the remaining instructions for use shall also be followed.

Remember! When using the device, protect children and other bystanders.

2.3. Personal safety

- a) Do not use the device when tired, ill or under the influence of alcohol, narcotics or medication which can significantly impair the ability to operate the device.
- b) The device is not designed to be handled by persons (including children) with limited mental and sensory functions or persons lacking relevant experience and/or knowledge unless they are supervised by a person responsible for their safety or they have received instruction on how to operate the device.
- c) The device can be handled only by physically fit persons who are capable of handling it, properly trained, familiar with this manual and trained within the scope of occupational health and safety.
- d) When working with the device, use common sense and stay alert. Temporary loss of concentration while using the device may lead to serious injuries.
- e) Use personal protective equipment as required for working with the device, specified in section 1 (Legend).
The use of correct and approved personal protective equipment reduces the risk of injury.
- f) To prevent the device from accidentally switching on, make sure the switch is on the OFF position before connecting to a power source.
- g) Do not wear loose clothing or jewellery. Keep hair, clothes and gloves away from moving parts. Loose clothing, jewellery or long hair may get caught in moving parts.
- h) The device is not a toy. Children must be supervised to ensure that they do not play with the device.

2.4. Safe device use

- a) Do not overload the device. Use the appropriate tools for the given task. A correctly-selected device will perform the task for which it was designed better and in a safer manner.
- b) Do not use the device if the ON/OFF switch does not function properly (does not switch the device on and off). Devices which cannot be switched on and off using the ON/OFF switch are hazardous, should not be operated and must be repaired.
- c) Disconnect the device from the power supply before commencement of adjustment, cleaning and maintenance. Such a preventive measure reduces the risk of accidental activation.
- d) When not in use, store in a safe place, away from children and people not familiar with the device who have not read the user manual. The device may pose a hazard in the hands of inexperienced users.
- e) Keep the device in perfect technical condition. Before each use check for general damage and especially check for cracked parts or elements and for any other conditions which may impact the safe operation of the device. If damage is discovered, hand over the device for repair before use.
- f) Device repair or maintenance should be carried out by qualified persons, only using original spare parts. This will ensure safe use.
- g) To ensure the operational integrity of the device, do not remove factory-fitted guards and do not loosen any screws.
- h) When transporting and handling the device between the warehouse and the destination, observe the occupational health and safety principles for manual transport operations which apply in the country where the device will be used.
- i) Do not move, adjust or rotate the device in the course of work.
- j) Do not leave this appliance unattended while it is in use.
- k) Clean the device regularly to prevent stubborn grime from accumulating.
- l) The device is not a toy. Cleaning and maintenance may not be carried out by children without supervision by an adult person.

- m) It is forbidden to interfere with the structure of the device in order to change its parameters or construction.
- n) Keep the device away from sources of fire and heat.
- o) It is forbidden to run the empty device! The minimum water level in the device is 50 mm above the build plate.

ATTENTION! Despite the safe design of the device and its protective features, and despite the use of additional elements protecting the operator, there is still a slight risk of accident or injury when using the device. Stay alert and use common sense when using the device.

3. Use guidelines

The device mediates the cooling of laboratory vessels with their contents. The device is a vessel filled with a suitable substance that can be heated in order to transfer heat through the walls of the laboratory vessel to the substance to be heated.

The product is intended for home use only!

The user is liable for any damage resulting from unintended use of the device.

3.1. Device description

SBS-TWB-2000	 <p>The image shows the SBS-TWB-2000 device with its lid closed. A green power switch is labeled '1'. The digital display shows '22.1' and 'E--1'. The control panel is labeled '2'.</p>
SBS-TWB-2002	 <p>The image shows the SBS-TWB-2002 device with its lid removed, revealing a heating plate. The green power switch is labeled '1'. The digital display shows '8888'. The control panel is labeled '2'. A red emergency stop button is labeled '3'. The device body is labeled '4'.</p>

1 - Main switch ON/OFF.

2 - Control panel.

3 - ON/OFF switch for pumping water out of the unit - during normal operation this switch must be turned OFF - OFF position (models SBS-TWB-2002, SBS-TWB-2004, SBS-TWB-2006 only).

4 - Water drain (not visible on the picture), (only models SBS-TWB-2002, SBS-TWB-2004, SBS-TWB-2006).

Control panel:

1 - Actual temperature display "PV"

2 - Set temperature display / time display "SV"

3 - Indicator lights: RUN (unit is running) / OUT (heating is on) / AT (AUTO function) / ST (set time) / ALM (alarm)

4 - SET button

5 - SHIFT function button

6 - Buttons for decreasing / increasing the entered values

3.2. Preparing for use

APPLIANCE LOCATION

The temperature of environment must not be higher than 40°C and the relative humidity should be less than 85%. Ensure good ventilation in the room in which the device is being used. There should be at least 10 cm distance between each side of the device and the wall or other objects. The device should always be used when positioned on an even, stable, clean, fireproof and dry surface, and be out of the reach of children and persons with limited mental and sensory functions. Position the device such that you always have access to the power plug. The power cord connected to the appliance must be properly grounded and correspond to the technical details on the product label.

Disassemble the device and all its components and clean them before the first use.

3.3. Device use

- Open the lid and add clean water or distilled water to the water tank. The water level must be higher than the heating element and the temperature sensor. NOTE: If an "Er-1" error appears on the display, add non-distilled water.
- Connect the appropriate power supply.
- The upper display of the unit (PV) shows the actual temperature and the lower display (SV) shows the set temperature or time.
- Press the SET button to proceed to temperature setting. Use the decrease/increase input buttons to set the desired temperature value. Confirm by pressing the SET button again.

Set the desired operating time:

- Press the SET button twice.
- Set the desired time value by means of the decreasing/increasing entry keys.
- Confirm by pressing the SET button.
- The device will start timing when the set temperature value is reached.

ATTENTION! If only the temperature value is set (time value = 0) the unit will operate in continuous mode.

The AT function is designed to eliminate the possibility of control errors based on parameter group settings (proportional, integral, differential):

- Hold down the SHIFT function key for few seconds. The "AT" indicator lamp will start flashing.
- When finished, the "AT" indicator light will turn off and the unit will return to normal operation.
- To interrupt the process, hold down the SHIFT function button again.

ATTENTION! For models SBS-TWB-2002, SBS-TWB-2004, SBS-TWB-2006 the ON/OFF switch "OUTLET" for draining water from the unit - during normal operation this switch must be turned off (OFF position). When finished working, pour out the water through the water drain - turn OUTLET switch to ON position.

3.4. Cleaning and maintenance

- a) Unplug the mains plug and allow the device to cool completely before each cleaning, adjustment or replacement of accessories, or if the device is not being used.
- b) Use only non-corrosive cleaners to clean the surface.
- c) After cleaning the device, all parts should be dried completely before using it again.
- d) Store the unit in a dry, cool place, free from moisture and direct exposure to sunlight.
- e) Do not spray the device with a water jet or submerge it in water.
- f) Do not allow water to get inside the device through vents in the housing of the device.
- g) Clean the vents with a brush and compressed air.
- h) The device must be regularly inspected to check its technical efficiency and spot any damage.
- i) Use a soft, damp cloth for cleaning.
- j) Do not use sharp and/or metal objects for cleaning (e.g. a wire brush or a metal spatula) because they may damage the surface material of the appliance.
- k) Do not clean the device with an acidic substance, agents of medical purposes, thinners, fuel, oils or other chemical substances because it may damage the device.

DISPOSING OF USED DEVICES:

Do not dispose of this device in municipal waste systems. Hand it over to an electric and electrical device recycling and collection point. Check the symbol on the product, instruction manual and packaging. The plastics used to construct the device can be recycled in accordance with their markings. By choosing to recycle you are making a significant contribution to the protection of our environment.

Contact local authorities for information on your local recycling facility.

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3.1. Description de l'appareil

SBS-TWB-2002	 <p>The image shows the SBS-TWB-2002 water bath with its lid removed. The control panel features a digital display showing '88.88', a green power button (1), a control panel (2), and a red emergency stop button (3). A drainage port (4) is visible on the right side.</p>
SBS-TWB-2004	 <p>The image shows the SBS-TWB-2004 water bath with its lid on. The control panel features a digital display showing '22.5', a green power button (1), a control panel (2), and a red emergency stop button (3). A drainage port (4) is visible on the right side.</p>
SBS-TWB-2006	 <p>The image shows the SBS-TWB-2006 water bath with its lid on. The control panel features a digital display showing '22.7', a green power button (1), a control panel (2), and a red emergency stop button (3). A drainage port (4) is visible on the right side.</p>

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3.1. Descripción del aparato

SBS-TWB-2002	 The image shows the SBS-TWB-2002 water bath with its lid removed. The lid is a flat metal plate with two circular heating elements. The main unit is a white rectangular box with a control panel on the front. The control panel features a digital display showing '88.88', a green power button (1), a red emergency stop button (3), and several smaller buttons for temperature and mode selection (2). A drain outlet (4) is visible on the right side of the unit.
SBS-TWB-2004	 The image shows the SBS-TWB-2004 water bath. It has a similar design to the 2002 model but with a lid that has four heating elements. The digital display shows '22.5'. The control panel (2) includes a green power button (1) and a red emergency stop button (3). A drain outlet (4) is on the right side.
SBS-TWB-2006	 The image shows the SBS-TWB-2006 water bath. It has a lid with four heating elements. The digital display shows '22.7'. The control panel (2) includes a green power button (1) and a red emergency stop button (3). A drain outlet (4) is on the right side.

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