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USER MANUAL

DIGITAL PRECISION SCALE

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LABORATORNÍ VÁHA

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BALANCE DE PRECISION

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BILANCIA DI PRECISIONE

MANUAL DE INSTRUCCIONES

BALANZA DE PRECISIÓN

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BRUGSANVISNING

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DE | EN | PL | CZ | FR | IT | ES | HU | DA | PT

SBS-LW-1200-MAX | SBS-LW-200-MAX
SBS-LW-300-MAX | SBS-LW-600-MAX

DE	Produktname:	Präzisionswaage
EN	Product name:	Digital Precision Scale
PL	Nazwa produktu:	Waga laboratoryjna
CZ	Název výrobku:	Laboratorní váha
FR	Nom du produit:	Balance de précision
IT	Nome del prodotto:	Bilancia di precisione
ES	Nombre del producto:	Balanza de precisión
HU	Termék neve:	Precíziós mérleg
DA	Produktnavn:	Præcisionsvægt
PT	Nome do produto:	Balança de Precisão Digital

DE	Modell:	SBS-LW-1200-MAX SBS-LW-200-MAX SBS-LW-300-MAX SBS-LW-600-MAX
EN	Product model:	
PL	Model produktu:	
CZ	Model výrobku:	
FR	Modèle:	
IT	Modello:	
ES	Modelo:	
HU	Modell:	
DA	Model:	
PT	Modelo:	

DE	Hersteller:	expondo Polska sp. z o.o. sp. k.
EN	Manufacturer:	
PL	Producent:	
CZ	Výrobce:	
FR	Fabricant:	
IT	Produttore:	
ES	Fabricante:	
HU	Termelő:	
DA	Producent:	
PT	Fabricante:	

DE	Anschrift des Herstellers:	ul. Nowy Kisielin – Innowacyjna 7, 66-002 Zielona Góra Poland, EU
EN	Manufacturer Address:	
PL	Adres producenta:	
CZ	Adresa výrobce:	
FR	Adresse du fabricant:	
IT	Indirizzo del produttore:	
ES	Dirección del fabricante:	
HU	A gyártó címe:	
DA	Producentens adresse:	
PT	Endereço do fabricante:	

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This User Manual has been translated for your convenience using machine translation. Reasonable efforts have been made to provide an accurate translation; however, no automated translation is perfect nor is it intended to replace human translators. The official User Manual is the English version. Any discrepancies or differences created in the translation are not binding and have no legal effect for compliance or enforcement purposes. If any questions arise related to the accuracy of the information contained in the User Manual, please refer to the English version of those contents which is the official version.

Technical data

Parameter description	Parameter value			
Product name	Digital precision scale			
Model	SBS-LW-1200-MAX	SBS-LW-200-MAX	SBS-LW-300-MAX	SBS-LW-600-MAX
Battery	7.4V DC / 2200mAh			
Power supply	12 V DC / 1 A			
Maximum load	1.2 kg	210 g	310 g	610 g
Minimum load	0.2 g	0.02 g	0.02 g	0.2 g
Breakdown	0.01 g	0.001 g	0.001 g	0.01 g

1. General Description

This manual is intended to assist you for safe and reliable use. The product is designed and manufactured strictly according to technical specifications using the latest technology and components and maintaining the highest quality standards.

CAREFULLY READ AND UNDERSTAND THIS MANUAL BEFORE STARTING THE WORK.

To ensure long and reliable operation of the device, make sure to operate and maintain it properly in accordance with the guidelines in this instruction manual. The technical data and specifications in this manual are up-to-date. The manufacturer reserves the right to make changes in order to improve the quality. Taking the technical progress and the possibility of reducing noise into account, the unit is designed and built in such a way so that risks resulting from noise emissions are reduced to the lowest possible level.

Explanation of symbols

	The product complies with applicable safety standards.
	Read the manual before use.
	Recyclable product.
	CAUTION! or WARNING! or REMEMBER! describing a given situation (general warning sign).
	CAUTION! Risk of electric shock!

CAUTION! The figures in this manual are illustrative only and may vary in some details from the actual appearance of the product.

2. Safety of use

CAUTION! Read all safety warnings and instructions. Failure to follow the warnings and instructions may result in electric shock, fire and/or serious injury or death.

The term "device" or "product" in the warnings and in the description of the instructions refers to digital precision scale.

Digital precision scale

2.1. Electrical safety

Applies to the battery charger:

- The plug of this device must fit into the outlet. Do not modify the plug in any way. Original plugs and matching outlets reduce the risk of electric shock.
- Avoid touching grounded parts, such as pipes, heaters, ovens, and refrigerators. There is an increased risk of electric shock if your body is grounded and touches the device while exposed to direct rain, wet pavement, or while working in a damp environment. If water enters the device, there is an increased risk of damage to the unit and electric shock.
- Do not touch the device with wet or damp hands.

- d) Do not use the cord in an unintended manner. Never use it to carry the device or to pull the plug out of the socket. Keep the cord away from heat sources, oil, sharp edges or moving parts. Damaged or tangled cords increase the risk of electric shock.
- e) When operating the product outdoors, use an extension cord designed for outdoor use. Using an extension cord designed for outdoor use reduces the risk of electric shock.
- f) If you cannot avoid using the product in a wet environment, use a residual current device (RCD) to connect the product to electrical mains. Using an RCD reduces the risk of electric shock.
- g) Do not use the device if the power cord is damaged or shows signs of wear. A damaged power cord should be replaced by a qualified electrician or the manufacturer's service department.
- h) To avoid electric shock, do not immerse the cable, plug, or the device itself in water or other liquid. Do not use the appliance on wet surfaces.
- i) When cleaning or using the appliance, never immerse it in water or other liquids.
- j) Do not allow the machine to get wet. Risk of electric shock!
- k) Use only the included power adapter!

2.2. Safety in the workplace

- a) Do not use the unit in a hazardous zone, for example in the presence of flammable liquids, gases or dust.
- b) If in doubt as to whether the product is working properly or found to be damaged, contact the manufacturer's technical service.
- c) Only the manufacturer's service department may repair the appliance. Do not attempt to repair the product on your own!
- d) In the event of ignition or a fire, use dry powder or CO2 extinguishers only to suppress the fire if the appliance is live.
- e) Keep this manual for future reference. If the product is to be handed over to a third party, hand it over with this user manual.
- f) Keep packaging components and small installation parts out of the reach of children.

Remember! Keep children and other bystanders safe while operating the appliance.

2.3. Personal safety

- a) The product is not intended to be used by individuals (including children) with reduced mental, sensory or intellectual capacity or lack of experience and/or knowledge, unless they are supervised by an individual responsible for their

safety or have been given instructions by the responsible individual on how to operate the product.

- b) The appliance is not a toy. Children should be watched to ensure that they do not play with the appliance.

2.4. Safe use of the product

- a) Keep unused appliance out of the reach of children and anyone unfamiliar with the appliance or this manual. Products are dangerous when used by inexperienced users.
- b) Keep the product in good working order. If damaged, have the device repaired before use.
- c) Keep the product out of the reach of children.
- d) Repairs and maintenance should be carried out by qualified personnel using only original spare parts. This will ensure the safety of use.
- e) To ensure the designed operational integrity of the device, do not remove factory-installed covers or loosen screws.
- f) Clean the device regularly to prevent permanent dirt build-up.
- g) The product is not a toy. Cleaning and maintenance must not be performed by children without adult supervision.
- h) Do not tamper with the device to alter its performance or design.
- i) Keep the unit away from sources of fire and heat.
- j) Do not overload the device.


CAUTION! Although the product has been designed to be safe, with adequate safeguards, and despite the additional safety features provided to the user, there is still a slight risk of accident or injury when handling the unit. Caution and common sense are advised when using the product.

3. Instructions for use


The device is designed to measure the weight of elements placed on the pan. Do not exceed the maximum load of the scale - it may damage it!

The user is responsible for any damage resulting from misuse.


3.1. Product overview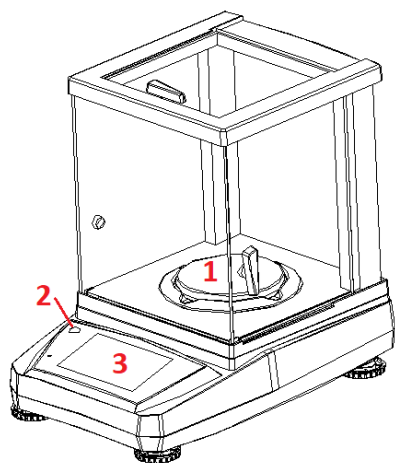

- 1 - pan
- 2 - level
- 3 - display

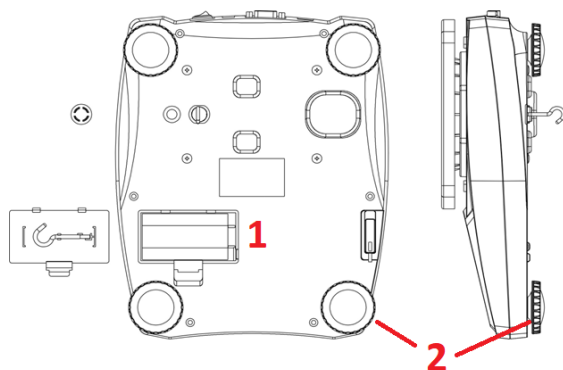- 1 - battery compartment
- 2 - leveling feet

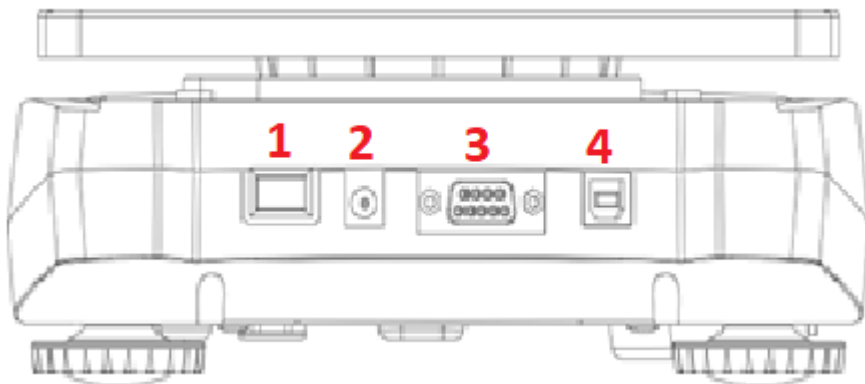- 1 - Switch
- 2 - Power socket
- 3 - RS232 socket
- 4 - USB socket

3.2. Preparation for operation

POSITIONING OF THE UNIT

The ambient temperature must not exceed 40°C and ambient humidity should not exceed 85%. Place the unit in a way that ensures good air circulation. Maintain a minimum clearance of 10 cm from any wall of the unit. Keep the unit away from any hot surfaces. Always operate the unit on a level, stable, clean, fireproof and dry surface and out of the reach of children and persons of impaired mental, sensory and intellectual functions. Place the unit in such a way that the mains plug can be reached at any time. Ensure that the power supply to the unit corresponds to that specified on the identification plate!

ASSEMBLY OF THE DEVICE

In order for the scale to function properly, it must be perfectly level. The scale is equipped with one level and adjustable feet for level control, with which it is possible to compensate for slight differences in height and/or any unevenness in the surface on which the scale is standing.

Adjust the feet so that the air bubble is exactly in the middle of the vial.

Place the scale horizontally and align the bubble inside the level.

For accurate measurements, re-level the scale carefully after each repositioning.

1


2

1 - the scale is properly leveled

2 - the scale is not properly leveled

3.3. Working with the device

3.3.1. Description of the display: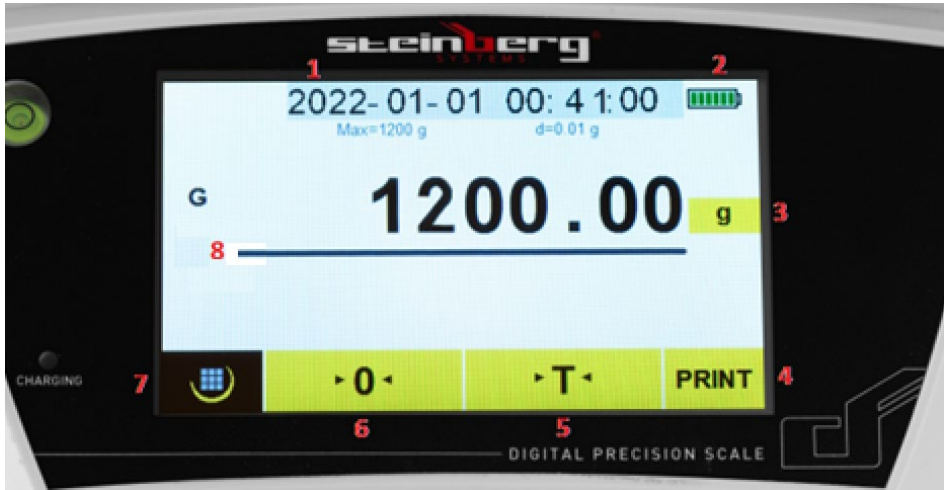

- 1 - Date and time
- 2 - Battery indicator
- 3 - Weighing unit
- 4 - Print button
- 5 - Tare button
- 6 - Reset button
- 7 - Menu button
- 8 - Weighing measurement range

The main switch of the scale is located on the back.

The zeroing function should only be used when there is nothing on the pan and the display does not show "0". The scale may not read "0" due to dust, dirt or moisture on the weighing pan or when a heavy object is dropped on the scale. The zeroing function essentially corrects the true zero on the scale and informs it that there is nothing on the pan.


The tare function is the weight of the package itself. It is the difference between the total weight of the goods with packaging and the weight of the goods without packaging. The tare is used to remove the weight of an item that should not be included in the overall weight measurement.

By pressing the **weighing units** button, you can set the desired unit:


<	Unit				
	g	oz	lb	dwt	
	ozt	ct	tl.T	tl.H	
	tl.J	GN	dr	MM	
	tola	gsm	T/A/R	T/M/R	

3.3.2. Scale functions: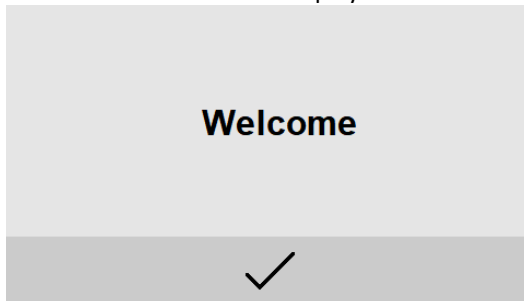

Turn on the scale and the display will show:

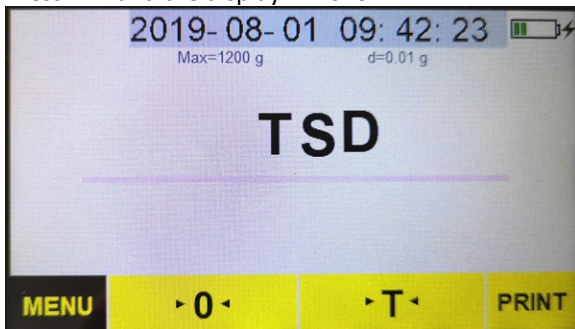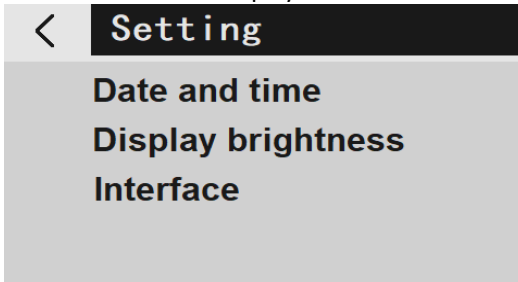
Press ✓ and the display will show:

Press the menu button **MENU** and the display will show:

 CAL  	Weighing	Counting
	Percentage	Density

Press  and the display will show:

3.3.2.1. Setting the date and time:

Press "Date and time" and the display will show:

20YY-MM-DD HH:QQ:SS			✓
1	2	3	←
4	5	6	
7	8	9	C
0			

After entering the date and time, press ✓.

3.3.2.2. Display brightness setting:


Press "Display brightness" and the display will show:

Press **CHANGE** to select "Medium", "Slight", "Bright" and press ✓.

3.3.2.3. Interface settings:

Press "Interface" and the display will show: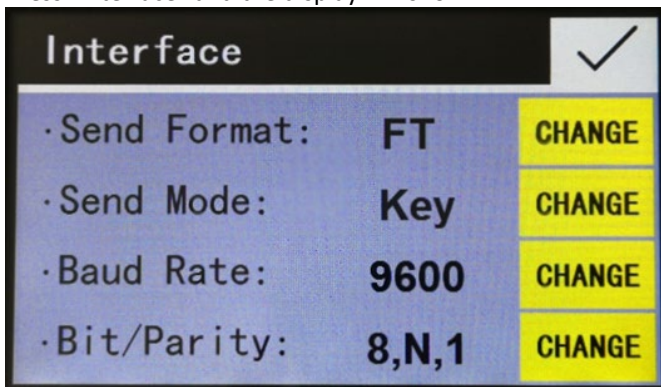

Press **CHANGE** to choose:

"Send Format:" PM/FT, default setting: PM

"Send Mode:" Key/Stb/Con/OFF, default setting: Stb

"OFF" means that the serial data output is turned off

"Key" indicates the manual mode by pressing the [PRINT] key

"Stb" means automatic printout when the scale is stable

"Con" stands for continuous printing

Baud Rate (transmission speed): 1200/2400/4800/9600, Default Setting: 9600


Bit/parity: 8, N, 1/7, O, 1/7, E, 1, default setting: 8, N, 1. (7, O, 1 means 7 data bits with odd parity, 7, E, 1 means 7 data bits with parity, 8, N, 1 means 8 data bits without parity)

Press ✓ to return to the setting mode.


Press **<** then **←** to return to the weighing mode.

3.3.2.4. Application options:


Turn on the scale and the display will show: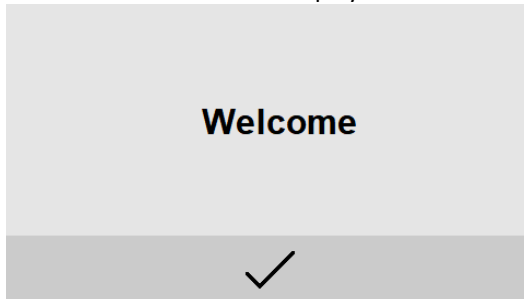

Press ✓ and the display will show:

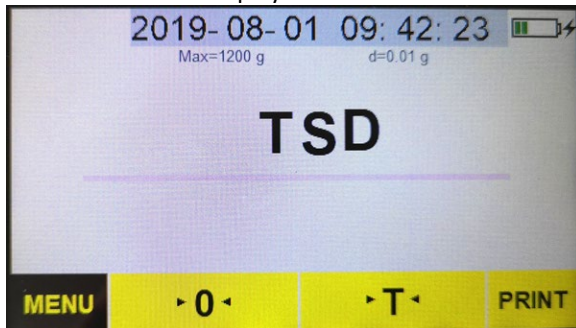Press **MENU** and the display will show:

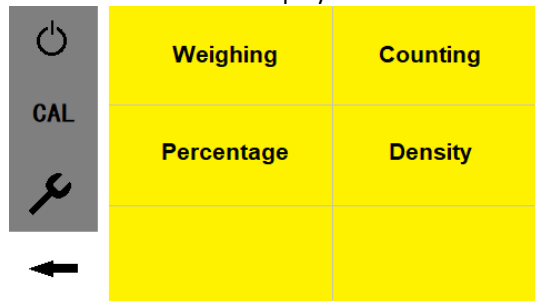
3.3.2.5. Counting function

Press "Counting" and the display will show:


Place the container on the weighing pan and press "T" to tare the scale. Press

Reference pcs and the display will show: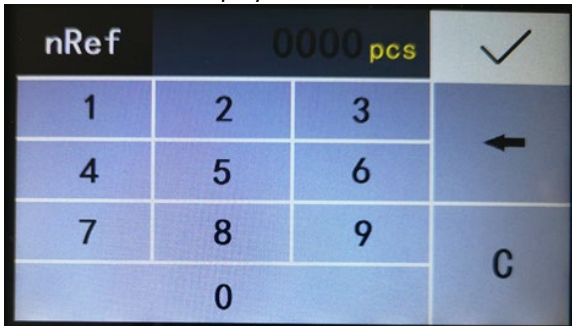

Enter the reference number of samples (example: 50 pieces), then press ✓ and the display will show:

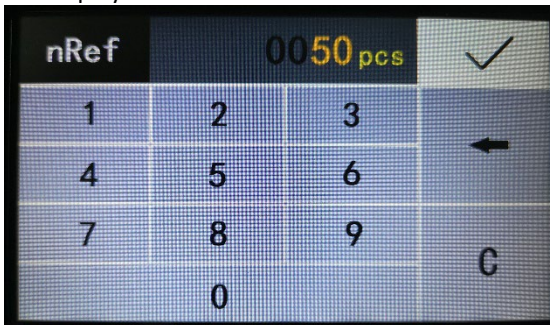
Put in the container the number of samples corresponding to the entered reference number, then press START and the display will show:

Put the total number of reference samples and the display will show:

Remove the samples, press END to exit the counting function and the display will show: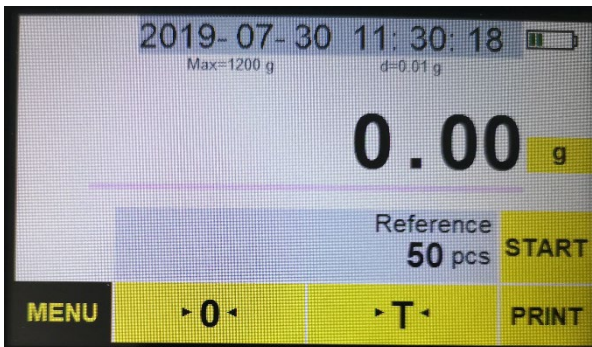

3.3.2.6. Percentage function


Press "Percentage" and the display will show:


Press Reference % and the display will show:


Enter the percentage of the reference sample (example: 10%):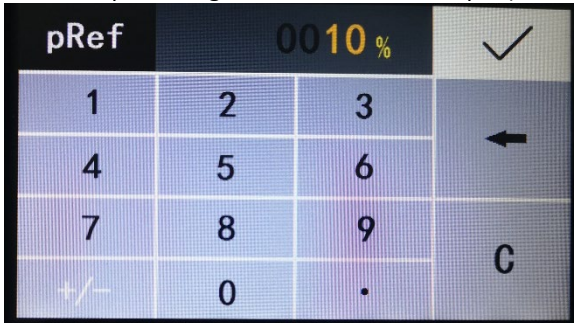


And press ✓ :

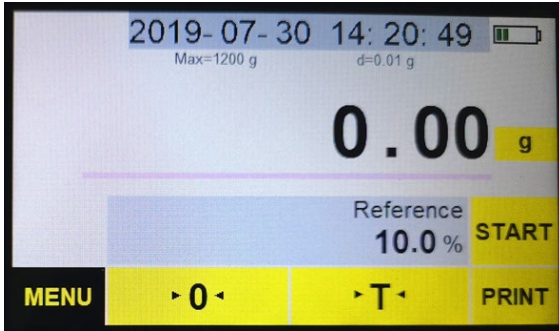Place the samples in the container on the pan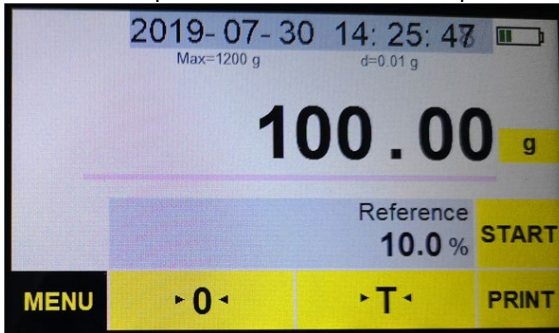


Press START and the display will show the percentage of the reference sample.

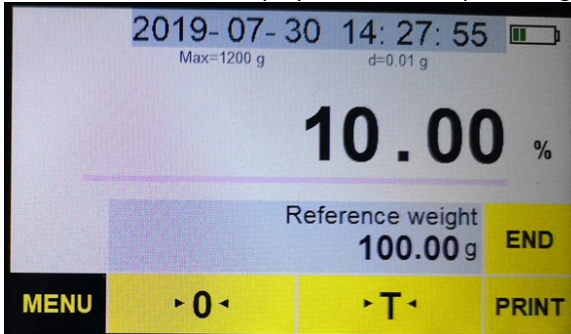Put the total number of reference samples into the container and the display will show the percentage of samples:


Remove the samples, press END to exit the percentage function: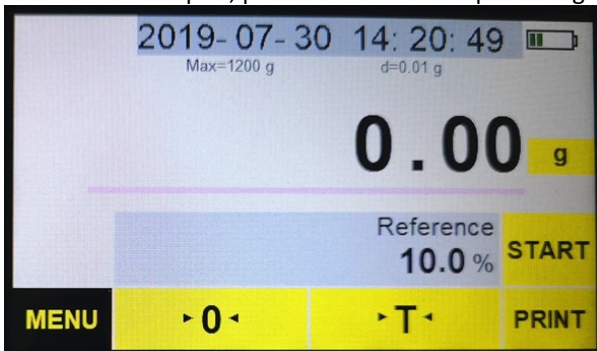


3.3.2.7. Density function

To perform this function, you need a density measuring tool - **this tool is not part of the scale kit - it is not included in the scope of delivery of the scale.**


Press "Density" and the display will show: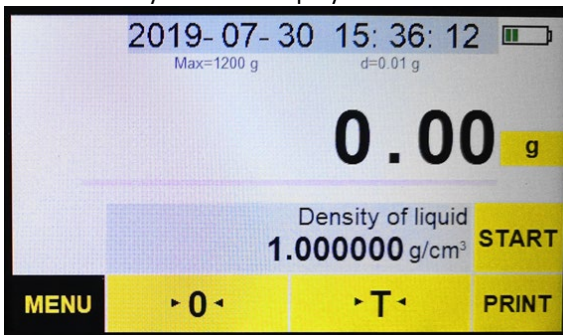


Press Density of liquid
g/cm³ and the display will show: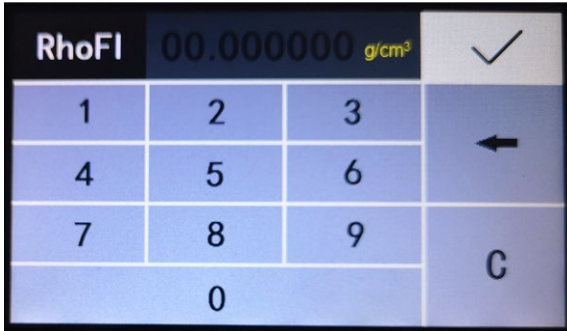

Enter the medium density of the liquid, then press ✓ :

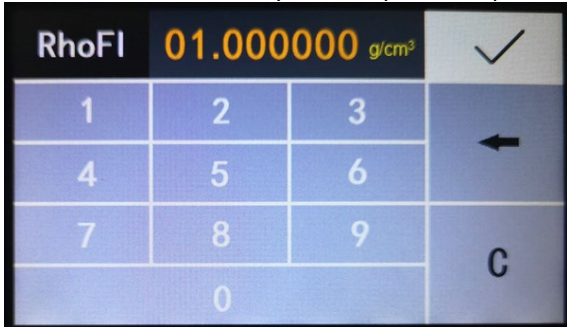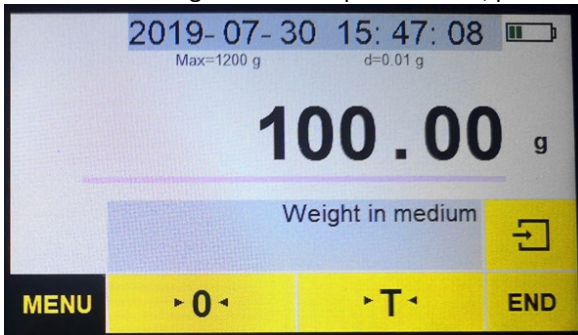
Remove the tare from the scale and set to zero, press START

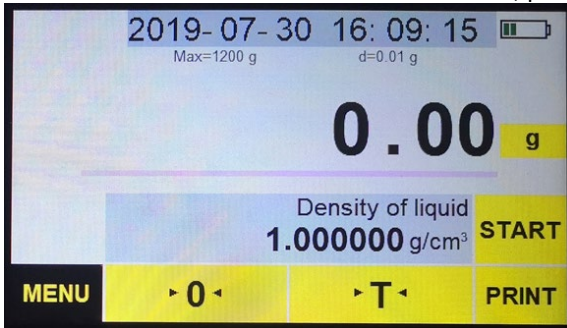Place the sample on the upper platform of the densimeter .


To check the weight of the sample in the air, press  and the display will show:

Press  and the display will show the density of the solid sample


Remove the load, press END to exit the density function: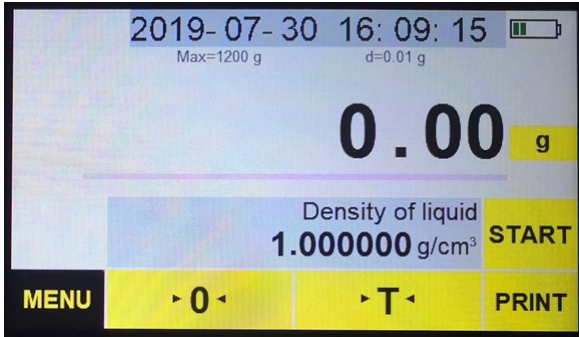


3.3.2.8. Calibration

Calibration is strongly recommended when the scale is initially installed or when the scale is moved far from its original location, or when environmental conditions change significantly. Before weighing and calibrating, it is always necessary to warm up the scale for 30 minutes.


Turn on the scale and the display will show: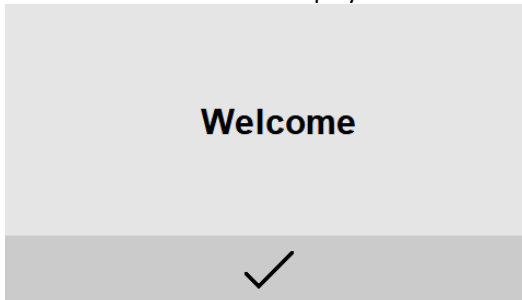


Press ✓ and the display will show:

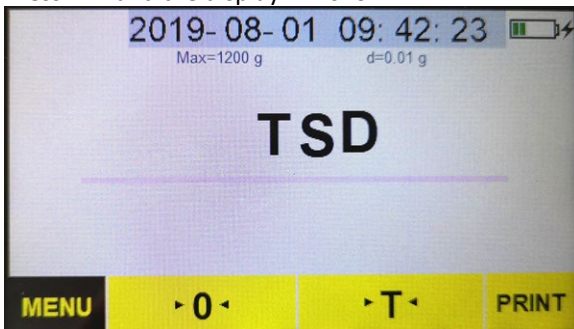
Press **MENU** and the display will show: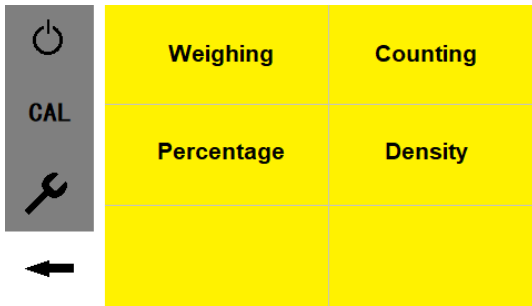


Press "CAL" and the display will show:

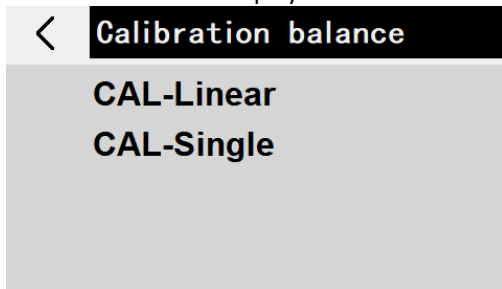
Press "CAL-Single" and the display will show: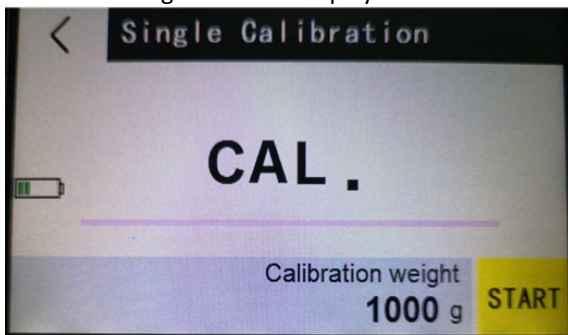


After a few seconds, the display will show a series of numbers which are referred to as A/D values: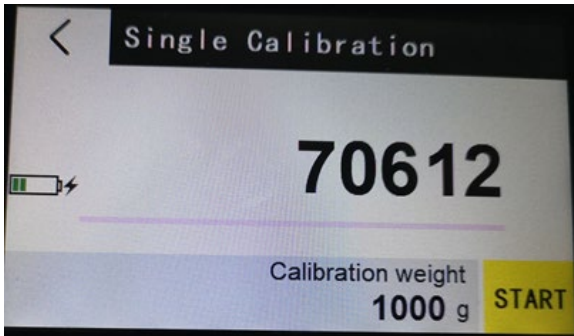


Press "Calibration weight" and the display will show: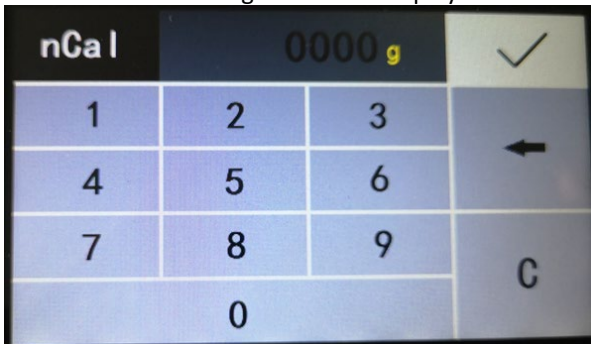


Enter the value of the calibration load using digits (the calibration load can be set according to the user's needs; it is recommended that the minimum calibration weight is at least 50% of the maximum load range of the scale)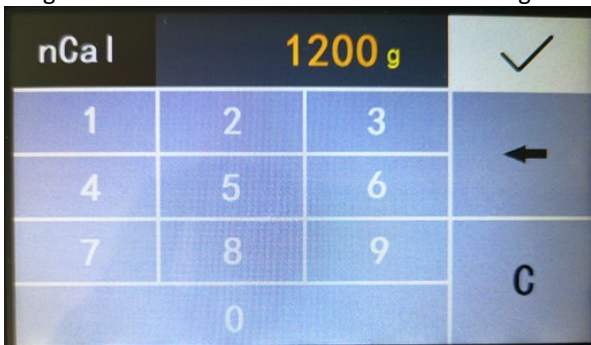

Press the ✓ button and the display will show: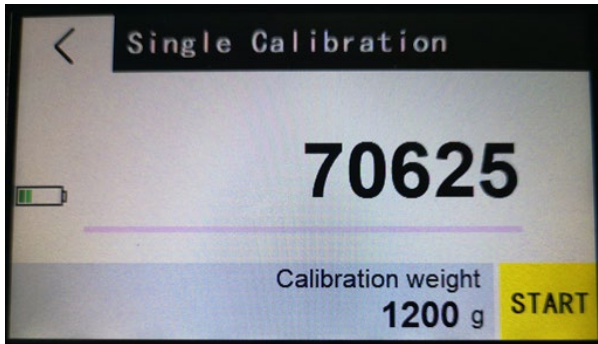

Press the START button and the display will show a flashing calibration value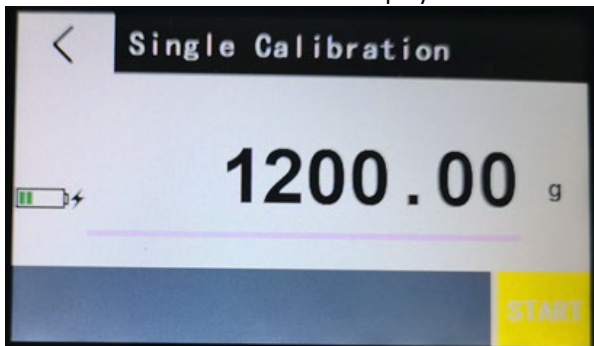

Place the appropriate test weight in the center of the pan.

Wait for a steady indicator and the display will return to weighing mode.


Calibration is complete

Place a known weight on the pan to check that it is accurate.

If the displayed weight does not match the value of the calibration weight, repeat the above steps.

3.3.2.9. Cooperation with a computer

Connect the scale to the computer with the RS232 cable: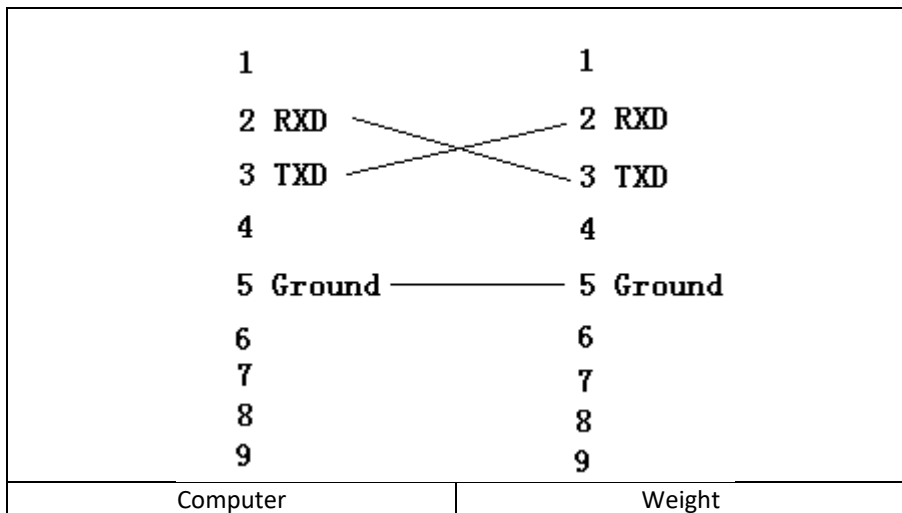

Turn on the scale. Run the "AccessPort.exe" program (to be downloaded from the expondo website).

In the menu, select "View" -> "Language" -> "English"


In the menu, select: "Tools" -> "Configuration" - "Display":

NOTE: Select the COM port on the computer.


Press OK.

View: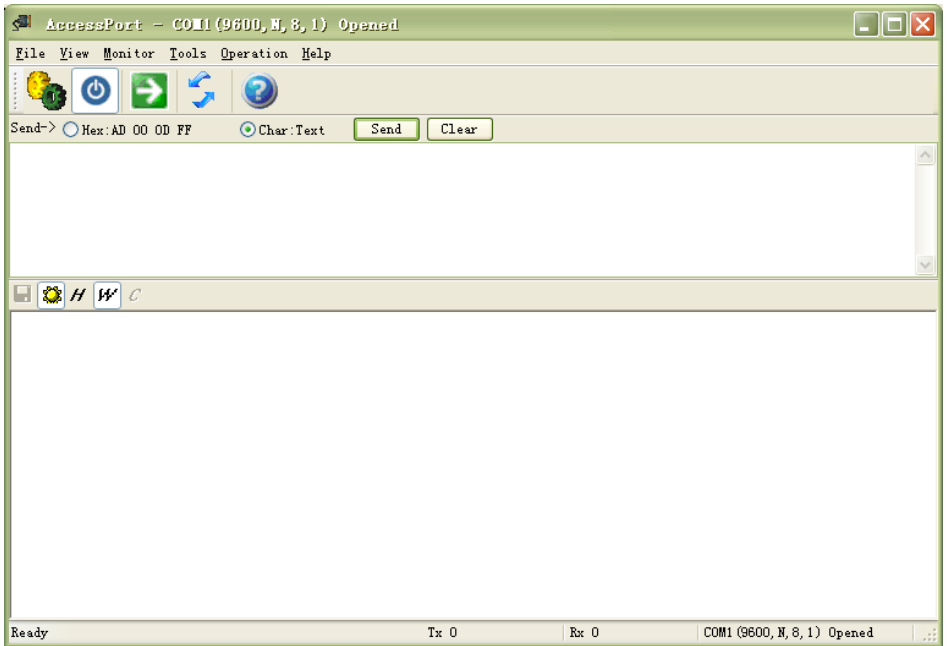

Press PRINT on the scale, the result will be shown on the screen:

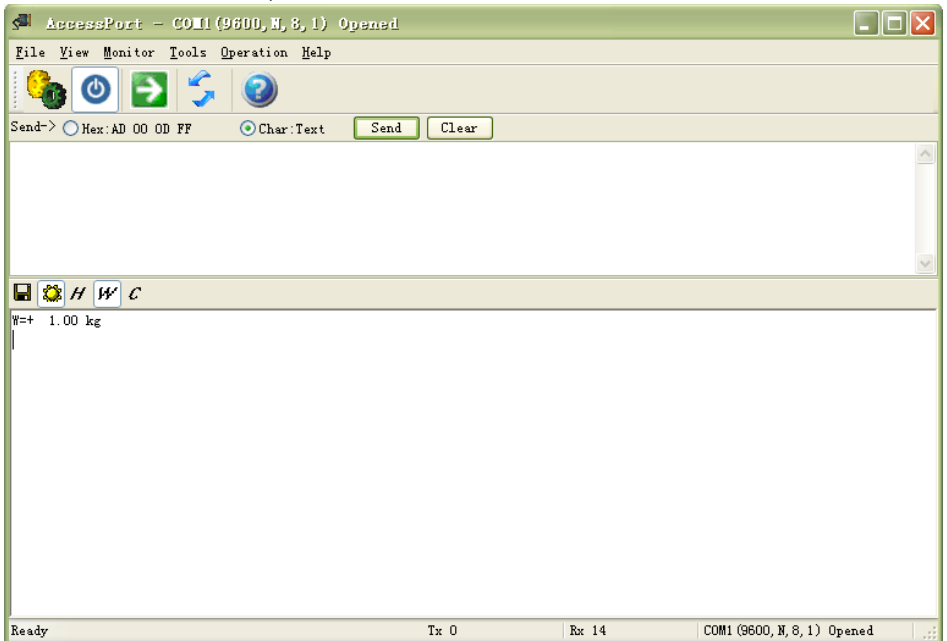3.4. Cleaning and maintenance

- a) Pull out the mains plug before each cleaning and when the unit is not in use.
- b) Use only non-corrosive cleaning agents for cleaning the surfaces.
- c) After each cleaning, all parts should be thoroughly dried before the device is reused.
- d) Store the unit in a dry and cool place protected from moisture and direct sunlight.
- e) Do not spray the unit with a stream of water or immerse it in water.
- f) Make sure that no water enters through the ventilation openings in the casing.
- g) Perform regular inspections of the unit checking technical fitness and any damages.
- h) Use a soft cloth for cleaning.
- i) Do not leave the batteries in the unit when it will not be used for a long time.
- j) Do not use sharp and/or metal objects (e.g. wire brush or metal spatulas) as they may damage the surface of the material from which the unit is made.
- k) Do not clean the unit with acidic substances, medical agents, thinners, fuel, oils or other chemicals. It may cause damage to the device.

INSTRUCTIONS FOR THE SAFE DISPOSAL OF BATTERIES.

A 7.4 V DC / 2200 mAh battery is installed in the device.

Remove used batteries from the unit in the same way you put them in.

Return the batteries to a unit responsible for their disposal.

DISPOSAL OF USED UNITS.

At the end of its useful life, this product should not be disposed of with normal household waste but should be taken to a collection point for the recycling of electrical and electronic equipment. This is indicated by the symbol on the product, operating instructions or packaging. The materials used in this appliance are recyclable according to their marking. By reuse, recycling or applying other forms of use of waste machines, you make a significant contribution to the protection of our environment.

Your local administration will provide you with information about the appropriate disposal point for used appliances.

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