USER MANUAL

Bedienungsanleitung
Instrukcja obsługi
Návod k použití
Manuel d'utilisation
Istruzioni per l'uso
Manual de instrucciones
Használati útmutató
Bruksanvisning
Käyttöohje
Gebruiksaanwijzing
Bruksanvisning
Instruções de utilização
Používateľská príručka
Ръководство за употреба
Οδηγίες χρήσης
Upute za uporabu
Naudojimo instrukcija
Manual de utilizare
Navodila za uporabo

MULTICHANNEL PIPETTE

DE	Produktname	Pipette
EN	Product name	Multichannel Pipette
PL	Nazwa produktu	Pipeta wielokanałowa
CZ	Název výrobku	Vícekanálová pipeta
FR	Nom du produit	Pipette multicanaux
IT	Nome del prodotto	Pipetta multicanale
ES	Nombre del producto	Pipeta multicanal
HU	Termék neve	Többcsatornás pipetta
DA	Produktnavn	Flerkanalpipette
FI	Tuotteen nimi	Monikanavainen pipetti
NL	Productnaam	Multikanaalpipet
NO	Produktnavn	Flerkanalpipette
SE	Produktnamn	Flervägspipett
PT	Nome do produto	Pipeta multicanal
SK	Názov produktu	Viackanálová pipeta
BG	Име на продукта	Многоканална пипета
EL	Όνομα προϊόντος	Πολυκάναλη πιπέτα
HR	Naziv proizvoda	Višekanalna pipeta
LT	Produktu pavadinimas	Daugiakanalė pipetė
RO	Numele produsului	Pipetă multicanal
SL	Ime izdelka	Večkanalna pipeta
DE Modell EN Product model PL Model produktu CZ Model výrobku FR Modèle IT Modello ES Modelo HU Modell DA Model FI Tuotteen malli NL Productmodel NO Produktmodell SE Produktmodell PT Modelo do produto SK Model BG Модел на продукт EL Μοντέλο προϊόντος HR Model proizvoda LT : Gaminio modelis RO : Model de produs SL : Model izdelka		SBS-LAB-117 SBS-LAB-118 SBS-LAB-119 SBS-LAB-120 SBS-LAB-121 SBS-LAB-122
DE Hersteller EN Manufacturer PL Producent CZ Výrobce FR Fabricant IT Produttore ES Fabricante HU Termelő DA Producent FI Valmistaja NL Producent NO Produsent SE Tillverkare PT Fabricante SK Výrobca BG Производител EL Κατασκευαστής HR Proizvođač LT Gamintojas RO Producător SL Proizvajalec		expondo Polska sp. z o.o. sp. k.
DE Anschrift des Herstellers EN Manufacturer Address PL Adres producenta CZ Adresa výrobce FR Adresse du fabricant IT Indirizzo del produttore ES Dirección del fabricante HU A gyártó címe DA Producentens adresse FI Valmistajan osoite NL Adres producent NO Produsentens adresse SE Tillverkarens adress PT Endereço do fabricante SK Adresa výrobcu BG Адрес на производителя EL : Διεύθυνση κατασκευαστή HR Adresa proizvođača LT Gamintojo adresas RO Adresa producătorului SL Naslov proizvajalca		ul. Nowy Kisielin – Innowacyjna 7, 66-002 Zielona Góra Poland, EU

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This User Manual has been translated for your convenience using machine translation. Reasonable efforts have been made to provide an accurate translation; however, no automated translation is perfect nor is it intended to replace human translators. The official User Manual is the English version. Any discrepancies or differences created in the translation are not binding and have no legal effect for compliance or enforcement purposes. If any questions arise related to the accuracy of the information contained in the User Manual, please refer to the English version of those contents which is the official version.


Technical data

Parameter description	Parameter value		
Product name	Multichannel Pipette		
Model	SBS-LAB-117	SBS-LAB-118	SBS-LAB-119
Capacity [μ l]	0.5-10	50-300	5-50
Number of pipettes [pcs]	12	12	12
Weight [kg]	0.2	0.2	0.2
Product name	Multichannel Pipette		
Model	SBS-LAB-120	SBS-LAB-121	SBS-LAB-122
Capacity [μ l]	0.5-10	50-300	5-50
Number of pipettes [pcs]	8	8	8
Weight [kg]	0.2	0.2	0.2

1. Product description

This pipette is a general-purpose instrument designed for sampling and dispensing precise liquid volumes. It operates on the air displacement principle and utilizes disposable tips for accuracy and cleanliness.

2. Product overview

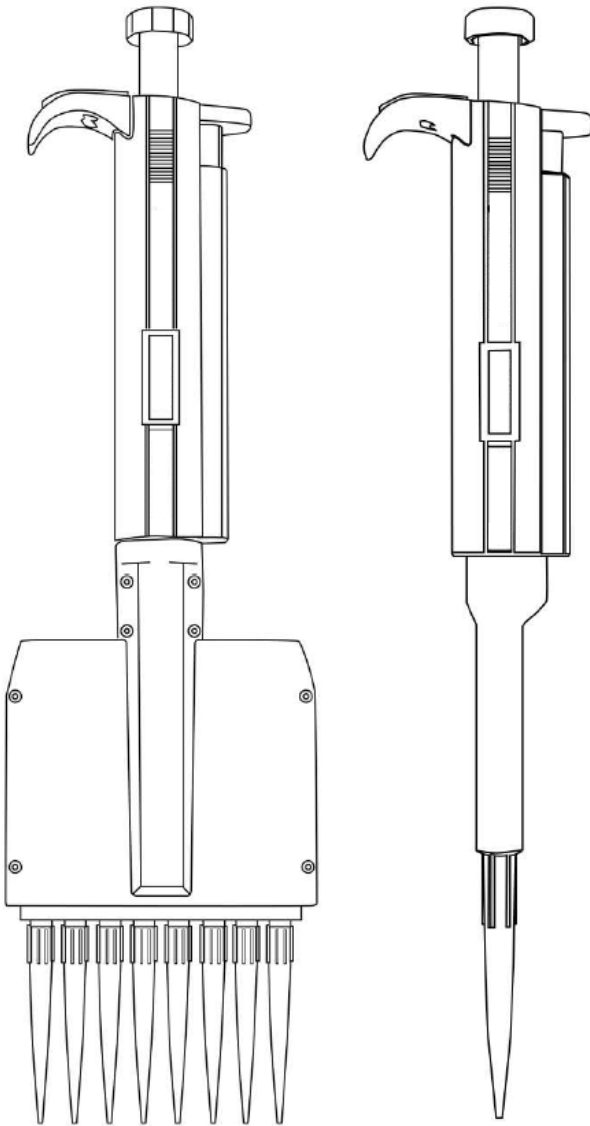3. How to use

Volume setting

The volume of the pipette is clearly shown through the handle grip window. The delivery volume (variable volume pipettes only) is set by turning the thumb button clockwise or anticlockwise. When setting the volume, please make sure that:

The desired delivery volume clicks into place. The digits are completely visible in the display window. The selected volume is within the pipette's specified range.

Note: Using excessive force to turn the push button outside the range may jam the mechanism and damage the pipette.

Sealing and ejecting tips

Before fitting a tip, make sure that the pipette tip cone is clean. Press the cone of the pipette firmly to ensure an airtight seal. The seal is tight when a visible sealing ring forms between the tip and the blank tip cone


Each pipette is equipped with a tip ejector to help eliminate safety hazards related to contamination. To properly eject the tip, press the ejector firmly downwards.

Removing the tip ejector

Please removing the tip ejector collar by following instructions

- Push the tip ejector down. (1)
- Push the opening tool pin between the ejector bar and ejector collar to release the locking mechanism. (2)
- Pull the ejector collar off. (3)
- Pushing the collar locking pin into the ejector bar hole. Keeping the tip ejector pushed down.
- Release the tip ejector and test that it runs smoothly.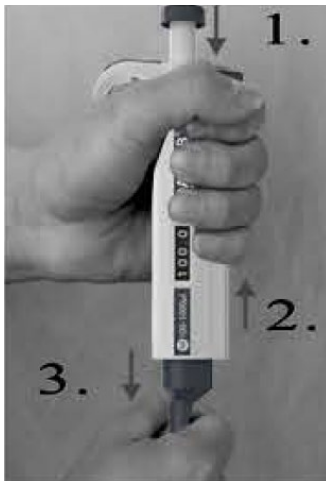

Pipetting techniques

Rotate the operating button to adjust volume setting within the permissible range;

Check the set volume in the display window.

Press the tip tightly on the tip cone

Forward Pipetting:

Press the operating button to the first stop point. **(B)**

Gently immerse the tip 2-3mm below the liquid surface. Slowly release the button to return to the starting position **(A)**. Carefully withdraw the tip from the liquid, touching it against the edge of the container to remove any excess liquid.

After a brief pause, press the button to the second stop point **(C)** (blowout) to fully empty the tip and ensure accurate liquid transfer. Release the button to the starting position **(A)**. If needed, replace the tip and continue pipetting. Press the ejector to remove and discard the tip.

Starting position **(A)**

First stop **(B)**

Second stop **(C)**

Reverse pipetting:

For viscous liquid, do reverse pipetting

Press the operating button to the second stop point (C) and immerse the pipette tip 2-3mm below the liquid surface.

Slowly release the button back to the starting position (A). Carefully withdraw the tip from the liquid, touching it against the edge of the container to remove any excess liquid.

Gently press the button to the first stop point (B) to dispense the liquid. Hold the operating button at the first stop (B) to ensure that any liquid remaining in the tip is not included in the delivery.

Discard the excess liquid along with the tip, or return the remaining liquid to the original container.

Starting position (A)

First stop (B)

Second stop (C)


4. Recalibration:

- place the calibration tool into the holes of the calibration adjustment lock (under the operating button)

- Turn the adjustment lock anticlockwise to decrease and clockwise to increase the volume.
- Repeat performance test procedure from step 1 until the pipetting results are correct.

1. Hold down the tip pipette.
2. Place the tooth of the opening tool between the tip ejector and the tip ejector collar to release the locking mechanism. **(D)**
3. Carefully release the tip ejector and remove the ejector collar.
4. Place the wrench end of the opening tool over the tip cone, turning it anticlockwise. Do not use any other tools. **(E)**
5. Wipe the piston, the O-ring and the tip cone with ethanol and a lint-free cloth

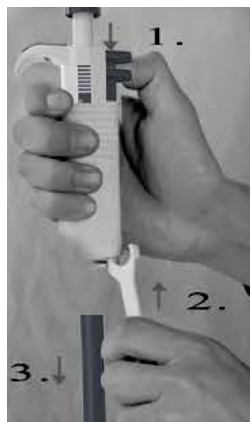**(D)**

Note: Models up to 10ul have a fixed O-ring locked inside the tip cone. Therefore, the O-ring cannot be accessed for maintenance.

6. Before replacing tip cone it is recommended to grease the piston slightly using the silicone grease provided

Note: Excessive use of grease may jam the piston.

7. After reassembling use the pipette (without liquid) several times to make
8. Check the pipette calibration

**(E)**

TROUBLESHOOTING

Problem	Possible cause	Solution
Droplets left inside the tip leakage or pipetted volume too small	Unsuitable tip Non-uniform wetting of the plastic	Use original tips Attach new tip
	Tip incorrectly attached Foreign particles between tip and tip cone Instrument contaminated	Attach firmly Use original tips Clean the tip cone, attach new tip
	Insufficient amount of grease on piston and O-ring	Clean and grease O-ring and piston, clean the tip cone Grease accordingly
	O-ring not correctly	Change the O-ring
	Positioned or damaged Incorrect operation	Follow instructions carefully
	Calibration altered or unsuitable for the liquid	Recalibrate according to instructions
	Instrument damage	Send for service
Push button jammed or moves erratically	Piston contaminated Penetration of solvent vapors	Clean and grease O-ring and piston, clean the tip cone
Pipette blocked; aspirated volume too small	Liquid has penetrated tip cone and dried	Clean and grease O-ring and piston, clean the tip cone
Tip ejector jammed or moves erratically	Tip cone and/or ejector collar contaminated	Clean the tip cone and the ejector collar

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