## USER MANUAL

Bedienungsanleitung  
Instrukcja obsługi  
Návod k použití  
Manuel d'utilisation  
Istruzioni per l'uso  
Manual de instrucciones  
Használati útmutató  
Brugsanvisning  
Käyttöohje  
Gebruiksaanwijzing  
Bruksanvisning  
Instruções de utilização  
Používateľská príručka

---

---

# OVERHEAD STIRRER

DE	Produktname	LABORRÜHRER
EN	Product name	OVERHEAD STIRRER
PL	Nazwa produktu	MIESZADŁO LABORATORYJNE
CZ	Název výrobku	LABORATORNÍ MÍCHADLO
FR	Nom du produit	AGITATEUR DE LABORATOIRE
IT	Nome del prodotto	AGITATORE MECCANICO
ES	Nombre del producto	AGITADOR MECÁNICO
HU	Termék neve	LABORATÓRIUMI KEVERŐ
DA	Produktnavn	ELEKTRISK OMRØRER LABORATORIE
FI	Tuotteen nimi	LABORATORIOSEKOITIN
NL	Productnaam	LABORATORIUMROERDER
NO	Produktnavn	LABORATORIERØRER
SE	Produktnamn	LABORATORIEOMRÖRARE
PT	Nome do produto	AGITADOR MECÂNICO
SK	Názov produktu	LABORATÓRNE MIEŠADLO
DE	Modell	SBS-LAB-126
EN	Product model	
PL	Model produktu	
CZ	Model výrobku	
FR	Modèle	
IT	Modello	
ES	Modelo	
HU	Modell	
DA	Model	
FI	Tuotteen malli	
NL	Productmodel	
NO	Produktmodell	
SE	Produktmodell	
PT	Modelo do produto	
SK	Model	
DE	Hersteller	expondo Polska sp. z o.o. sp. k.
EN	Manufacturer	
PL	Producent	
CZ	Výrobce	
FR	Fabricant	
IT	Produttore	
ES	Fabricante	
HU	Termelő	
DA	Producent	
FI	Valmistaja	
NL	Producent	
NO	Produsent	
SE	Tillverkare	
PT	Fabricante	
SK	Výrobca	
DE	Anschrift des Herstellers	ul. Nowy Kisielin – Innowacyjna 7, 66-002 Zielona Góra   Poland, EU
EN	Manufacturer Address	
PL	Adres producenta	
CZ	Adresa výrobce	
FR	Adresse du fabricant	
IT	Indirizzo del produttore	
ES	Dirección del fabricante	
HU	A gyártó címe	
DA	Producentens adresse	
FI	Valmistajan osoite	
NL	Adres producent	
NO	Produsentens adresse	
SE	Tillverkarens adress	
PT	Endereço do fabricante	
SK	Adresa výrobcu	

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This User Manual has been translated for your convenience using machine translation. Reasonable efforts have been made to provide an accurate translation; however, no automated translation is perfect nor is it intended to replace human translators. The official User Manual is the English version. Any discrepancies or differences created in the translation are not binding and have no legal effect for compliance or enforcement purposes. If any questions arise related to the accuracy of the information contained in the User Manual, please refer to the English version of those contents which is the official version.

## Technical data

Parameter description	Parameter value
Product name	Overhead Stirrer
Model	SBS-LAB-126
Rated voltage [V~] / frequency [Hz]	230/50
Rated power [W]	60
Max. rotation speed [rev/min]	50 – 1200
Dimensions [width x depth x height; mm]	300 x 200 x 650
Weight [kg]	7

### 1. General description

The user manual is designed to assist in the safe and trouble-free use of the device. The product is designed and manufactured in accordance with strict technical guidelines, using state-of-the-art technologies and components. Additionally, it is produced in compliance with the most stringent quality standards.

**DO NOT USE THE DEVICE UNLESS YOU HAVE THOROUGHLY READ AND UNDERSTOOD THIS USER MANUAL.**

To increase the product life of the device and to ensure trouble-free operation, use it in accordance with this user manual and regularly perform maintenance tasks. The technical data and specifications in this user manual are up to date. The manufacturer reserves the right to make changes associated with quality improvement. The device is designed to reduce noise emission risks to a minimum, taking into account technological progress and noise reduction opportunities.

#### Legend

The product satisfies the relevant safety standards.

Read instructions before use.

The product must be recycled.

**WARNING! or CAUTION! or REMEMBER!**

General warning sign

Wear protective goggles when working with harmful substances.

Wear protective gloves when working with harmful substances.

Wear a protective apron when working with harmful substances.

**ATTENTION! Fire hazard!**  
Applies when working with flammable materials.

**WARNING! Danger of poisoning!**  
Applies when working with toxic substances.

**ATTENTION! Electric shock warning!**

Only use indoors.

**PLEASE NOTE! Drawings in this manual are for illustration purposes only and in some details may differ from the actual product.**

## 2. Usage safety

**ATTENTION!** Read all safety warnings and all instructions. Failure to follow the warnings and instructions may result in electric shock, fire and/or serious injury or even death.

The terms "device" or "product" are used in the warnings and instructions to refer to:  
Overhead Stirrer

### 2.1. Electrical safety

- a) The plug must fit the socket. Do not modify the plug in any way. Using original plugs and matching sockets reduces the risk of electric shock.
- b) Avoid touching earthed elements such as pipes, heaters, boilers and refrigerators. There is an increased risk of electric shock if the earthed device is exposed to rain, comes into direct contact with a wet surface or is operating in a damp environment. Water getting into the device increases the risk of damage to the device and of electric shock.
- c) Do not touch the device with wet or damp hands.
- d) Use the cable only for its designated use. Never use it to carry the device or to pull the plug out of a socket. Keep the cable away from heat sources, oil, sharp edges or moving parts. Damaged or tangled cables increase the risk of electric shock.
- e) If using the device in a damp environment cannot be avoided, a residual current device (RCD) should be applied. The use of an RCD reduces the risk of electric shock.
- f) Do not use the device if the power cord is damaged or shows obvious signs of wear. A damaged power cord should be replaced by a qualified electrician or the manufacturer's service centre.
- g) To avoid electric shock, do not immerse the cord, plug or device in water or other liquids. Do not use the device on wet surfaces.

### 2.2. Safety in the workplace

- a) Make sure the workplace is clean and well lit.
- b) If there are any doubts as to the correct operation of the device, contact the manufacturer's support service.
- c) Only the manufacturer's service point may repair the device. Do not attempt any repairs independently!
- d) In case of fire, use a powder or carbon dioxide (CO<sub>2</sub>) fire extinguisher (one intended for use on live electrical devices) to put it out.
- e) Children or unauthorised persons are forbidden to enter a work station.
- f) Use the device in a well-ventilated space.
- g) Regularly inspect the condition of the safety labels. If the labels are illegible, they must be replaced.
- h) Please keep this manual available for future reference. If this device is passed on to a third party, the manual must be passed on with it.

- i) Keep the device away from children and animals.

## 2.3. Personal safety

- a) Do not use the device when tired, ill or under the influence of alcohol, narcotics or medication which can significantly impair the ability to operate the device.
- b) The machine may be operated by physically fit persons who are able to handle the machine, are properly trained, who have reviewed this operating manual and have received training in occupational health and safety.
- c) The machine is not designed to be handled by persons (including children) with limited mental and sensory functions or persons lacking relevant experience and/ or knowledge unless they are supervised by a person responsible for their safety or they have received instruction on how to operate the machine.
- d) When working with the device, use common sense and stay alert.
- e) Use personal protective equipment as required for working with the device, specified in section 1 (Legend).  
The use of correct and approved personal protective equipment reduces the risk of injury if the substance spills or sparkles.
- f) The device is not a toy. Children must be supervised to ensure that they do not play with the device.

## 2.4. Safe device use

- a) Disconnect the device from the power supply before commencement of adjustment, cleaning and maintenance. Such a preventive measure reduces the risk of accidental activation.
- b) Keep the device out of the reach of children.
- c) Device repair or maintenance should be carried out by qualified persons, only using original spare parts. This will ensure safe use.
- d) To ensure the operational integrity of the device, do not remove factory-fitted guards and do not loosen any screws.
- e) Do not touch articulated parts or accessories unless the device has been disconnected from the power source.
- f) Do not move, adjust or rotate the device in the course of work.
- g) Do not leave this appliance unattended while it is in use.
- h) Keep the device away from sources of fire and heat.
- i) It is forbidden to interfere with the structure of the device in order to change its parameters or construction.
- j) Do not overload the device.
- k) Observe if the substance doesn't spill out of the vessel during operation.
- l) Pathogenic materials must be stirred only in closed containers.
- m) When mixing viscous liquids reduce the rotational speed and mix a small volume of liquid only.
- n) Follow biosafety regulations and check the resistance of materials when working with chemical, corrosive, infectious, pathogenic and biological materials.

**ATTENTION! Despite the safe design of the device and its protective features, and despite the use of additional elements protecting the operator, there is still a slight risk of accident or injury when using the device. Stay alert and use common sense when using the device.**

## 3. Use guidelines

The unit is designed for use in laboratories, workshops, etc. It is used for mixing liquids in laboratory vessels. The product is outside the scope of the MDR regulations (EU 2017/745) for medical devices.

**The user is liable for any damage resulting from unintended use of the device.**

### 3.1. Device description

- 1 - Clamp
- 2 - Stand
- 3 - Display
- 4 - Knob for on/off and time adjustment
- 5 - Base

## 3.2. Preparing for use

### APPLIANCE LOCATION

The temperature of environment must not be higher than 40°C and the relative humidity should be less than 85%. Ensure good ventilation in the room in which the device is being used. There should be at least 10 cm distance between each side of the device and the wall or other objects. The device should always be used when positioned on an even, stable, clean, fireproof and dry surface, and be out of the reach of children and persons with limited mental and sensory functions. Position the device such that you always have access to the power plug. The power cord connected to the appliance must be properly grounded and correspond to the technical details on the product label.

## 3.3. Device use

Connect the power supply.

A single press of the knob toggles between speed mode and time mode setting.

Turn the dial to set the desired parameters after selecting the mode (time --: -- continuous mode).

The appliance starts when the knob is pressed for 2 seconds.

The appliance stops when the knob is pressed once during use.

During use, only the speed can be adjusted, but the time cannot be adjusted.

## 3.4. Cleaning and maintenance

- a) Unplug the mains plug and allow the device to cool completely before each cleaning, adjustment or replacement of accessories, or if the device is not being used.
  - Wait for the rotating elements to stop.
- b) Use only non-corrosive cleaners to clean the surface.
- c) After cleaning the device, all parts should be dried completely before using it again.
- d) Store the unit in a dry, cool place, free from moisture and direct exposure to sunlight.

- e) Do not spray the device with a water jet or submerge it in water.
- f) Do not allow water to get inside the device through vents in the housing of the device.
- g) Clean the vents with a brush and compressed air.
- h) The device must be regularly inspected to check its technical efficiency and spot any damage.
- i) Use a soft, damp cloth for cleaning.
- j) Do not use sharp and/or metal objects for cleaning (e.g. a wire brush or a metal spatula) because they may damage the surface material of the appliance.
- k) Do not clean the device with an acidic substance, agents of medical purposes, thinners, fuel, oils or other chemical substances because it may damage the device.

**DISPOSING OF USED DEVICES:**

Do not dispose of this device in municipal waste systems. Hand it over to an electric and electrical device recycling and collection point. Check the symbol on the product, instruction manual and packaging. The plastics used to construct the device can be recycled in accordance with their markings. By choosing to recycle you are making a significant contribution to the protection of our environment. Contact local authorities for information on your local recycling facility.

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