USER MANUAL

Bedienungsanleitung
Instrukcja obsługi
Návod k použití
Manuel d'utilisation
Istruzioni per l'uso
Manual de instrucciones
Használati útmutató
Bruksanvisning
Käyttöohje
Gebruiksaanwijzing
Bruksanvisning
Instruções de utilização
Používateľská príručka
Ръководство за употреба
Οδηγίες χρήσης
Urpute za uporabu
Naudojimo instrukcija
Manual de utilizare
Navodila za uporabo

VORTEX MIXER SET

DE	Produktname	Vortex-Mischer Set
EN	Product name	VORTEX MIXER SET
PL	Nazwa produktu	Mikser wirowy
CZ	Název výrobku	Sada vortexového mixéru
FR	Nom du produit	Ensemble mélangeur vortex
IT	Nome del prodotto	Set mixer vortex
ES	Nombre del producto	Conjunto de mezclador vortex
HU	Termék neve	Vortex keverő készlet
DA	Produktnavn	Vortex mixer sæt
FI	Tuotteen nimi	Vortex-sekoitinsetti
NL	Productnaam	Vortexmixer set
NO	Produktnavn	Vortex mikser sett
SE	Produktnamn	Vortex mixer set
PT	Nome do produto	Conjunto de misturador vórtice
SK	Názov produktu	Sada vortexového miešača
BG	Име на продукта	Комплект вихрового міксера
EL	Όνομα προϊόντος	Σετ αναμεικτήρα δίνης
HR	Naziv proizvoda	Vortex mikser set
LT	Produkto pavadinimas	Vortex maišyklės rinkinys
RO	Numele produsului	Set mixer vortex
SL	Ime izdelka	Set mešalnika za vrtince
DE Modell EN Product model PL Model produktu CZ Model výrobku FR Modèle IT Modello ES Modelo HU Modell DA Model FI Tuotteen malli NL Productmodel NO Produktmodell SE Produktmodell PT Modelo do produto SK Model BG Модел на продукт EL Μοντέλο προϊόντος HR Model proizvoda LT : Gaminio modelis RO : Model de produs SL : Model izdelka		SBS-LAB-136
DE Hersteller EN Manufacturer PL Producent CZ Výrobce FR Fabricant IT Produttore ES Fabricante HU Termelő DA Producent FI Valmistaja NL Producent NO Produsent SE Tillverkare PT Fabricante SK Výrobca BG Производител EL Κατασκευαστής HR Proizvođač LT Gamintojas RO Producător SL Proizvajalec		expondo Polska sp. z o.o. sp. k.
DE Anschrift des Herstellers EN Manufacturer Address PL Adres producenta CZ Adresa výrobce FR Adresse du fabricant IT Indirizzo del produttore ES Dirección del fabricante HU A gyártó címe DA Producentens adresse FI Valmistajan osoite NL Adres producent NO Produsentens adresse SE Tillverkarens adress PT Endereço do fabricante SK Adresa výrobcu BG Адрес на производителя EL : Διεύθυνση κατασκευαστή HR Adresa proizvođača LT Gamintojo adresas RO Adresa producătorului SL Naslov proizvajalca		ul. Nowy Kisielin – Innowacyjna 7, 66-002 Zielona Góra Poland, EU

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This User Manual has been translated for your convenience using machine translation. Reasonable efforts have been made to provide an accurate translation; however, no automated translation is perfect nor is it intended to replace human translators. The official User Manual is the English version. Any discrepancies or differences created in the translation are not binding and have no legal effect for compliance or enforcement purposes. If any questions arise related to the accuracy of the information contained in the User Manual, please refer to the English version of those contents which is the official version.


FOR YOUR SAFETY AND CORRECT OPERATION, PLEASE CAREFULLY READ THIS INSTRUCTION BEFORE USING IT.

NOTE: All of the information reported herein is based on data available at the moment of printing. The factory reserves the right to modify its own products at any moment without notice and incur any sanction. So, it is suggested to always verify possible updates.

1. TECHNICAL DATA

Parameter description	Parameter value
Product name	Vortex Mixer
Model	SBS-LAB-136
Rated voltage [V~] / Frequency [Hz]	230 / 50
Rated power [W]	20
Touch [rpm]	0 – 3000
Run [rpm]	100 – 3000
IP protection class	IP21

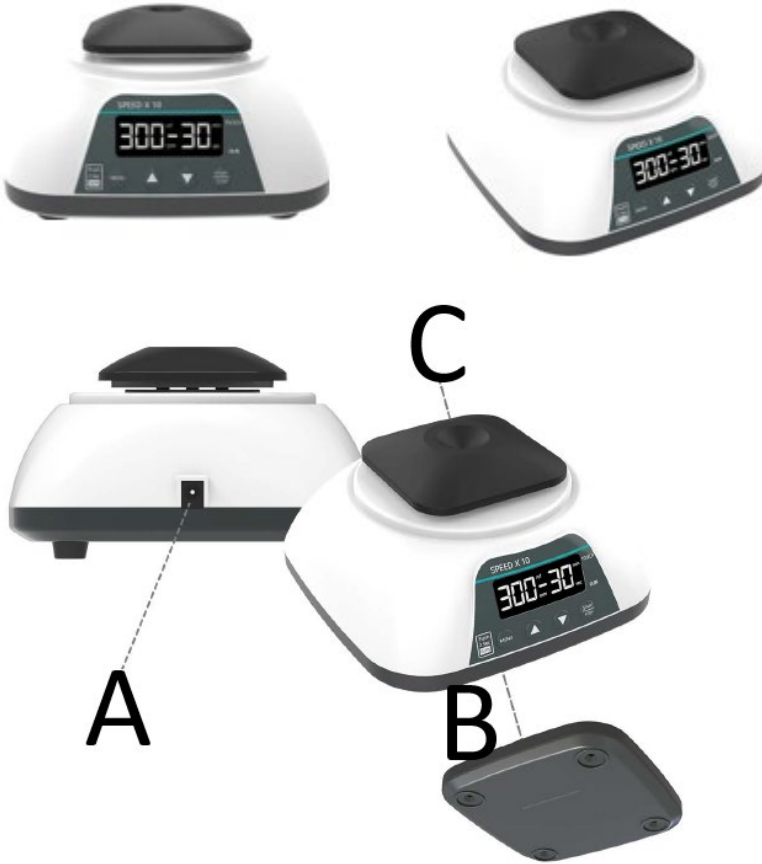2. PRODUCT OVERVIEW

- A DC Power Outlet
- B Zinc aluminum alloy base
- C Module

3. STEPS FOR USAGE

- Plug in
- Non-digital display model: use switch to select mode, use knob to adjust the number of revolutions
- Digital display model: long press for 2 seconds to switch the mode, press the button to adjust the speed
- In press-run mode, you need to press to start

4. Safety

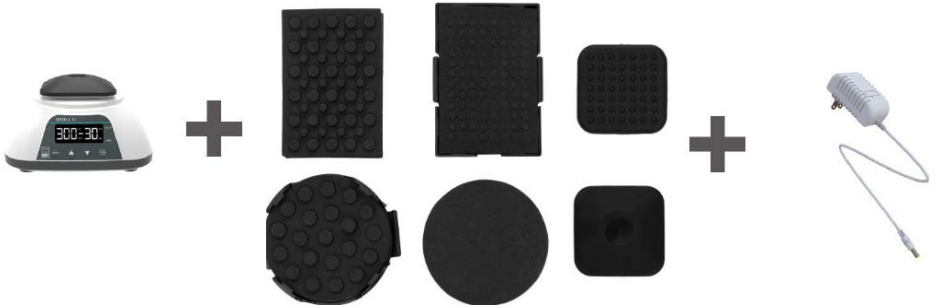
It is strictly forbidden to unplug and plug the power connector and toggle the power button when the hands are wet with liquid!

It is strictly forbidden to unplug the power plug when the device is on!

It is strictly forbidden to maintain and clean the instrument when it is powered on!

It is strictly forbidden to install the instrument on an uneven, swaying and vibrating work surface!

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