## USER MANUAL

Bedienungsanleitung  
Instrukcja obsługi  
Návod k použití  
Manuel d'utilisation  
Istruzioni per l'uso  
Manual de instrucciones  
Használati útmutató  
Bruksanvisning  
Käyttöohje  
Gebruiksaanwijzing  
Bruksanvisning  
Instruções de utilização  
Používateľská príručka  
Ръководство за употреба  
Οδηγίες χρήσης  
Upute za uporabu  
Naudojimo instrukcija  
Manual de utilizare  
Navodila za uporabo

---

---

## MAGNETIC HOTPLATE STIRRER

<b>DE</b>	Produktname	Magnetrührer
<b>EN</b>	Product name	Magnetic Hotplate Stirrer
<b>PL</b>	Nazwa produktu	Mieszadło magnetyczne z podgrzewaniem
<b>CZ</b>	Název výrobku	Magnetické míchačky s ohřevem
<b>FR</b>	Nom du produit	Agitateur magnétique chauffant
<b>IT</b>	Nome del prodotto	Agitatore magnetico riscaldante
<b>ES</b>	Nombre del producto	Agitador magnético con placa calefactora
<b>HU</b>	Termék neve	Mágneses fűtőlap keverő
<b>DA</b>	Produktnavn	Magnetisk varmeplade omrører
<b>FI</b>	Tuotteen nimi	Magneettinen kuumalevyn sekoitin
<b>NL</b>	Productnaam	Magnetische verwarmingsplaatmenger
<b>NO</b>	Produktnavn	Magnetisk varmeplateomrører
<b>SE</b>	Produktnamn	Magnetisk varmplatta omrörare
<b>PT</b>	Nome do produto	Agitador magnético com placa de aquecimento
<b>SK</b>	Názov produktu	Magnetická miešačka s ohrevom
<b>BG</b>	Име на продукта	Магнитна бъркалка с нагревателна плоча
<b>EL</b>	Όνομα προϊόντος	Μαγνητικός αναδευτήρας με θερμαινόμενη πλάκα
<b>HR</b>	Naziv proizvoda	Magnetna mešalna plošča s segrevanjem
<b>LT</b>	Produkto pavadinimas	Magnetinė kaitinimo plokštelė su maišykle
<b>RO</b>	Numele produsului	Agitator magnetic cu plită de încălzire
<b>SL</b>	Ime izdelka	Magnetni mešalnik s grelno ploščo
<b>DE</b> Modell   <b>EN</b> Product model   <b>PL</b> Model produktu   <b>CZ</b> Model výrobku   <b>FR</b> Modèle   <b>IT</b> Modello   <b>ES</b> Modelo   <b>HU</b> Modell   <b>DA</b> Model   <b>FI</b> Tuotteen malli   <b>NL</b> Productmodel   <b>NO</b> Produktmodell   <b>SE</b> Produktmodell   <b>PT</b> Modelo do produto   <b>SK</b> Model   <b>BG</b> Модел на продукт   <b>EL</b> Μοντέλο προϊόντος   <b>HR</b> Model proizvoda   <b>LT</b> : Gaminio modelis   <b>RO</b> : Model de produs   <b>SL</b> : Model izdelka		SBS-LAB-137
<b>DE</b> Hersteller   <b>EN</b> Manufacturer   <b>PL</b> Producent   <b>CZ</b> Výrobce   <b>FR</b> Fabricant   <b>IT</b> Produttore   <b>ES</b> Fabricante   <b>HU</b> Termelő   <b>DA</b> Producent   <b>FI</b> Valmistaja   <b>NL</b> Producent   <b>NO</b> Produsent   <b>SE</b> Tillverkare   <b>PT</b> Fabricante   <b>SK</b> Výrobca   <b>BG</b> Производител   <b>EL</b> Κατασκευαστής   <b>HR</b> Proizvođač   <b>LT</b> Gamintojas   <b>RO</b> Producător   <b>SL</b> Proizvajalec		expondo Polska sp. z o.o. sp. k.
<b>DE</b> Anschrift des Herstellers   <b>EN</b> Manufacturer Address   <b>PL</b> Adres producenta   <b>CZ</b> Adresa výrobce   <b>FR</b> Adresse du fabricant   <b>IT</b> Indirizzo del produttore   <b>ES</b> Dirección del fabricante   <b>HU</b> A gyártó címe   <b>DA</b> Producentens adresse   <b>FI</b> Valmistajan osoite   <b>NL</b> Adres producent   <b>NO</b> Produsentens adresse   <b>SE</b> Tillverkarens adress   <b>PT</b> Endereço do fabricante   <b>SK</b> Adresa výrobcu   <b>BG</b> Адрес на производителя   <b>EL</b> : Διεύθυνση κατασκευαστή   <b>HR</b> Adresa proizvođača   <b>LT</b> Gamintojo adresas   <b>RO</b> Adresa producătorului   <b>SL</b> Naslov proizvajalca		ul. Nowy Kisielin – Innowacyjna 7, 66-002 Zielona Góra   Poland, EU

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This User Manual has been translated for your convenience using machine translation. Reasonable efforts have been made to provide an accurate translation; however, no automated translation is perfect nor is it intended to replace human translators. The official User Manual is the English version. Any discrepancies or differences created in the translation are not binding and have no legal effect for compliance or enforcement purposes. If any questions arise related to the accuracy of the information contained in the User Manual, please refer to the English version of those contents which is the official version.

## Technical data

Parameter description	Parameter value
Product name	Hotplate Magnetic Stirrer
Model	SBS-LAB-137
Rated voltage [V~] / Frequency [Hz]	230 / 50
Rated power [W]	230
Working plate [mm]	φ 135
Plate temperature [°C]	280
Rotation speed [rpm]	50-1000
Dimensions [Width x Depth x Height; mm]	225 x 155 x 410
Weight [kg]	1.5

### Product brief

The heating magnetic stirrer adopts international advanced design concept and manufacturing technology. It has the advantages of smooth and beautiful shape, compact and firm structure, high versatility and simple operability. It is a rare instrument for scientific research and experiments. This heating magnetic stirrer is mainly used to stir and heat low viscosity liquid or solid-liquid mixture.

### Instrument using environment

- > Indoor use
- >Altitude: ≤2000m
- >The working temperature range of the instrument is +5 °C~+40 °C
- > The applicable temperature range of the instrument is ≤80%
- > There is no vibration and airflow that affects performance around.

## Product introduction

1. Bracket
2. Heating plate
3. Cross clamp
4. Temperature probe
5. Operation panel
6. Speed control knob: Rotate the knob to set the speed. Press the knob lightly to start/pause.
7. Temperature control knob: Rotate the knob to set the temperature. Press the knob lightly to start/pause.

## Correct usage

### Application

> It is mainly used for stirring and heating liquid or solid-liquid mixture with low viscosity.

### Usage

> It is widely used as an experimental tool in colleges and universities, environmental protection, health, epidemic prevention, chemical industry, tap water, medical treatment and other units.

### Steps for usage

> Turn on the power

> Rotate the left knob to adjust the temperature and the right knob to adjust the speed.

> Press any knob to start.

(Note: This model without timing function)

### Safety Precautions

> It is strictly forbidden to unplug and plug the power connector and toggle the power button when the hands are wet with liquid!

> It is strictly forbidden to unplug the power plug when the instrument is live!

> It is strictly forbidden to maintain and clean the instrument when it is powered on!

> It is strictly forbidden to install the instrument on an uneven, swaying and vibrating work surface!

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