## USER MANUAL

Bedienungsanleitung  
Instrukcja obsługi  
Návod k použití  
Manuel d'utilisation  
Istruzioni per l'uso  
Manual de instrucciones  
Használati útmutató  
Bruksanvisning  
Käyttöohje  
Gebruiksaanwijzing  
Bruksanvisning  
Instruções de utilização  
Používateľská príručka  
Ръководство за употреба  
Οδηγίες χρήσης  
Upute za uporabu  
Naudojimo instrukcija  
Manual de utilizare  
Navodila za uporabo

---

---

## MULTI-TUBE VORTEX MIXER

<b>DE</b>	Produktname	Vortex Mixer Multitube
<b>EN</b>	Product name	Multi-Tube Vortex Mixer
<b>PL</b>	Nazwa produktu	Mieszadło wirowe wielorurowe
<b>CZ</b>	Název výrobku	Vícezkumavkový vortex mixer
<b>FR</b>	Nom du produit	Mélangeur vortex multi-tube
<b>IT</b>	Nome del prodotto	Miscelatore vortex multi-tubo
<b>ES</b>	Nombre del producto	Mezclador vortex multi-tubo
<b>HU</b>	Termék neve	Többcsöves vortex keverő
<b>DA</b>	Produktnavn	Multi-tube vortex mixer
<b>FI</b>	Tuotteen nimi	Moniputkinen vortex-mikseri
<b>NL</b>	Productnaam	Multi-buis vortex mixer
<b>NO</b>	Produktnavn	Multi-rør vortex mikser
<b>SE</b>	Produktnamn	Multi-rör vortexmixer
<b>PT</b>	Nome do produto	Misturador vortex multi-tubos
<b>SK</b>	Názov produktu	Viaczkumavkový vortex miešač
<b>BG</b>	Име на продукта	Вихров миксер с много тръби
<b>EL</b>	Όνομα προϊόντος	Πολυσωλήνας vortex αναδευτήρας
<b>HR</b>	Naziv proizvoda	Vortex mikser za više epruveta
<b>LT</b>	Produkto pavadinimas	Daugiavamzdis „Vortex“ maišytuvas
<b>RO</b>	Numele produsului	Mixer vortex multi-tuburi
<b>SL</b>	Ime izdelka	Multi-cevni vortex mešalnik
<b>DE</b> Modell   <b>EN</b> Product model   <b>PL</b> Model produktu   <b>CZ</b> Model výrobku   <b>FR</b> Modèle   <b>IT</b> Modello   <b>ES</b> Modelo   <b>HU</b> Modell   <b>DA</b> Model   <b>FI</b> Tuotteen malli   <b>NL</b> Productmodel   <b>NO</b> Produktmodell   <b>SE</b> Produktmodell   <b>PT</b> Modelo do produto   <b>SK</b> Model   <b>BG</b> Модел на продукт   <b>EL</b> Μοντέλο προϊόντος   <b>HR</b> Model proizvoda   <b>LT</b> : Gaminio modelis   <b>RO</b> : Model de produs   <b>SL</b> : Model izdelka		SBS-LAB-138
<b>DE</b> Hersteller   <b>EN</b> Manufacturer   <b>PL</b> Producent   <b>CZ</b> Výrobce   <b>FR</b> Fabricant   <b>IT</b> Produttore   <b>ES</b> Fabricante   <b>HU</b> Termelő   <b>DA</b> Producent   <b>FI</b> Valmistaja   <b>NL</b> Producent   <b>NO</b> Produsent   <b>SE</b> Tillverkare   <b>PT</b> Fabricante   <b>SK</b> Výrobca   <b>BG</b> Производител   <b>EL</b> Κατασκευαστής   <b>HR</b> Proizvođač   <b>LT</b> Gamintojas   <b>RO</b> Producător   <b>SL</b> Proizvajalec		expondo Polska sp. z o.o. sp. k.
<b>DE</b> Anschrift des Herstellers   <b>EN</b> Manufacturer Address   <b>PL</b> Adres producenta   <b>CZ</b> Adresa výrobce   <b>FR</b> Adresse du fabricant   <b>IT</b> Indirizzo del produttore   <b>ES</b> Dirección del fabricante   <b>HU</b> A gyártó címe   <b>DA</b> Producentens adresse   <b>FI</b> Valmistajan osoite   <b>NL</b> Adres producent   <b>NO</b> Produsentens adresse   <b>SE</b> Tillverkarens adress   <b>PT</b> Endereço do fabricante   <b>SK</b> Adresa výrobcu   <b>BG</b> Адрес на производителя   <b>EL</b> : Διεύθυνση κατασκευαστή   <b>HR</b> Adresa proizvođača   <b>LT</b> Gamintojo adresas   <b>RO</b> Adresa producătorului   <b>SL</b> Naslov proizvajalca		ul. Nowy Kisielin – Innowacyjna 7, 66-002 Zielona Góra   Poland, EU

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This User Manual has been translated for your convenience using machine translation. Reasonable efforts have been made to provide an accurate translation; however, no automated translation is perfect nor is it intended to replace human translators. The official User Manual is the English version. Any discrepancies or differences created in the translation are not binding and have no legal effect for compliance or enforcement purposes. If any questions arise related to the accuracy of the information contained in the User Manual, please refer to the English version of those contents which is the official version.

## Technical data

Parameter description	Parameter value
Product name	Multi-Tube Vortex Mixer
Model	SBS-LAB-138
Rated voltage [V~] / Frequency [Hz]	230 / 50
Rated power [W]	100
Timer duration	1min – 24h
Ambient temperature [°C]	5-40
Rotation speed [rpm]	500-3000
Dimensions [Width x Depth x Height; mm]	445 x 282 x 330
Weight [kg]	14.85

### Product brief

The multi-tube vortex mixer uses brushless moto and microcomputer control technology.

The unique knob operation mode is simple and easy to use, and by changing different tube fixation sponges, a variety of commonly used tubes can be mixed and cultured. It is suitable for the fields of biotechnology, microbiology and medica analysis.

### Instrument using environment

- > Indoor use
- > Altitude: ≤2000m
- > The working temperature range of the instrument is +5 °C~+40 °C
- > The applicable temperature range of the instrument is ≤80%
- > There is no vibration and airflow that affects performance around.

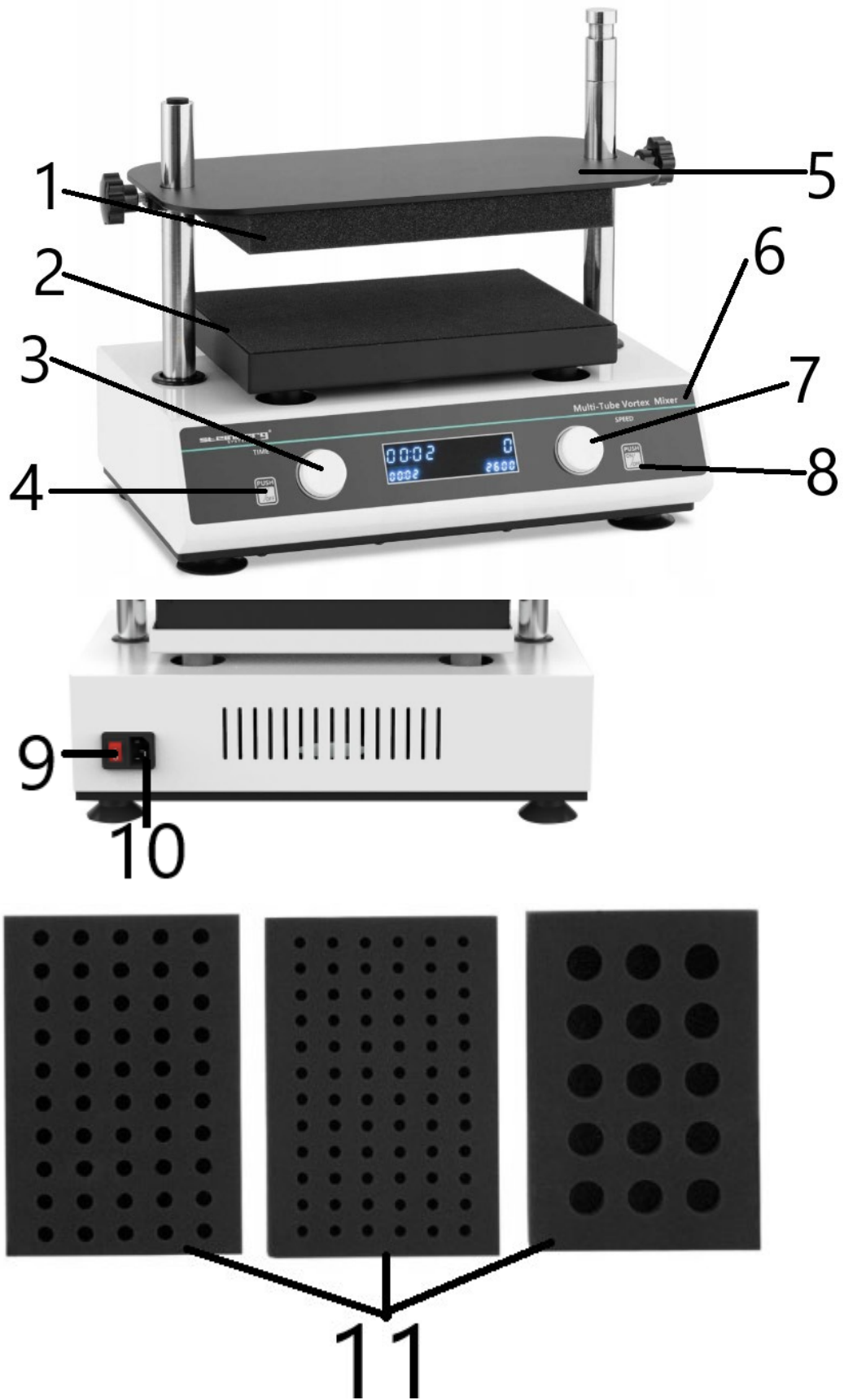
### **Important Safety Operation Information**

- Please make sure to understand the use of the instrument and read the instruction manual carefully before operating the instrument safely.
- Please read the safety instructions and instructions below carefully to understand all the precautions to avoid possible accidental damage to the instrument operation.

### **Safety**

- The following basic safety precautions must be observed during all operation, maintenance, and repair of this instrument. Failure to comply with these measures or other warnings in this manual may affect the intended scope of use of the protection box provided by the instrument.
- Please read the manual carefully and train the operators before operation.
- Operators should not open the maintenance instrument by themselves, which will cause you to lose the warranty qualification, may also lead to electric shock and other dangers. If you need to repair, please contact us first.
- Before connecting the power supply, ensure that the voltage of the power supply is the same as the voltage required by the instrument. And ensure that the rated load of the power socket is not less than the requirements of the instrument.
- If the power cable is damaged, it must be replaced. Replace the power cable with a power cable of the same type and specifications.
- Do not press anything on the power line when the instrument is in use. Do not put the power cord in the place where people walk.
- Be sure to hold the plug in hand when plugging in and out the power cord. When inserting the plug, ensure that the plug is fully inserted into the socket. Do not pull the power cord when pulling out the plug.
- • The instrument should be placed in a place with low humidity, less dust and away from water sources and avoid direct sunlight and strong light. Indoor ventilation should be maintained, without corrosive gas or strong magnetic interference, and away from heating, furnace, and other heat sources. Do not place the instrument in a damp or dusty area.
- • When stopping work, the power should be turned off. When the instrument is not used for a long time, the power plug should be unplugged, and the instrument should be covered with a soft cloth or plastic paper to prevent dust from entering.
- • In the following cases, the power plug of the instrument should be immediately removed from the power socket and contact the supplier:
  - There is liquid spilled into the instrument.
  - The instrument is exposed to rain or water.
  - The instrument does not work properly, especially with abnormal sound or smell.
  - The instrument falls or the shell is damaged.
  - The function of the instrument has been significantly changed.

## Product introduction

- 1 Foam mat of upper tray
- 2 Vortex shaking platform
- 3 Time adjustment knob
- 4 Start/Stop button
- 5 Fixed tray above test tube
- 6 Operation panel
- 7 Speed adjustment knob
- 8 Touch running button
- 9 Power switch
- 10 Power interface and fuse
- 11 Foam test tube holder

## Operation guide

### Speed and timing setting

A. **Turn on the power switch**, the screen will appear as shown in the picture, and the instrument will enter the initialization, accompanied by a "beep" prompt sound.

B. **The system enters the parameter configuration mode.** B. The system enters the parameter configuration mode. The right side of the screen displays the currently set speed, and the left side displays the set countdown time and flashes a number. The time separator on the left side does not flash.

C. **At this time, press the knob button** to move the flashing position to the left and select the number to be adjusted. Turn the knob clockwise or counterclockwise to adjust the value.

D. **When the value is less than 1 minute, '00:00' is displayed.** Timing function is disabled, enabling the start counting function. The system must be stopped manually by the user.

### Run, stop function

A. **After the parameter setting is complete, press the X button** and the system will play a 'Beep' prompt tone. At this time, the system starts to run.

B. **When the speed is adjusted to the preset value**, the system starts the countdown, the system sends out two 'Beep' prompts, and the time separator in the window starts to blink.

C. **When the speed is adjusted to the preset value**, the system starts the countdown, the system sends out two 'Beep' prompts, and the time separator in the right window starts to blink.

- After the countdown, the system stops running. The system plays a 'Beep' prompt to return to the configuration mode.

D. **Long press the X button to force the operation to stop.**

**Touch running mode**

A. **When you hold down the button**, the system will play a 'Beep' prompt tone, and the system will start to run.

- After the speed is adjusted to the preset value, the system will play a 'Beep' prompt tone twice, and the system will time and display. At this time, the time separator in the right window will blink. Release the button and the system stops running mode.

**NOTICE:**

- When the time parameter is set to '00:00', the instrument continues to run in countdown mode.
- After the end of the timing, the instrument will stop automatically and give a 'Beep' prompt tone and enter the configuration mode automatically.
- The value adjustment of the left window is controlled by the left knob, and the value adjustment of the right window is controlled by the right knob.
- When using the timing function, the maximum time is 99 hours and 69 minutes, and the system continues to run after the maximum time, but the timing stops.
- Non-professionals should not handle faults with power on.

## TROUBLESHOOTING

Problem	Possible cause	Solution
The display does not light up after turning	The power is not connected.	Check the power supply
	The switch is damaged.	Change the switch
There is no vortex in the liquid after operation	The speed setting is not reasonable.	Reset the speed
	The upper tray is pressed too loose.	Adjust the EVA Foam position and press again
Large instrument vibration	Asymmetry of mixing position	Place the mixture symmetrically
The displayed speed does not match the actual speed.	Controller failure.	Contact the supplier
The screen on the left shows 8888.	Motor drive circuit failure.	Contact the supplier
The knob does not work.	Broken button.	Contact the supplier
Excessive noise	The handwheel is not tightened.	Tighten the left and right handwheels
	The parts are loose, etc.	Contact the supplier

**Additional Notes:**

A. **Check the power cord connection** to ensure that the external power supply meets the requirements.

B. **Restart the instrument** after replacing the fuse. If the problem is still not resolved, please contact the supplier.

C. **Restart the instrument**; if the problem still exists, please contact the supplier.

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