## USER MANUAL

Bedienungsanleitung  
Instrukcja obsługi  
Návod k použití  
Manuel d'utilisation  
Istruzioni per l'uso  
Manual de instrucciones  
Használati útmutató  
Bruksanvisning  
Käyttöohje  
Gebruiksaanwijzing  
Bruksanvisning  
Instruções de utilização  
Používateľská príručka  
Ръководство за употреба  
Οδηγίες χρήσης  
Upute za uporabu  
Naudojimo instrukcija  
Manual de utilizare  
Navodila za uporabo

---

---

## ROLLER MIXER

<b>DE</b>	Produktname	Rollenmischer
<b>EN</b>	Product name	Roller mixer
<b>PL</b>	Nazwa produktu	Mieszadło laboratoryjne
<b>CZ</b>	Název výrobku	Válečkový míchač
<b>FR</b>	Nom du produit	Mélangeur à rouleaux
<b>IT</b>	Nome del prodotto	Miscelatore a rulli
<b>ES</b>	Nombre del producto	Mezclador de rodillos
<b>HU</b>	Termék neve	Görgős keverő
<b>DA</b>	Produktnavn	Rulleblander
<b>FI</b>	Tuotteen nimi	Rullasekoitin
<b>NL</b>	Productnaam	Rollenmenger
<b>NO</b>	Produktnavn	Rullemikser
<b>SE</b>	Produktnamn	Rullblandare
<b>PT</b>	Nome do produto	Misturador de rolos
<b>SK</b>	Názov produktu	Valčekový miešač
<b>BG</b>	Име на продукта	Валцов смесител
<b>EL</b>	Όνομα προϊόντος	Μίκτης κυλίνδρων
<b>HR</b>	Naziv proizvoda	Mješač valjaka
<b>LT</b>	Produktu pavadinimas	Volelių maišytuvas
<b>RO</b>	Numele produsului	Mixer cu role
<b>SL</b>	Ime izdelka	Valjčni mešalnik
<b>DE</b> Modell   <b>EN</b> Product model   <b>PL</b> Model produktu   <b>CZ</b> Model výrobku   <b>FR</b> Modèle   <b>IT</b> Modello   <b>ES</b> Modelo   <b>HU</b> Modell   <b>DA</b> Model   <b>FI</b> Tuotteen malli   <b>NL</b> Productmodel   <b>NO</b> Produktmodell   <b>SE</b> Produktmodell   <b>PT</b> Modelo do produto   <b>SK</b> Model   <b>BG</b> Модел на продукт   <b>EL</b> Μοντέλο προϊόντος   <b>HR</b> Model proizvoda   <b>LT</b> : Gaminio modelis   <b>RO</b> : Model de produs   <b>SL</b> : Model izdelka		SBS-LAB-140
<b>DE</b> Hersteller   <b>EN</b> Manufacturer   <b>PL</b> Producent   <b>CZ</b> Výrobce   <b>FR</b> Fabricant   <b>IT</b> Produttore   <b>ES</b> Fabricante   <b>HU</b> Termelő   <b>DA</b> Producent   <b>FI</b> Valmistaja   <b>NL</b> Producent   <b>NO</b> Produsent   <b>SE</b> Tillverkare   <b>PT</b> Fabricante   <b>SK</b> Výrobca   <b>BG</b> Производител   <b>EL</b> Κατασκευαστής   <b>HR</b> Proizvođač   <b>LT</b> Gamintojas   <b>RO</b> Producător   <b>SL</b> Proizvajalec		expondo Polska sp. z o.o. sp. k.
<b>DE</b> Anschrift des Herstellers   <b>EN</b> Manufacturer Address   <b>PL</b> Adres producenta   <b>CZ</b> Adresa výrobce   <b>FR</b> Adresse du fabricant   <b>IT</b> Indirizzo del produttore   <b>ES</b> Dirección del fabricante   <b>HU</b> A gyártó címe   <b>DA</b> Producentens adresse   <b>FI</b> Valmistajan osoite   <b>NL</b> Adres producent   <b>NO</b> Produsentens adresse   <b>SE</b> Tillverkarens adress   <b>PT</b> Endereço do fabricante   <b>SK</b> Adresa výrobcu   <b>BG</b> Адрес на производителя   <b>EL</b> : Διεύθυνση κατασκευαστή   <b>HR</b> Adresa proizvođača   <b>LT</b> Gamintojo adresas   <b>RO</b> Adresa producătorului   <b>SL</b> Naslov proizvajalca		ul. Nowy Kisielin – Innowacyjna 7, 66-002 Zielona Góra   Poland, EU

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**DE**

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This User Manual has been translated for your convenience using machine translation. Reasonable efforts have been made to provide an accurate translation; however, no automated translation is perfect nor is it intended to replace human translators. The official User Manual is the English version. Any discrepancies or differences created in the translation are not binding and have no legal effect for compliance or enforcement purposes. If any questions arise related to the accuracy of the information contained in the User Manual, please refer to the English version of those contents which is the official version.

## Technical data

Parameter description	Parameter value
Product name	Roller mixer
Model	SBS-LAB-140
Rated voltage [V~] / Frequency [Hz]	230 / 50
Rated power [W]	25
Rotation speed [rpm]	20-80
Number of rollers	6
Dimensions [Width x Depth x Height; mm]	445 x 260 x 103
Weight [kg]	2.45

### Instrument using environment

- > Indoor use
- > Altitude:  $\leq 2000\text{m}$
- > The instrument is suitable for ambient temperature range of  $+5^{\circ}\text{C} \sim +40^{\circ}\text{C}$
- > Instrument operation suitable for relative humidity range of  $\leq 80\%$
- > No vibrations and airflow affecting performance are present
- > There are no conductive dust, explosive gases and corrosive gases in the surrounding air
- > Keep a safety distance of 10cm between the front and back left and right when the instrument is working

## Product brief

This mixer is designed and manufactured according to international advanced design concepts and technologies.

The mixer is highly efficient and gentle and is suitable for different experiments such as blood analysis and sample mixing.

## Product overview

## Product introduction

1. Operating panel
2. Rollers (made of silicone)
3. Switch

## Correct usage

### Application

> This instrument is **suitable for a variety of mixing operations** in different applications, including:

- Electron microscopy analysis
- Blood cell suspensions
- Homogeneous solid blood and plasma mixing
- Tissue specimen processing
- Washing precipitates
- Disintegration sheet analysis
- Extraction processes
- Dissolution of slow reactions
- Rotary sealing operations

### Usage

> Widely used in scientific research institutions, medical schools, health care and other units, molecular biology, virology, microbiology, pathology, immunology and other laboratories.

### Steps for usage

- > Plug in the power supply and turn on the switch.
- > Place the sample tube/reagent bottle on the roller.
- > MENU key to switch time (min) /speed (rpm), up and down keys to adjust parameters.
- > Press the START/STOP keys to start/stop the device.

## Safety Precautions

- > It is strictly forbidden to unplug the power connector and toggle the switching power button when your hands are stained with liquid!
- > It is strictly forbidden to unplug the power plug when the instrument is live!
- > It is strictly forbidden to maintain and scrub the instrument when it is powered on!
- > It is strictly forbidden to install the instrument on an uneven, shaky and shaky work surface!

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