USER MANUAL

Bedienungsanleitung
Instrukcja obsługi
Návod k použití
Manuel d'utilisation
Istruzioni per l'uso
Manual de instrucciones
Használati útmutató
Bruksanvisning
Käyttöohje
Gebruiksaanwijzing
Bruksanvisning
Instruções de utilização
Používateľská príručka
Ръководство за употреба
Οδηγίες χρήσης
Upute za uporabu
Naudojimo instrukcija
Manual de utilizare
Navodila za uporabo

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DE	Produktname	Zentrifuge
EN	Product name	Mini Centrifuge
PL	Nazwa produktu	Mała wirówka laboratoryjna
CZ	Název výrobku	Mini centrifuga
FR	Nom du produit	Mini-centrifuge
IT	Nome del prodotto	Mini centrifuga
ES	Nombre del producto	Mini centrifugadora
HU	Termék neve	Mini centrifuga
DA	Produktnavn	Mini-centrifuge
FI	Tuotteen nimi	Minisentrifugi
NL	Productnaam	Mini-centrifuge
NO	Produktnavn	Mini-sentrifuge
SE	Produktnamn	Mini-centrifug
PT	Nome do produto	Mini centrifuga
SK	Názov produktu	Mini centrifúga
BG	Име на продукта	Мини центрофуга
EL	Όνομα προϊόντος	Μίνι φυγόκεντρος
HR	Naziv proizvoda	Mini centrifuga
LT	Produkto pavadinimas	Mini centrifuga
RO	Numele produsului	Mini centrifugă
SL	Ime izdelka	Mini centrifuga
DE Modell EN Product model PL Model produktu CZ Model výrobku FR Modèle IT Modello ES Modelo HU Modell DA Model FI Tuotteen malli NL Productmodel NO Produktmodell SE Produktmodell PT Modelo do produto SK Model BG Модел на продукт EL Μοντέλο προϊόντος HR Model proizvoda LT : Gaminio modelis RO : Model de produs SL : Model izdelka		SBS-LAB-141
DE Hersteller EN Manufacturer PL Producent CZ Výrobce FR Fabricant IT Produttore ES Fabricante HU Termelő DA Producent FI Valmistaja NL Producent NO Produsent SE Tillverkare PT Fabricante SK Výrobca BG Производител EL Κατασκευαστής HR Proizvođač LT Gamintojas RO Producător SL Proizvajalec		expondo Polska sp. z o.o. sp. k.
DE Anschrift des Herstellers EN Manufacturer Address PL Adres producenta CZ Adresa výrobce FR Adresse du fabricant IT Indirizzo del produttore ES Dirección del fabricante HU A gyártó címe DA Producentens adresse FI Valmistajan osoite NL Adres producent NO Produsentens adresse SE Tillverkarens adress PT Endereço do fabricante SK Adresa výrobcu BG Адрес на производителя EL : Διεύθυνση κατασκευαστή HR Adresa proizvođača LT Gamintojo adresas RO Adresa producătorului SL Naslov proizvajalca		ul. Nowy Kisielin – Innowacyjna 7, 66-002 Zielona Góra Poland, EU

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This User Manual has been translated for your convenience using machine translation. Reasonable efforts have been made to provide an accurate translation; however, no automated translation is perfect nor is it intended to replace human translators. The official User Manual is the English version. Any discrepancies or differences created in the translation are not binding and have no legal effect for compliance or enforcement purposes. If any questions arise related to the accuracy of the information contained in the User Manual, please refer to the English version of those contents which is the official version.

Technical data

Parameter description	Parameter value
Product name	Mini centrifuge
Model	SBS-LAB-141
Rated voltage [V~] / frequency [Hz]	100-240~/50
Rated power [W]	60
Max capacity	Centrifuge tube 15ml*8 Blood tube 5/10ml*8
Speed range [rpm]	500-4000
Increment [rpm]	100
Timing range [s]	30-3590
Max RCF (xg)	1967
Semidiameter of rotor [cm]	11
Max operating altitude [m]	52000
Allowable ambient temperature [°C]	5-40
Dimensions [width x depth x height; cm]	29 x 40 x 20
Weight [kg]	8

1. General description

The user manual is designed to assist in the safe and trouble-free use of the device. The product is designed and manufactured in accordance with strict technical guidelines, using state-of-the-art technologies and components. Additionally, it is produced in compliance with the most stringent quality standards.

DO NOT USE THE DEVICE UNLESS YOU HAVE THOROUGHLY READ AND UNDERSTOOD THIS USER MANUAL.

To increase the product life of the device and to ensure trouble-free operation, use it in accordance with this user manual and regularly perform maintenance tasks. The technical data and specifications in this user manual are up to date. The manufacturer reserves the right to make changes associated with quality improvement. The device is designed to reduce noise emission risks to a minimum, taking into account technological progress and noise reduction opportunities.

Legend

	The product satisfies the relevant safety standards.
	Read instructions before use.
	The product must be recycled.
	WARNING! or CAUTION! or REMEMBER! Applicable to the given situation. (general warning sign)

	Wear protective goggles.
	Wear protective gloves.
	Wear protective clothing.
	Wear a protective apron.
	Read the instructions.
	ATTENTION! Electric shock warning!
	WARNING! Toxic substances, danger of poisoning!
	Only use indoors.

PLEASE NOTE! Drawings in this manual are for illustration purposes only and in some details may differ from the actual product.

2. Usage safety

ATTENTION! Read all safety warnings and all instructions. Failure to follow the warnings and instructions may result in electric shock, fire and/or serious injury or even death.

The terms "device" or "product" are used in the warnings and instructions to refer to:
Mini centrifuge

2.1. Electrical safety

- The plug must fit the socket. Do not modify the plug in any way. Using original plugs and matching sockets reduces the risk of electric shock.
- Avoid touching earthed elements such as pipes, heaters, boilers and refrigerators. There is an increased risk of electric shock if the earthed device is exposed to rain, comes into direct contact with a wet surface or is operating in a damp environment. Water getting into the device increases the risk of damage to the device and of electric shock.
- Do not touch the device with wet or damp hands.
- Use the cable only for its designated use. Never use it to carry the device or to pull the plug out of a socket. Keep the cable away from heat sources, oil, sharp edges or moving parts. Damaged or tangled cables increase the risk of electric shock.
- Do not use in very humid environments or in the direct vicinity of water tanks.
- Before the first use, please check whether the main voltage type and current comply with the indicated data on the type plate.

2.2. Safety in the workplace

- Make sure the workplace is clean and well lit. A messy or poorly lit workplace may lead to accidents. Try to think ahead, observe what is going on and use common sense when working with the device.

- b) If you are unsure about whether the product is operating correctly or if you find damage, please contact the manufacturer's service centre.
- c) Only the manufacturer's service centre may make repairs to the device. Do not attempt to make repairs yourself!
- d) In case of fire, use a powder or carbon dioxide (CO₂) fire extinguisher (one intended for use on live electrical devices) to put it out.
- e) Children or unauthorised persons are forbidden to enter a work station. A distraction may result in loss of control over the device.
- f) Use the device in a well-ventilated space.
- g) Regularly inspect the condition of the safety labels. If the labels are illegible, they must be replaced.
- h) Please keep this manual available for future reference. If this device is passed on to a third party, the manual must be passed on with it.
- i) Keep packaging elements and small assembly parts in a place not available to children.
- j) Keep the device away from children and animals.
- k) If this device is used together with another equipment, the remaining instructions for use shall also be followed.

Remember! When using the device, protect children and other bystanders.

2.3. Personal safety

- a) Do not use the device when tired, ill or under the influence of alcohol, narcotics or medication which can significantly impair the ability to operate the device.
- b) The device can be handled only by physically fit persons who are capable of handling it, properly trained, familiar with this manual and trained within the scope of occupational health and safety.
- c) When working with the device, use common sense and stay alert. Temporary loss of concentration while using the device may lead to serious injuries.
- d) Use personal protective equipment as required for working with the device, specified in section 1 "Legend". The use of correct and approved personal protective equipment reduces the risk of injury.
- e) To prevent the device from accidentally switching on, make sure the switch is on the OFF position before connecting to a power source.
- f) Do not overestimate your abilities. When using the device, keep your balance and remain stable at all times. This will ensure better control over the device in unexpected situations.
- g) The device is not a toy. Children must be supervised to ensure that they do not play with the device.
- h) Do not put your hands or other items inside the device while it is in use!

2.4. Safe device use

- a) Do not use the device if the "ON/OFF" switch does not function properly (does not switch the device on and off). Devices which cannot be switched on and off using the "ON/OFF" switch are hazardous, should not be operated and must be repaired.
- b) Make sure the plug is disconnected from the socket before attempting any adjustments, accessory replacements or before putting the device aside. Such precautions will reduce the risk of accidentally activating the device.
- c) When not in use, store in a safe place, away from children and people not familiar with the device who have not read the user manual. The device may pose a hazard in the hands of inexperienced users.
- d) Keep the device in perfect technical condition. Before each use check for general damage, especially check moving components for cracked parts or elements, and for any other conditions which may impact the safe operation of the device. If damage is discovered, hand over the device for repair before use.
- e) Keep the device out of the reach of children.
- f) Device repair or maintenance should be carried out by qualified persons, only using original spare parts. This will ensure safe use.
- g) To ensure the operational integrity of the device, do not remove factory-fitted guards and do not loosen any screws.
- h) Do not move, adjust or rotate the device in the course of work.

- i) Do not leave this appliance unattended while it is in use.
- j) Clean the device regularly to prevent stubborn grime from accumulating.
- k) The device is not a toy. Cleaning and maintenance may not be carried out by children without supervision by an adult person.
- l) It is forbidden to interfere with the structure of the device in order to change its parameters or construction.
- m) Keep the device away from sources of fire and heat.
- n) Do not overload the device.
- o) Do not cover the ventilation openings!
- p) Pathogenic materials may only be mixed in closed vessels.
- q) Do not cover the device while it is in operation.
- r) Keep the device away from strong magnetic fields.
- s) When mixing viscous liquids, it is recommended to reduce the rotation speed and mix a small volume of liquid.
- t) Follow biological safety regulations and check the vessel's resistance to given substances when working with chemicals, corrosive, infectious, pathogenic and biological substances.

ATTENTION! Despite the safe design of the device and its protective features, and despite the use of additional elements protecting the operator, there is still a slight risk of accident or injury when using the device. Stay alert and use common sense when using the device.

3. Use guidelines

The product separates samples in various tube sizes and microplates, with adjustable time, speed, and temperature settings.

The user is liable for any damage resulting from unintended use of the device.

3.1. Device description

Number	Description
1	Dual E-lock
2	Rotor
3	External circulation heat emission hole
4	Emergency switch

3.2. Device use

1. Plug the product in and turn the switch on.
2. Press the OPEN button to open the lid. Insert the casing and the sample, ensuring the sample is balanced and placed symmetrically.
3. Close the lid. Press the MENU button to adjust the time, speed, and safe temperature.
4. Press START/STOP to start.
5. Press the P button to memorize multiple groups of modes and recall the common mode to run.

Note:

- Make sure there is no vibration and airflow that affects the performance of the device.
- Keep a safety distance of 10cm before and after device use.

3.3. Package content

Part	Quantity	Unit
Centrifuge	1	Pc
Power cord	1	Pc
Adapters for 2/5/10 ml test tubes	8	Pc
Holder for 15ml test tubes	8	Pc

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