## USER MANUAL

Bedienungsanleitung  
Instrukcja obsługi  
Návod k použití  
Manuel d'utilisation  
Istruzioni per l'uso  
Manual de instrucciones  
Használati útmutató  
Bruksanvisning  
Käyttöohje  
Gebruiksaanwijzing  
Bruksanvisning  
Instruções de utilização  
Používateľská príručka  
Ръководство за употреба  
Οδηγίες χρήσης  
Upute za uporabu  
Naudojimo instrukcija  
Manual de utilizare  
Navodila za uporabo

---

---

## HOT PLATE

<b>DE</b>	Produktname	Wasserbad mit Magnetrührer
<b>EN</b>	Product name	Water bath with magnetic stirrer
<b>PL</b>	Nazwa produktu	łaźnia wodna z mieszadłem magnetycznym
<b>CZ</b>	Název výrobku	Vodní lázeň s magnetickým míchadlem
<b>FR</b>	Nom du produit	Bain-marie avec agitateur magnétique
<b>IT</b>	Nome del prodotto	Bagno d'acqua con agitatore magnetico
<b>ES</b>	Nombre del producto	Baño de agua con agitador magnético
<b>HU</b>	Termék neve	Vízfürdő mágneses keverővel
<b>DA</b>	Produktnavn	Vandbad med magnetomrører
<b>FI</b>	Tuotteen nimi	Vesikylpy magneettisekoittimella
<b>NL</b>	Productnaam	Waterbad met magnetische roerder
<b>NO</b>	Produktnavn	Vannbad med magnetisk omrører
<b>SE</b>	Produktnamn	Vattenbad med magnetomrörare
<b>PT</b>	Nome do produto	Banho-maria com agitador magnético
<b>SK</b>	Názov produktu	Vodný kúpeľ s magnetickým miešačom
<b>BG</b>	Име на продукта	Водна баня с магнитна бъркалка
<b>EL</b>	Όνομα προϊόντος	Λουτρό νερού με μαγνητικό αναδευτήρα
<b>HR</b>	Naziv proizvoda	Vodena kupelj s magnetnim miješalom
<b>LT</b>	Produkto pavadinimas	Vandens vonia su magnetine maišykle
<b>RO</b>	Numele produsului	Baie de apă cu agitator magnetic
<b>SL</b>	Ime izdelka	Vodna kopel z magnetnim mešalom
<b>DE</b> Modell   <b>EN</b> Product model   <b>PL</b> Model produktu   <b>CZ</b> Model výrobku   <b>FR</b> Modèle   <b>IT</b> Modello   <b>ES</b> Modelo   <b>HU</b> Modell   <b>DA</b> Model   <b>FI</b> Tuotteen malli   <b>NL</b> Productmodel   <b>NO</b> Produktmodell   <b>SE</b> Produktmodell   <b>PT</b> Modelo do produto   <b>SK</b> Model   <b>BG</b> Модел на продукт   <b>EL</b> Μοντέλο προϊόντος   <b>HR</b> Model proizvoda   <b>LT</b> : Gaminio modelis   <b>RO</b> : Model de produs   <b>SL</b> : Model izdelka		SBS-LAB-142
<b>DE</b> Hersteller   <b>EN</b> Manufacturer   <b>PL</b> Producent   <b>CZ</b> Výrobce   <b>FR</b> Fabricant   <b>IT</b> Produttore   <b>ES</b> Fabricante   <b>HU</b> Termelő   <b>DA</b> Producent   <b>FI</b> Valmistaja   <b>NL</b> Producent   <b>NO</b> Produsent   <b>SE</b> Tillverkare   <b>PT</b> Fabricante   <b>SK</b> Výrobca   <b>BG</b> Производител   <b>EL</b> Κατασκευαστής   <b>HR</b> Proizvođač   <b>LT</b> Gamintojas   <b>RO</b> Producător   <b>SL</b> Proizvajalec		expondo Polska sp. z o.o. sp. k.
<b>DE</b> Anschrift des Herstellers   <b>EN</b> Manufacturer Address   <b>PL</b> Adres producenta   <b>CZ</b> Adresa výrobce   <b>FR</b> Adresse du fabricant   <b>IT</b> Indirizzo del produttore   <b>ES</b> Dirección del fabricante   <b>HU</b> A gyártó címe   <b>DA</b> Producentens adresse   <b>FI</b> Valmistajan osoite   <b>NL</b> Adres producent   <b>NO</b> Produsentens adresse   <b>SE</b> Tillverkarens adress   <b>PT</b> Endereço do fabricante   <b>SK</b> Adresa výrobcu   <b>BG</b> Адрес на производителя   <b>EL</b> : Διεύθυνση κατασκευαστή   <b>HR</b> Adresa proizvođača   <b>LT</b> Gamintojo adresas   <b>RO</b> Adresa producătorului   <b>SL</b> Naslov proizvajalca		ul. Nowy Kisielin – Innowacyjna 7, 66-002 Zielona Góra   Poland, EU

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This User Manual has been translated for your convenience using machine translation. Reasonable efforts have been made to provide an accurate translation; however, no automated translation is perfect nor is it intended to replace human translators. The official User Manual is the English version. Any discrepancies or differences created in the translation are not binding and have no legal effect for compliance or enforcement purposes. If any questions arise related to the accuracy of the information contained in the User Manual, please refer to the English version of those contents which is the official version.

## Technical data

Parameter description	Parameter value
Product name	Water bath with magnetic stirrer
Model	SBS-LAB-142
Rated voltage [V~] / Frequency [Hz]	230 / 50
Rated power [W]	800
Speed range [rpm]	100 - 1200
Temperature range [°C]	0 – 100
Dimensions [Width x Depth x Height; mm]	265 x 260 x 330
Weight [kg]	2.75

### Instrument Usage Environment:

- **Indoor use only**
- **Altitude:** Must be  $\leq 2000$  meters above sea level
- **Working temperature range:**  $+5^{\circ}\text{C}$  to  $+40^{\circ}\text{C}$
- **Applicable humidity range:**  $\leq 80\%$
- Ensure there is **no vibration** or **airflow** around that could affect the instrument's performance.

# EN

## Product brief

The **Water Bath with Magnetic Stirrer** utilizes advanced international design concepts and manufacturing technology. It boasts a sleek, aesthetically pleasing appearance with a compact, firm structure that ensures high versatility and ease of operation. This makes it a valuable, high-quality instrument for scientific research experiments.

### Key Features:

- Smooth and elegant design
- Compact and durable structure
- Highly versatile
- Easy to operate

### Applications:

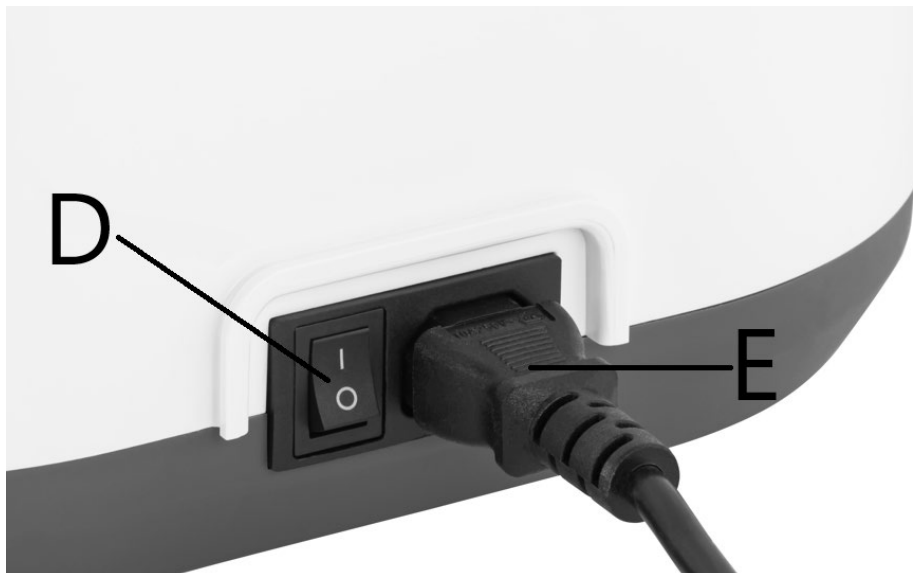
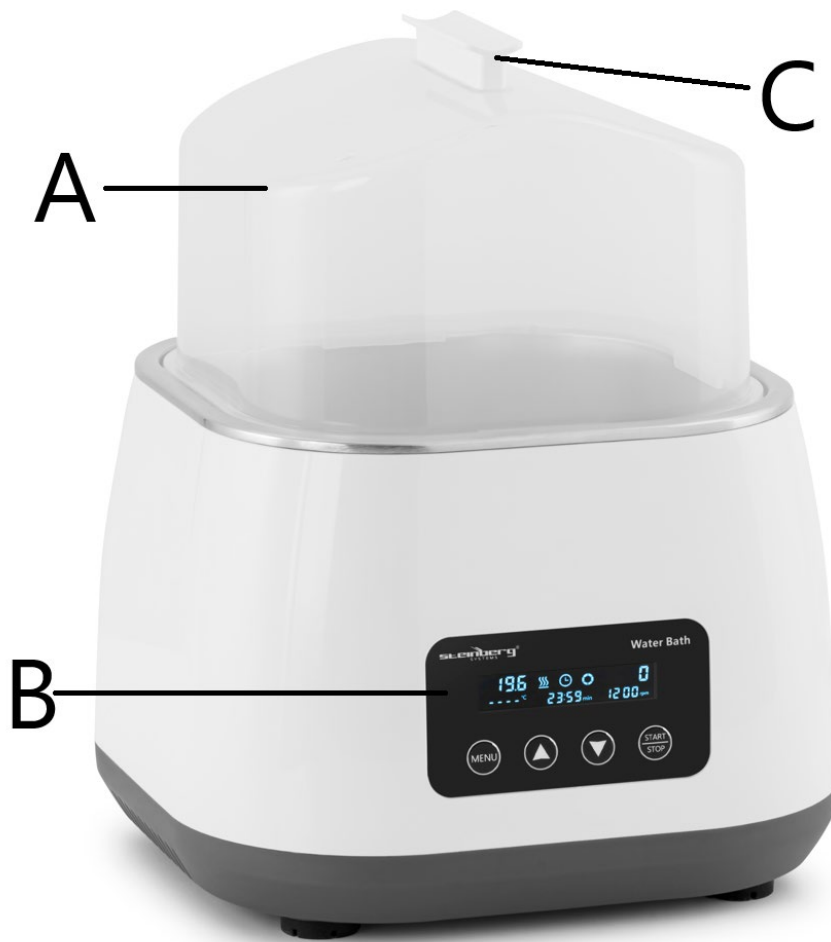
This instrument is widely employed in various laboratory and research settings for tasks such as:

- Drying
- Concentration
- Distillation
- Impregnation of chemical reagents
- Impregnation of medical and biological products

### Ideal for Use In:

- Biology
- Genetics
- Aquatic products
- Environmental protection
- Medical and sanitation fields
- Biochemical laboratories
- Analysis rooms
- Educational institutions
- Scientific research

Product introduction

- A. Transparent tall cover
- B. Operation panel
- C. Safety handle
- D. Power switch
- E. Power cord

## Correct usage

### Application:

The instrument is primarily used for:

- **Drying**
- **Concentrating**
- **Distilling**
- **Impregnating chemical reagents**
- **Impregnating medical products**
- **Heating and stirring biological products**

### Usage:

It is widely employed in various fields, including:

- **Universities**
- **Environmental protection**
- **Health and epidemic prevention**
- **Chemical industry**
- **Tap water testing**
- **Medical research and experimental tools**

## Steps for usage

### 1. Prepare the Water Bath:

- Add clean water to the water bath.
- Place the stirrer bar in the water bath and set it on the steaming rack.

### 2. Power On:

- Plug in the water bath and turn it on.

### 3. Place the Container:

- Insert the thermostatic container with the stirrer bar into the water bath and close the lid.

### 4. Set Time and Temperature:

- Use the **MENU key** to toggle between time and temperature settings.
- Adjust the parameters using the **up or down** buttons.

## EN

### 5. Start the Water Bath:

- Press the **START/STOP** key to start or stop the water bath's heating function.

### 6. After Use:

- Remove the container from the water bath.
- Turn off the power and unplug the device.
- Pour out the water inside the container and ensure it is clean and dry.

### Safety Precautions:

1. **Do not unplug or plug the power connector** or toggle the power button when your hands are wet with liquid.
2. **Do not unplug the power plug** while the instrument is still live or powered on.
3. **Do not perform maintenance or cleaning** of the instrument while it is powered on.
4. **Do not install the instrument** on an uneven, swaying, or vibrating work surface.

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