USER MANUAL

Bedienungsanleitung
Instrukcja obsługi
Návod k použití
Manuel d'utilisation
Istruzioni per l'uso
Manual de instrucciones
Használati útmutató
Bruksanvisning
Käyttöohje
Gebruiksaanwijzing
Bruksanvisning
Instruções de utilização
Používateľská príručka
Ръководство за употреба
Οδηγίες χρήσης
Upute za uporabu
Naudojimo instrukcija
Manual de utilizare
Navodila za uporabo

PRECISION BALANCE

DE	Produktname	Präzisionswaage
EN	Product name	Precision balance
PL	Nazwa produktu	Precyzyjna waga
CZ	Název výrobku	Přesné vyvážení
FR	Nom du produit	Balance de précision
IT	Nome del prodotto	Bilancia di precisione
ES	Nombre del producto	Balanza de precisión
HU	Termék neve	Precíziós mérleg
DA	Produktnavn	Præcisionsvægt
FI	Tuotteen nimi	Tarkka tasapaino
NL	Productnaam	Precisiebalans
NO	Produktnavn	Presisjonsbalanse
SE	Produktnamn	Precisionsbalans
PT	Nome do produto	Balança de precisão
SK	Názov produktu	Presné vyváženie
BG	Име на продукта	Прецизен баланс
EL	Όνομα προϊόντος	Ισορροπία ακριβείας
HR	Naziv proizvoda	Precizna vaga
LT	Produkto pavadinimas	Tikslus balansas
RO	Numele produsului	Balanta de precizie
SL	Ime izdelka	Natančna tehtnica
DE Modell EN Product model PL Model produktu CZ Model výrobku FR Modèle IT Modello ES Modelo HU Modell DA Model FI Tuotteen malli NL Productmodel NO Produktmodell SE Produktmodell PT Modelo do produto SK Model BG Модел на продукт EL Μοντέλο προϊόντος HR Model proizvoda LT : Gaminio modelis RO : Model de produs SL : Model izdelka		SBS-LW-1030 SBS-LW-1040 SBS-LW-1050 SBS-LW-1060 SBS-LW-1070
DE Hersteller EN Manufacturer PL Producent CZ Výrobce FR Fabricant IT Produttore ES Fabricante HU Termelő DA Producent FI Valmistaja NL Producent NO Produsent SE Tillverkare PT Fabricante SK Výrobca BG Производител EL Κατασκευαστής HR Proizvođač LT Gamintojas RO Producător SL Proizvajalec		expondo Polska sp. z o.o. sp. k.
DE Anschrift des Herstellers EN Manufacturer Address PL Adres producenta CZ Adresa výrobce FR Adresse du fabricant IT Indirizzo del produttore ES Dirección del fabricante HU A gyártó címe DA Producentens adresse FI Valmistajan osoite NL Adres producent NO Produsentens adresse SE Tillverkarens adress PT Endereço do fabricante SK Adresa výrobcu BG Адрес на производителя EL : Διεύθυνση κατασκευαστή HR Adresa proizvođača LT Gamintojo adresas RO Adresa producătorului SL Naslov proizvajalca		ul. Nowy Kisielin – Innowacyjna 7, 66-002 Zielona Góra Poland, EU

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This User Manual has been translated using machine translation. We have made every effort to ensure the translation is accurate, but please note that automated translations are not perfect and are not meant to replace human translators. The official version of the User Manual is in English. Any differences between the translated version and the original English are not legally binding. If you have any questions about the accuracy of the translation, please refer to the English version, which is the official reference. More language versions are available upon request via info@expondo.com.

Technical data

Parameter description	Parameter value				
Product name	Precision balance				
Model	SBS-LW-1030	SBS-LW-1040	SBS-LW-1050	SBS-LW-1060	SBS-LW-1070
Max load [kg]	3	3	6	10	20
Min load [g]	0.2	2			
Dimensions (Width x Length x Height) (mm)	205 x 300 x 95	210 x 315 x 95	205 x 300 x 95	210 x 310 x 95	205 x 300 x 95
Weight [kg]	1.6	1.65			
Division [g]	0.01	0.1			
Power adapter [V/A]	12 / 1.0				
Battery	6 x AA Batteries				
Protection class	II				

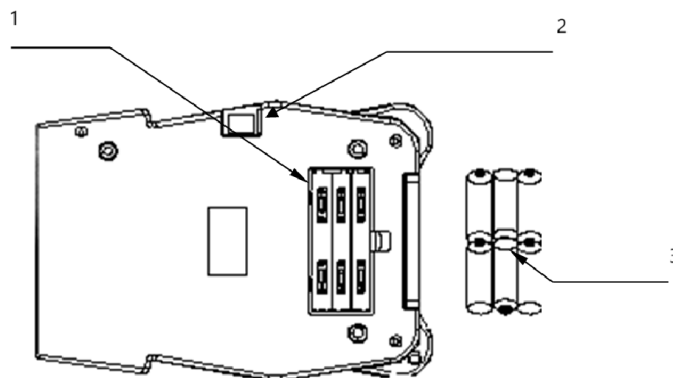
Purpose

The product is used to measure the weight of objects with high accuracy and precision, typically within a milligram or even finer increments. Its purpose is essential in environments where precise measurements are crucial, such as in laboratories, research facilities, pharmaceuticals, and industrial applications.

Usage

Preparation

Batteries

- 1- Battery compartment
- 2- Switch power
- 3- 6 x AA Batteties

Setting up the scale/balance

CAUTION

- Avoid installing the scale/balance in direct sunlight, as it may lead to discoloration or malfunction.
- Ensure the scale/balance is set up on a sturdy weighing table so that it remains level.
- The scale/balance will not provide accurate measurements if it is not properly leveled.
- For accurate weighing, always place the scale/balance on a firm, flat, and level surface.

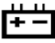
Weighing

Before Weighing

Whenever possible, allow the balance to warm up for 10 minutes after turning it on. This ensures the balance will function properly and provide accurate measurements.

NOTE

- Error messages
 - Err-0: Overload
 - Err-Z: Exceed zero track range when turning on the power
 - Err- S: The scale is unstable when turning on the power
 - : Low battery

Weighing Procedures

1. Turn on the Balance

Press  to power on the balance. When the power is turned on, all display segments will briefly appear on the LCD, and after a few seconds, "0" will be displayed.

2. Select the Weighing Unit

Press  to choose the desired weighing unit (e.g., "g", "ct", etc.). The selected unit will be shown alongside the weight value.

3. Start Weighing

• Without a container

- 1) Ensure the display reads "0". If not, press  to reset it to "0".
- 2) Place the object on the weighing platform.
- 3) Once the reading stabilizes, the stability indicator will appear.

• With a container

- 1) Place the empty container on the platform.
- 2) Wait for the stability indicator to appear, then press  to reset the display to "0".
- 3) Add the objects to be weighed into the container. When the reading becomes stable, the stability indicator will appear.

Calibration

When to Calibrate

Calibration may be necessary when the scale/balance is first installed or moved a significant distance, as the

gravitational force affecting the weight may vary between locations. Additionally, over time and with regular use, mechanical deviations can occur, making recalibration necessary.

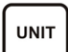
How to Calibrate

1. Prepare for Calibration

Before entering calibration mode, locate the switch key on the bottom of the scale and turn it on.

Then, press  to power on the scale and allow it to warm up for 1 minute.

2. Enter Calibration Mode

- 1) Press  to power on the scale.
- 2) Once the display shows flashing digits followed by "-----", press . The display will show "CAL". Press  again, and the display will show "X0000" (X is the flashing digit).
- 3) Use the  key to move the flashing digit to the right, and press  or  to increase the value. (X represents the calibration weight, which should be set based on your specific calibration requirements).

3. Calibration with the Weight

- 1) After selecting the appropriate calibration weight, place the weight on the platform.
- 2) Wait for the stability indicator to appear, then press .
- 3) The display will return to weighing mode, indicating that calibration is complete.

4. Verify Calibration

- 1) Turn off the scale by pressing , then turn it back on.
- 2) Place the calibration weight on the platform to verify the accuracy of the weighing.
- 3) If the weight is not accurate, repeat steps 2-4.

Counting Procedures

1. Turn on the Scale

Press  to power on the scale. Wait until "0" appears on the display.

2. Start the Count Procedure

If necessary, press  to reset the display to "0".

3. Place Samples on the Pan

Place a predetermined number of sample items (either 10, 20, 50, or 100 pieces) on the pan. The

weight of the samples will be displayed. Press  to enter PCS (piece counting) mode. Then press , and the display will show "P=XX pcs". Press  again to select the sample size (10, 20, 50, or 100).

4. Confirm the Sample Size

Press  to confirm. The display will show "XX pcs", indicating the sample size you selected. The scale will store the sample size, and you can now remove the samples if you want the display to return to "0".

5. Return to Weighing Mode

Press  again to return to the standard weighing mode.

Function Settings

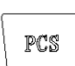
- Press  to turn on the scale. The display will show flashing digits, followed by "-----". Press , and the display will show "b xxx BAU" (baud rate). Use the  key to select from 1200, 2400, 4800, or 9600 (default: 9600). Press  to enter Parity mode. The display will show "P xxx PAR". Use  to select from odd, even, or none:

- Odd: 7 data bits with odd parity
- Even: 7 data bits with even parity
- None: 8 data bits without parity (default: none)

Press  to enter Print mode. The display will show "n xxx Mod". Use  to choose from:

- OFF: Serial data output disabled
- KEY: Data prints on-demand when  is pressed (default)
- STB: Data prints automatically when the scale stabilizes
- Con: Continuous printing Press  again to return to weighing mode.

- Press  to turn on the scale. The display will show flashing digits followed by "-----". Press  to enter unit selection mode. The display will show "ON X" (where X stands for units like g, oz, lb, dwt, etc.). Press  to toggle between ON (activate the selected unit) or OFF (deactivate the unit). Press  to select and confirm the unit mode, then press  again to return to weighing mode.

- Switch the power to "ON" and press  to turn on the scale. The display will show "F-E-9 to 0" followed by "-----". Press  to enter Zero Track settings. The display will show "Zrd x.x d". Use

Press **UNIT** to select the Zero Track value (0.0d to 2.0d). Press **ZERO** to enter Zero Track Time settings, displayed as "Zrt x.x sec". Use **UNIT** to select the time (0.5 to 4.0 seconds). Press **ZERO** to set Zero value, displayed as "rzd x d". Use **UNIT** to choose between 1d, 3d, 5d, or 8d.

Press **ZERO** to enter Filter Range (FIL) settings. Use **UNIT** to select the filter range (0-3). Press **ZERO** to enter the zero range adjustment for power-on. Use **UNIT** to select percentage values (PZR 1, PZR 2, PZR 3, etc.). Press **ZERO** to set the Zero range, then use **UNIT** to select percentage values for zero range (KZR 3, 4, 5, 10).

Press **ZERO** to enter TARE range settings. Use **UNIT** to select percentage (KTR 50 or KTR 100). Press **ZERO** to enter overload protection settings. Use **UNIT** to select overload values (0, OVR-2, OVR-5, OVR-10).

Press **ZERO** to set the beep function. Use **UNIT** to toggle between "bz ON" (activate) or "bz OFF" (deactivate). Press **ZERO** to unite the TARE and ZERO functions. Use **UNIT** to switch between "ZT ON" (activated) and "ZT OFF" (deactivated).

Press **ZERO** to enter Auto-Off settings. The display will show "A xxx". Use **UNIT** to select the auto-off time (3, 5, 10 minutes, or OFF). Press **ZERO** to adjust the backlight. The display will show "L X". Use **UNIT** to select:

- OFF: Disable backlight
- ON: Backlight always on
- Au: Auto backlight mode

4. Press **ZERO** to return to weighing mode.

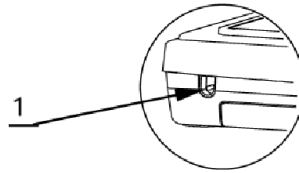
Disposing of Used Devices

Do not dispose of this device in municipal waste systems. Hand it over to an electric and electrical device recycling and collection point. Check the symbol on the product, instruction manual and packaging. The plastics used to construct the device can be recycled in accordance with their markings. By choosing to recycle you are making a significant contribution to the protection of our environment.

Contact local authorities for information on your local recycling facility.

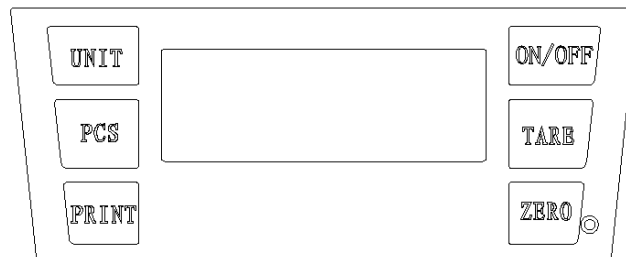
Parts

Back side

1- AC adaptor socket

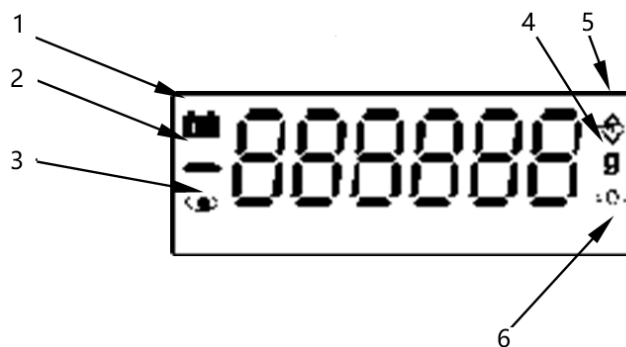
Top

Keypad Functions

1.  : Turns the power on or off.
2.  Resets the scale to zero; the zero-setting range is 4% of the total capacity.
3.  Deducts the weight of the container from the total weight.
4.  Prints the data currently displayed.
5.  Toggles between different weighing units.
6.  Activates the counting function.

Display

- 1- Low battery indicator
- 2- Negative value indicator
- 3- STABLE indicator (when the reading is stable)
 - 4- Weighing unit
 - 5- TARE indicator
 - 6- ZERO indicator

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