USER MANUAL

Bedienungsanleitung
Instrukcja obsługi
Návod k použití
Manuel d'utilisation
Istruzioni per l'uso
Manual de instrucciones
Használati útmutató
Bruksanvisning
Käyttöohje
Gebruiksaanwijzing
Bruksanvisning
Instruções de utilização
Používateľská príručka
Ръководство за употреба
Οδηγίες χρήσης
Upute za uporabu
Naudojimo instrukcija
Manual de utilizare
Navodila za uporabo

MICROSCOPE

DE	Produktname	Mikroskop
EN	Product name	Microscope
PL	Nazwa produktu	Mikroskop
CZ	Název výrobku	Mikroskop
FR	Nom du produit	Microscope
IT	Nome del prodotto	Microscopio
ES	Nombre del producto	Microscopio
HU	Termék neve	Mikroszkóp
DA	Produktnavn	Mikroskop
FI	Tuotteen nimi	Mikroskooppi
NL	Productnaam	Microscoop
NO	Produktnavn	Mikroskop
SE	Produktnamn	Mikroskop
PT	Nome do produto	Microscópio
SK	Názov produktu	Mikroskop
BG	Име на продукта	Микроскоп
EL	Όνομα προϊόντος	Μικροσκόπιο
HR	Naziv proizvoda	Mikroskop
LT	Produkto pavadinimas	Mikroskopas
RO	Numele produsului	Microscop
SL	Ime izdelka	mikroskop
DE Modell EN Product model PL Model produktu CZ Model výrobku FR Modèle IT Modello ES Modelo HU Modell DA Model FI Tuotteen malli NL Productmodel NO Produktmodell SE Produktmodell PT Modelo do produto SK Model BG Модел на продукт EL Μοντέλο προϊόντος HR Model proizvoda LT : Gaminio modelis RO : Model de produs SL : Model izdelka		SBS-MK-8
DE Hersteller EN Manufacturer PL Producent CZ Výrobce FR Fabricant IT Produttore ES Fabricante HU Termelő DA Producent FI Valmistaja NL Producent NO Produsent SE Tillverkare PT Fabricante SK Výrobca BG Производител EL Κατασκευαστής HR Proizvođač LT Gamintojas RO Producător SL Proizvajalec		expondo Polska sp. z o.o. sp. k.
DE Anschrift des Herstellers EN Manufacturer Address PL Adres producenta CZ Adresa výrobce FR Adresse du fabricant IT Indirizzo del produttore ES Dirección del fabricante HU A gyártó címe DA Producentens adresse FI Valmistajan osoite NL Adres producent NO Produsentens adresse SE Tillverkarens adress PT Endereço do fabricante SK Adresa výrobcu BG Адрес на производителя EL : Διεύθυνση κατασκευαστή HR Adresa proizvođača LT Gamintojo adresas RO Adresa producătorului SL Naslov proizvajalca		ul. Nowy Kisielin – Innowacyjna 7, 66-002 Zielona Góra Poland, EU

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This User Manual has been translated using machine translation. We have made every effort to ensure the translation is accurate, but please note that automated translations are not perfect and are not meant to replace human translators. The official version of the User Manual is in English. Any differences between the translated version and the original English are not legally binding. If you have any questions about the accuracy of the translation, please refer to the English version, which is the official reference. More language versions are available upon request via info@expondo.com.

Technical data

Parameter description	Parameter value
Product name	Microscope
Model	SBS-MK-8
Voltage [V]	5
Magnification ratio	1600X
Dimensions (Width x Length x Height) (mm)	230 x 180 x 290
Weight [kg]	1.05
Protection class	II

Purpose

The product is used to magnify small objects or details that are too small to be seen clearly with the naked eye. By enlarging these objects, the product allows users to observe fine structures and details, making it an essential tool in scientific research, biology, medicine, material science, and various other fields.

Product overview

This digital microscope features continuous zoom and a wide range of viewing options, offering magnification levels from 1x to 1600x. It ensures you can always find the ideal magnification for any observation need.

Installation

Works with Windows (Windows 10 Example)

1. **Check for Windows Camera:** Ensure your PC has the "Windows Camera" app installed. If not, download it from the Windows/Microsoft Store or search for it on Google.
2. **Power On the Microscope:** Turn on the microscope and connect it to your PC with the USB cable.
3. **Select PC Camera Mode:** On the microscope, choose "PC Camera" when prompted (options include "Mass Storage," "PC Camera," and "REC-mode"), and confirm.
4. **Open Windows Camera:** Start the "Windows Camera" app. Your PC should detect the microscope and display the live feed.

NOTE

If the camera feed does not appear, unplug the USB cable and reconnect it.

Works with Mac/iMac (Using Photo Booth)

1. **Check for Photo Booth:** Ensure your Mac has "Photo Booth" installed. If not, download it from the App Store.
2. **Power On the Microscope:** Turn on the microscope and connect it to your Mac with the USB cable (a USB hub may be needed).
3. **Select PC Camera Mode:** On the microscope, choose "PC Camera" when prompted (options include "Mass Storage," "PC Camera," and "Charging-mode"), and confirm.
4. **Open Photo Booth:**
 - Go to **Finder > Applications > Photo Booth**.
 - Select **GENERAL-UVC#2 (Default)** from the menu bar.

Usage

: Press and hold for 2 seconds to turn the device on or off.

: Access the function settings.

Screensaver Options:

- Off
- 30 seconds
- 1 minute
- 2 minutes (optional)

Automatic Shutdown Options:

- Off
- 3 minutes

- 5 minutes
- 10 minutes

Hz Light Source:

- 50 Hz
- 60 Hz

Language Options: 16 languages available.

Date Adjustment:

Set the year, month, day, hour, minute, and second.

In Video Mode

- **Resolution:**
 - 1080 FHD
 - 720P
 - VGA
- **Circulating camera:**
 - Off
 - 3 minutes
 - 5 minutes
 - 10 minutes

NOTE When the memory card is full, the oldest video data will be overwritten with new recordings.

- **HDR (High Dynamic Range):** On/Off

NOTE When HDR is enabled, video recording starts automatically if motion is detected under the lens, even without manual operation.

- **Exposure Compensation:**
 - 7 levels available for adjustment.
- **Date Label:** Display/Close
 - *Date and Time can be displayed or hidden on the screen.*

In Photo Mode

- **Photograph Delay:**
 - Single
 - 2 seconds
 - 5 seconds
 - 10 seconds

- **Photo Pixels:**
 - 1.3 M
 - 2 M
 - 5 M
 - 8 M
 - 10 M
 - 12 M
- **Continuous Shooting:**
 - 3 consecutive shots
- **Photo Quality:**
 - Quality
 - Standard
 - Compression
- **Sharpness:**
 - Strong
 - Standard
 - Soft
- **Color Mode:**
 - Natural
 - Black and White
 - Nostalgia
- **ISO:**
 - Automatic
 - 100
 - 200
 - 400
- **Exposure Compensation:**
 - 7 levels available for adjustment
- **Photo Assist:**
 - Anti-hand shock: On/Off
- **Date Label:**
 - On/Off
- **Storage Options:**

- View Files: Up/Down
- Delete
- Protection

: Mode key for switching between modes.

- Camera Mode
- Photo Mode
- Playback Mode

: Up

: Down

: OK

For Photos or Videos

1. Power Supply (DC Interface): Micro USB port.
2. Flash Card Slot: MicroSD card slot (supports 1-128GB, Class 10 or higher).
3. Light Brightness Regulator: Adjusts brightness.
4. Reset Key: System restore function. If the device becomes unresponsive, insert a needle or similar object into this hole to restart the machine.
5. LED Lamps: 8 high-brightness LEDs with a lifespan of up to 100,000 hours.

Disposing of Used Devices

Do not dispose of this device in municipal waste systems. Hand it over to an electric and electrical device recycling and collection point. Check the symbol on the product, instruction manual and packaging. The plastics used to construct the device can be recycled in accordance with their markings. By choosing to recycle you are making a significant contribution to the protection of our environment.

Contact local authorities for information on your local recycling facility.

Parts

- 1- Switch Machine (short press for two seconds to turn on/off)
- 2- Function Keys (screen saver, auto power off, light source, language, date, format, default settings)
- 3- Mode Button (Camera/Photo/Playback Mode)
- 4- OK Key(Photo/Video, On)
- 5- Up Arrow
- 6- Down Arrow
- 7- Object Expansion Adjustment Key
- 8- Bright 8 LED Lights
- 9- Power DC Interface
- 10- REST Key, System Restore Button (If the machine is unresponsive, use a sharp needle to press the button inside this hole to restart the machine)
- 11- Flash Card Socket
- 12- Light Modulator

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