USER MANUAL

Bedienungsanleitung
Instrukcja obsługi
Návod k použití
Manuel d'utilisation
Istruzioni per l'uso
Manual de instrucciones
Használati útmutató
Brugsanvisning
Käyttöohje
Gebruiksaanwijzing
Bruksanvisning
Instruções de utilização
Používateľská príručka
Ръководство за употреба
Οδηγίες χρήσης
Upute za uporabu
Naudojimo instrukcija
Manual de utilizare
Navodila za uporabo

OVERHEAD STIRRER

DE	Produktname	Laborrührer
EN	Product name	Overhead stirrer
PL	Nazwa produktu	Mieszadło laboratoryjne
CZ	Název výrobku	Horní míchadlo
FR	Nom du produit	Agitateur suspendu
IT	Nome del prodotto	Agitatore aereo
ES	Nombre del producto	Agitador de varilla
HU	Termék neve	Keverőfejes keverő
DA	Produktnavn	Elektrisk omrører laboratorie
FI	Tuotteen nimi	Yläpuolinen sekoitin
NL	Productnaam	Bovenste roerder
NO	Produktnavn	Overhead rører
SE	Produktnamn	Overhead omrörare
PT	Nome do produto	Agitador de teto
SK	Názov produktu	Horné miešadlo
BG	Име на продукта	Горна бъркалка
EL	Όνομα προϊόντος	Αναδευτήρας από πάνω
HR	Naziv proizvoda	Gornja mješalica
LT	Produktu pavadinimas	Viršutinė maišyklė
RO	Numele produsului	Agitator deasupra capului
SL	Ime izdelka	Zgornji mešalnik
DE Modell EN Product model PL Model produktu CZ Model výrobku FR Modèle IT Modello ES Modelo HU Modell DA Model FI Tuotteen malli NL Productmodel NO Produktmodell SE Produktmodell PT Modelo do produto SK Model BG Модел на продукт EL Μοντέλο προϊόντος HR Model proizvoda LT : Gaminio modelis RO : Model de produs SL : Model izdelka		SBS-MR-2020 SBS-MR-2040 SBS-MR-2060 SBS-MR-1830-FR
DE Hersteller EN Manufacturer PL Producent CZ Výrobce FR Fabricant IT Produttore ES Fabricante HU Termelő DA Producent FI Valmistaja NL Producent NO Produsent SE Tillverkare PT Fabricante SK Výrobca BG Производител EL Κατασκευαστής HR Proizvođač LT Gamintojas RO Producător SL Proizvajalec		expondo Polska sp. z o.o. sp. k.
DE Anschrift des Herstellers EN Manufacturer Address PL Adres producenta CZ Adresa výrobce FR Adresse du fabricant IT Indirizzo del produttore ES Dirección del fabricante HU A gyártó címe DA Producentens adresse FI Valmistajan osoite NL Adres producent NO Produsentens adresse SE Tillverkarens adress PT Endereço do fabricante SK Adresa výrobcu BG Адрес на производителя EL : Διεύθυνση κατασκευαστή HR Adresa proizvođača LT Gamintojo adresas RO Adresa producătorului SL Naslov proizvajalca		ul. Nowy Kisielin – Innowacyjna 7, 66-002 Zielona Góra Poland, EU

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This User Manual has been translated using machine translation. We have made every effort to ensure the translation is accurate, but please note that automated translations are not perfect and are not meant to replace human translators. The official version of the User Manual is in English. Any differences between the translated version and the original English are not legally binding. If you have any questions about the accuracy of the translation, please refer to the English version, which is the official reference.

Technical data

Parameter description	Parameter value			
Product name	Overhead stirrer			
Model	SBS-MR-2020	SBS-MR-2040	SBS-MR-2060	SBS-MR-1830-FR
Rated voltage [V~] / frequency [Hz]	100-240 / 50-60			
Rated power [W]	75	120	200	150
Max mixing (H2O) [L]	20	40	60	30
Speed range [rpm]	100-2000	100-2000	100-2000	50-1800
Protection class	III			
Dimensions [width * length * height; mm]	200*420*600	200*462*595	200*365*600	200*410*600
Weight [kg]	7.55	8.3	8.4	8.4

Description

The product is designed for mixing and stirring low, medium, and high-viscosity liquids in laboratory settings. The main body is constructed from high-quality engineering plastics, ensuring stable operation. The output end features a variable speed booster mechanism that increases output torque. The stirring speed can be adjusted as needed for the experiment, and the digital display provides clear, real-time speed information for accurate data collection.

The user is liable for any damage resulting from unintended use of the device.

Installation

1. Working Environment Requirements:

- 1) The workspace should be clean and dry.
- 2) Place the instrument on a horizontal, stable surface.
- 3) Minimize vibration in the workspace.
- 4) Keep the instrument away from magnetic objects and sources.
- 5) Do not use the instrument in explosive environments.
- 6) Avoid using the instrument in areas with high humidity or dust for extended periods.

2. Equipment Installation:

- 1) After unpacking and removing all packaging materials, place the base plate on a level and flat surface.
- 2) Attach the vertical rod to the base plate and tighten it securely. Insert the sub-rod into the main rod and tighten it as well.

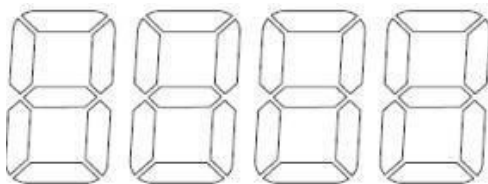
- 3) Put the bottom plate on a stable place, install the column in the hole of the bottom plate, and tighten it with a wrench to prevent shaking.
- 4) Slide the small-hole cross-fixing frame onto the vertical pole and tighten it at the bottom. Screw the large-hole cross-fixing frame onto the pole and install the universal clamp on the lower cross-fixing frame, tightening it securely. Insert the rod into the upper cross-fixing frame and tighten it.
- 5) One hand holds the main shaft of the stirring paddle, while holding the lower part of the chuck and turning the upper part of the chuck with one hand, it will automatically get stuck during rotation.
- 6) Insert the stirring rod into the drill chuck and slowly tighten it to ensure proper alignment and prevent off-center positioning.
- 7) Plug the power cord into the power socket, connect the power source to the back of the instrument, and the whole machine is installed.

Operation

SBS-MR-2020 & SBS-MR-2040 & SBS-MR-2060

1. Basic version

1) Operation interface


- Knob: Long press to turn the instrument on or off. A short press controls the start, pause, and speed adjustment. Rotate clockwise to increase the speed and counterclockwise to decrease it.

2) Operation instructions

- Place the sample container on the bottom plate. Loosen the motor's cross-fixing frame and adjust the motor's height so the stirring blade extends into the container, ensuring the blade is approximately 1 cm from the bottom. Tighten the motor's fixing frame and adjust the container's position to centre the stirring paddle, securing it with the universal clamp.
- Press and hold the knob to power on the device. Turn the knob clockwise to increase the speed or counterclockwise to decrease it.
- Press the knob to start the operation, and the speed will gradually increase. To pause, press the knob again, and press it once more to resume.
- Once the stirring is complete, press the knob to stop the operation, or press and hold the knob to turn off the device. Wait for the stirring blade to stop before removing the sample.

2. Updated version

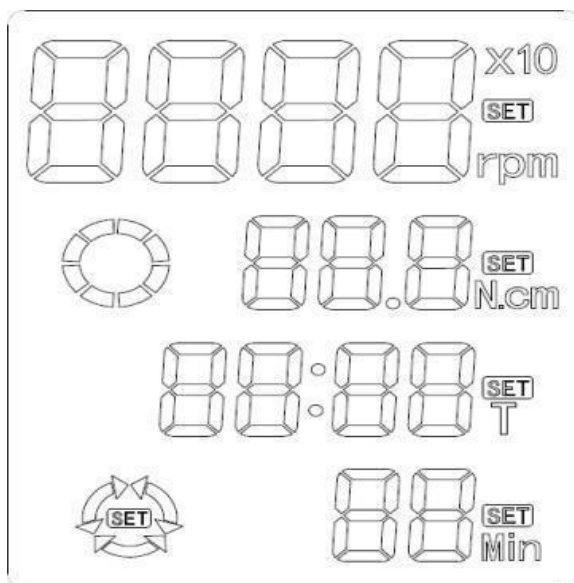1) Operation interface

- **SET Key:** Used to select rotation speed and set the timer. Press and hold to enter programming mode.
- **ON/OFF Key:** Short press to power the device on or off.
- **Knob Encoder:** Rotate to adjust speed and set the timer. Turning clockwise increases speed, while counterclockwise decreases it. Short press to start or pause operation and long press to stop the device.

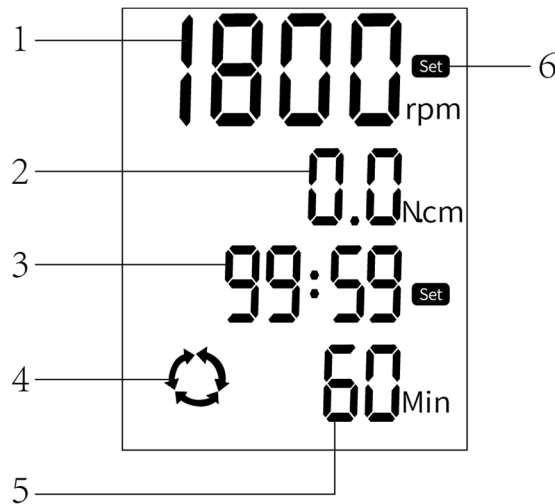
2) Operation instructions

- Place the sample container on the bottom plate. Loosen the motor's cross-fixing frame and adjust the motor height so the stirring blade extends into the container, positioning the blade approximately 1 cm from the container's bottom. Tighten the motor's cross-fixing frame, adjust the container's position to center the stirring paddle, and secure it with the universal clamp.
- **Conventional Mode:**
 - a) **Speed Setting:** Short press the ON/OFF button to power on the device. Short press the SET button to enter speed adjustment mode and use the knob to adjust the speed. Turning the knob clockwise increases the speed, while counterclockwise decreases it.
 - b) **Timer Setting:** After setting the speed, press the SET button again to enter the timer adjustment mode. Rotate the knob to set the timing duration—clockwise to increase and counterclockwise to decrease. The minimum time is 1 minute, and setting the timer to 0 allows the device to run continuously.
 - c) **Start/Pause/Stop:** Short press the knob to start operation, press again to pause, and long press to stop and return to the settings interface.
- **Program Control Mode:**
 - a) **Entering Program Mode:** Short press the ON/OFF button to turn on the power. Long press the SET button to enter multi-stage programming mode, displaying "Program 1" by default. Adjust the speed using the knob, turning clockwise to increase and counterclockwise to decrease. Press SET again to set the timer, using the knob to adjust time, turning clockwise to increase and counterclockwise to decrease.
 - b) **Additional Programs:** Short press the SET button to move to "Program 2" and refer to the instructions from [a)] for setting methods.

- c) **Program-Controlled Operation:** While in the "Program 2" interface, short press the knob to start running from "Program 1" through "Program 2." In the "Program 3" interface, short press the knob to start running from "Program 1" through "Program 3," and so on. Up to 6 programs can be set.
- d) **Shut Down:** After running, short press the ON/OFF button to turn off the device.

SBS-MR-1830-FR

1. Operation interface

- 1- Speed value display
- 2- Torque value display
- 3- Timing time display
- 4- Forward and reverse tips
- 5- Intermittent forward and reverse time display
- 6- Setting icon display

2. Button function

	<p>Set key Press this key to switch between speed setting and timing time setting.</p>
	<p>Forward and reverse setting key Press this key to set forward, reverse, forward and reverse operation mode.</p>
	<p>On/off key Press this key to turn on/off the stirring function.</p>

3. Operation instructions

- a) Press the "SET" key to enter speed adjustment mode. Use the knob to adjust the speed: turning it clockwise decreases the speed, while turning it counterclockwise increases the speed.
- b) Once the speed is set, press the "SET" key again to enter the timer setting mode. Rotate the knob to adjust the timer: turning it clockwise decreases the time, while counterclockwise increases it. If the timer is set to 0, the device will run continuously.

- c) Press the "Switch" key to enter the forward and reverse mode settings. Use the "Switch" button to choose between forward, reverse, or forward and reverse running modes. Adjust the interval between forward and reverse by turning the knob: clockwise decreases the interval, while counterclockwise increases it.
- d) Short press the "ON/OFF" key to start the instrument. Press it again to pause operation. Long press the "ON/OFF" key to return to the main settings menu.

Maintenance

1. Avoid using hard objects to clean the machine to prevent surface damage.
2. Unplug the power cord if the machine is not in use for an extended period. Store it in a cool, dry place.
3. Before restarting after long inactivity, check the machine's functions to ensure proper operation.

Troubleshooting

SBS-MR-2020 & SBS-MR-2040 & SBS-MR-2060


Problem	Possible cause	Solution
Unable to boot	<ol style="list-style-type: none"> 1. Loose power cord. 2. The power supply is damaged. 	<ol style="list-style-type: none"> 1. Re-plug the power cord. 2. Replace the power supply.
No mixing	<ol style="list-style-type: none"> 1. Motor damage. 2. Instrument damage. 	<ol style="list-style-type: none"> 1. Replace the motor 2. Replace the meter.

SBS-MR-1830-FR

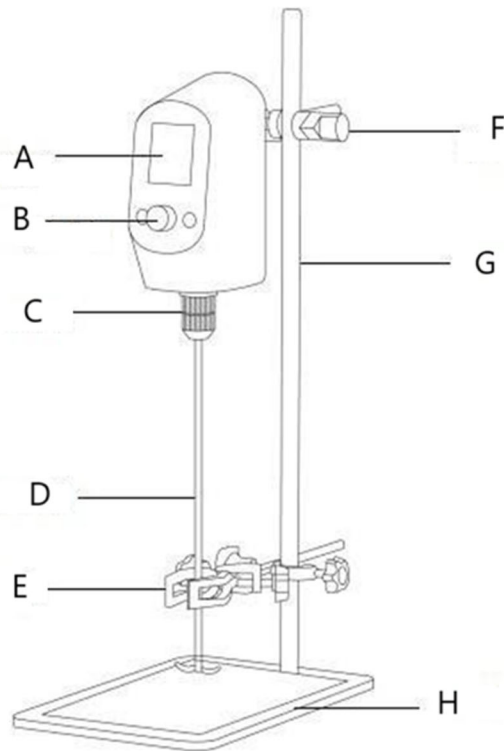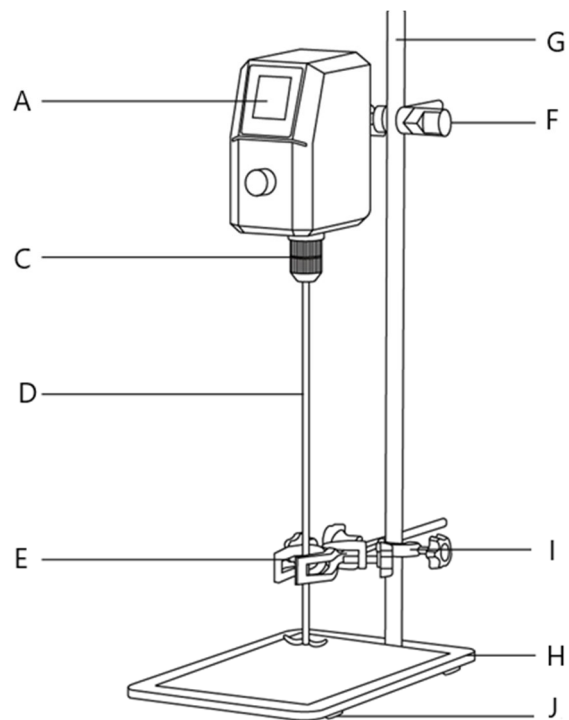
Problem	Possible cause	Solution
The speed cannot reach the set value.	The set speed does not match the current medium viscosity	Restart after turning down the speed.
Can't boot normally.	The power cord is not connected tightly.	Check whether the power cord is tightly connected or connected to other power outlets.
ERR1 is displayed in the value setting area.	The plug wire falls off.	Check the internal cable of the instrument and re-plug.

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