USER MANUAL

Bedienungsanleitung
Instrukcja obsługi
Návod k použití
Manuel d'utilisation
Istruzioni per l'uso
Manual de instrucciones
Használati útmutató
Bruksanvisning
Käyttöohje
Gebruiksaanwijzing
Bruksanvisning
Instruções de utilização
Používateľská príručka
Ръководство за употреба
Οδηγίες χρήσης
Upute za uporabu
Naudojimo instrukcija
Manual de utilizare
Navodila za uporabo

MINI VORTEX MIXER

DE	Produktname	Mini-Wirbelmischer
EN	Product name	Mini vortex mixer
PL	Nazwa produktu	Miniaturowy mikser wirowy
CZ	Název výrobku	Mini vortexový mixér
FR	Nom du produit	Mini mélangeur vortex
IT	Nome del prodotto	Mini miscelatore vortex
ES	Nombre del producto	Mini mezclador de vórtice
HU	Termék neve	Mini örvénykeverő
DA	Produktnavn	Mini vortex-mixer
FI	Tuotteen nimi	Mini vortex-sekoitin
NL	Productnaam	Mini-vortexmixer
NO	Produktnavn	Mini vortex mikser
SE	Produktnamn	Mini vortexblandare
PT	Nome do produto	Mini misturador vortex
SK	Názov produktu	Mini vortexový mixér
BG	Име на продукта	Мини вихров миксер
EL	Όνομα προϊόντος	Μίνι μίξερ vortex
HR	Naziv proizvoda	Mini vortex miješalica
LT	Produkto pavadinimas	Mini sūkurinis maišytuvas
RO	Numele produsului	Mini mixer vortex
SL	Ime izdelka	Mini vrtnični mešalnik
DE Modell EN Product model PL Model produktu CZ Model výrobku FR Modèle IT Modello ES Modelo HU Modell DA Model FI Tuotteen malli NL Productmodel NO Produktmodell SE Produktmodell PT Modelo do produto SK Model BG Модел на продукт EL Μοντέλο προϊόντος HR Model proizvoda LT : Gaminio modelis RO : Model de produs SL : Model izdelka		SBS-LAB-151 SBS-LAB-152
DE Hersteller EN Manufacturer PL Producent CZ Výrobce FR Fabricant IT Produttore ES Fabricante HU Termelő DA Producent FI Valmistaja NL Producent NO Produsent SE Tillverkare PT Fabricante SK Výrobca BG Производител EL Κατασκευαστής HR Proizvođač LT Gamintojas RO Producător SL Proizvajalec		expondo Polska sp. z o.o. sp. k.
DE Anschrift des Herstellers EN Manufacturer Address PL Adres producenta CZ Adresa výrobce FR Adresse du fabricant IT Indirizzo del produttore ES Dirección del fabricante HU A gyártó címe DA Producentens adresse FI Valmistajan osoite NL Adres producent NO Produsentens adresse SE Tillverkarens adress PT Endereço do fabricante SK Adresa výrobcu BG Адрес на производителя EL : Διεύθυνση κατασκευαστή HR Adresa proizvođača LT Gamintojo adresas RO Adresa producătorului SL Naslov proizvajalca		ul. Nowy Kisielin – Innowacyjna 7, 66-002 Zielona Góra Poland, EU

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This User Manual has been translated using machine translation. We have made every effort to ensure the translation is accurate, but please note that automated translations are not perfect and are not meant to replace human translators. The official version of the User Manual is in English. Any differences between the translated version and the original English are not legally binding. If you have any questions about the accuracy of the translation, please refer to the English version, which is the official reference. More language versions are available upon request via info@expondo.com.

Technical data

Parameter description	Parameter value	
Product name	Mini vortex mixer	
Model	SBS-LAB-151	SBS-LAB-152
Voltage [V] / Frequency [Hz]	100-240 / 50-60	
Output voltage [V]	12	15
Output power [W]	7.2	10
IP class	IP42	IP41
Dimensions (Width x Length x Height) [mm]	98x98x82	170x155x82
Weight [kg]	1.05	1.3
Protection class	II	

Purpose

The product is a compact, laboratory instrument designed for efficiently mixing small volumes of liquids in test tubes or vials. It generates a rapid vortex motion to thoroughly blend samples, making it ideal for applications such as preparing reagents, mixing solutions, or homogenizing samples in research, clinical, and educational labs.

The user is liable for any damage resulting from unintended use of the product.

Correct Use

The device is designed for shaking test tubes. Gently push the test tube into the soft attachment to start shaking. Shaking stops as soon as you lift out the test tube.

Unpacking

Please unpack the device carefully and check it for damage. It is important that any transport damage is detected when the device is unpacked. As supplied, the package contains:

- The device,
- A plug-in power supply unit,
- An instruction manual.

Commissioning

Connect the device to a socket.

Maintenance and Cleaning

NOTE

- Wear the proper protective gloves during cleaning of the devices.
- Electrical devices may not be placed in the cleansing agent for the purpose of cleaning.
- Before using other cleaning agents than the recommended ones, make sure this method does not destroy the instrument.

The device is maintenance-free. It is subject only to the natural wear and tear of components and their statistical failure rate. Use only cleansing agents which have been approved to clean the devices.

Use the corresponding clean agent based on the material in the following table:

Material	Cleaning Agent
Dyes	Isopropyl alcohol
Constructions materials	Water-containing tenside Isopropyl alcohol
Cosmetics	Water-containing tenside/Isopropyl alcohol
Foodstuffs	Water-containing tenside
Fuels	Water-containing tenside

Disposing of Used Devices

Do not dispose of this device in municipal waste systems. Hand it over to an electric and electrical device recycling and collection point. Check the symbol on the product, instruction manual and packaging. The plastics used to construct the device can be recycled in accordance with their markings. By choosing to recycle you are making a significant contribution to the protection of our environment.

Contact local authorities for information on your local recycling facility.

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