ULTRASONIC CLEANER

PROCLEAN 3.2S

PROCLEAN 4.5S

PROCLEAN 6.5S

PROCLEAN 10.0S

PROCLEAN 30.0S

USER MANUAL

BEDIENUNGSANLEITUNG
INSTRUKCJA OBSŁUGI
NÁVOD K POUŽITÍ
MANUEL D'UTILISATION
ISTRUZIONI PER L'USO
MANUAL DE INSTRUCCIONES

DE	3
EN	7
PL	11
CZ	15
FR	19
IT	23
ES	27

PRODUKTNAME	ULTRASCHALLREINIGER
PRODUCT NAME	ULTRASONIC CLEANER
NAZWA PRODUKTU	MYJKA ULTRADŹWIĘKOWA
NÁZEV VÝROBKU	ULTRAZVUKOVÁ ČISTIČKA
NOM DU PRODUIT	NETTOYEUR À ULTRASONS
NOME DEL PRODOTTO	PULITORE ULTRASUONI
NOMBRE DEL PRODUCTO	LIMPIADOR ULTRASÓNICO
MODELL	
PRODUCT MODEL	PROCLEAN 3.2S
MODEL PRODUKTU	PROCLEAN 4.5S
MODEL VÝROBKU	PROCLEAN 6.5S
MODÈLE	PROCLEAN 10.0S
MODELLO	PROCLEAN 30.0S
MODELO	
IMPORTEUR	
IMPORTER	
IMPORTER	
DOVOZCE	EXPONDO POLSKA SP. Z O.O. SP. K.
IMPORTATEUR	
IMPORTATORE	
IMPORTADOR	
ADRESSE VON IMPORTEUR	
IMPORTER ADDRESS	
ADRES IMPORTERA	
ADRESA DOVOZCE	UL. NOWY KISIELIN-INNOWACYJNA 7, 66-002 ZIELONA GÓRA POLAND, EU
ADRESSE DE L'IMPORTATEUR	
INDIRIZZO DELL'IMPORTATORE	
DIRECCIÓN DEL IMPORTADOR	

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Parameter - Beschreibung	Parameter - Wert		
Produktname	Ultraschallreiniger		
Modell	PROCLEAN 3.2S	PROCLEAN 4.5S	PROCLEAN 6.5S
Kapazität [L]	3,2	4,5	6,5
Innenmaße des Behälters [mm]	100 ×140 ×240	100 ×155 ×300	150 ×155 ×300
Nennspannung [V~] / Frequenz [Hz]	230/50		
Schallfrequenz [kHz]	40		
Ultraschallstärke [W]	120	120	180
Ablasshahn	Nein		Ja
Temperatur [°C]	0-80		

Produktname	Ultraschallreiniger	
Modell	PROCLEAN 10.0S	PROCLEAN 30.0S
Kapazität [L]	10	30
Innenmaße des Behälters [mm]	150 ×240 ×300	200 ×300 ×500
Nennspannung [V~] / Frequenz [Hz]	230/50	
Schallfrequenz [kHz]	40	
Ultraschallstärke [W]	240	600
Ablasshahn	Ja	
Temperatur [°C]	0-80	

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Parameter description	Parameter value	
Product name	Ultrasonic Cleaner	
Model	PROCLEAN 10.0S	PROCLEAN 30.0S
Capacity [L]	10	30
Internal dimensions of tank [mm]	150 ×240 ×300	200 ×300 ×500
Rated voltage [V~] / Frequency [Hz]	230/50	
Sound frequency [kHz]	40	
Ultrasonic power [W]	240	600
Drain tap	Yes	
Temperature [°C]	0-80	

1. GENERAL DESCRIPTION

The user manual is designed to assist in the safe and trouble-free use of the device. The product is designed and manufactured in accordance with strict technical guidelines, using state-of-the-art technologies and components. Additionally, it is produced in compliance with the most stringent quality standards.

DO NOT USE THE DEVICE UNLESS YOU HAVE THOROUGHLY READ AND UNDERSTOOD THIS USER MANUAL

To increase the product life of the device and to ensure trouble-free operation, use it in accordance with this user manual and regularly perform maintenance tasks. The technical data and specifications in this user manual are up to date. The manufacturer reserves the right to make changes associated with quality improvement. The device

is designed to reduce noise emission risks to a minimum, taking into account technological progress and noise reduction opportunities.

LEGEND

	The product satisfies the relevant safety standards.
	Read instructions before use.
	The product must be recycled.
	WARNING! or CAUTION! or REMEMBER! Applicable to the given situation (general warning sign).
	Use ear protection. Exposure to loud noise may result in hearing loss.
	ATTENTION! Electric shock warning!
	ATTENTION! Hot surface, risk of burns!

PLEASE NOTE! Drawings in this manual are for illustration purposes only and in some details may differ from the actual product.

The original operation manual is written in German. Other language versions are translations from the German.

2. USAGE SAFETY

ATTENTION! Read all safety warnings and all instructions. Failure to follow the warnings and instructions may result in electric shock, fire and/or serious injury or even death.

The terms "device" or "product" are used in the warnings and instructions to refer to "Ultrasonic Cleaner". Do not use in very humid environments or in the direct vicinity of water tanks. Prevent the device from getting wet. Risk of electric shock! Do not put your hands or other items inside the device while it is in use! Do not cover the ventilation openings!

2.1. ELECTRICAL SAFETY

- a) The plug must fit the socket. Do not modify the plug in any way. Using original plugs and matching sockets reduces the risk of electric shock.
- b) Avoid touching earthed elements such as pipes, heaters, boilers and refrigerators. There is an increased risk of electric shock if the earthed device is exposed to rain, comes into direct contact with a wet surface or is operating in a damp environment. Water getting into the device increases the risk of damage to the device and of electric shock.
- c) Do not touch the device with wet or damp hands.
- d) Use the cable only for its designated use. Never use it to carry the device or to pull the plug out of a socket. Keep the cable away from heat sources, oil, sharp edges or moving parts. Damaged or tangled cables increase the risk of electric shock.
- e) If using the device in a damp environment cannot be avoided, a residual current device (RCD) should be applied. The use of an RCD reduces the risk of electric shock.
- f) Do not use the device if the power cord is damaged or shows obvious signs of wear. A damaged power cord should be replaced by a qualified electrician or the manufacturer's service centre.
- g) To avoid electric shock, do not immerse the cord, plug or device in water or other liquids. Do not use the device on wet surfaces.

2.2. SAFETY IN THE WORKPLACE

- Make sure the workplace is clean and well lit. A messy or poorly lit workplace may lead to accidents. Try to think ahead, observe what is going on and use common sense when working with the device.
- If you discover damage or irregular operation, immediately switch the device off and report it to a supervisor without delay.
- If there are any doubts as to the correct operation of the device, contact the manufacturer's support service.
- Only the manufacturer's service point may repair the device. Do not attempt any repairs independently!
- In case of fire, use a powder or carbon dioxide (CO₂) fire extinguisher (one intended for use on live electrical devices) to put it out.
- Children or unauthorised persons are forbidden to enter a work station. (A distraction may result in loss of control over the device).
- Use the device in a well-ventilated space.
- Regularly inspect the condition of the safety labels. If the labels are illegible, they must be replaced.
- Please keep this manual available for future reference. If this device is passed on to a third party, the manual must be passed on with it.
- Keep packaging elements and small assembly parts in a place not available to children.
- Keep the device away from children and animals.

REMEMBER! When using the device, protect children and other bystanders.

2.3. PERSONAL SAFETY

- Do not use the device when tired, ill or under the influence of alcohol, narcotics or medication which can significantly impair the ability to operate the device.
- The device can be handled only by physically fit persons who are capable of handling it, properly trained, familiar with this manual and trained within the scope of occupational health and safety.
- When working with the device, use common sense and stay alert. Temporary loss of concentration while using the device may lead to serious injuries.
- To prevent the device from accidentally switching on, make sure the switch is on the OFF position before connecting to a power source.
- Do not overestimate your abilities. When using the device, keep your balance and remain stable at all times. This will ensure better control over the device in unexpected situations.
- The device is not a toy. Children must be supervised to ensure that they do not play with the device.

2.4. SAFE DEVICE USE

- Do not overload the device. Use the appropriate tools for the given task. A correctly-selected device will perform the task for which it was designed better and in a safer manner.
- Do not use the device if the ON/OFF switch does not function properly (does not switch the device on and off). Devices which cannot be switched on and off using the ON/OFF switch are hazardous, should not be operated and must be repaired.
- Disconnect the device from the power supply before commencement of adjustment, cleaning and maintenance. Such a preventive measure reduces the risk of accidental activation.

- When not in use, store in a safe place, away from children and people not familiar with the device who have not read the user manual. The device may pose a hazard in the hands of inexperienced users.
- Keep the device in perfect technical condition. Before each use check for general damage and especially check for cracked parts or elements and for any other conditions which may impact the safe operation of the device. If damage is discovered, hand over the device for repair before use.
- Keep the device out of the reach of children.
- Device repair or maintenance should be carried out by qualified persons, only using original spare parts. This will ensure safe use.
- To ensure the operational integrity of the device, do not remove factory-fitted guards and do not loosen any screws.
- When transporting and handling the device between the warehouse and the destination, observe the occupational health and safety principles for manual transport operations which apply in the country where the device will be used.
- Do not move, adjust or rotate the device in the course of work.
- Do not leave this appliance unattended while it is in use.
- Clean the device regularly to prevent stubborn grime from accumulating.
- The device is not a toy. Cleaning and maintenance may not be carried out by children without supervision by an adult person.
- Before adjusting, replacing accessories, or storing the appliance, unplug the appliance from the power source. This will decrease the possibility of accidentally switching it on.
- Do not run the device when empty.
- It is forbidden to interfere with the structure of the device in order to change its parameters or construction.
- Do not dive parts of the body or parts of living things (animals, plants) into the basin, especially when the unit is in operation.
- Avoid the contact with eyes or skin of the detergents used.
- Do not inhale the detergent.
- Do not use the device if it's empty. The tank of the device must be filled with minimum level of the cleaning liquid in accordance with the below table:
- Do not fill the tank with water above 50°C.
- Do not place items that have been in contact with acids, chloride ions and other aggressive substances in the basin.
- Do not use flammable liquids or other chemicals containing chloride ions, or leading to their release (such as some disinfectants, detergents or dishwashing liquids).
- During the heating of the cleaning liquid, be sure to stir the liquid every 15 minutes so as to avoid overheating and boiling.
- During ultrasonic cleaning, you will hear the sound of cavitation. When using the device for a long time, please wear hearing protection within a radius of 2 m around the device.
- The cleaning fluid heats up when it is used long. Watch the temperature if you clean objects sensitive to heat.

ATTENTION! Despite the safe design of the device and its protective features, and despite the use of additional elements protecting the operator, there is still a slight risk of accident or injury when using the device. Stay alert and use common sense when using the device.

3. USE GUIDELINES

The ultrasonic cleaner is a device used to remove operational dirt from surfaces, tools and functional parts of machines.

The user is liable for any damage resulting from unintended use of the device.

3.1. DEVICE DESCRIPTION

- Cover Handle
- Control panel (varies depending on model)
- Handle
- Power socket and fuse socket
- Portable rack
- Tub
- Drain plug

3.2. PREPARING FOR USE

APPLIANCE LOCATION

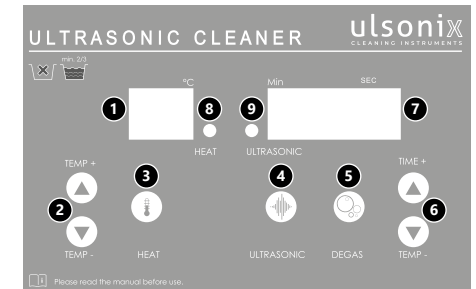
Temperatura otoczenia nie może przekraczać 40°C The temperature of environment must not be higher than 40°C and the relative humidity should be less than 85%. Ensure good ventilation in the room in which the device is being used. There should be at least 10 cm distance between each side of the device and the wall or other objects. The device should always be used when positioned on an even, stable, clean, fireproof and dry surface, and be out of the reach of children and persons with limited mental and sensory functions. Position the device such that you always have access to the power plug. The power cord connected to the appliance must be properly grounded and correspond to the technical details on the product label.

ASSEMBLING THE APPLIANCE

- Unpack the device.
- Check that the device is undamaged and that all components are included in the set. If it is damaged or any elements are missing, contact the manufacturer.
- Place the portable rack into the tub

- Connect the power cable to the device socket.
- Place the lid on top of the device to cover the tub

3.3. DEVICE USE

- Temperature indicator
- Temperature adjustment buttons
- „HEAT”
- „ULTRASONIC”
- Degas button
- Operation time adjustment buttons
- Minute indicator
- Indicator for heating in process
- Indicator for ultrasonic cleaning in process

HOW TO USE THE DEVICE

- Fill the tank with cleaning solution; Connect the power cord to the power source; When the cleaner has been started, the temperature indicator will display the current ambient temperature.
- Press the operation time adjustment button to set the washing cleaner operation time, pressing the button once changes the value by one minute, pressing and holding the button changes the value by ten minutes. To end the cleaner operation, press and hold the "4" button.
PLEASE NOTE: the cleaning cycle should not exceed one hour.
- Press the temperature control button to set the cleaner temperature, pressing the button once will change the value by one degree Celsius, pressing and holding the button will change the value by ten degrees.
- Press the "HEAT" or "4" button to start both functions at once. After pressing the "4" button again, the ultrasonic function and "3" function will be turned off; in order to continue of heating, press the "3" button.
- To start the degassing function press the "5" button
- when you have finished work, empty the tub.

3.4. CLEANING AND MAINTENANCE

- Unplug the mains plug and allow the device to cool completely before each cleaning, adjustment or replacement of accessories, or if the device is not being used.
- Use only non-corrosive cleaners to clean the surface.
- After cleaning the device, all parts should be dried completely before using it again.
- Do not spray the device with a water jet or submerge it in water.
- Do not allow water to get inside the device through vents in the housing of the device.

- f) The device must be regularly inspected to check its technical efficiency and spot any damage.
 g) Use a soft cloth for cleaning.
 h) Do not use sharp and/or metal objects for cleaning (e.g. a wire brush or a metal spatula) because they may damage the surface material of the appliance.

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NAMEPLATE TRANSLATIONS

ulsonix
CLEANING INSTRUMENTS

1 PRODUCT NAME ULTRASONIC CLEANER
2 MODEL
3 ULTRASONIC POWER
4 HEATING POWER
5 TANK DIMENSIONS
6 TANK CAPACITY
7 PRODUCTION YEAR
8 SERIAL NO.

10 ELECTRIC SHOCK | DON'T OPEN

11 ATTENTION: THIS MACHINE MUST BE GROUNDED!
12 BEWARE: TO MINIMIZE THE RISK OF ELECTRIC SHOCK, PLEASE DON'T REMOVE ANY SCREWS OR THE HOUSING! PLEASE DON'T REPLACE ANY PARTS YOURSELF, TO BE REPLACED ONLY BY A CERTIFIED SPECIALIST! TO PREVENT THE RISK OF THE FIRE AND ELECTRIC SHOCK, PLEASE DO NOT EXPOSE THE MACHINE TO RAIN OR MOISTURE!

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For the disposal of the device please consider and act according to the national and local rules and regulations.

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