USER MANUAL

Bedienungsanleitung
Instrukcja obsługi
Návod k použití
Manuel d'utilisation
Istruzioni per l'uso Manual
de instrucciones Használati
útmutató Brugsanvisning
Käyttöohje
Gebruiksaanwijzing
Bruksanvisning Instruções
de utilização Používateľská
príručka Ръководство за
употреба Οδηγίες χρήσης
Upute za uporabu
Naudojimo instrukcija
Manual de utilizare
Navodila za uporabo

P R E C I S I O N S C A L E

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For products shipped in October 2025 or later

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Do not place the unit close to windows, doors, or fans that may cause unstable results due to air currents.

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4.3.4. TARE

This button is used to calibrate the scale. Place an empty container on the weighing plate until the balance indicates a stable value. Then press TARE . Your balance now displays "0". To exit the tare function, remove the container and press TARE . The scale will go back to normal.

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4.3.6. Energy saving function

The scale turns off automatically after 300 seconds.

4.3.7. First use

Press the ON / OFF button to turn on the scale. After turning on the scale, the display will first show "000000" to "999999", and then "0.0g". Let the scale acclimate for 15 minutes to achieve optimum precision results.

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4.3.8. Overload

When the weight of the object exceeds the capacity of the scale, it displays "OverLd". Remove immediately the weight of the scale to prevent damage.

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4.3.10. Calibration

- 1) Press and hold the MODE button for 4 seconds to switch to the calibration mode. The total capacity of the scale is displayed.
- 2) Place the corresponding standard weight on the scale until the word "PASS" is displayed.
- 3) Your scale is then calibrated.

4.3.11. Optional weight calibration

- When the scale displays the total capacity value, press the "Tare" to enter the self-selected weight value, and press the "TARE" again to move the number.
- Press PCS to select the number.
- Press "TARE" to confirm.
- Place the corresponding weight on the scale until the word "PASS" is displayed.
- Your scale is then calibrated.

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For products shipped in September 2025 or earlier

This User Manual has been translated using machine translation. We have made every effort to ensure the translation is accurate, but please note that automated translations are not perfect and are not meant to replace human translators. The official version of the User Manual is in English. Any differences between the translated version and the original English are not legally binding. If you have any questions about the accuracy of the translation, please refer to the English version, which is the official reference. More language versions are available upon request via info@expondo.com.

1. Technical data

Table 1: Technical data of the product

Parameter description	Parameter value	
Product name	Precision scale	
Model	SBS-LW-7500A	SBS-LW-10000A
Max. load [g]	7500	10000
Min. load [g]	0.5	
Division [g]	0.1	
Standard (e)	e=10d	
Battery	4 x AA	
Dimensions [width x depth x height; mm]	170 x 260 x 80	
Weight [kg]	1.5	

2. General description

The user manual is designed to assist in the safe and trouble-free use of the device. The product is designed and manufactured in accordance with strict technical guidelines, using state-of-the-art technologies and components. Additionally, it is produced in compliance with the most stringent quality standards.

DO NOT USE THE DEVICE UNLESS YOU HAVE THOROUGHLY READ AND UNDERSTOOD THIS USER MANUAL.

To increase the product life of the device and to ensure trouble-free operation, use it in accordance with this user manual and regularly perform maintenance tasks. The technical data and specifications in this user manual are up to date. The manufacturer reserves the right to make changes associated with quality improvement. The device is designed to reduce noise emission risks to a minimum, taking into account technological progress and noise reduction opportunities.

2.1. Legend

Icon	Description
------	-------------

	The product satisfies the relevant safety standards.
	Read instructions before use.
	The product must be recycled.
	WARNING! or CAUTION! or REMEMBER! Applicable to the given situation. (general warning sign)
	ATTENTION! Electric shock warning!

PLEASE NOTE! DRAWINGS IN THIS MANUAL ARE FOR ILLUSTRATION PURPOSES ONLY AND IN SOME DETAILS MAY DIFFER FROM THE ACTUAL PRODUCT.

3. Usage safety

ATTENTION! READ ALL SAFETY WARNINGS AND ALL INSTRUCTIONS. FAILURE TO FOLLOW THE WARNINGS AND INSTRUCTIONS MAY RESULT IN ELECTRIC SHOCK, FIRE AND/OR SERIOUS INJURY OR EVEN DEATH.

The terms "device" or "product" are used in the warnings and instructions to refer to:

Precision scale

3.1. Electrical safety

- a) Do not touch the device with wet or damp hands.
- b) **ATTENTION! DANGER TO LIFE!** While cleaning, never immerse the device in water or other liquids.
- c) Do not use in very humid environments or in the direct vicinity of water tanks.
- d) Prevent the device from getting wet. Risk of electric shock!

3.2. Safety in the workplace

- a) Make sure the workplace is clean and well lit. A messy or poorly lit workplace may lead to accidents. Try to think ahead, observe what is going on and use common sense when working with the device.

- b) If you discover damage or irregular operation, immediately switch the device off and report it to a supervisor without delay.
- c) If you are unsure about whether the product is operating correctly or if you find damage, please contact the manufacturer's service centre.
- d) Only the manufacturer's service centre may make repairs to the product. Do not attempt to make repairs yourself!
- e) In case of fire, use a powder or carbon dioxide (CO₂) fire extinguisher (one intended for use on live electrical devices) to put it out.
- f) Regularly inspect the condition of the safety labels. If the labels are illegible, they must be replaced.
- g) Please keep this manual available for future reference. If this device is passed on to a third party, the manual must be passed on with it.
- h) Keep packaging elements and small assembly parts in a place not available to children.

REMEMBER! WHEN USING THE DEVICE, PROTECT CHILDREN AND OTHER BYSTANDERS.

3.3. Personal safety

- a) Do not use the device when tired, ill or under the influence of alcohol, narcotics or medication which can significantly impair the ability to operate the device.
- b) The device is not designed to be handled by persons (including children) with limited mental and sensory functions or persons lacking relevant experience and/or knowledge unless they are supervised by a person responsible for their safety or they have received instruction on how to operate the device.

3.4. Safe device use

- a) Do not overload the device. Use the appropriate tools for the given task. A correctly-selected device will perform the task for which it was designed better and in a safer manner.
- b) Keep the device in perfect technical condition. Before each use check for general damage, especially check moving components for cracked parts or elements, and for any other conditions which may impact the safe operation of the device. If damage is discovered, hand over the device for repair before use.
- c) Device repair or maintenance should be carried out by qualified persons, only using original spare parts. This will ensure safe use.
- d) To ensure the operational integrity of the device, do not remove factory-fitted guards and do not loosen any screws.

- e) When transporting and handling the device between the warehouse and the destination, observe the occupational health and safety principles for manual transport operations which apply in the country where the device will be used.
- f) Clean the device regularly to prevent stubborn grime from accumulating.
- g) Ensure that the weight placed on the scales is stable.
- h) Check that all screws, rods, and brackets are secure.
- i) Do not weigh in case of extreme humidity and rain.
- j) Avoid working at high ambient temperatures.

ATTENTION! DESPITE THE SAFE DESIGN OF THE DEVICE AND ITS PROTECTIVE FEATURES, AND DESPITE THE USE OF ADDITIONAL ELEMENTS PROTECTING THE OPERATOR, THERE IS STILL A SLIGHT RISK OF ACCIDENT OR INJURY WHEN USING THE DEVICE. STAY ALERT AND USE COMMON SENSE WHEN USING THE DEVICE.


4. Use guidelines

Thank you for choosing our product. Each scale has a range of options and features that are ideal both for use in laboratory, classroom, or for private use.

We hope you enjoy working with our balance!

The user is liable for any damage resulting from unintended use of the device.

4.1. Product description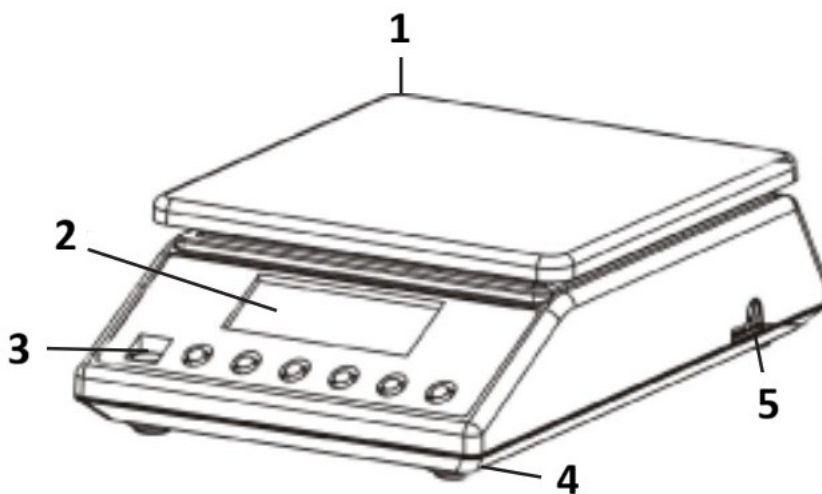

1) Weighing plate of stainless steel

2) Display

- 3) Level bubble
- 4) Plastic feet
- 5) Power cable plug

4.2. Preparing for use

4.2.1. Inserting the batteries

Remove the battery box cover and insert the batteries (AA)

4.2.2. Placement of the device and precautionary measures


To keep your balance in good condition, we recommend you follow these tips:


Avoid extreme temperatures. Do not place the unit in direct sunlight or near the air conditioning.

Make sure the scale is placed on a stable table and is not subject to vibrations.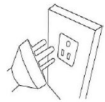

Avoid unstable power sources. Do not use close to large electrical appliances such as welding machines or large motors. Do not leave batteries in the device if you do not use it for a long time.

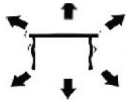Avoid vibrations. Do not place your scale near vibrating heavy machinery.

Avoid places where moisture can lead to condensation. Avoid direct contact with water.

Do not place the unit close to windows, doors or fans that may cause unstable results due to air currents.

Keep your machine clean. Do not store products on the scale when it is not used.

4.3. Device use

4.3.1. ON / OFF

- 1) Press the ON / OFF button to start the unit.
- 2) Please allow the scale to warm up for at least 5 minutes for more accurate weighing results.
- 3) Hold the ON / OFF button pressed for 3 seconds to turn off the scale.

4.3.2. UNIT

The unit change is possible by pressing the MODE button. Then press up to reach the desired unit.

There are 13 units: g, ct, lb, oz, dr, GN, ozt, dwt, MM, tl.J, tl.T, tl.H, t

4.3.3. PCS - Counting

Counting mode: Press PCS to enter the counting function. The scales switch from the weighing mode to the counting mode. By pressing this button, you can switch from one mode to another.

4.3.4. TARE

This button is used to calibrate the scale. Place an empty container on the weighing plate until the balance indicates a stable value. Then press ZERO. Your balance now displays "0". To exit the tare function, remove the container and press ZERO. The scale will back to normal.

4.3.5. Green backlight of the display

The backlight of the display turns off automatically if there is no load on the scale. The scale goes into standby mode. During weighing, the backlight of the display is turned on.

4.3.6. Energy saving function

The scale turns off automatically after the time scheduled (2, 5, or 8 minutes of inactivity). This function is turned off by factory default settings. To turn it on, a full calibration procedure must be performed in accordance with the instructions below.

Calibration requires standard weights 5 kg and 7.5 kg / or 10 kg, depending on the scale model:

- 1) Turn off the scale.
- 2) Press TARE and ON / OFF simultaneously for about 5 seconds. The display will show SET and then OFF (FOO).
- 3) Press ON / OFF – display will show 0(F01).
- 4) Press ON / OFF – display will show 1.0(F02).
- 5) Press ON / OFF – display will show 0.5(F03).
- 6) Press ON / OFF – display will show 1.0(F04).
- 7) Press ON / OFF – display will show X(F05) – where X can be 0 (AUTO-OFF is off), 2 (scale will turn off after 2 minutes of inactivity), 5 (scale will turn off after 5 minutes of inactivity), or 8 (scale will turn off after 8 minutes of inactivity). X can be changed by pressing the MODE button.

- 8) Press ON / OFF – display will show 1(F06).
- 9) Press ON / OFF – display will show 4(F07).
- 10) Press ON / OFF – display will show 75000(CAP) or 100000(CAP) depending on the scale model.
- 11) Press ON / OFF – display will show 1(div).
- 12) Press ON / OFF – display will show 0.1(dEC).
- 13) Press ON / OFF – display will show on(LCP).
- 14) Press ON / OFF – display will show 5000.0(WE1).
- 15) Press ON / OFF – display will show 7500.0(WE2) or 10000.0 (WE2) depending on the scale model.
- 16) Press ON / OFF – display will show 30.....(SCL).
- 17) Press ON / OFF – display will show LoAd and blinking 50000 (WE1).
- 18) Put on the scale standard weight 5 kg – display will show some numbers and then the word “PASS” and the weight value.
- 19) Then the display will show LoAd and blinking 75000 (WE2) or LoAd and blinking 100000 (WE2) depending on the scale model.
- 20) Put on the scale standard weight 7.5 kg or 10 kg depending on the scale model – display will show some numbers and then the word “PASS” and the weight value.
- 21) End of calibration.

4.3.7. First use

Press the ON / OFF button to turn on the scale. After turning on the scale, the display will first show “000000” to “999999”, and then “0.0g”. Let the scale acclimate for 15 minutes to achieve optimum precision results.

NOTICE!

- Use the scale only as explained above in the technical details.
- The unit should be levelled using the built-in spirit level.
- Use the device on a stable and flat surface. Protect the device from shocks, humidity, radiation, and magnetic fields.
- If the screen does not show a stable result, make sure the support of the platform is not damaged.

4.3.8. Overload

When the weight of the object exceeds the capacity of the scale, it displays “HHHH”. Remove immediately the weight of the scale to prevent damage.

4.3.9. Counting function

Put a number of countable parts on the scale (10, 20, 30, 40, 50, 60, 70, 80, 90, 100) and then press PCS. You will pass in counting mode. Press the PCS button again to select the number corresponding to the number of pieces set. Then press the ON / OFF button to confirm. Your balance is now in counting mode. To return to weighing mode, simply press the MODE button.

4.3.10. Calibration (requires standard weight 5kg)

22) Press and hold the MODE button for 4 seconds to switch to the calibration mode. The total capacity of the scale is displayed.

23) Place the corresponding standard weight on the scale until the word "PASS" is displayed.

24) Your scale is then calibrated.

4.3.11. Standards requirements

- For scales of 7500g: 5 kg and 7.5kg
- For scales of 10000g: 5 kg and 10kg

4.4. Cleaning and maintenance

4.4.1. General instructions

- a) Use only non-corrosive cleaners to clean the surface.
- b) After cleaning the device, all parts should be dried completely before using it again.
- c) Store the unit in a dry, cool place, free from moisture and direct exposure to sunlight.
- d) Do not spray the device with a water jet or submerge it in water.
- e) Do not allow water to get inside the device through vents in the housing of the device.
- f) The device must be regularly inspected to check its technical efficiency and spot any damage.
- g) Use a soft cloth for cleaning.
- h) Do not leave the battery in the device if it will not be used for a longer period of time.
- i) Do not use sharp and/or metal objects for cleaning (e.g. a wire brush or a metal spatula) because they may damage the surface material of the appliance.
- j) Do not clean the device with an acidic substance, agents of medical purposes, thinners, fuel, oils or other chemical substances because it may damage the device.

4.4.2. Safe removal of batteries and rechargeable batteries

Batteries are installed in the devices.

Remove used batteries from the device using the same procedure by which you installed them.

Recycle batteries with the appropriate organisation or company.

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