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LABORZENTRIFUGE

## **USER MANUAL**

ELECTRIC CENTRIFUGE

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LABORATORNÍ CENTRIFUGA

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CENTRIFUGA DA LABORATORIO

## **MANUAL DE INSTRUCCIONES**

CENTRIFUGADORA DE LABORATORIO

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**SBS-LZ-4000/20-6 | SBS-LZ-4000/20-12 | SBS-MZ-4000/6-1.5**

## INHALT | CONTENU | CONTENT | CONTENUTO | CONTENIDO | TREŠĆ | OBSAH

▪ Deutsch	3
▪ English	7
▪ Polski	11
▪ Česky	15
▪ Français	19
▪ Italiano	23
▪ Español	27

NAZWA PRODUKTU	WIRÓWKA LABORATORYJNA
PRODUKTNAME	LABORZENTRIFUGE
PRODUCT NAME	ELECTRIC CENTRIFUGE
NOM DU PRODUIT	CENTRIFUGEUSE DE LABORATOIRE
NOME DEL PRODOTTO	CENTRIFUGA DA LABORATORIO
NOMBRE DEL PRODUCTO	CENTRIFUGADORA DE LABORATORIO
NÁZEV VÝROBKU	LABORATORNÍ CENTRIFUGA
MODEL PRODUKTU	
MODELL	
PRODUCT MODEL	SBS-LZ-4000/20-6
MODÈLE	SBS-LZ-4000/20-12
MODELLO	SBS-MZ-4000/6-1.5
MODELO	
MODEL VÝROBKU	
PRODUCENT	
MANUFACTURER	
PRODUCENT	
FABRICANT	EXPONDO POLSKA SP. Z O.O. SP. K.
PRODUTTORE	
FABRICANTE	
VÝROBCE	
ADRES PRODUCENTA	
ANSCHRIFT DES HERSTELLERS	
MANUFACTURER ADDRESS	
ADRESSE DU FABRICANT	UL. NOWY KISIELIN – INNOWACYJNA 7, 66-002 ZIELONA GÓRA   POLAND, EU
INDIRIZZO DEL PRODUTTORE	
INDIRIZZO DEL PRODUTTORE	
ADRESA VÝROBCE	

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### SYMBOLE

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## USER MANUAL

### SYMBOLS

	The operation manual must be reviewed.
	Never dispose of electrical equipment together with household waste.
	This machine conforms to the CE declarations.
	Attention! Risk of accident.
	Safety goggles must be worn.
	Attention! Wear protective gloves.

 **PLEASE NOTE!** Drawings in this manual are for illustration purposes only and in some details may differ from the actual machine.

The original operation manual is in German. Other language versions are translations from German.

## I. USER MANUAL

### SAFETY INSTRUCTIONS

#### GENERAL SAFETY INFORMATION FOR THE USE OF ELECTRICAL DEVICES

To avoid any risk of injury from fire or electric shock, please follow the basic safety instructions when using this device. Please read the instructions carefully and make sure that you have understood them well. Keep the manual near the equipment to be able to read it at any time. Always use current sources connected to the ground that provide the correct voltage (indicated on the label on the device). If you have any doubts, let an electrician check that the connection is properly grounded. Never use a defective power cable. Do not open the unit in damp or wet environment, or if your hands or body are damp or wet. Protect the unit from solar radiation. Use the device in a protected place to avoid damaging the equipment or putting at risk others. Make sure the device is able to cool and avoid placing it too close to other devices that produce heat. Before cleaning the machine, disconnect it from the power source. Use a soft damp cloth for cleaning. Avoid using detergents and make sure that no liquid enters the unit. No internal element of this device needs to be maintained by the user. Opening the device without our approval leads to a loss of warranty!

### II. SAFETY GUIDELINES

1. Please read and understand these instructions carefully before using the device! Please observe the safety guidelines carefully in order to prevent damages through improper use!
2. Please keep this manual available for future reference. If this device is passed on to third parties, the manual must be passed on with it.
3. Only use this device indoors and for its intended use.
4. We do not offer any warranty for damages resulting from improper use or incorrect operation.
5. Before using this device for the first time, please check whether the main voltage type and current comply with the indicated data on the type plate.
6. This device is not intended to be used by persons (including children) with limited physical, sensory or mental capabilities or lack of experience and/or knowledge, unless they are supervised by a person responsible for their safety or have received instruction as to how the device is used.
7. **ELECTRIC SHOCK HAZARD!** Do not try to repair the unit yourself. In case of device failures, repairs must be done by qualified experts.
8. Please check the main plug and power cable regularly. If the power cable for this device is damaged, it has to be replaced by the manufacturer or its customer service or another qualified person in order to prevent hazards.
9. Prevent damages to the power cable by avoiding squeezing or bending it or rubbing it on sharp edges. Also keep the power cable away from hot surfaces and open flames.
10. **ATTENTION! DANGER TO LIFE!** While cleaning, never immerse the device in water or other liquids.
11. Under no circumstances should the housing be opened.
12. Do not place your hands or legs near the rotating parts.

### III. TECHNISCHE DATEN

Product name	ELECTRIC CENTRIFUGE		
Model	SBS-LZ-4000/20-6	SBS-LZ-4000/20-12	SBS-MZ-4000/6-1.5
Power supply [V]/Frequency [Hz]	230~/50		
Rated power [W]	110	180	15
Rotation speed [r/min]	3200	3000	3500
Time range [min]	0-30	0-30	-
Ambient temperature [°C]	0-30	0-30	≤10
Maximum number of containers	6	12	6 / 16
Container volume [ml]	20	20	0,5 - 1,5
Max. Relative Centrifugal Force [xg]	1200	1150	700

### IV. SCOPE OF THE APPLICATION

The laboratory centrifuge is designed for general use in industrial laboratories and on other sites, such as schools or another educational establishments for scientific purposes.

**ATTENTION! It is not a medical device and should not be used for medical research and tests in hospitals and other health facilities.**

**The user is liable for any damage resulting from non-intended use of the device!**

### V. BEFORE THE FIRST USE

Upon receipt of the goods, check the packaging for defects and open the packaging if there are no damages. If the packaging is damaged, please contact your transport company and distributor within 3 days, and document the damages as detailed as possible. Do not turn the package upside down! When transporting the package, please ensure that it is kept horizontal and stable.

### VI. DISPOSING OF THE PACKAGING

Please keep all packaging materials (cardboard, plastic tapes and Styrofoam), so that in case of a problem, the device can be sent back to the service centre in accurate condition!

### VII. INSTALLING THE APPLIANCE

#### POSITIONING THE MACHINE

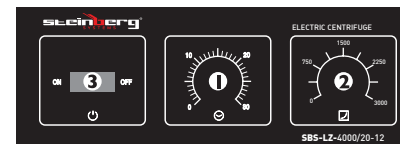
The machine should be positioned on a dry and stable surface, vertically so as to prevent movement during operation. The machine should be positioned on a flat, level surface, with sufficient load capacity for the machine and its contents. Ensure unobstructed access to the electricity plug at all times. Make sure the power source reflects the parameters on the rating plate! Prior to first use remove and wash all elements and wash the entire machine.

### VIII. MACHINE OPERATION – BASICS

Device description

1. Rotor
2. Protective cover
3. Housing

Device control panel description

1. Time adjustment knob
2. Speed adjustment knob
3. ON/OFF switch

### DEVICE OPERATION

1. Set both switches to "0".
2. Connect the power cable to a power socket.
3. Then insert the containers symmetrically into the device rotor.

Use special adaptors to secure smaller containers in the rotor slots.

4. For the SBS-MZ-4000/6-1.5 model there is an option to change the round rotor for a rectangular one equipped with more slots for containers.
5. Set the time and revs using the knobs (applicable to SBS-LZ-4000/20-6, SBS-LZ-4000/20-12), close the protective lid and activate the device.

### ROTOR CHANGE (APPLIES TO SBS-MZ-4000/6-1.5)

Use an Allen key to loosen the mounting bolt.

Attach a rectangular rotor using the mounting bolt.

Then lift the entire rotor.

Then tighten the bolt so that the rotor does not move out of position when in use.

### CENTRIFUGAL FORCE CALCULATIONS

$$RCF = 1.118 * 10^{-5} * R * (RPM)^2$$

RCF – Centrifugal force [g]

R – radius (distance from the axis of revolution) [cm]

RPM – Rotational speed [rev/min]

### IX. TRANSPORTATION AND STORAGE

During transport, the machine should be protected from shaking, crashing and turning upside down. Store it in a properly ventilated place with dry air and without any corrosive gas.

### X. CLEANING AND MAINTENANCE

- Disconnect the machine from the electricity when cleaning it.
- Use cleaner without corrosive substances to clean surface.

- After cleaning, all parts should be able to dry before the machine is to be reused.
- Store the unit in a dry, cool place, free from moisture.
- Never spray the device with water jets.

#### XI. REGULAR CONTROL OF THE DEVICE

Check regularly whether the device is damaged. If it should be damaged, please stop using it immediately and contact your customer service to solve the problem.

What to do in case of a problem?

Please contact your customer service and prepare the following information:

- Invoice number and serial number (the latter is to be found on the technical plate on the device).
- If relevant, a picture of the damaged, broken or defective part.
- It will be easier for your customer service clerk to determine the source of the problem if you give a detailed and precise description of the matter. The more detailed your information, the better the customer service will be able to solve your problem rapidly and efficiently!

**CAUTION:** Never open the device without the authorization of your customer service. This can lead to a loss of warranty!

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