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USER MANUAL

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1. GENERAL DESCRIPTION

The user manual is designed to assist in the safe and trouble-free use of the device. The product is designed and manufactured in accordance with strict technical guidelines, using state-of-the-art technologies and components. Additionally, it is produced in compliance with the most stringent quality standards.

DO NOT USE THE DEVICE UNLESS YOU HAVE THOROUGHLY READ AND UNDERSTOOD THIS USER MANUAL

To increase the product life of the device and to ensure trouble-free operation, use it in accordance with this user manual and regularly perform maintenance tasks. The technical data and specifications in this user manual are up to date. The manufacturer reserves the right to make changes associated with quality improvement. The device is designed to reduce noise emission risks to a minimum, taking into account technological progress and noise reduction opportunities.

LEGEND

	The product satisfies the relevant safety standards.
	Read instructions before use.
	The product must be recycled.

	WARNING! or CAUTION! or REMEMBER! Applicable to the given situation (general warning sign).
	ATTENTION! Electric shock warning!
	ATTENTION! Hot surface, risk of burns!
	Connect the device to a grounded power supply.
	Do not use the device to experiment with flammable, explosive, poisonous and highly caustic substances.

PLEASE NOTE! Drawings in this manual are for illustration purposes only and in some details may differ from the actual product.

The original operation manual is written in German. Other language versions are translations from the German.

2. USAGE SAFETY

ATTENTION! Read all safety warnings and all instructions. Failure to follow the warnings and instructions may result in electric shock, fire and/or serious injury or even death.

The terms "device" or "product" are used in the warnings and instructions to refer to "Digital & Magnetic Stirring Heating Mantles, Magnetic Stirring Heating Mantles". Do not use in very humid environments or in the direct vicinity of water tanks. Prevent the device from getting wet. Risk of electric shock! Do not cover the ventilation openings!

2.1. ELECTRICAL SAFETY

- a) The plug must fit the socket. Do not modify the plug in any way. Using original plugs and matching sockets reduces the risk of electric shock.
- b) Avoid touching earthed elements such as pipes, heaters, boilers and refrigerators. There is an increased risk of electric shock if the earthed device is exposed to rain, comes into direct contact with a wet surface or is operating in a damp environment. Water getting into the device increases the risk of damage to the device and of electric shock.
- c) Do not touch the device with wet or damp hands.
- d) Use the cable only for its designated use. Never use it to carry the device or to pull the plug out of a socket. Keep the cable away from heat sources, oil, sharp edges or moving parts. Damaged or tangled cables increase the risk of electric shock.
- e) If using the device in a damp environment cannot be avoided, a residual current device (RCD) should be applied. The use of an RCD reduces the risk of electric shock.
- f) Do not use the device if the power cord is damaged or shows obvious signs of wear. A damaged power cord should be replaced by a qualified electrician or the manufacturer's service centre.
- g) To avoid electric shock, do not immerse the cord, plug or device in water or other liquids. Do not use the device on wet surfaces.

2.2. SAFETY IN THE WORKPLACE

- a) Make sure the workplace is clean and well lit. A messy or poorly lit workplace may lead to accidents. Try to think ahead, observe what is going on and use common sense when working with the device.
- b) Do not use the device in a potentially explosive environment, for example in the presence of flammable liquids, gases or dust. The device generates sparks which may ignite dust or fumes.

- c) If you discover damage or irregular operation, immediately switch the device off and report it to a supervisor without delay.
- d) If there are any doubts as to the correct operation of the device, contact the manufacturer's support service.
- e) Only the manufacturer's service point may repair the device. Do not attempt any repairs independently!
- f) In case of fire, use a powder or carbon dioxide (CO₂) fire extinguisher (one intended for use on live electrical devices) to put it out.
- g) Please keep this manual available for future reference. If this device is passed on to a third party, the manual must be passed on with it.
- h) Keep packaging elements and small assembly parts in a place not available to children.
- i) Keep the device away from children and animals.

REMEMBER! When using the device, protect children and other bystanders.

2.3. PERSONAL SAFETY

- a) Do not use the device when tired, ill or under the influence of alcohol, narcotics or medication which can significantly impair the ability to operate the device.
- b) The device is not designed to be handled by persons (including children) with limited mental and sensory functions or persons lacking relevant experience and/or knowledge unless they are supervised by a person responsible for their safety or they have received instruction on how to operate the device.
- c) When working with the device, wear personal protective equipment that prevents:
 - splashing or contact with evaporating liquid,
 - contact with toxic or flammable gases.
- d) To prevent accidental starting, make sure the knobs are in off position before connecting to the power source.
- e) The device is not a toy. Children must be supervised to ensure that they do not play with the device.

2.4. SAFE DEVICE USE

- a) Do not use the device if the knobs or control buttons do not function properly. Devices that cannot be controlled by the controls are dangerous, may not be operated and must be repaired.
- b) Disconnect the device from the power supply before commencement of adjustment, cleaning and maintenance. Such a preventive measure reduces the risk of accidental activation.
- c) When not in use, store in a safe place, away from children and people not familiar with the device who have not read the user manual. The device may pose a hazard in the hands of inexperienced users.
- d) Keep the device in perfect technical condition. Before each use check for general damage and especially check for cracked parts or elements and for any other conditions which may impact the safe operation of the device. If damage is discovered, hand over the device for repair before use.
- e) Keep the device out of the reach of children.
- f) Device repair or maintenance should be carried out by qualified persons, only using original spare parts. This will ensure safe use.
- g) To ensure the operational integrity of the device, do not remove factory-fitted guards and do not loosen any screws.

- h) Do not move, adjust or rotate the device in the course of work.
- i) Do not leave this appliance unattended while it is in use.
- j) Clean the device regularly to prevent stubborn grime from accumulating.
- k) The device is not a toy. Cleaning and maintenance may not be carried out by children without supervision by an adult person.
- l) It is forbidden to interfere with the structure of the device in order to change its parameters or construction.
- m) Keep the device away from sources of fire and heat.
- n) **WARNING!** When the heating function is active, the heating mantle heats up to a high temperature. Do not touch hot parts of the device as this may cause burns.
- o) When working with flammable materials, pay special attention that the set heating temperature is not too high.
- p) Make sure that the liquid container can withstand the set heating temperature. A melted container may cause a fire.
- q) Never heat an empty container.
- r) Control the amount of liquid in containers. The liquid may evaporate.
- s) Spilling the flammable contents of the container onto the heating elements of the device may cause a fire.
- t) The set heating temperature must be at least 25°C lower than the ignition temperature of the stirred substance.
- u) Special attention must be paid to the hazards arising from:
 - flammable materials and substances with low boiling point,
 - overfilled containers,
 - containers made of brittle material or of low melting point.
- v) Pathogenic materials must be stirred only in closed containers.
- w) Do not cover the device while it is in operation.
- x) Make sure that the power cord does not touch any heated parts of the device.
- y) Keep the device away from strong magnetic fields.

ATTENTION! Despite the safe design of the device and its protective features, and despite the use of additional elements protecting the operator, there is still a slight risk of accident or injury when using the device. Stay alert and use common sense when using the device.

3. USE GUIDELINES

The device is designed for use in laboratories, schools, universities, and similar institutions. It is used for non-contact stirring and for heating liquids in laboratory vessels. **The user is liable for any damage resulting from unintended use of the device.**

3.1. DEVICE DESCRIPTION

SBS-LHM-1000/DH

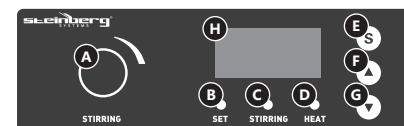
SBS-LHM-1000/H, SBS-LHM-500/H

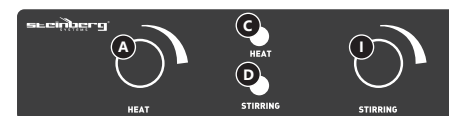
1. control panel
2. chamber with heating mantle
3. probe
4. probe clamp
5. the element for adjusting the height and extension of the clamp
6. probe support rod
7. probe cable

CONTROL PANEL SBS-LHM-1000/DH

CONTROL PANEL SBS-LHM-1000/H, SBS-LHM-500/H

- A. heating temperature control knob
- B. heating temperature control indicator light
- C. active stirring indicator light
- D. active heating indicator light
- E. button for activating the function of controlling the heating temperature and displaying the heating temperature
- F. increasing the heating temperature
- G. decreasing the heating temperature
- H. display of liquid temperature or the set heating temperature
- I. stirring speed control knob

3.2. PREPARING FOR USE

AMBIENT CONDITIONS

The device should be used in a working environment with the following parameters:

Ambient temperature	5–40°C
Ambient humidity	≤90%

APPLIANCE LOCATION

Ensure good ventilation in the room in which the device is being used. There should be at least 10 cm distance between each side of the device and the wall or other objects. The device should always be used when positioned on an even, stable, clean, fireproof and dry surface, and be out of the reach of children and persons with limited mental and sensory functions. Position the device such that you always have access to the power plug. The power cord connected to the appliance must be properly grounded and correspond to the technical details on the product label.

ASSEMBLING THE DEVICE

SBS-LHM-1000/DH

Before starting to use the device, install the temperature probe correctly.

- a) Screw the probe support rod elements together.
- b) Position the rod in the opening on the housing of the device. The spline in the opening should fit into the groove at the bottom of the rod.
- c) Fit the adjusting element and tighten the clamping knob.
- d) Place the probe clamp into the adjusting element and tighten the clamping knob.
- e) Place the probe in the clamp and make sure it is correctly clamped. The probe should be oriented with the cable upwards.
- f) Connect the probe cable to the suitable socket in the housing of the device.
- g) The device is ready to use.

3.3. DEVICE USE

- a) Pour the liquid into the container and insert the magnetic stirrer into the liquid.
- b) Place the container with the liquid inside the heating mantle chamber.
For the SBS-LHM-1000/DH model, insert the end of temperature probe into the liquid to a depth of at least 20 mm.
When using a metal container, do not place temperature probe on the bottom of the container. Placing the temperature probe on the bottom of the container may result in an incorrect measurement reading. The end of the measuring probe must be at least 5 mm from the bottom of the container; a distance of 10 mm is the best.
- c) Plug the power cord into the device and into the electrical outlet.
- d) Activate stirring with the knob [I]. When the stirring is active, the stirring indicator lamp lights up.
CAUTION! Do not set the stirring speed to high from the start. Increase the stirring speed gradually.

Depending on the model:

SBS-LHM-1000/DH

- e) Set the heating temperature by following these steps:
 - activate the heating temperature setting mode by pressing the [E] button,

- increase or decrease the set heating temperature using the [F] or [G] buttons.

SBS-LHM-1000/H, SBS-LHM-500/H

- Activate heating (if required) with the knob [A]. When the heating is on, the active heating indicator lamp lights up.
CAUTION! Do not activate heating under an empty container.
- After finishing work with the device, switch off stirring and heating and then disconnect the device from power supply.

3.4. CLEANING AND MAINTENANCE

- Unplug the mains plug and allow the device to cool completely before each cleaning, adjustment or replacement of accessories, or if the device is not being used.
- Use only non-corrosive cleaners to clean the surface.
- After cleaning the device, all parts should be dried completely before using it again.
- Store the unit in a dry, cool place, free from moisture and direct exposure to sunlight.
- Do not spray the device with a water jet or submerge it in water.
- Do not allow water to get inside the device through vents in the housing of the device.
- Clean the vents with a brush and compressed air.
- The device must be regularly inspected to check its technical efficiency and spot any damage.
- Use a soft cloth for cleaning.

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