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PH WERT TESTER

USER MANUAL

PH METER

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MIERNIK PH

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PH-METRE

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MANUAL DE INSTRUCCIONES

MEDIDOR DE PH

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PH MÉRŐ

BRUGSANVISNING

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DE | EN | PL | CZ | FR | IT | ES | HU | DA

SBS-PH-100 | SBS-PH-300

DE	Produktname:	PH-WERT TESTER
EN	Product name:	PH METER
PL	Nazwa produktu:	MIERNIK PH
CZ	Název výrobku	MĚŘIČ HODNOTY PH
FR	Nom du produit:	PH-METRE
IT	Nome del prodotto:	MISURATORE-PH
ES	Nombre del producto:	MEDIDOR DE PH
HU	Termék neve	PH MÉRŐ
DA	Produktnavn	PH-VÆRDI-TESTER
DE	Modell:	SBS-PH-100 SBS-PH-300
EN	Product model:	
PL	Model produktu:	
CZ	Model výrobku	
FR	Modèle:	
IT	Modello:	
ES	Modelo:	
HU	Modell	
DA	Model	
DE	Hersteller	expondo Polska sp. z o.o. sp. k.
EN	Manufacturer	
PL	Producent	
CZ	Výrobce	
FR	Fabricant	
IT	Produttore	
ES	Fabricante	
HU	Termelő	
DA	Producent	
DE	Anschrift des Herstellers	ul. Nowy Kisielin – Innowacyjna 7, 66-002 Zielona Góra Poland, EU
EN	Manufacturer Address	
PL	Adres producenta	
CZ	Adresa výrobce	
FR	Adresse du fabricant	
IT	Indirizzo del produttore	
ES	Dirección del fabricante	
HU	A gyártó címe	
DA	Producentens adresse	

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1. General Description

The instruction manual is intended to assist in safe and reliable use. The product is designed and manufactured strictly according to technical specifications using the latest technology and components and maintaining the highest quality standards.

**PLEASE CAREFULLY READ
AND UNDERSTAND THIS INSTRUCTION MANUAL BEFORE
OPERATION,**

To ensure long and reliable operation of the unit, make sure to operate and maintain it properly in accordance with the guidelines in this instruction manual. The technical data and specifications contained in this instruction manual are up to date. The manufacturer reserves the right to make changes in order to improve the quality.

Explanation of symbols

	The product complies with applicable safety standards.
	Please read the instructions before use.
	Recyclable product.

CAUTION! The illustrations in this instruction manual are for reference only and may differ from the actual product in some details.

The original instruction manual is in the German language version. Other language versions are translations from German.

2. Safety of use

CAUTION! Read all safety warnings and all instructions. Failure to follow warnings and instructions could result in serious injury or even death.

The term "appliance" or "product" in the warnings and in the description of the instructions refers to the pH METER

2.1. Safety in the workplace

- a) Keep the work area tidy and well lit. Disorder or poor lighting can lead to accidents. Be foresighted, watch what you are doing and use common sense when using the unit.
- b) If you have any doubts as to whether the unit is working properly or if it is damaged, contact the manufacturer's service department.
- c) Only the manufacturer's service department may repair the unit. Do not carry out repairs yourself!
- d) Keep these instructions for use for future reference. If the unit is to be passed on to a third party, the operating instructions must also be handed over together with the unit.
- e) Keep the packaging and small assembly parts out of the reach of children.
- f) Keep the unit away from children and animals.
- g) When using this unit together with other units, also follow the other instructions for use.

**Please
note!**

Keep children and other bystanders safe while operating the unit.

2.2. Personal safety

- a) The unit is not intended to be used by persons (including children) with reduced mental, sensory or intellectual functions or persons who lack experience and/or knowledge unless they are supervised or have been instructed by a person responsible for their safety on how to operate the unit.
- b) The unit is not a toy. Children should be watched to ensure that they do not play with the unit.

2.3. Safe use of the unit

- a) Keep unused equipment out of the reach of children and out of the reach of anyone unfamiliar with the unit or this instruction manual. These units is dangerous in the hands of inexperienced users.
- b) Keep the unit in good working condition. Check before each use for general damage or damage to moving parts (cracks in parts and components or any other condition that may affect the safe operation of the unit). If damaged, have the unit repaired before use.

-
- c) Keep the unit out of the reach of children.
 - d) To ensure the designed operational integrity of the unit, do not remove factory-installed covers or loosen screws.
 - e) The unit is not a toy. Cleaning and maintenance must not be performed by children without adult supervision.
 - f) Do not tamper with the unit to alter its performance or design.
 - g) Keep the unit away from sources of fire and heat.

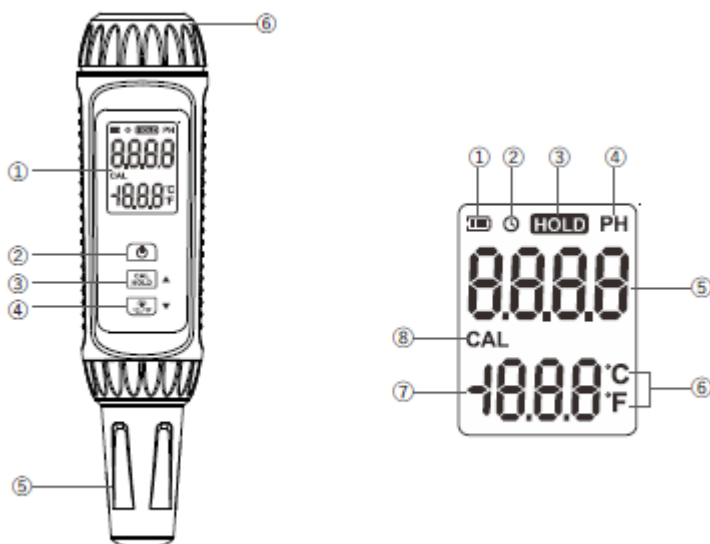
CAUTION! Although the product has been designed to be safe, with adequate safeguards, and despite the additional safety features provided to the user, there is still a slight risk of accident or injury when handling the unit. You are advised to use caution and common sense when using this product.

3. Rules of use

The device is designed to measure solution pH and solution temperature.

The user is responsible for any damage resulting from misuse.

3.1. Description

1. LCD display board
2. ON/ OFF button
3. UP button (HOLD/calibration function)
4. DOWN button (backlight/change unit)
5. Probe cover
6. Battery cover

1. Battery status indicator
2. Symbol for shut-off function after a certain time
3. HOLD function icon
4. pH display icon
5. Display area for pH value
6. Temperature unit
7. Temperature display area
8. Calibration icon

3.2. Working with the unit

TAKING THE MEASUREMENT

1. Before each measurement, wash the probe with distilled water and wipe it dry with a cotton cloth.
2. Start the instrument by pressing ON/OFF button (2).
3. Dip the probe tip into the solution to be tested.
4. Read the value on the display.

-
5. Clean the probe tip as in step 1, press and hold ON/OFF button to turn off the instrument.
 6. Pour the probe humidification fluid into the cover and close the cover.

BACKLIGHT ON THE SCREEN

To turn on the screen backlight, press the DOWN button after turning on the unit. Pressing the button again will turn the screen backlight off.

CAUTION! When the pH of the substance being measured exceeds the alarm level (pH>12.5 or pH<3.5) the screen will illuminate red.

CALIBRATION

The device should be calibrated if:

- it is being used for the first time,
- the probe has been replaced with a new one,
- the device has not been used for a long time,
- after measuring in a very strong solution,
- if the instrument indicates a measurement of pH>12 or pH<2,
- a fluoride containing solution was measured and the instrument indicated pH<7.

3-POINT CALIBRATION PROCEDURE:

1. Start the instrument and leave it in the measurement state.
2. Hold down the UP button until the CAL symbol appears on the display to enter the calibration state.
3. Once in calibration mode, the temperature display zone will display 4.00. At this point, wash the probe with distilled water and wipe with a cotton cloth to prepare it for measurement.
Then insert the probe into the pH 4.00 solution and wait 3-5 minutes.
When the pH is stable, briefly press the ON/OFF button to proceed to the next calibration point.
4. Repeat the probe cleaning.
5. Insert the probe into the pH 6.86 solution and wait 3-5 minutes. When pH is stable, briefly press the ON/OFF button again to go to the next calibration point.
6. Repeat the cleaning process.
7. Insert the probe into the pH 9.18 solution and wait 3-5 minutes. When the pH is stable, briefly press the ON/OFF button again to complete the calibration.
8. If:
 - calibration was successful- the display will show PASS,
 - calibration failed- the display will show ERR indicating that the pH of the calibration solution is not within the calibration point and the instrument

needs to be recalibrated.

CAUTION! If the pH of the solution used for calibration differs from the default value it is possible to adjust these values during calibration using the UP and DOWN buttons.

CHANGING THE TEMPERATURE UNIT

To change the unit of temperature measurement, start the instrument and press and hold the DOWN button (4) in the measurement state.

HOLD MODE

This function is used to freeze the value on the display. To activate this function press UP (3) button during measurement.

The display will show the HOLD symbol. Press the UP button again briefly to cancel the hold.

SWITCH OFF FUNCTION AFTER A SET TIME

To activate this function, switch on the device and briefly press the ON/OFF button while in measurement mode. In this way the device will automatically switch off after 15 minutes of inactivity. This function is indicated by the clock symbol on the display.

3.3. Cleaning and maintenance

- a) Store the unit in a dry and cool place protected from moisture and direct sunlight.
- b) Perform regular inspections of the unit checking technical fitness and any damages.
- c) Do not leave the batteries in the unit when it will not be used for a long time.
- d) Avoid touching the probe directly with your hand, this can damage it!
- e) Always put the probe cover back on when the device will not be used for a long time.
- f) The probe should be kept moist, therefore a small amount of special liquid should be placed in the probe cover after each use.

The use of 3.0 mol/l potassium chloride solution is recommended. Do not use tap water or demineralized water for this purpose!

-
- g) After each use the probe should be cleaned by rinsing it with demineralized water and wiping with a cotton cloth. In order to thoroughly clean the probe from contaminants it is recommended to regularly wash the probe with 0.1 mol/l diluted hydrochloric acid.
 - h) A new probe should be immersed in 3 mol/l potassium chloride solution for 24 hours.

INSTRUCTIONS FOR THE SAFE DISPOSAL OF BATTERIES.

The units are equipped with 1.5V batteries.

To replace the batteries, unscrew the battery cover (6), place the batteries in the compartment according to the polarity markings and screw the cover back on.

Return the batteries to a unit responsible for their disposal.

DISPOSAL OF USED UNITS.

At the end of its useful life, this product should not be disposed of with normal household waste but should be taken to a collection point for the recycling of electrical and electronic equipment. This is indicated by the symbol on the product, operating instructions or packaging. The materials used in this unit are recyclable according to their marking. You will be making an important contribution to protecting our environment by reusing, recycling or otherwise disposing of used units.

Your local administration will provide you with information about the appropriate disposal point for used units.

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